
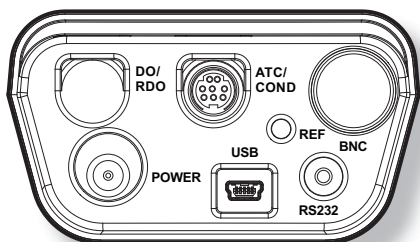


## Preparation

### Power Source

1. Power adapter (sold separately)
  - a. Select the appropriate wall socket plug plate.
  - b. Slide off the clear plastic cover from the plug plate.
  - c. Slide the plug plate into the groove on the back of the power adapter.
  - d. Connect the power adapter to the meter and power outlet.
2. Batteries (included with meter)
  - a. Select four AA alkaline batteries.
  - b. Confirm that the meter is powered off.
  - c. Remove the battery compartment cover – loosen the screws holding the battery cover, release the top portion of the battery cover from the meter (use a coin or your finger) and release the bottom portion of the battery cover.
  - d. Orientate the batteries as shown in the battery compartment housing and insert batteries.
  - e. Replace the battery compartment cover and screws.

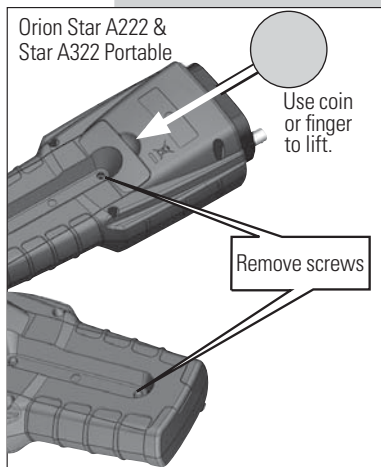


### Electrodes and Other Connections

1. Prepare the conductivity cell and any other applicable electrodes according to the directions in the electrode user guide.
2. Connect the appropriate items as labeled on the meter and as shown in the figure above:

For additional information on meter setup and operation, refer to the reference guide. The reference guide is on the included CD and available at

[www.thermoscientific.com/water](http://www.thermoscientific.com/water).

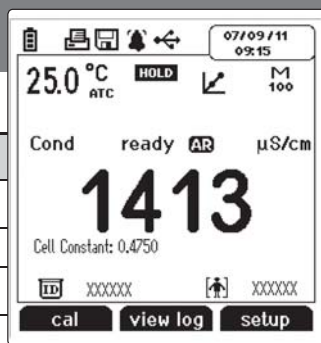

## Display Information

Display Icon	Description
	Shown when the meter is running on AC power.
	Shown when the meter has batteries installed.
	Indicates data is being sent to a computer or printer.
	Indicates data is being sent to the data log.
	Shown when an alarm is set and the alarm value is reached.
	Indicates the meter is set to be interfaced with a printer or computer via the RS232 port.
	Indicates the meter is set to be interfaced with a printer or computer via the USB port.
	Displays the time and date entered in the setup menu.
	Displays the current temperature based on the temperature probe reading or entered temperature value. Shows the origin of the temperature as MAN (entered temperature) or ATC (temperature probe).
	Shown when  is pressed and the displayed measurement is frozen.
	Indicates a calibration was successfully completed.
	Indicates a method is in use and the number of the method being used.
<b>Cond</b>	Indicates the type of measurement and determines the type of calibration that will be performed.
<b>ready</b>	Specifies the stability of the electrode as <b>stabilizing</b> or <b>ready</b> .
	Shown when the meter is in AUTO-READ mode. The  icon will blink while the reading is stabilizing and stop blinking when the reading is stable and the measurement is locked on the display.
<b>1413 μS/cm</b>	Displays the measurement value based on the last saved calibration and current electrode reading. Units are shown to the right of the value.
<b>Cell Constant: 0.4750</b>	Shows the cell constant in use from the last saved calibration.
XXXXXX	Shows the operator assigned sample ID number. <i>Star A322 meter only.</i>
XXXXXX	Shows the operator assigned user ID number. <i>Star A322 meter only.</i>
	Displays the action that will be performed when <b>f1</b> is pressed.
	Displays the action that will be performed when <b>f2</b> is pressed.
	Displays the action that will be performed when <b>f3</b> is pressed.

## Keypad Display Information



Press the f1, f2 and f3 function keys to perform the action shown above each key on the display.



Press to turn the meter on.

When the meter is on, press and quickly release to turn the display backlight on or off or hold down to turn the meter off.



In the measurement mode, press to take a measurement.

In the setup, calibration and other menus, press to escape the current menu and return to the measurement mode.



In the measurement mode, press to enter the setup menu.

In the setup, calibration and other menus, press to scroll up through a list of options.



In the continuous measurement mode, press to freeze the displayed measurement and press again to unfreeze the measurement.

In the setup, calibration and other menus, press to scroll left through a list of options.



In the measurement mode, press to change the displayed measurement mode. Options are Cond (conductivity), TDS (total dissolved solids), Salinity and Res (resistivity).









In the setup, calibration and other menus, press to scroll right through a list of options.



In the measurement mode, press to log or print a measurement.

In the setup, calibration and other menus, press to scroll down through a list of options.

## Keypad




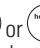
1. Press  to power the meter on. When the meter is on, press and quickly release  to turn the backlight on or off or press and hold  for about three seconds to power the meter off.
2. Press  to exit any meter function and return to the measurement mode.
3. The **f1**, **f2**, and **f3** function keys perform a variety of meter operations. The menu-specific operation is shown above each key. For example, press **f1** in the measurement mode to start a calibration.
4. The , ,  or  keys are used as navigation keys (up, right, down, left) when selecting from a fixed list or grid of meter options. In the measurement mode, these keys are used to access the setup menu, change the measurement mode, manually log or print a measurement and hold (freeze) a displayed measurement.

## Conductivity Calibration




One to five conductivity standards can be used for calibration. Always use fresh standards and select standards that are near the sample conductivity. Prepare the conductivity cell according to the instructions in the conductivity cell use guide. Connect the conductivity cell and any other electrodes to be used to the meter. Power on the meter and set the measurement mode to conductivity.




**Note:** For an automatic calibration, the nominal cell constant of the conductivity cell must be entered in the setup menu before the calibration is performed and Thermo Scientific Orion 100  $\mu\text{S}/\text{cm}$ , 1413  $\mu\text{S}/\text{cm}$  and/or 12.9  $\text{mS}/\text{cm}$  conductivity standards must be used.

### Automatic and Direct Calibration














1. In the measurement mode, press **f1 (cal)**.
2. Rinse the conductivity cell and any other electrodes in use with distilled water, blot dry with a lint-free tissue and place into the standard.
3. When the conductivity cell and standard are ready, press **f3 (start)**.
4. Wait for the conductivity value on the meter to stabilize and stop flashing and perform one of the following actions:
  - a. Press **f2 (accept)** to accept the displayed conductivity value.
  - b. Press **f3 (edit)** to access the numeric entry screen and edit the conductivity standard value.
    - i. Press , ,  or  to highlight a number or decimal point, press **f3 (enter)** to select the highlighted item and repeat until the standard value at the measured temperature is shown.
    - ii. Press **f2 (done)** to exit the numeric entry screen.
    - iii. Press **f2 (accept)** to accept the entered conductivity value.
5. Press **f2 (next)** to proceed to the next standard and repeat steps 2 through 4 or press **f3 (cal done)** to save and end the calibration. If five standards are used, the calibration will save and end once the fifth conductivity standard value is accepted.
6. The meter will display the calibration summary including the average calculated cell constant. Press **f1 (meas)** to export the data to the calibration log or press **f2 (print)** to export the data to the calibration log and a printer or computer. The meter will automatically proceed to the measurement mode.

## Measurement

Press  while taking a measurement in the continuous measurement mode to freeze the display and press  a second time to unfreeze the display and continue the measurement. Press  while taking a measurement to manually export the measurement to the data log, if the data log is enabled in the setup menu.




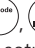










1. Ensure the measurement mode is set correctly. To set the measurement, press  to scroll through the conductivity, TDS, salinity and resistivity measurement modes.
2. Rinse the conductivity cell electrode and any other electrodes in use with distilled water, blot dry with a lint-free tissue and place into the sample.
3. Start the measurement and wait for it to stabilize.
  - a. If the meter is in **AUTO-READ** mode (default setting), press  to start the measurement. When the **AR** icon stops flashing, record the conductivity, TDS, salinity or resistivity and temperature of the sample. Press  again to start a new measurement.
  - b. If the meter is in continuous mode, the meter will immediately start taking a measurement and update the display whenever the measurement changes. Wait for the display to show **ready** and record the conductivity, TDS, salinity or resistivity and temperature of the sample.
  - c. *Star A322 meter only* – If the meter is in timed mode, the meter will log measurements at the preselected time interval, regardless of the measurement stability. The meter will update the display whenever the measurement changes, so the conductivity, TDS, salinity or resistivity and temperature of the sample can be recorded when the display shows **ready**.
4. Remove the conductivity cell from the sample, rinse with distilled water, blot dry and place into the next sample.
5. Repeat steps 3 and 4 for all samples.
6. When all samples have been measured, store the conductivity cell according to its user guide.

## Read Type Selection

1. In the measurement mode, press .
2. Press    or  to highlight **COND Channel** and press **f3 (select)**.
3. Press  or  to highlight **Mode and Settings** and press **f3 (select)**.
4. Press  or  to highlight **Read Type** and press **f3 (select)**.
5. Press  or  to highlight **Auto**, **Continuous** or **Timed** (*Star A322 meter only*) and press **f3 (select)**.
  - a. *Star A322 meter only* – If **Timed** is selected and the time interval needs to be changed – highlight **Timed**; press  to highlight hours (HH), minutes (MM) or seconds (SS); press **f3 (edit)** to access the numeric entry screen; use the numeric entry screen to change the values and press **f1 (back)** when the time interval is correct.
6. Press  to return to the measurement mode.
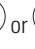
## Setup Menu
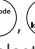
### Navigating the Setup Menu















1. In the measurement mode, press  to enter the main setup menu.
2. Press , ,  or  to scroll through the main setup menu options. Press **f3 (select)** to select a main setup menu option.
3. Press  or  to scroll through setup submenu options. Press **f3 (select)** to select a setup submenu option.
4. Perform the appropriate actions to set the desired parameter in the setup submenus.
  - a. To select a value from a list of options, press  or  to highlight the desired value and press **f3 (select)** to set the value.
  - b. To enter a numeric value, use the numeric entry screen.
    - i. Select the value to be entered by pressing **f3 (select)** or **f3 (edit)**. The numeric entry screen will popup on the display.
    - ii. Press , ,  or  to highlight a number, decimal place or negative sign; press **f3 (enter)** to select the highlighted item and repeat until the desired value is shown on the top of the numeric entry screen.
    - iii. Press **f2 (done)** to save the value and exit the numeric entry screen.
5. Press **f1 (back)** and then  key to return to the measurement mode at any time.
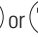
### Setup Menu Overview

COND Channel	Settings	Log View	Diagnostics
<p><b>Method</b></p> <hr/> <p><b>Mode &amp; Settings</b></p> <ul style="list-style-type: none"> <li>• Measure Mode</li> <li>• Read Type</li> <li>• Cell K</li> <li>• Stability</li> <li>• Averaging</li> <li>• Ref Temp</li> <li>• Temp Comp.</li> <li>• Temp Coeff</li> <li>• Alarm Settings</li> <li>• Sample ID</li> <li><small>*Star A322 meter only</small></li> </ul> <hr/> <p><b>Temperature</b></p> <ul style="list-style-type: none"> <li>• Manual Temp Value</li> <li>• Temperature Unit</li> <li>• Temperature Calibration</li> </ul>	<ul style="list-style-type: none"> <li>• Export Data</li> <li>• Data Log</li> <li>• Date / Time</li> <li>• Language</li> <li>• Key Press Beep</li> <li>• Alarm Beep</li> <li>• Contrast</li> <li>• Auto Shut Off</li> <li>• User ID</li> <li><small>*Star A322 meter only</small></li> </ul>	<ul style="list-style-type: none"> <li>• Data Log</li> <li>• Calibration Log</li> </ul>	<ul style="list-style-type: none"> <li>• Meter Self Test</li> <li>• Factory Reset</li> <li>• About Meter</li> </ul>













## Nominal Cell Constant Entry

The nominal cell constant value is used during an automatic conductivity calibration and allows the meter to determine which Thermo Scientific Orion conductivity standard is being used for the calibration.

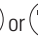1. In the measurement mode, press .
2. Press , ,  or  to highlight *COND Channel* and press **f3 (select)**.
3. Press  or  to highlight *Mode and Settings* and press **f3 (select)**.
4. Press  or  to highlight *Cell K* and press **f3 (select)**.
5. Press **f3 (select)** to access the numeric entry screen and enter the nominal cell constant of the conductivity cell.
  - a. Press , ,  or  to highlight a number or decimal point, press **f3 (enter)** to select the highlighted item and repeat until the nominal cell constant value is shown.
  - b. Press **f2 (done)** to exit the numeric entry screen.
6. Press **f1 (back)** and then press  to return to the measurement mode.













## Reference Temperature Selection

The reference temperature is the temperature that all conductivity measurements will be reported at by the meter if temperature compensation is enabled.

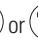













1. In the measurement mode, press .
2. Press , ,  or  to highlight *COND Channel* and press **f3 (select)**.
3. Press  or  to highlight *Mode and Settings* and press **f3 (select)**.
4. Press  or  to highlight *Ref Temp.* and press **f3 (select)**.
5. Press  or  to highlight **5 °C (Star A322 meter only), 10 °C (Star A322 meter only), 15 °C, 20 °C or 25 °C** and press **f3 (select)**.
6. Press  to return to the measurement mode.

## Temperature Compensation Selection











The temperature compensation can be turned off or set to Linear, nLFn (non-linear ultra pure non-degassed water), nLFu (non-linear ultra pure degassed water) or EP (temperature compensation off and warning is displayed if conductivity values are outside EP requirements for ultra pure water) and is used to report all conductivity measurements at the selected reference temperature.

1. In the measurement mode, press .
2. Press , ,  or  to highlight *COND Channel* and press **f3 (select)**.
3. Press  or  to highlight *Mode and Settings* and press **f3 (select)**.
4. Press  or  to highlight *Temp. Comp.* and press **f3 (select)**.
5. Press  or  to highlight **Off, Linear, nLFn, nLFu (Star A322 meter only) or EP (Star A322 meter only)** and press **f3 (select)**.
6. Press  to return to the measurement mode.

## Viewing the Calibration Log

1. In the measurement mode, press .
2. Press , ,  or  to highlight **View Log** and press **f3 (select)**.
3. Press  or  to highlight **Calibration Log** and press **f2 (accept)**.
4. Press  or  to highlight **Conductivity – Channel** and press **f2 (select)**.
5. Press  or  to highlight **Conductivity, Resistivity, TDS, Salinity or Temperature** as the calibration type and press **f2 (select)**.
6. The meter will display a list of calibrations for the selected channel and calibration type. The list shows the sequential number of the calibration and the date and time it was saved (**07/01/2011 12:45**).
7. To view the calibration data, press  or  to highlight a specific calibration and press **f2 (select)**. Press **f2 (print)** to print the calibration or press **f1 (back)** to return to the list of calibrations.
8. Press  to return to the measurement mode.

## Viewing the Data Log

1. In the measurement mode, press .
2. Press , ,  or  to highlight **View Log** and press **f3 (select)**.
3. Press  or  to highlight **Data Log** and press **(accept)**.
4. The meter will display a list of the data points. The list shows the sequential number of the data point and the date and time the data point was saved (**07/01/2011 12:45**).
5. To view the measurement information for an individual data point, press  or  to highlight the data point and press **f2 (select)**. Press **f2 (print)** to print the data point or press **f1 (back)** to return to the list of data points.
6. Press  to return to the measurement mode.