


## Quick-Action Spigot Use and Care

**Warning:**  
**Do not use spigot as a handle!**

- Spigot spout fits 5/8-in. I.D. tubing. Adapter included takes tubing with 1/4- to 5/16-in I.D.


### To attach spigot:

1. Screw lock nut of spigot onto boss of carboy until tight.
2. Push spigot body firmly over nipple of threaded boss to seat O-ring. You will feel the Spigot body snap into place. Spigot body is seated properly when nut turns easily again.
3. Point spout towards bottom of container and turn lock nut until tight again.

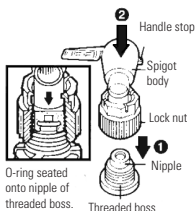

### To remove and disassemble spigot for cleaning:

1. Unscrew lock nut from from container. Nut and spigot housing are permanently attached; do not separate them.
2. Unscrew spigot handle counter clockwise from spigot body lifting handle up over handle-stop.

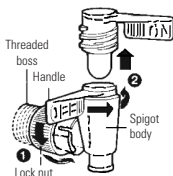

### To clean spigot:

1. **General:** Wash in a mild detergent; rinse with tap water; repeat rinse with distilled water.
2. **Dishwasher:** Use plastics cycle; avoid excessive abrasion by covering metal projection; weigh down spigot; use hot water. Detergent usually not required.

### To reassemble spigot:

1. Insert handle into spigot body with handle facing away (180°) from handle-stop.
2. Turn handle clockwise, lifting handle gently over handle-stop. Turn until handle stops. Spigot should be in "full off" position.

