Alexa Fluor[®] Reference Standards

Pub. no. MAN0009935

by *life* technologies

molecular probes[®]

Store at ≤-20°C, protected from light.

Label	Cat. no.	Ex/Em max*	Quantity
Alexa Fluor® 350 carboxylic acid	A33076	346/442	5 mg
Alexa Fluor® 488 carboxylic acid	A33077	495/519	5 mg
Alexa Fluor® 532 carboxylic acid	A33078	531/554	5 mg
Alexa Fluor® 546 carboxylic acid	A33079	556/575	5 mg
Alexa Fluor® 555 carboxylic acid	A33080	555/565	5 mg
Alexa Fluor® 568 carboxylic acid	A33081	578/603	5 mg
Alexa Fluor [®] 594 carboxylic acid	A33082	590/617	5 mg
Alexa Fluor® 633 carboxylic acid	A33083	632/647	5 mg
Alexa Fluor® 647 carboxylic acid	A33084	650/668	5 mg
Alexa Fluor® 750 carboxylic acid	A33085	749/775	5 mg

* Approximate excitation and emission maxima, in nm.



For Research Use Only. Not for use in diagnostic procedures.

Guidelines for use

- Dissolve the Alexa Fluor[®] carboxylic acid compound in aqueous buffer (e.g. PBS), or high-quality, anhydrous dimethylsulfoxide (DMSO) or dimethylformamide (DMF) at a final concentration of 1–10 mg/mL. Once reconstitued, this solution is stable for 1 year, when stored at 2–8°C, protected from light.
- Prepare working solutions immediately before use, and then discard the working solutions after use.
- Because the actual fluorescence units can vary as a function of several factors including temperature and instrument settings, it is important to rigorously control these variables. For example, dilute the stock solution with the same buffer as that used with the analyte of interest and use exactly the same instrument settings.
- A 100-fold dilution of the standard solution is recommended as a starting point before measuring fluorescence.

Limited Product Warranty

Life Technologies Corporation and/or its affiliate(s) warrant their products as set forth in the Life Technologies' General Terms and Conditions of Sale found on Life Technologies' website at www.lifetechnologies.com/termsandconditions. If you have any questions, please contact Life Technologies at www.lifetechnologies.com/support.

BIGLANGE. LET TECHNOLOGIES CORPORTION AND/ORITIS AFFLUATES) DISCLAM ALL WARRANTES WITH RESPECT TO THE DOLLMENT, DPRESSED OR MAREED, MICLIDNE UNIT DI LIMITE DI THE DISC OF MERCINALILUIT, FINTESS CAR A BARTICLLAR PURCHES, OR NAN-AFFRICAN AND MAREET, THE DISCLAM ALL MAREET, A CONTRACT, TOTAL MARANTI, ON MARA

© 2014 Life Technologies Corporation. All rights reserved. The trademarks mentioned herein are the property of Life Technologies Corporation and/or its affiliate(s) or their respective owners.

For support visit lifetechnologies.com/support

lifetechnologies.com

13 January 2014