

BackDrop® Background Suppressor for Microplate Assays

Cat. nos. B10511, B10512

Pub. no. MAN0010825

Revision A.0

Product	Cat. no.	Amount	Storage
BackDrop® Background Suppressor for Microplate Assays	B10511	5 mL	• 2–8°C • Do not freeze
	B10512	50 mL	

Quick Protocol



1. Prepare a 1X Microplate Assay Reagent by diluting the BackDrop[®] Background Suppressor reagent 1:10 in your microplate assay reagent of choice.



2. Remove media from cells.



3. Add 100 μ L per well into 96-well or 25 μ L per well into a 384-well plate of the 1X Microplate Assay Reagent containing the BackDrop[®] Background Suppressor reagent.



4. Add drug or compound library to the wells and proceed with analysis.

Limited Product Warranty

Life Technologies Corporation and/or its affiliate[s] warrant their products as set forth in the Life Technologies' General Terms and Conditions of Sale found on Life Technologies' website at www.lifetechnologies.com/termsandconditions. If you have any questions, please contact Life Technologies at www.lifetechnologies.com/support.

DISCLAIMER - LIFE TECHNOLOGIES CORPORATION AND/OR ITS AFFILIATE[S] DISCLAIM ALL WARRANTIES WITH RESPECT TO THIS DOCUMENT, EXPRESSED OR IMPLIED, INCLUDING BUT NOT LIMITED TO THOSE OF MERCHANTABILITY, FITNESS FOR A PARTICULAR PURPOSE, OR NON-INFRINGEMENT. TO THE EXTENT ALLOWED BY LAW, IN NO EVENT SHALL LIFE TECHNOLOGIES AND/OR ITS AFFILIATE[S] BE LIABLE, WHETHER IN CONTRACT, TORT, WARRANTY, OR UNDER ANY STATUTE OR ON ANY OTHER BASIS FOR SPECIAL, INCIDENTAL, INDIRECT, PUNITIVE, MULTIPLE OR CONSEQUENTIAL DAMAGES IN CONNECTION WITH OR ARISING FROM THIS DOCUMENT, INCLUDING BUT NOT LIMITED TO THE USE THEREOF.

©2014 Thermo Fisher Scientific Inc. All rights reserved. All trademarks are the property of Thermo Fisher Scientific and its subsidiaries unless otherwise specified.

For support visit lifetechnologies.com/support
lifetechnologies.com