MitoProbe[™] TMRM Kit for Flow Cytometry

Catalog No. M20036

Pub. No. MAN0017633

Rev. B.0

Product information

Presence of a mitochondrial membrane potential ($\Delta \Psi m$), the electrochemical gradient between the interior and exterior of the mitochondria, is a sign of healthy cells. As a cationic and lipophilic fluorescent dye, TMRM can be used to detect changes in the membrane potential by tracking the redistribution of the dye. A healthy cell will display a robust fluorescent signal; however, the fluorescence diminishes when the mitochondrial membrane depolarizes due to changes such as induction of apoptosis.

Table 1. Contents and storage

Product	Amount	Concentration	Storage*
TMRM (Component A)	100 µL	20 µM solution in DMSO	• Store in freezer (–5°C to
CCCP (Carbonyl cyanide 3-chlorophenylhydrazone) (Component B)	125 µL	50 mM solution in DMSO	–30°C) • Desiccate • Protect from light
* When stored as directed, the product is stable for at least 6 months from the date receipt. Approximate Ex/Em maxima: 561 nm/585 nm			

Materials required, but not
provided• Cells of interest• Phosphate-buffered saline (PBS) or similar protein-free buffer• Culture media containing protein such as FBS or BSA• Flow cytometer

Caution No data are available addressing the mutagenicity or toxicity of the TMRM reagent. Handle the DMSO and DMSO containing reagents with caution because DMSO is known to facilitate the entry of organic molecules into tissues. Always wear suitable protective clothing, gloves and eye/face protection when handling this reagent. Dispose of the reagents in compliance with all pertaining local regulations.

Storage and handlingUpon receipt, store the kit components desiccated at ≤-20°C until required for use.
When stored as directed, TMRM and CCCP reagents are stable for at least 6 months.
Allow the products to warm to room temperature before opening the vials. Stock
solution can be frozen after use, but should be aliquoted to avoid repeated freezing and
thawing.




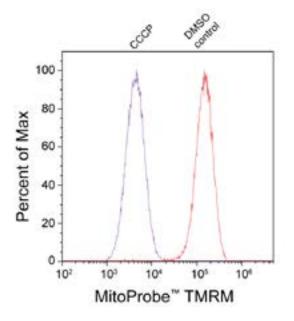


Figure 1. MitoProbe[™] TMRM detects changes in mitochondrial membrane potential. Jurkat cells, a human T-lymphocyte cell line, were treated with DMSO (control) or 50 µM CCCP to induce mitochondrial membrane depolarization. Cells were subsequently stained with MitoProbe[™] TMRM reagent and analyzed on an Attune[™] NxT Flow Cytometer using a 561-nm laser for excitation and 585/16BP-nm filter for emission.

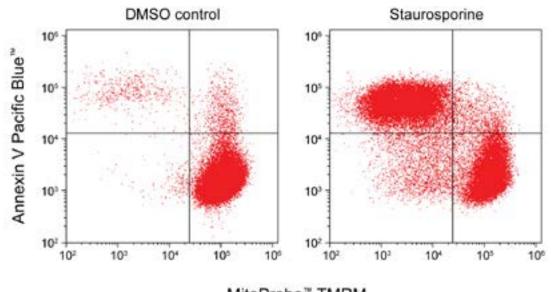

MitoProbe[™] TMRM

Figure 2. Detection of mitochondrial membrane potential changes in apoptotic cells. Jurkat cells, a human T-lymphocyte cell line, were treated with DMSO (control) or 500 nM staurosporine for 2 hours. Cells were subsequently stained with 20 nM MitoProbe[™] TMRM reagent for 30 minutes at 37°C, followed by a wash and additional stain with Annexin V Pacific Blue[™] reagent. Data were acquired on an Attune[™] NxT Flow Cytometer using a 561-nm laser and a 585/16-nm emission filter for the MitoProbe[™] TMRM reagent and a 405-nm laser and 440/50-nm emission filter for the Annexin V Pacific Blue[™] reagent. Staurosporine induced apoptosis, resulting in a mixed population of cells containing a population of healthy MitoProbe[™] TMRM-positive cells as well as a population of apoptotic, Annexin V Pacific Blue[™]-positive, MitoProbe[™] TMRM-low cells.

Prepare reagent	toProbe [™] TMRM reagent is provided as a 1000X stock solution at a concentration o M in DMSO. To use it, simply add 1 μL of 20 μM TMRM reagent to 1 mL of a 10 ⁶ cells.	of
Stain cultured cells with TMRM reagent	fore beginning your experiment, ensure that the vial of CCCP has equilibrated to om temperature.	
	r each sample, resuspend the cells in cell culture medium or PBS at approximately $^{10^6}$ cells/mL.	
	r the control sample, add 1 μ L of 50 mM CCCP to the cells, then incubate for ninutes at 37°C, 5% CO ₂ .	
	r experimental samples, add 1 μ L from the 20 μ M stock TMRM reagent solution 0 nM final concentration) and incubate for 30 minutes at 37°C, 5% CO ₂ .	
	<i>tional:</i> Wash the cells once in 1 mL of culture medium or PBS, then resuspend the ls in 500 µL of PBS (or other suitable buffer).	
	alyze the cells on a flow cytometer with 561-nm excitation using emission filters propriate for R-phycoerythrin.	
	te: TMRM is optimally excited at 561 nm; however, it can also be excited by the 2-nm or 488-nm laser lines. In each case, the PE equivalent bandpass filter should b ed to detect TMRM.	e

Ordering information

Cat. No.	Product	Unit size
M20036	MitoProbe [™] TMRM Kit for Flow Cytometry	100 tests

Documentation and support

Customer and Technical Support

Visit thermofisher.com/support for the latest in services and support, including:

- Worldwide contact telephone numbers
- Product support, including:
 - Product FAQs
 - Software, patches, and updates
 - Training for many applications and instruments
- · Order and web support
- Product documentation, including:
 - User guides, manuals, and protocols
 - Certificates of Analysis
- Safety Data Sheets (SDSs; also known as MSDSs)

Note: For SDSs for reagents and chemicals from other manufacturers, contact the manufacturer.

Limited Product Warranty

Life Technologies Corporation and/or its affiliate(s) warrant their products as set forth in the Life Technologies' General Terms and Conditions of Sale found on Life Technologies' website at www.thermofisher.com/us/en/home/global/terms-and-conditions.html. If you have any questions, please contact Life Technologies at thermofisher.com/support.

The information in this guide is subject to change without notice.

For Research Use Only. Not for use in diagnostic procedures.

DISCLAIMER: TO THE EXTENT ALLOWED BY LAW, LIFE TECHNOLOGIES AND/OR ITS AFFILIATE(S) WILL NOT BE LIABLE FOR SPECIAL, INCIDENTAL, INDIRECT, PUNITIVE, MULTIPLE OR CONSEQUENTIAL DAMAGES IN CONNECTION WITH OR ARISING FROM THIS DOCUMENT, INCLUDING YOUR USE OF IT.

Revision history: Pub. No. MAN0017633

Revision	Date	Description
B.0	21 June 2018	Product name correction in the Methods section
A.0	16 February 2018	New user guide

Important Licensing Information: These products may be covered by one or more Limited Use Label Licenses. By use of these products, you accept the terms and conditions of all applicable Limited Use Label Licenses.

Manufacturer: Life Technologies Corporation | 29851 Willow Creek Road | Eugene, OR 97402

Trademarks: All trademarks are the property of Thermo Fisher Scientific and its subsidiaries unless otherwise specified.

©2018 Thermo Fisher Scientific Inc. All rights reserved.

