



PRODUCT INFORMATION

**Glycogen,
RNA grade**

Inert co-precipitant of nucleic acids. Derived from oysters.

Pub. No. MAN0013149

Rev. Date 09 May 2017 (Rev. C.00)

#R0551

Lot: _ Expiry Date: _

Store at -20 °C

Ordering Information

Component	#R0551
Glycogen, RNA grade, Concentration: 20 mg/mL aqueous solution.	2 x 0.1 mL

www.thermofisher.com

For Research Use Only. Not for use in diagnostic procedures.

Description

Glycogen is a highly purified polysaccharide derived from oysters. It is an inert carrier, free of host DNA/RNA. Glycogen is insoluble in ethanol solution; in the presence of salts it forms a precipitate that traps the target nucleic acids. During centrifugation, a visible pellet is formed, which greatly facilitates handling of nucleic acids. Glycogen quantitatively precipitates nucleic acids from diluted solutions with a higher efficiency than that of tRNA, linear polyacrylamide or sonicated DNA (1-4). Glycogen molecules are highly branched structures composed of thousands of glucose molecules bonded to each other. The molecular weight of the largest individual glycogen molecule containing about 50,000 glucose molecules appears to be 8 million.



Protocol for RNA Precipitation from Diluted Solutions

1. Add 1/10 volume of 3 M sodium acetate (or 2 M sodium chloride, or 5 M ammonium acetate) to RNA solution.
Use DEPC-treated tubes and tips when working with RNA.
2. Add glycogen to a final 0.05-1 $\mu\text{g}/\mu\text{L}$ concentration.
Use up to 1 μL of glycogen per 20 μL of the solution.
For precipitation of oligonucleotides, do not use higher than 1 $\mu\text{g}/\mu\text{L}$ final glycogen concentration.
3. Add 1 volume of isopropanol (or 2.5 volumes of ethanol) to the solution. Mix gently but thoroughly.
Use ethanol for < 200 bp fragments.
4. Incubate the mixture at -20 °C for 60 min, or at -70 °C for 30 min.
Longer incubation time and lower temperature provide better recovery of nucleic acids.
5. Centrifuge the mixture for 10-15 min at 10,000 rpm.
6. Discard the supernatant.
7. Rinse the pellet with cold 70% ethanol.
8. Air-dry the pellet.
Avoid over-drying the pellet, as it then takes more time to dissolve.
9. Dissolve RNA in DEPC-treated Water (#R0601).

Note

- 5 μg of Glycogen (0.25 μL) forms a visible pellet.
- Optimal for recovery of oligonucleotides (>8 bases) and low amounts (>20 pg) of nucleic acids from diluted solutions.
Up to a final concentration of 8 $\mu\text{g}/\mu\text{L}$, glycogen does not interfere with most downstream applications, e.g. cDNA synthesis.
- Up to a final concentration of 0.4 $\mu\text{g}/\mu\text{L}$, glycogen does not affect *in vitro* transfection of eukaryotic cells.
- Does not interfere with gel electrophoresis of nucleic acids.
- Does not interfere with spectrophotometrical determination of nucleic acids concentration ($A_{260-280}$ measurements).

(continued on back page)

CERTIFICATE OF ANALYSIS

Nucleic Acids Precipitation Assay

RNA was precipitated with glycogen and ethanol. 85 ± 15 % of RNA was recovered.

Nicking Activity Assay

No detectable degradation was observed after incubation of supercoiled plasmid DNA with Glycogen.

Labeled Oligonucleotide (LO) Assay

No detectable degradation after incubation of single-stranded or double-stranded radiolabeled oligonucleotides with Glycogen.

Ribonuclease Assay

No detectable degradation was observed after incubation of [³H]-RNA with Glycogen.

Protease Assay

No detectable degradation of protease substrate after incubation of FTC-casein with Glycogen.

Quality authorized by:

 Jurgita Zilinskiene

References

1. Tracy, S., Improved rapid methodology for the isolation of nucleic acids from agarose gels, Prep. Biochem., 11, 251-268, 1981.
2. Helms, C., A new method for purifying lambda DNA from phage lysates, DNA, 4, 39-49, 1985.
3. Hengen, P. N., Methods and reagents – Carriers for precipitating nucleic acids, TIBS, 21, 224-225, 1996.
4. Sambrook, J., Russell, D.W., Molecular Cloning: A Laboratory Manual, the Third edition, Cold Spring Harbor Laboratory Press, Cold Spring Harbor, New York, A8. 12-8.13, 2001.

LIMITED USE LABEL LICENSE: Internal Research and Development Use Only.

The purchase of this product conveys to the buyer the limited, non-exclusive, non-transferable right (without the right to resell, repackage, or further sublicense) to use this product for internal research and development purposes. No other license is granted to the buyer whether expressly, by implication, by estoppel or otherwise. In particular, the purchase of the product does not include or carry any right or license to use, develop, or otherwise exploit this product commercially and no rights are conveyed to the buyer to use the product or components of the product for purposes including but not limited to provision of services to a third party, generation of commercial databases or clinical diagnostics. This product is sold pursuant to authorization from Thermo Fisher Scientific and Thermo Fisher Scientific reserves all other rights. For information on purchasing a license for uses other than internal research and development purposes, please contact outlicensing@lifetech.com or Out Licensing, Life Technologies Inc., 5781 Van Allen Way, Carlsbad, California 92008.

PRODUCT USE LIMITATION

This product is developed, designed and sold exclusively for research purposes and in vitro use only. The product was not tested for use in diagnostics or for drug development, nor is it suitable for administration to humans or animals. Please refer to www.thermofisher.com for Material Safety Data Sheet of the product.

© 2017 Thermo Fisher Scientific Inc. All rights reserved. All trademarks are the property of Thermo Fisher Scientific Inc. and its subsidiaries.