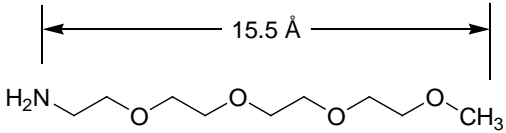
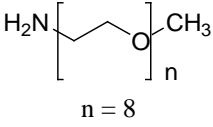
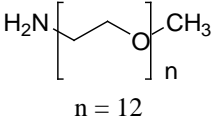
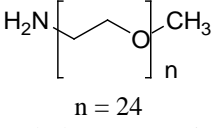


MA(PEG)_n Reagents

Methyl-amine PEGylation reagents

2074.2

Number	Description	
26110 26111	MA(PEG)₄ , 100mg MA(PEG)₄ , 1g Form: Colorless to pale yellow liquid Molecular Weight: 207.27 Spacer Arm: 15.5Å	 <p>Methyl-PEG₄-Amine</p>
26112 26113	MA(PEG)₈ , 100mg MA(PEG)₈ , 1g Form: Colorless to pale yellow liquid at ambient or solid at -20°C Molecular Weight: 383.48 Spacer Arm: 29.7Å	 <p>Methyl-PEG₈-Amine</p>
26114 26115	MA(PEG)₁₂ , 100mg MA(PEG)₁₂ , 1g Form: Colorless to pale yellow liquid at ambient or solid at -20°C Molecular Weight: 559.69 Spacer Arm: 43.9Å	 <p>Methyl-PEG₁₂-Amine</p>
26116 26117	MA(PEG)₂₄ , 100mg MA(PEG)₂₄ , 1g Form: White to pale yellow waxy solid Molecular Weight: 1088.32 Spacer Arm: 86.1Å	 <p>Methyl-PEG₂₄-Amine</p>

Storage: Upon receipt store product at -20°C protected from moisture. Product is shipped at ambient temperature.

Introduction

The methyl-PEG_n-amine (MA[PEG]_n) PEGylation reagents are methyl ether-terminated PEG amines that are used for modifying proteins or surfaces such as beads, nanoparticles and self-assembled monolayers. Modification of proteins adds polyethylene glycol (PEG) spacers, which impart increased water solubility, reduced immunogenicity of the labeled molecule and enhanced *in vivo* stability in solution.¹ Functionalization of solid surfaces, such as quantum dots, self-assembled monolayers and nanoparticles, with polyethylene glycol spacers significantly reduces nonspecific protein binding.²⁻⁷ MA(PEG)_n reagents used with CA(PEG)_n reagents in surface modification can form a hydrophilic “lawn” of methyl ether-terminated PEGs with periodic exposed carboxy-terminated PEGs. The exposed carboxy groups can be coupled to affinity ligands using the carbodiimide coupling reaction with EDC and sulfo-NHS.

Typical PEGylation reagents contain heterogeneous mixtures of different PEG chain lengths; however, our PEGylation reagents are homogenous compounds of defined molecular weight and spacer length, providing precision in optimizing modification applications.

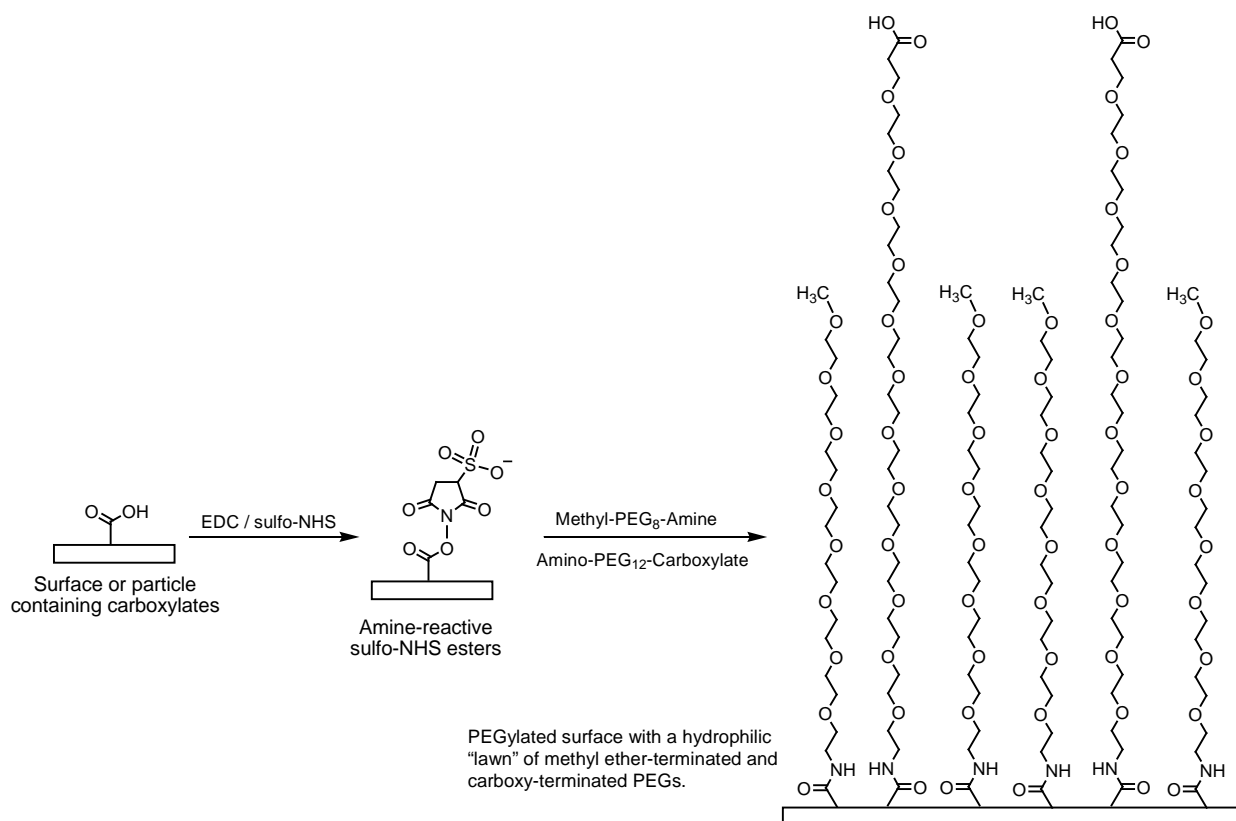


Figure 1. Two-step modification of surfaces with CA(PEG)_n-MA(PEG)_n reagents.

Important Product Information

- The MA(PEG)_n reagents are low-melting solids that are difficult to weigh and dispense. To facilitate handling, make a stock solution by dissolving the reagent in dry water-miscible solvent (e.g., DMF or DMSO).
- Store unused stock solution in a moisture-free condition (e.g., capped under and inert gas such as argon or nitrogen) at -20°C. Equilibrate reagent vial to room temperature before opening to avoid moisture condensation inside the container. To minimize air exposure, cap the stock solution with a septum and use a syringe to remove the solution.
- Use the MA(PEG)_n reagents (see Related Thermo Scientific Products) in combination with CA(PEG)_n reagents to modify surfaces and minimize nonspecific binding.
- Use non-amine-containing buffers at pH 7-9 such as PBS (20mM sodium phosphate, 150mM NaCl; pH 7.4) (Product No. 28372); 20mM HEPES; 100mM carbonate/bicarbonate; or 50mM borate. Do not use buffers that contain primary amines, such as Tris or glycine, which compete with acylation.
- The CA(PEG)_n-to-MA(PEG)_n ratio and the reagent mixture-to-surface carboxylic acid molar ratio in the reaction affects the number of carboxy groups modified on the surface and the number of new carboxylic acid residues available for further modification. Optimize these ratios to obtain the modification level needed for the specific application.

Procedure for Coupling CA(PEG)_n-MA(PEG)_n Mixtures to Carboxylated Surfaces

The following protocol, adapted from a procedure described by Grabarek and Gergely⁸ is a two-step coupling reaction using EDC and NHS or Sulfo-NHS. The CA(PEG)_n-MA(PEG)_n mixture is coupled to a carboxylated surface without exposing the carboxylic acids on CA(PEG)_n to EDC. The activation reaction requires quenching with a thiol-containing compound.

The activation reaction with EDC and Sulfo-NHS is most efficient at pH 4.5-7.2; however, the reaction of Sulfo-NHS-activated molecules with primary amines is most efficient at pH 7-8. For best results, perform the first reaction in MES buffer (or other non-amine, non-carboxy buffer) at pH 5-6, then raise the pH to 7.2-7.5 with phosphate buffer (or other non-amine buffer) immediately before reacting with the CA(PEG)_n-MA(PEG)_n mixture. Use DTT to quench the activation reaction. The conjugation reaction is quenched using hydroxylamine, Tris or glycine.

Materials Required

- Water-miscible organic solvent (molecular sieve-treated) such as dimethylsulfoxide (DMSO, Product No. 20684) or dimethylformamide (DMF, Product No. 20673) for preparing the reagent stock solution
- Small-volume, non-coring syringes for dispensing the reagent stock solution while minimizing exposure to air
- EDC (Product No. 77149)
- Activation Buffer: MES-buffered saline (0.1M MES, 0.5M NaCl; pH 6.0 or 0.1M MES, 0.9% NaCl; pH 4.7; Product No. 28390)
- Conjugation Buffer: Phosphate-buffered saline, PBS (20mM sodium phosphate, 0.15M NaCl; pH 7.2, Product No. 28372)
- NHS or Sulfo-NHS (Product No. 24500 and 24510, respectively)
- Dithiothreitol (DTT; Product No. 20290 or 20291)
- Hydroxylamine • HCl (Product No. 26103)

Procedure

1. Equilibrate EDC, NHS or sulfo-NHS, CA(PEG)_n, and MA(PEG)_n to room temperature before opening bottles.
2. Prepare CA(PEG)_n and MA(PEG)_n stock solutions by dissolving 100mg of each reagent (~100µL) in the desired amount of dry water-miscible solvent (e.g., DMF or DMSO).
3. Cap, store and handle stock solutions as directed in the Important Product Information Section.
4. Add appropriate amounts of EDC and NHS or sulfo-NHS to the appropriate amount of carboxylated surface in Activation Buffer and react for 15 minutes at room temperature.
5. Add DTT to quench the EDC.

Note: For surfaces that can be easily washed, the quenching step can be skipped and the surface washed with Coupling Buffer to remove any remaining EDC and NHS.

6. Add the CA(PEG)_n-MA(PEG)_n mixture prepared in Conjugation Buffer to the activated surface and react for 2 hours at room temperature.
7. To quench the reaction, add hydroxylamine or another amine-containing buffer. Hydroxylamine hydrolyzes non-reacted NHS on the solid surface and results in hydroxamate formation. Other quenching methods involve adding Tris, lysine, glycine or ethanolamine; however, these primary amine-containing compounds modify carboxyls.

Note: The newly introduced carboxy groups can be further modified by repeating Steps 4 and 5.

8. Add the desired amine-containing substrate, prepared in Coupling Buffer, to the activated surface and react for 2 hours at room temperature.
9. Quench the reaction as described in Step 7.

Related Thermo Scientific Products

26120	CA(PEG) ₄ , 100mg
26121	CA(PEG) ₄ , 1g
26122	CA(PEG) ₈ , 100mg
26123	CA(PEG) ₈ , 1g
26124	CA(PEG) ₁₂ , 100mg
26125	CA(PEG) ₁₂ , 1g
26126	CA(PEG) ₂₄ , 100mg
26127	CA(PEG) ₂₄ , 1g
28390	BupH™ MES Buffered Saline, 10 packs
28372	BupH Phosphate Buffered Saline, 40 packs
77149	EDC, 10mg
24500	NHS (N-hydroxy succinimide), 25g
24510	Sulfo-NHS (sulfo N-hydroxy succinimide), 500mg
20290	DTT, 5g
20291	DTT, No-Weigh™ Format, 48 tubes × 7.7mg
26103	Hydroxylamine, 25g

Cited References

1. Morar, A.S., *et al.* (2006). PEGylation of proteins: A structural approach. *BioPharm Int* April 34-46.
2. Prime, K.L. and Whitesides, G.M. (1991). Self-assembled organic monolayers: model systems for studying absorption of proteins at surfaces. *Science* **252**:1164.
3. Bentzen, E.L., *et al.* (2005). Surface modification to reduce non-specific binding of quantum dots in live cell assays. *Bioconjugate Chem* **16**:1488-94.
4. Lin, P-C., *et al.* (2006). Ethylene glycol-protected magnetic nanoparticles for a multiplexed immunoassay in human plasma. *Small* **2(4)**:485-9.
5. Zheng, M., *et al.* (2003). Ethylene glycol monolayer protected nanoparticles for eliminating nonspecific binding with biological molecules. *J Am Chem Soc* **125**:7790-1.
6. Verma, A. and Rotello, V.M. (2005). Surface recognition of biomacromolecules using nanoparticle receptors. *Chem Commun* **3**:303-12.
7. Kidambi, S., *et al.* (2004). Selective depositions on polyelectrolyte multilayers: self-assembled monolayers of m-dPEG acid as molecular template. *J Am Chem Soc* **126**:4697-03.
8. Grabarek, Z. and Gergely, J. (1990). Zero-length crosslinking procedure with the use of active esters. *Anal Biochem* **185**:131-5.

Products are warranted to operate or perform substantially in conformance with published Product specifications in effect at the time of sale, as set forth in the Product documentation, specifications and/or accompanying package inserts ("Documentation"). No claim of suitability for use in applications regulated by FDA is made. The warranty provided herein is valid only when used by properly trained individuals. Unless otherwise stated in the Documentation, this warranty is limited to one year from date of shipment when the Product is subjected to normal, proper and intended usage. This warranty does not extend to anyone other than Buyer. Any model or sample furnished to Buyer is merely illustrative of the general type and quality of goods and does not represent that any Product will conform to such model or sample.

NO OTHER WARRANTIES, EXPRESS OR IMPLIED, ARE GRANTED, INCLUDING WITHOUT LIMITATION, IMPLIED WARRANTIES OF MERCHANTABILITY, FITNESS FOR ANY PARTICULAR PURPOSE, OR NON INFRINGEMENT. BUYER'S EXCLUSIVE REMEDY FOR NON-CONFORMING PRODUCTS DURING THE WARRANTY PERIOD IS LIMITED TO REPAIR, REPLACEMENT OF OR REFUND FOR THE NON-CONFORMING PRODUCT(S) AT SELLER'S SOLE OPTION. THERE IS NO OBLIGATION TO REPAIR, REPLACE OR REFUND FOR PRODUCTS AS THE RESULT OF (I) ACCIDENT, DISASTER OR EVENT OF FORCE MAJEURE, (II) MISUSE, FAULT OR NEGLIGENCE OF OR BY BUYER, (III) USE OF THE PRODUCTS IN A MANNER FOR WHICH THEY WERE NOT DESIGNED, OR (IV) IMPROPER STORAGE AND HANDLING OF THE PRODUCTS.

Unless otherwise expressly stated on the Product or in the documentation accompanying the Product, the Product is intended for research only and is not to be used for any other purpose, including without limitation, unauthorized commercial uses, in vitro diagnostic uses, ex vivo or in vivo therapeutic uses, or any type of consumption by or application to humans or animals.

Current product instructions are available at www.thermoscientific.com/pierce. For a faxed copy, call 800-874-3723 or contact your local distributor.

© 2012 Thermo Fisher Scientific Inc. All rights reserved. Unless otherwise indicated, all trademarks are property of Thermo Fisher Scientific Inc. and its subsidiaries. Printed in the USA.