


Superspeed Rotors

Thermo Scientific A21-24x15c Rotor



ROTOR SPECIFICATIONS

Type	Fixed Angle
Material	Aluminum
Capacity	24 x 15 mL
Maximum Speed	21,500 rpm
Tube Dimensions Ø x L	121.5 x 16.5 mm
Angle	42/30°
K Factor	379
Net Weight	8.7 kg /19.1 lbs
Warranty	10 years
Certifications	

	RCF	RADIUS
Maximum	60,049 x g (Outer)	122 mm (Outer)
	63,049 x g (Inner)	122 mm (Inner)
Minimum	31,525 x g (Outer)	61 mm (Outer)
	31,525 x g (Inner)	44 mm (Inner)

ROTOR AND ACCESSORIES ORDERING INFORMATION

		THERMO SCIENTIFIC	
		SORVALL LYNX 6000	SORVALL LYNX 4000
A21-24x15c	Max. Speed	21,500 rpm	19,500 rpm
	Max. RCF	63,049 x g	51,865 x g

A21-24x15c Rotor Package includes:

Description	Cat. No.
A21-24x15c Rotor	75003004
Vacuum Grease	76003500
Lubricant	75003786
Instruction Manual	50136234
Retaining Ring Pliers	65614

Accessories:

Description	Cat. No.
Rotor Stand	75003711
Replacement O-Ring	75007003

 Biocontainment certification by Public Health England, Porton Down, UK.

LABWARE ORDERING INFORMATION

Tube Vol. (mL)	Vol. Fill (mL)	Description	Cat. No.	Qty. /Pkg.	Tubes /Rotor	Max. Speed (rpm)	Max. RCF ¹ (x g)	Dims ØxL (mm)	REQUIRED TUBE CLOSURES, ADAPTERS, TOOLS & ACCESSORIES			
									Type	Cat. No.	Qty. /Pkg.	Description
15.0	15.0	Nunc™ PP Conical Disposable Tube ¹	339651	500	24	11,600	18,500	–	Closure	Included	–	–
15.0	15.0	Greiner™ PP Conical Disposable Tube ¹	–	–	24	21,500	62,532	–	Closure	Included	–	–
15.0	15.0	Greiner Conical Disposable Tube ¹	–	–	24	21,500	62,532	–	Closure	Included	–	–
15.0	15.0	Falcon™ PP Conical Disposable Tube ¹	–	–	24	21,500	62,532	–	Closure	Included	–	–
15.0	15.0	Corning™ PP Conical Disposable Tube ¹	–	–	24	21,500	62,532	–	Closure	Included	–	–
15.0	15.0	Corning PET Conical Disposable Tube ¹	–	–	24	21,500	62,532	–	Closure	Included	–	–
15.0	15.0	Sarstedt™ PP Conical Disposable Tube ¹	–	–	24	21,500	62,532	–	Closure	Included	–	–

¹ Listed Max Speeds for Conical Disposable tubes may be higher than OEM tube specifications and may result in crazing and other cosmetic damage to the tube. Do not use these tubes more than once.

Refer to manufacturer's recommendations for product performance information.

© 2012, 2013, 2014, 2021 Thermo Fisher Scientific Inc. All rights reserved. Greiner is a registered trademark of Greiner. Falcon and Corning are registered trademarks of Corning. Sarstedt is a registered trademark of Sarstedt, Walter DBA. Amicon is a registered trademark of Millipore Corp. All other trademarks are the property of Thermo Fisher Scientific Inc. and its subsidiaries. Specifications, terms and pricing are subject to change. Not all products are available in all countries. Please consult your local sales representative for details.

DSCFGLYNXA2124X15C COL34223 0721

ThermoFisher
SCIENTIFIC