



PRODUCT INFORMATION

T4 DNA Polymerase

Pub. No. MAN0012013

Rev. Date 13 June 2016 (B.00)

Lot: _

Expiry Date: _

Components	#EP0061	#EP0062
	100 U	500 U
Concentration	5 U/ μ L	5 U/ μ L
5X Reaction Buffer	0.35 mL	2 \times 1 mL

Store at -20 °C

www.thermofisher.com

For Research Use Only. Not for use in diagnostic procedures.

Description

T4 DNA Polymerase, a template-dependent DNA polymerase, catalyzes 5'→3' synthesis from primed single-stranded DNA. The enzyme has a 3'→5' exonuclease activity, but lacks 5'→3' exonuclease activity.

Applications

- Blunting of DNA ends: fill-in 5'-overhangs or/and removal of 3'-overhangs (1, 2), see protocol on back page.
- Blunting of PCR products with 3'-dA overhangs (6).
- Synthesis of labeled DNA probes by the replacement reaction (3).
- Oligonucleotide-directed site-specific mutagenesis (4).
- Ligation-independent cloning of PCR products (5).

Source

E.coli cells with a cloned gene 43 of bacteriophage T4.

Molecular Weight

104 kDa monomer.

Definition of Activity Unit

One unit of the enzyme catalyzes the incorporation of 10 nmol of deoxyribonucleotides into a polynucleotide fraction in 30 min at 37 °C.

Storage Buffer

The enzyme is supplied in: 20 mM potassium phosphate (pH 7.5), 200 mM KCl, 2 mM DTT, and 50% (v/v) glycerol.

5X Reaction Buffer

335 mM Tris-HCl (pH 8.8 at 25 °C), 33 mM MgCl₂, 5 mM DTT, 84 mM (NH₄)₂SO₄.

Inhibition and Inactivation

- Inhibitors: metal chelators, nucleotide analogs 2(*p-n*-butylanilino)-dATP, N²-(*p-n*-butylphenyl)-dGTP), SH-blocking compounds (7).
- Inactivated by heating at 75 °C for 10 min.

Note

- The 3'→5' exonuclease activity of T4 DNA Polymerase is stronger on single-stranded DNA than on double-stranded DNA and greater (more than 200 times) than that of DNA Polymerase I, *E.coli* (1).
- Activity in Thermo Scientific buffers, % (in comparison to activity in assay buffer)

Buffers	Activity, %
for restriction enzymes: Thermo Scientific™ FastDigest™, FastDigest Green, O, R, 1X Thermo Scientific™ Tango™, 2X Tango, BamHI, EcoRI, Ecl136II, KpnI, PaeI, SacI B, G	100 75-100
for PCR buffers: <i>Taq</i> buffer with KCl and <i>Pfu</i> buffer <i>Taq</i> buffer with (NH ₄) ₂ SO ₄	50 100
RT buffers	100

CERTIFICATE OF ANALYSIS

Endodeoxyribonuclease Assay

No detectable degradation was observed after incubation of supercoiled plasmid DNA with T4 DNA Polymerase.

Quality authorized by:

 Jurgita Zilinskiene

(continued on back page)

Protocol for blunting of 5'- or 3'-overhangs

1. Prepare the following reaction mixture:

5X reaction buffer	4 μ L
Linear DNA or PCR product	1 μ g
dNTP Mix, 2 mM each (#R0241)	1 μ L (0.1 mM final concentration)
T4 DNA Polymerase	0.2 μ L (1 U)
Water, nuclease-free (#R0581)	to 20 μ L

2. Mix thoroughly, spin briefly and incubate at 11 °C for 20 min or at room temperature for 5 min.
3. Stop the reaction by heating at 75 °C for 10 min.

References

1. Sambrook, J., Russell D.W., Molecular Cloning: A Laboratory Manual, the third edition, Cold Spring Harbor Laboratory Press, Cold Spring Harbor, New York, 2001.
2. Ausubel, F.M., et al., Current Protocols in Molecular Biology, vol. 1, John Wiley & Sons, Inc., Brooklyn, New York, 3.5.11-3.5.12, 1994-2004.
3. Challberg, M.D., Englund, P.T., Specific labeling of 3'-termini with T4 DNA polymerase, Methods Enzymol., 65, 39-43, 1980.
4. Kunkel, I.A., et al., Rapid and efficient site-specific mutagenesis without phenotypic selection, Methods Enzymol., 154, 367-382, 1987.
5. Haun, R.S., et al., Rapid, reliable ligation-independent cloning of PCR products using modified plasmid vectors, BioTechniques, 13, 515-518, 1992.
6. Wang, K., et al., A simple method using T4 DNA polymerase to clone polymerase chain reaction products, BioTechniques, 17, 236-238, 1994.
7. Eun, H-M., Enzymology Primer for Recombinant DNA Technology, Academic Press, Inc., 1996

LIMITED USE LABEL LICENSE: Internal Research and Development Use Only.

The purchase of this product conveys to the buyer the limited, non-exclusive, non-transferable right (without the right to resell, repackage, or further sublicense) to use this product for internal research and development purposes. No other license is granted to the buyer whether expressly, by implication, by estoppel or otherwise. In particular, the purchase of the product does not include or carry any right or license to use, develop, or otherwise exploit this product commercially and no rights are conveyed to the buyer to use the product or components of the product for purposes including but not limited to provision of services to a third party, generation of commercial databases or clinical diagnostics. This product is sold pursuant to authorization from Thermo Fisher Scientific and Thermo Fisher Scientific reserves all other rights. For information on purchasing a license for uses other than internal research and development purposes, please contact outlicensing@lifetech.com or Out Licensing, Life Technologies Inc., 5791 Van Allen Way, Carlsbad, California 92008.

PRODUCT USE LIMITATION

This product is developed, designed and sold exclusively *for research purposes and in vitro use only*. The product was not tested for use in diagnostics or for drug development, nor is it suitable for administration to humans or animals.

Please refer to www.thermofisher.com for Material Safety Data Sheet of the product.

© 2016 Thermo Fisher Scientific Inc. All rights reserved. All trademarks are the property of Thermo Fisher Scientific Inc. and its subsidiaries.