



Publication Part No. LC5688.PPS Publication No. MAN0009621 Rev B.0


 Package contents	Catalog Numbers LC5688	Size: 250 µL
---	----------------------------------	------------------------

 Storage conditions	<ul style="list-style-type: none"> Store at –30°C to –10°C. Avoid repeated freezing and thawing. 	
---	--	--

 Required materials	<ul style="list-style-type: none"> Polyacrylamide gel(s) Electrophoresis apparatus for protein analysis Appropriate sample buffer (when silver staining) 	
---	---	--


 Selection guide	Unstained Protein Standards Go online to view related products.	
--	---	--

- HiMark™ Unstained Protein Standard allows you to accurately determine molecular weight of high molecular weight proteins.

 Product description	<ul style="list-style-type: none"> This standard consists of 9 protein bands ranging in molecular weight from 40–500 kDa. 	
--	--	--

- Storage buffer: 250mM Tris-HCl (pH 8.5), 0.5 mM EDTA, 50 mM DTT, 10% glycerol, 2% LDS, 0.2 mM Coomassie G-250, and 0.175 mM Phenol Red

- Use with NuPAGE® 3–8% or 7% Tris-Acetate Gels and Tris-Acetate SDS Buffer System under denaturing conditions. Click here to view our online [Protein Gel Selection Guide](#).

 Important guidelines	<ul style="list-style-type: none"> Visualize bands using Coomassie or silver staining techniques. 	
---	--	--

- DO NOT apply heat >37°C or add a reducing agent. HiMark™ Unstained Standard is supplied ready-to-use.

 Online resources	Visit our product page for additional information and protocols. For support, visit www.lifetechnologies.com/support .	
---	--	--



For Research Use Only. Not for use in diagnostic procedures.

18 March 2014

Protocol

- Thaw the standard at room temperature.
- Vortex gently to ensure the solution is homogeneous.

Note: If there is precipitate in the standard, thaw for 10–15 minutes at room temperature and vortex. If necessary, warm at 30°C (**do not** exceed 37°C).

- (*Silver staining only*) Dilute the standard 1:20 in 1X Sample Buffer.
- Load the standard on the gel (see table for recommended volumes).

Gel Type	Volume
Mini gel (1.0-mm thick)	5 µL
Mini gel (1.5-mm thick)	7 µL

Limited product warranty

Life Technologies Corporation and/or its affiliate(s) warrant their products as set forth in the Life Technologies' General Terms and Conditions of Sale found on Life Technologies' website at www.lifetechnologies.com/termsandconditions. If you have any questions, please contact Life Technologies at www.lifetechnologies.com/support.

Important licensing information

This product may be covered by one or more Limited Use Label Licenses. By use of this product, you accept the terms and conditions of all applicable Limited Use Label Licenses.

Disclaimer

LIFE TECHNOLOGIES CORPORATION AND/OR ITS AFFILIATE(S) DISCLAIM ALL WARRANTIES WITH RESPECT TO THIS DOCUMENT, EXPRESSED OR IMPLIED, INCLUDING BUT NOT LIMITED TO THOSE OF MERCHANTABILITY, FITNESS FOR A PARTICULAR PURPOSE, OR NON-INFRINGEMENT, TO THE EXTENT ALLOWED BY LAW. IN NO EVENT SHALL LIFE TECHNOLOGIES AND/OR ITS AFFILIATE(S) BE LIABLE, WHETHER IN CONTRACT, TORT, WARRANTY, OR UNDER ANY STATUTE OR ON ANY OTHER BASIS FOR SPECIAL, INCIDENTAL, INDIRECT, PUNITIVE, MULTIPLE OR CONSEQUENTIAL DAMAGES IN CONNECTION WITH OR ARISING FROM THIS DOCUMENT, INCLUDING BUT NOT LIMITED TO THE USE THEREOF.

© 2014 Thermo Fisher Scientific Inc. All rights reserved. All trademarks are the property of Thermo Fisher Scientific and its subsidiaries unless otherwise specified.

kDa

500
290
240
160
116
97
66
55
40



5 µL of HiMark™ Unstained Protein Standard on a NuPAGE® 3–8% Tris-Acetate Gel with Tris-Acetate SDS Running Buffer.

Stain: SimplyBlue™ SafeStain.

life
technologies™

For support, visit www.lifetechnologies.com/support.