NovaBright[™] Secreted Placental Alkaline Phosphatase (SEAP) Enzyme Reporter Gene Chemiluminescent Detection System 2.0

Catalog nos. N10577, N10578

Table 1. Contents and storage information.

Material	Composition	N10577	N10578	Storage	Stability				
NovaBright [™] Secreted Placental Alkaline Phosphatase (SEAP) Enzyme Reporter Gene Chemiluminescent Detection Kit 2.0 2–6°C Components									
Assay buffer (Component A)	Contains a proprietary mixture of non-placental alkaline phosphatase inhibitors	14 mL	60 mL	• 2–6°C • Protect from	When stored as directed, this kit is stable for 1 year.				
Reaction buffer (Component B)	Contains CSPD® substrate and Emerald-III™ luminescence enhancer	14 mL	60 mL	light • DO NOT FREEZE					
NovaBright [™] Secreted Placental Alkaline Phosphatase (SEAP) Enzyme Reporter Gene Chemiluminescent Detection Kit 2.0 ≤-20°C Components									
Control enzyme	0.3 ng/μL purified human placental alkaline phosphatase in 150 mM Tris (pH 7.8), 50 mM NaCl, 50% glycerol	50 μL	50 μL	• ≤-20°C	When stored as directed, this kit is stable for 1 year.				
Number of assays: Sufficient material is supplied for 192 (Cat. no. N10577) or 960 (Cat. no. N10578) microplate assays based on the									

Introduction

protocol below.

The NovaBright[™] Secreted Placental Alkaline Phosphatase (SEAP) Enzyme Reporter Gene Chemiluminescent Detection System 2.0 is a chemiluminescent reporter gene assay system designed for the rapid and sensitive detection of secreted placental alkaline phosphatase (SEAP) in cell culture media. SEAP is a reporter protein that is secreted into the cell culture medium and detected by testing aliquots of the medium, leaving the cells intact for further experimentation.^{1,2} SEAP is a truncated form of human placental alkaline phosphatase (PLAP). Detection of non-secreted placental alkaline phosphatase is also possible (see PLAP (Non-Secreted) Detection Assay Performed Directly in the Cell Culture Microplate).

Description of the System	The next-generation NovaBright [™] Secreted Placental Alkaline Phosphatase (SEAP)
	Enzyme Reporter Gene Chemiluminescent Detection System 2.0 incorporates CSPD*
	chemiluminescent substrate and the next generation Emerald-III [™] luminescence enhancer
	for high sensitivity and wide dynamic range. This new detection assay provides higher signal
	intensity and higher assay signal/noise performance than the original assay system. All the
	reagents in the new assay system are provided ready-to-use without the need for any reagent
	preparation steps, which allows fewer assay steps and shorter assay duration. The original
	NovaBright [™] Secreted Placental Alkaline Phosphatase (SEAP) Enzyme Reporter Gene
	Chemiluminescent Detection System ^{3,4} has been used for the detection of secreted placental
	alkaline phosphatase reporter enzyme in cell culture media ^{5,6} and for the quantitation of non-
	secreted placental alkaline phosphatase in cell and tissue extracts. ^{7,8}

The NovaBright[™] Secreted Placental Alkaline Phosphatase (SEAP) Enzyme Reporter Gene Chemiluminescent Detection 2.0 assay is simple and rapid. Depending on the expression system, secreted placental alkaline phosphatase can be measured in as little as 6 hours after cell transfection. Cell culture medium is incubated first with an assay buffer that contains differential inhibitors of non-placental alkaline phosphatase (serum and endogenous cellular alkaline phosphatases) and then with the reaction buffer containing CSPD[®] substrate and Emerald-III[™] light emission enhancer until maximum light signal is reached in approximately 20 minutes. The light emission kinetics provide a persistent glow signal, with light emission half-life of approximately 2.5 hours, enabling measurement over a wide time interval. Light emission is measured in a luminometer without the need for automated injection capability.

Chemiluminescent reporter assays for secreted placental alkaline phosphatase may be performed in cells that have endogenous non-placental alkaline phosphatase activity. Endogenous non-placental enzyme activity is significantly reduced with a combination of heat inactivation and differential inhibitors that do not significantly inhibit the transfected placental alkaline phosphatase. It is important to determine the level of endogenous enzyme in the medium of non-transfected cells to establish an assay background. Certain cell lines such as HeLa and others derived from cervical cancers may express placental alkaline phosphatase, which may produce high assay backgrounds when shed into the medium.⁹ Therefore, we do **not** recommend the use of secreted alkaline phosphatase as a reporter system in these cell lines.

ApplicationsThe NovaBright™ Secreted Placental Alkaline Phosphatase (SEAP) Enzyme Reporter Gene
Chemiluminescent Detection System has been used widely for the following applications.
The new version 2.0 system is compatible with all of these applications.

- Reporter gene assays measuring gene expression in established cell lines¹⁰ and in transfected primary cells,^{11,12} including as a gene knockdown/RNA interference read-out.¹³
- Viral functional assays, including viral gene expression assays,^{14,15} viral replication,^{16,17} viral fusogenicity,¹⁸ virus neutralization and viral-mediated cell-cell fusion,¹⁹ and viral infectivity.²⁰
- Assaying serum samples from transgenic, transfected, or viral vector-infected animals to measure SEAP levels in mouse,²¹ rat,²² marmoset,²³ monkey,²⁴ and pig sera,²⁵ and in chicken egg allantoic fluid.²⁶
- Sensitive detection of mouse SEAP protein (mSEAP),²¹ which has been developed for improved SEAP protein stability in transgenic mice.
- Measuring SEAP as a functional reporter for receptor-ligand binding assays with a SEAP-ligand chimera,²⁷ protease-mediated secretion,²⁸ and for secretion pathway activity,^{29,30} including as a functional assay to measure effects of siRNA-mediated protein knockdown on specific protein secretion pathways.³¹
- Measuring non-placental alkaline phosphatase as a biomarker.¹⁰

Materials Required but Not Provided	 Mammalian cells in adherent or suspension culture and culture medium SEAP reporter expression vector 96-well luminometer microplates (solid white) Microplate luminometer (single-mode or multi-mode) <i>Optional</i>: 96-well tissue culture-treated luminometer microplates (white with clear bottom)
Caution	Components A and B are irritating to eyes and skin. In case of contact with eyes, rinse immediately with plenty of water and seek medical advice. Wear suitable protective clothing and eye/face protection.
General Guidelines	 Read the entire Experimental Protocols section before proceeding. Perform all assays at room temperature, unless otherwise indicated. The NovaBright[∞] Secreted Placental Alkaline Phosphatase (SEAP) Enzyme Reporter Gene Chemiluminescent Detection System 2.0 assay system is not provided with SEAP reporter expression vectors. These are available through various commercial suppliers for cell culture transfection as well as <i>in vivo</i> delivery and expression
	 We recommend using a single-mode luminometer or a multi-mode detection instrument set for luminescence measurement to measure the light emission from 96- or 384-well microplates. If using a multi-mode detection instrument, make sure that wavelength filters are not in use. Taking luminescence measurements with wavelength filters in place is not necessary in the absence of excitation light and because it reduces signal intensity, we do not recommended it. The assay can also be performed in a tube luminometer or a 384-well microplate format, with reagent volumes scaled as appropriate. If sample dilution is desired, we suggest using sterile water or standard Tris-saline solutions available in the laboratory.
Controls	 Recommended positive and negative controls are listed below. Positive Control The control enzyme supplied with the kit, purified human placental alkaline phosphatase (hPLAP), provides a positive control for the assay reagents, as well as a means to determine the range of detection of the luminometer instrumentation, if desired. The control enzyme dilution curve is not intended (or accurate) for absolute quantitation of reporter enzyme concentrations, as the specific activity of the control hPLAP enzyme and the reporter enzyme may differ significantly. Additional positive controls can include the use of control SEAP constructs that provide constitutive expression of reporter enzyme as a positive control for cell transfection. The control hPLAP enzyme is approximately 0.3 ng/μL (0.75 U/mL) in 150 mM Tris (pH 7.8), 50 mM NaCl, 50% glycerol. Generate an enzyme dilution curve by serially diluting the control enzyme in the desired cell culture medium. Detection curve should range from approximately 1 pg/mL to 10 ng/mL. Determine the total accent background by accenting a volume of culture medium from
	Determine the total assay background by assaying a volume of culture medium from mock-transfected cells that is equivalent to that from transfected cells. In experiments involving induction of reporter expression, assay uninduced cells to determine the total

involving induction of reporter expression, assay uninduced cells to determine the total assay background. Experimental results are typically expressed as a ratio of signal from experimental cells to assay background (signal/background) or as results from induced cells normalized to results obtained with uninduced cells.

SEAP Detection Assay Using a Separate Assay Microplate		This procedure is for detecting secreted placental alkaline phosphatase (SEAP) in 25 μL of conditioned cell culture medium removed from the cell culture microplate.
	1.1	Equilibrate the Assay Buffer (Component A, 50 $\mu L/well$) and the Reaction Buffer (Component B, 50 $\mu L/well$) to room temperature.
	1.2	Add 25 μ L of cell culture medium (or control enzyme) to assay microplate wells.
	1.3	Add 50 μL of Assay Buffer per well, and incubate the plate at 65°C for 5 minutes.
	1.4	Add 50 μ L of Reaction Buffer per well, and incubate the plate for 20 minutes.
	1.5	Place microplate in a luminometer and measure for 0.1–1 second per well.
SEAP Detection Assay Performed Directly in the Cell Culture Microplate		This procedure is for detecting secreted placental alkaline phosphatase (SEAP) in 100 μ L (or less) of conditioned cell culture medium directly in the cell culture microplate.
		Note: The microplate must be white with a clear bottom or solid white.
:	2.1	Equilibrate the Assay Buffer (Component A, 50 μ L/well) and the Reaction Buffer (Component B, 50 μ L/well) to room temperature.
:	2.2	Add 50 μ L of Assay Buffer directly to each well of a culture microplate containing 100 μ L of cells and conditioned cell culture medium, and incubate at 65°C for 5 minutes.
:	2.3	Add 50 μ L of Reaction Buffer per well, and incubate for 20 minutes.
	2.4	Place the microplate in a luminometer and measure chemiluminescence for 0.1–1 second per well.
PLAP (Non-Secreted) Detection Assay Performed Directly in the Cell Culture Microplate	3.1 3.2 3.3 3.4	This procedure is for adherent cells expressing non-secreted placental alkaline phosphatase (PLAP), cultured in 96-well tissue culture-treated luminometer plates. If the culture medium is removed from the wells before performing the assay, heat inactivation is not necessary with this protocol. We recommend removing the culture medium before the assay, which increases the assay sensitivity ~20-fold, because the serum AP is responsible for the majority of the assay background. Remove the culture medium from the assay wells. Add 50 µL of Assay Buffer (Component A) per well, and incubate for 5 minutes at ambient temperature. Add 50 µL of Reaction Buffer (Component B) per well, and incubate for 20 minutes. Place the microplate in a luminometer and measure chemiluminescence for 0.1–1 second per well.

References

Gene 66, 1 (1988); 2. Meth Enzymol 216, 362 (1992); 3. BioTechniques 17, 172 (1994); 4. Clin Chem 42, 1542 (1996); 5. J Endocrin 193, 421 (2007);
 J Biol Chem 271, 7019 (1996); 7. Mol Cell Biol 15, 4272 (1995); 8. J Biol Chem 272, 791 (1997); 9. Int J Cancer 27, 637 (1981); 10. J Biol Chem 280, 25111 (2005); 11. Proc Natl Acad Sci USA 98, 3756 (2001); 12. J Neurosci 23, 4420 (2003); 13. Cell Research 15, 111 (2005); 14. J Virol 75, 3391 (2001); 15. J Biol Chem 281, 8205 (2006); 16. Antimicrob Agents Chemother 48, 2876 (2006); 17. J Gen Virol 87, 2297 (2006); 18. J Virol 77, 6823 (2003); 19. J Virol 76, 2075 (2002); 20. J Gen Virol 80, 1241 (1999); 21. J Gene Med 5, 773 (2003); 22. J Gene Med 6, 111 (2004); 23. J Virol 72, 6770 (1998); 24. Human Gene Therapy 13, 1611 (2002); 25. DNA Cell Biol 22, 807 (2003); 26. J Gen Virol 84, 781 (2003); 27. Mol Endocrinol 18, 150 (2004); 28. Mol Cell 2, 505 (1998); 29. J Exp Med 199, 1201 (2004); 30. Mol Biol Cell 19, 722 (2008); 31. J Cell Biol 179, 1123 (2007).

Product List Current prices may be obtained from our website or from our Customer Service Department.

Cat. no.	Product Name	Unit Size
N10577	NovaBright™ Secreted Placental Alkaline Phosphatase (SEAP) Enzyme Reporter Gene Chemiluminescent Detection Kit 2.0	
	192 assays	1 kit
N10578	NovaBright™ Secreted Placental Alkaline Phosphatase (SEAP) Enzyme Reporter Gene Chemiluminescent Detection Kit 2.0	
	960 assays	1 kit

Contact Information

Corporate Headquarters

5791 Van Allen Way Carlsbad, CA 92008 Phone: (760) 603-7200 Fax: (760) 602-6500

European Headquarters

Inchinnan Business Park. 3 Fountain Drive Paisley PA4 9RF, UK Phone: +44 (0) 141 814 6100 Fax: +44 (0) 141 814 6117 Email: euroinfo@invitrogen.com Technical Services: eurotech@invitrogen.com

Japanese Headquarters

LÖOP-X Bldg. 6F. 3-9-15, Kaigan Minato-ku, Tokyo 108-0022 Phone: 81 3 5730 6509 Fax: 81 3 5730 6519 Email: jpinfo@invitrogen.com

For country-specific contact information, visit www.invitrogen.com.

Limited Use Label License No. 223: Labeling and Detection Technology

The purchase of this product conveys to the buyer the non-transferable right to use the purchased amount of the product and components of the product in research conducted by the buyer (whether the buyer is an academic or for-profit entity). The buyer cannot sell or otherwise transfer (a) this product (b) its components or (c) materials made using this product or its components to a third party or otherwise use this product or its components or materials made using this product or its components for Commercial Purposes. The buyer may transfer information or materials made through the use of this product to a scientific collaborator, provided that such transfer is not for any Commercial Purpose, and that such collaborator agrees in writing (a) to not transfer such materials to any third party, and (b) to use such transferred materials and/or information solely for research and not for Commercial Purposes. Commercial Purposes means any activity by a party for consideration and may include, but is not limited to: (1) use of the product or its components in manufacturing; (2) use of the product or its components to provide a service, information, or data; (3) use of the product or its components for therapeutic, diagnostic or prophylactic purposes; or (4) resale of the product or its components, whether or not such product or its components are resold for use in research. Invitrogen Corporation will not assert a claim against the buyer of infringement of the above patents based upon the manufacture, use or sale of a therapeutic, clinical diagnostic, vaccine or prophylactic product developed in research by the buyer in which this product or its components was employed, provided that neither this product nor any of its components was used in the manufacture of such product. If the purchaser is not willing to accept the limitations of this limited use statement, Invitrogen is willing to accept return of the product with a full refund. For information on purchasing a license to this product for purposes other than research, contact Molecular Probes, Inc., Business Development, 29851 Willow Creek Road, Eugene, OR 97402, Tel: (541) 465-8300. Fax: (541) 335-0354.

Limited Warranty

Invitrogen (a part of Life Technologies Corporation) is committed to providing our customers with high-quality goods and services. Our goal is to ensure that every customer is 100% satisfied with our products and our service. If you should have any questions or concerns about an Invitrogen product or service, contact our Technical Support Representatives. All Invitrogen products are warranted to perform according to specifications stated on the certificate of analysis. The Company will replace, free of charge, any product that does not meet those specifications. This warranty limits the Company's liability to only the price of the product. No warranty is granted for products beyond their listed expiration date. No warranty is applicable unless all product components are stored in accordance with instructions. The Company reserves the right to select the method(s) used to analyze a product unless the Company agrees to a specified method in writing prior to acceptance of the order. Invitrogen makes every effort to ensure the accuracy of its publications, but realizes that the occasional typographical or other error is inevitable. Therefore the Company makes no warranty of any kind regarding the contents of any publications or documentation. If you discover an error in any of our publications, please report it to our Technical Support Representatives. Life Technologies Corporation shall have no responsibility or any special, incidental, indirect or consequential loss or damage whatsoever. The above limited warranty is sole and exclusive. No other warranty is made, whether expressed or implied, including any warranty of merchantability or fitness for a particular purpose.

The trademarks mentioned herein are the property of Life Technologies Corporation or their respective owners.

Copyright 2009, Life Technologies Corporation. All rights reserved. For research use only. Not intended for any animal or human therapeutic or diagnostic use.