# Thermo Scientific DiluFlux Pro-3 & DiluFlux Pro-5 Automated Gravimetric Dilutors

## Version 1.2

- May 2016

EN

FR

**Thermo**  
SCIENTIFIC

# Table Of Contents

<b>Chapter 1</b>	1. Introduction . . . . .	5
<b>Chapter 2</b>	2. Safety Precautions . . . . .	7
<b>Chapter 3</b>	3. Setting-Up . . . . .	9
	3.1 Unpacking . . . . .	9
	3.1.1 Unpacking DiluFlux Pro-3 dilutor . . . . .	9
	3.1.2 Unpacking DiluFlux Pro-5 dilutor . . . . .	10
	3.2 Unit description . . . . .	11
	3.2.1 Description of the DiluFlux Pro-3 dilutor . . . . .	11
	3.2.2 Description of the DiluFlux Pro-5 dilutor . . . . .	12
	3.3 Setting up . . . . .	13
	3.3.1 Fixed arm (DiluFlux Pro-3 dilutor only) . . . . .	13
	3.3.2 BagOpen™ bag opener . . . . .	14
	3.3.3 Booster kit (optional) . . . . .	16
	3.3.4 Distribution assembly . . . . .	18
	3.3.5 Adjusting the pump and insertion of the tubings . . . . .	19
<b>Chapter 4</b>	4. How to use . . . . .	21
	4.1 Safety precautions . . . . .	21
	4.2 Keyboard . . . . .	21
	4.3 Dilution . . . . .	22
	4.4 Distribution . . . . .	24
	4.5 Calibration . . . . .	25
	4.5.1 How to do a calibration . . . . .	25
	4.5.2 Metrological data . . . . .	25
	4.6 Traceability . . . . .	26
	4.6.1 Printer . . . . .	26
	4.6.1.1 Loading roll paper / label roll . . . . .	26
	4.6.1.2 How to use . . . . .	27
	4.6.1.3 Loading transfer film . . . . .	28
	4.6.2 Monitoring software . . . . .	29
	4.6.2.1 Download . . . . .	29
	4.6.2.2 Use of software . . . . .	30
	4.6.2.3 Stickers . . . . .	32
	4.7 Error messages . . . . .	33
<b>Chapter 5</b>	5. Accessories (Optional) . . . . .	35
<b>Chapter 6</b>	6. Maintenance . . . . .	41
	6.1 Pump heads maintenance . . . . .	42
	6.2 DiluFlux Pro-5 dilutor: cleaning of robotic arm . . . . .	43
	6.3 DiluFlux Pro-5 dilutor: nozzle holder cleaning . . . . .	43
<b>Chapter 7</b>	7. Troubleshooting . . . . .	45
<b>Chapter 8</b>	8. Appendix . . . . .	49
	8.1 Technical specifications . . . . .	49

# 1. Introduction

Microbiological analyses are becoming increasingly demanding: sample-taking, diluting and blending are essential factors ensuring the quality and reproducibility of the results.

The Thermo Scientific™ DiluFlux™ Pro range of Automated Gravimetric Dilutors will answer your sample dilution needs. Easy-to-use, with a removable drip tray to make cleaning easy, it is an accurate device in compliance with the latest international standards: ISO 7218, ISO 6887-1, FDA-BAM.

The DiluFlux Pro instrument has been designed to facilitate work in a microbiological safety cabinet: it has the lowest height of any dilutor on the market.

Thanks to its 3 position adjustable arm and its robotic arm, the DiluFlux Pro dilutor enables you to have free space to manipulate the samples above the BagOpen™ bag holder.

# 2

## 2. Safety Precautions

To reduce fire risks, electric shocks or injuries during the use of the DiluFlux Pro dilutor, please read the safety precautions and carefully keep them next to the unit.

The DiluFlux Pro dilutor must be placed on a horizontal, rigid, stable surface, to prevent it from falling on the user.

Power should be connected with an undamaged power cord provided with a ground terminal.

If the DiluFlux Pro dilutor is powered by an extension cord or a portable device with electrical outlets, make sure that these devices are not exposed to water splashes or moisture.

Do not install the DiluFlux Pro dilutor in a very humid place. This may cause an electric shock or short circuit.

Do not put your finger in the pump as they may be pinched.

If liquid gets inside the DiluFlux Pro dilutor through the ventilation cut-outs at the back of the unit, you must:

- Avoid any contact with the unit
- Turn the power off
- Contact Thermo Fisher Scientific Technical Support

Only qualified technicians, duly authorized by Thermo Fisher Scientific Technical Support, can open the device.

Do not use or store flammable liquids (kerosene, gasoline, etc.) in your DiluFlux Pro dilutor or nearby. These substances may evaporate and present a risk of fire or explosion.

Please carefully follow all these operating and maintenance instructions.

# 3

## 3. Setting-Up

### 3.1 Unpacking

Please refer to the following diagram showing how to unpack the unit.

Retain the packaging (foam blocks & cardboard box) in case you have to move or ship the unit for maintenance purposes.

#### 3.1.1 Unpacking DiluFlux Pro-3 dilutor

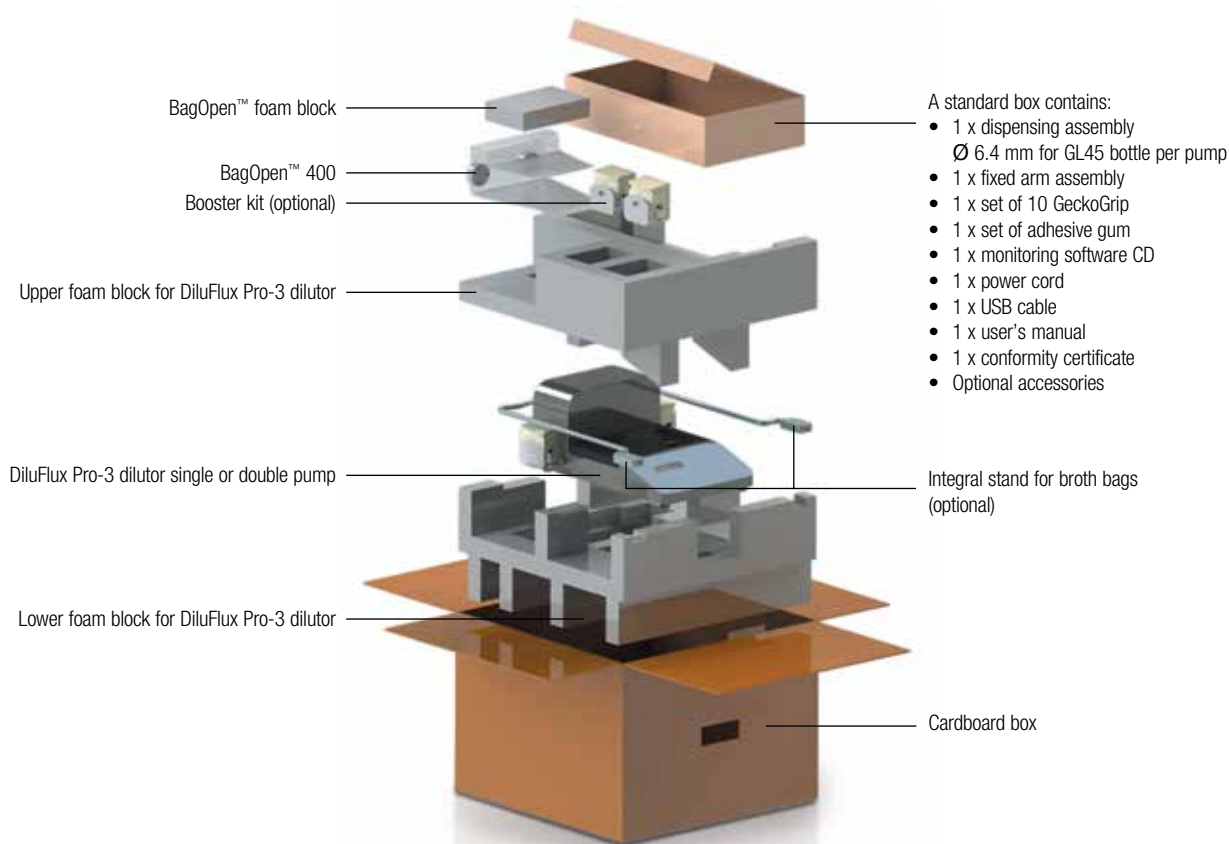
Figure 1

### 3.1.2 Unpacking DiluFlux Pro-5 dilutor

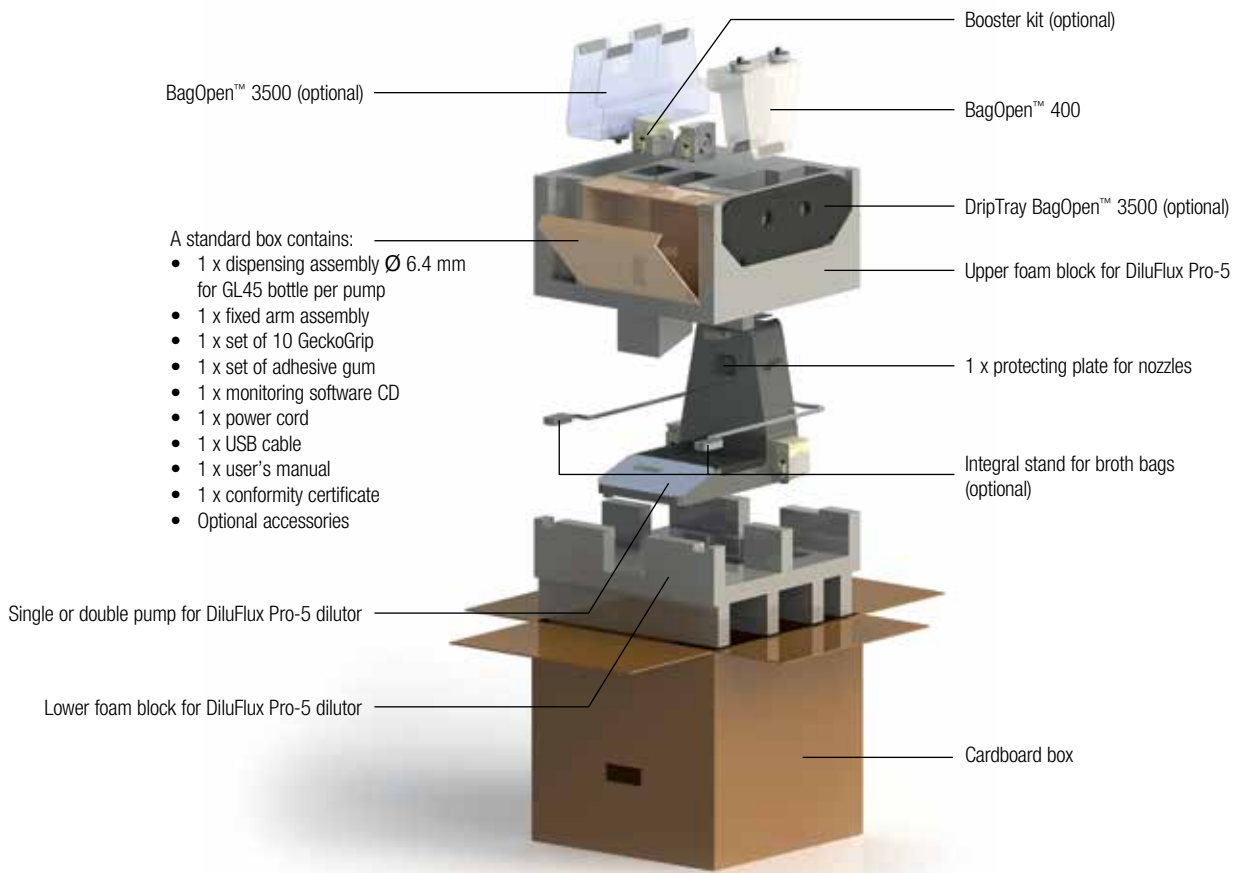
Figure 2

## 3.2 Unit description

### 3.2.1 Description of the DiluFlux Pro-3 dilutor

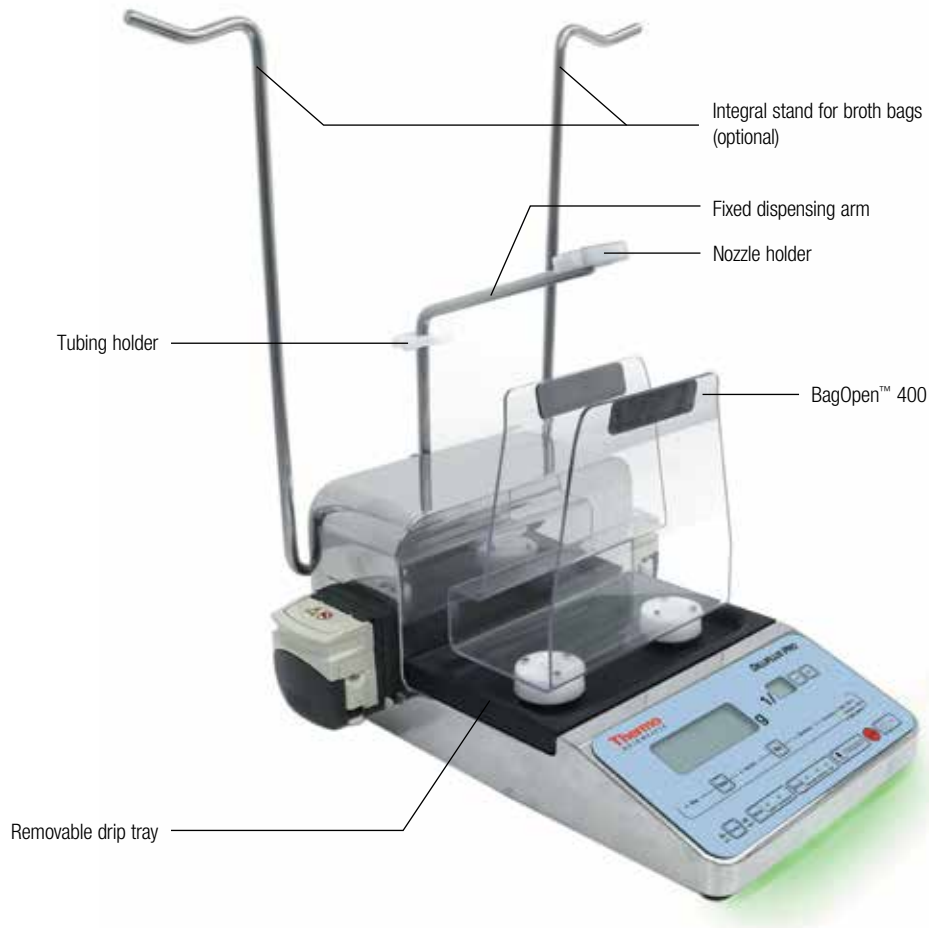
Figure 3

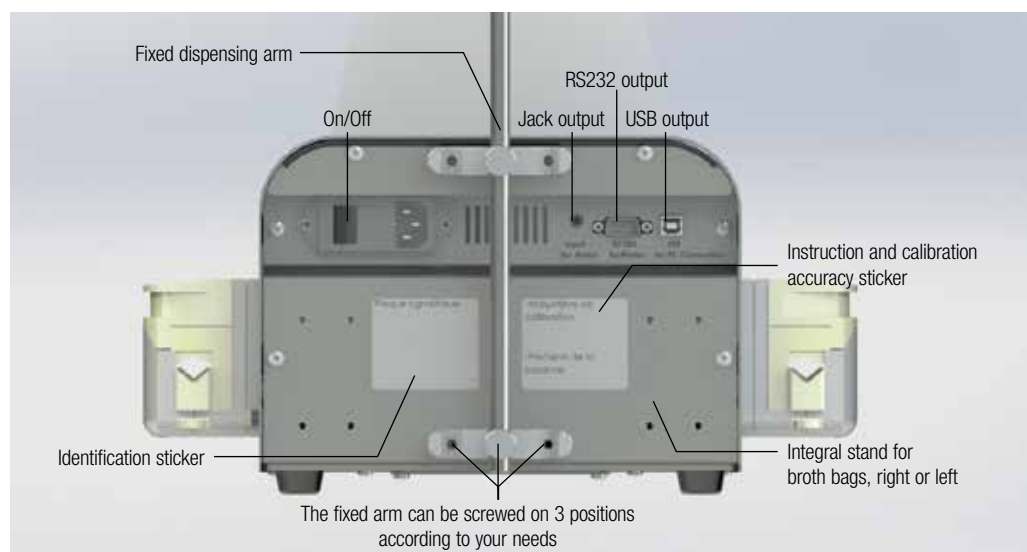
Figure 4—Back of the unit

- RS232 output: to connect a printer
- USB output: to connect a computer
- Jack output: to connect a foot pedal

### 3.2.2 Description of the DiluFlux Pro-5 dilutor

Figure 5

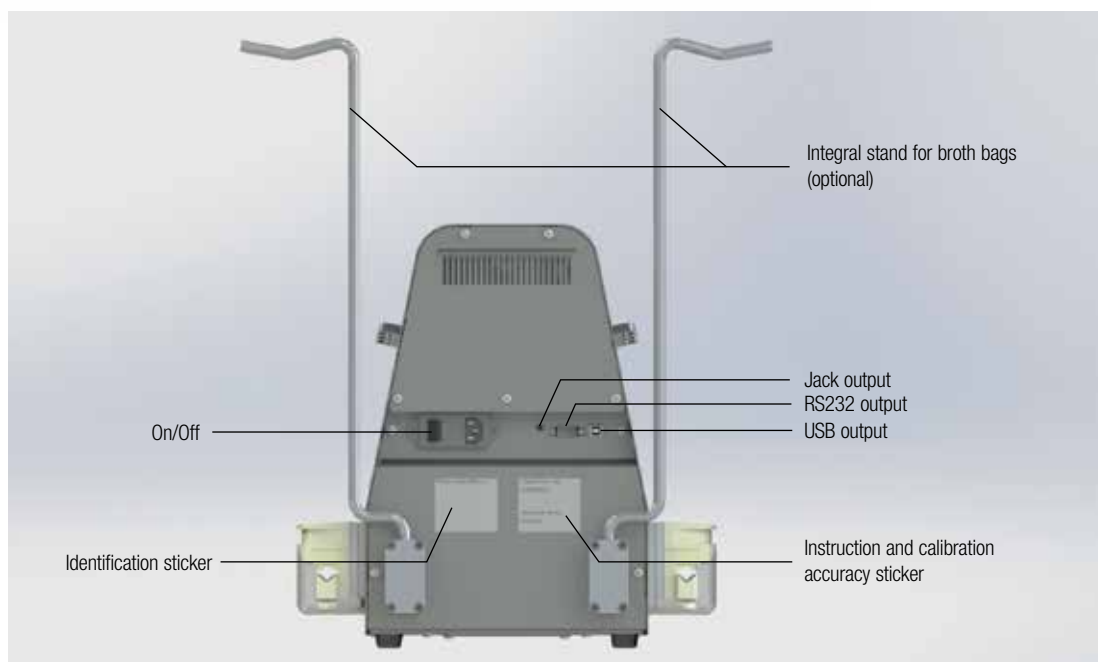
Figure 6—Back of the unit

- RS232 output: to connect a printer
- USB output: to connect a computer
- Jack output: to connect a foot pedal


## 3.3 Setting up

For best operation of your Diluflux Pro dilutor, you should:

- Install your DiluFlux Pro dilutor on a flat, stable, horizontal surface, which is important for the accuracy of the weighing platform. Also, keep it away from any vibrating devices.
- If you use a bottle as a diluent container, place it on the same horizontal surface, level or lower than the Diluflux Pro dilutor, to prevent the water from flowing after the pump stops. The bottle may be placed on the floor to save space on the bench.

### 3.3.1 Fixed arm (DiluFlux Pro-3 dilutor only)


Install the fixed arm at the back of the unit using the two screws provided with the arm. Choose between the three possible positions: a, b or c.

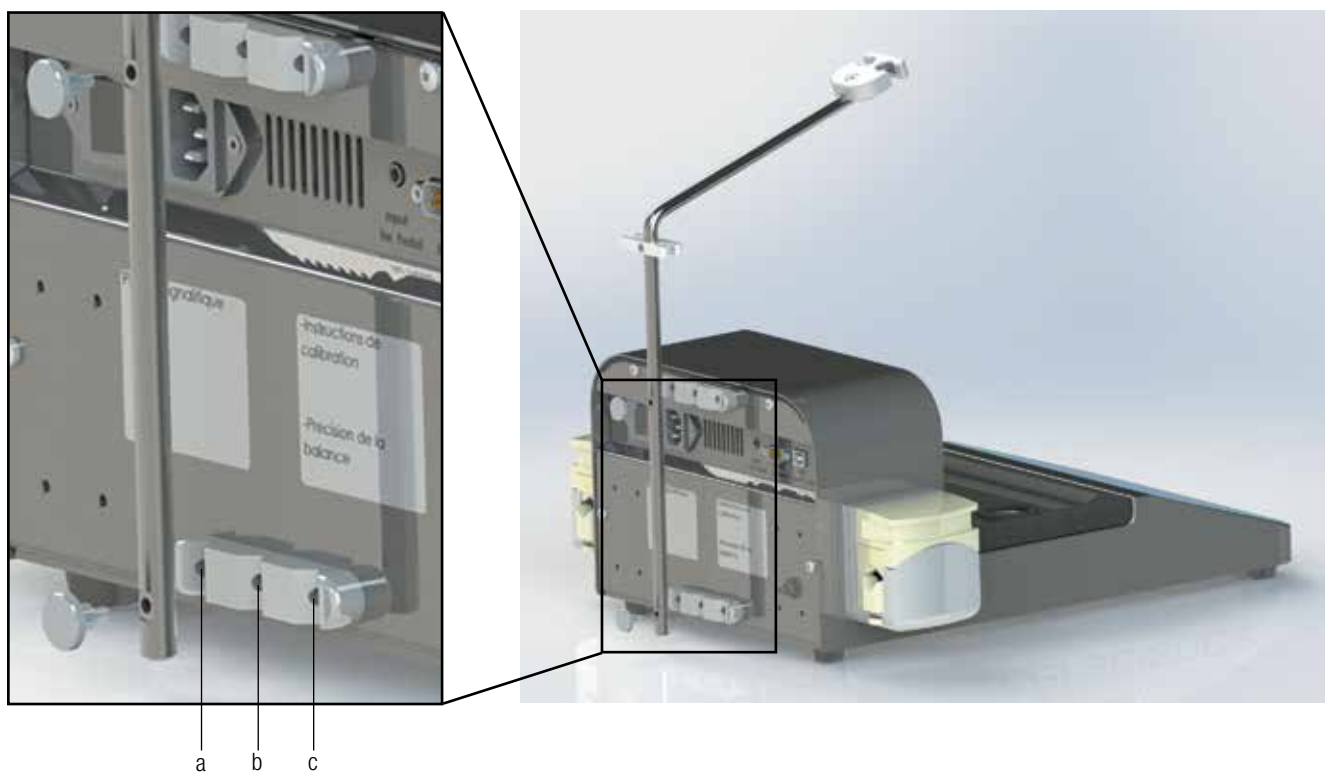
**Figure 7**—Screwing on the fixed arm

### 3.3.2 BagOpen™ bag opener

Clip the BagOpen™ bag holder into the two holes provided for this purpose. For DiluFlux Pro-5 dilutor only, also clip the magnetized plate for nozzle protection.

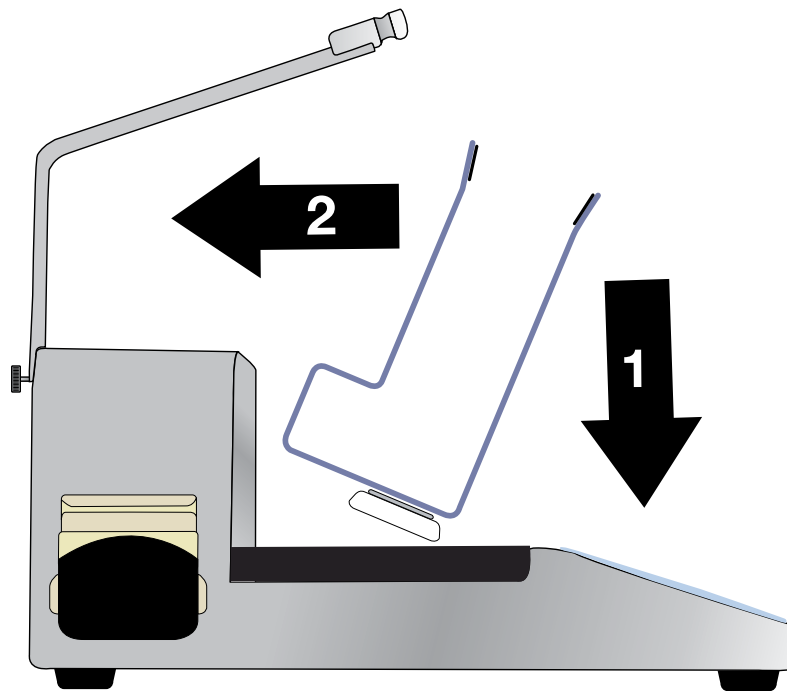
**Figure 8** –Setting-up of the BagOpen™ for DiluFlux Pro-3 dilutor

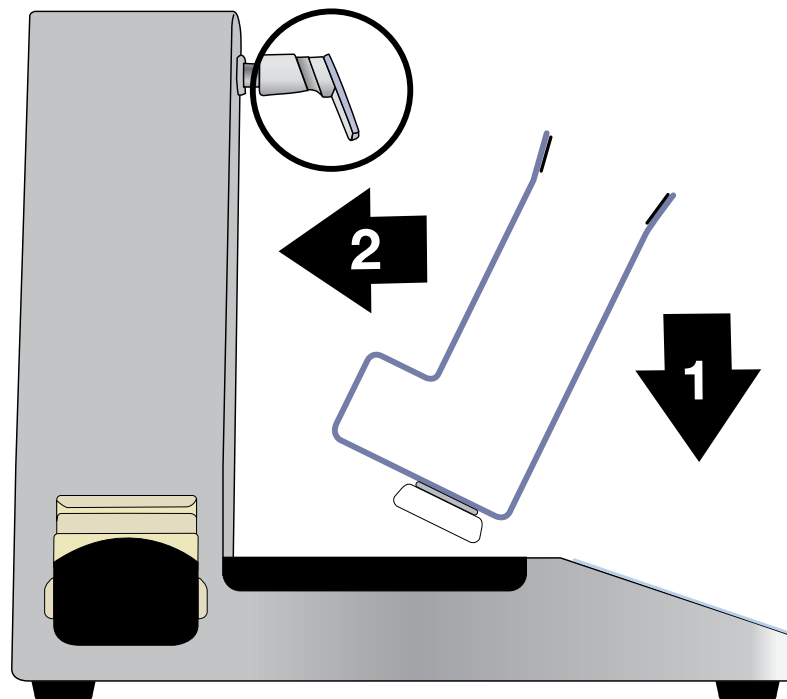
**Figure 9** –Setting-up of the BagOpen™ and the protecting plate for nozzles for the DiluFlux Pro-5 dilutor

**IMPORTANT:** The best way to remove the BagOpen™ is to tilt it towards you before lifting it up (see Figure 10 and Figure 11).

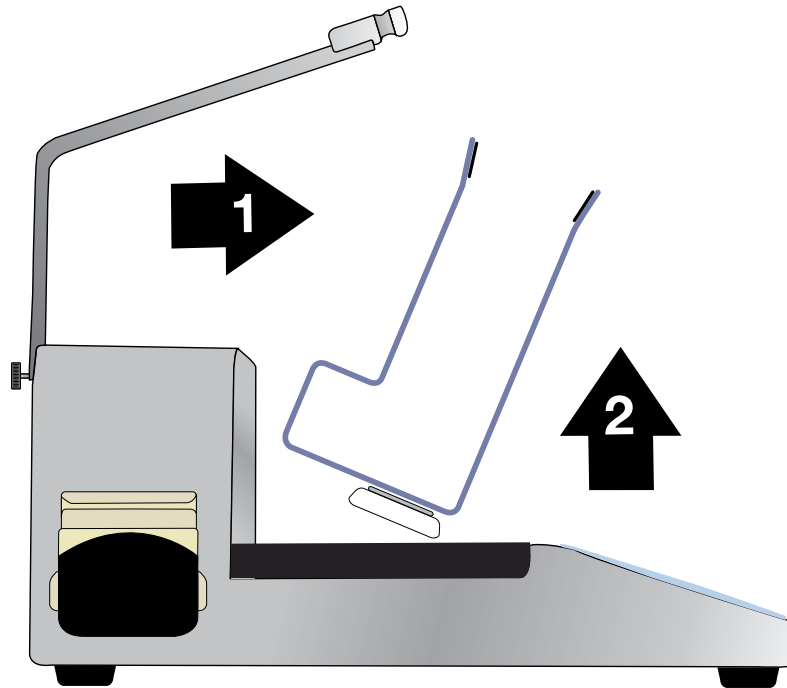**Figure 10** –Removal of the BagOpen™ for the DiluFlux Pro-3 dilutor

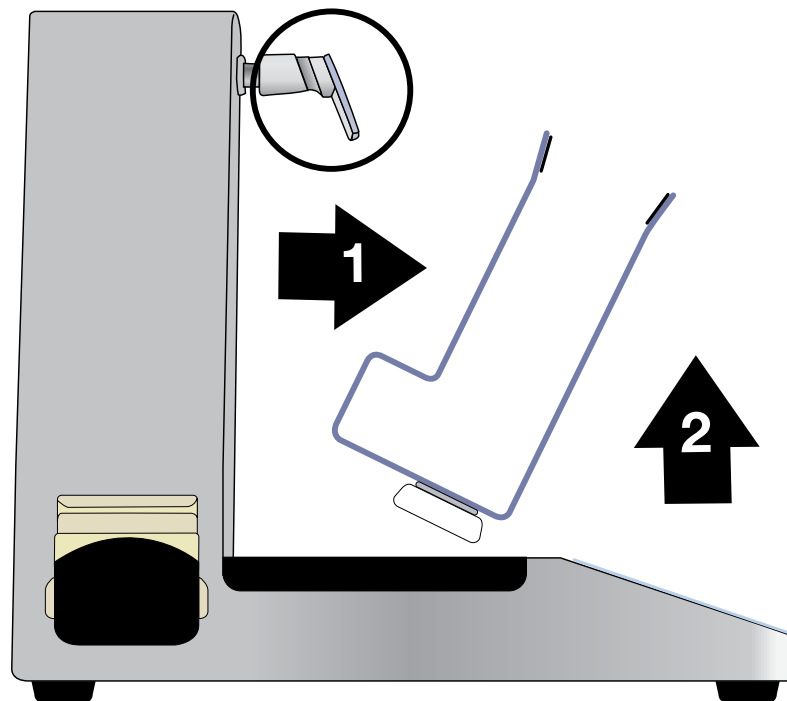**Figure 11** –Removal of the BagOpen™ and the protecting plate for nozzles for the DiluFlux Pro-5 dilutor

### 3.3.3 Booster kit (optional)

The booster kit comprises a pump and a pump extension.

To insert the pump extension between the pump and the pump holder:

1. Remove the tube from the pump.
2. Remove the pump: press the tab (see Figure 12) and at the same time turn it in the direction of the arrow. Release the pump.

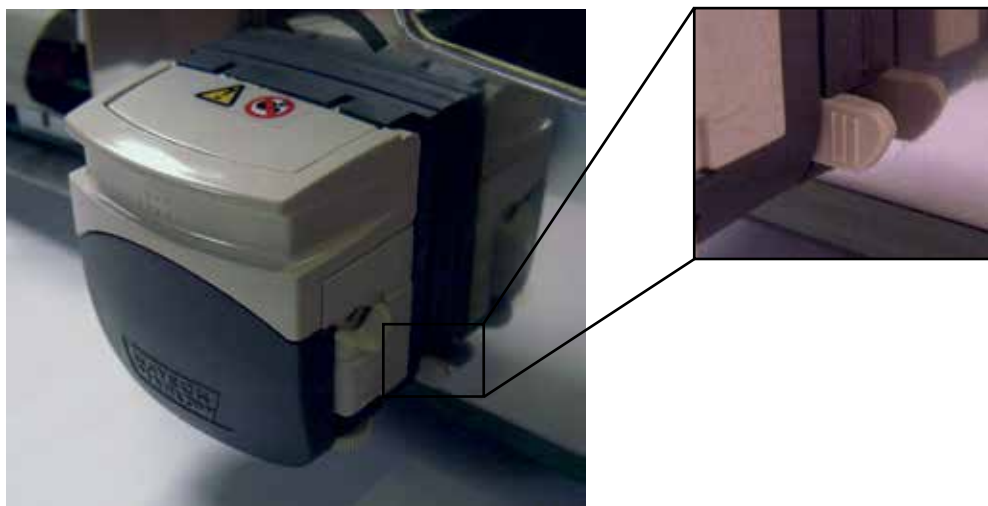
Figure 12

3. Clip the pump extension (see Figure 13) on to the pump and then open the two covers (see Figure 14):

Figure 13

Figure 14

4. Make sure that the rollers of the main pump and pump extension are NOT aligned but staggered in relation to each other:

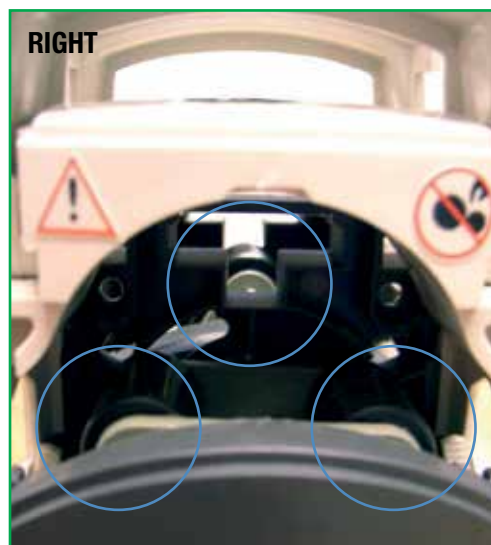
Figure 15

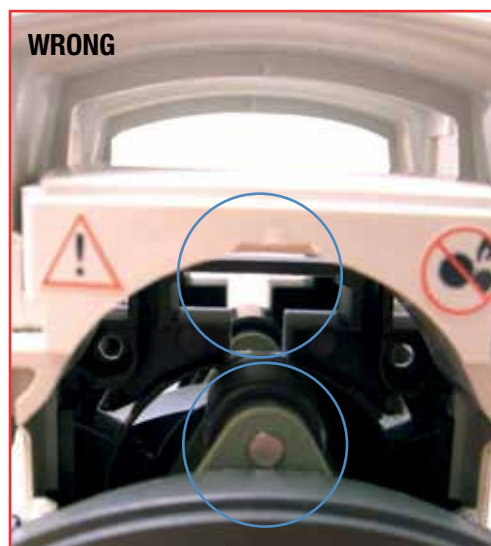
Figure 16

If WRONG, separate the pump from the pump extension by pushing the tab (see point 2 above). Move the roller wheels slightly to one side and clip the pump and the pump extension together again.

5. Replace the booster kit on the pump holder.

#### Assembly tubings for kit booster:

Remove the pump tubing from the current assembly and replace it with the double tubing for kit booster.

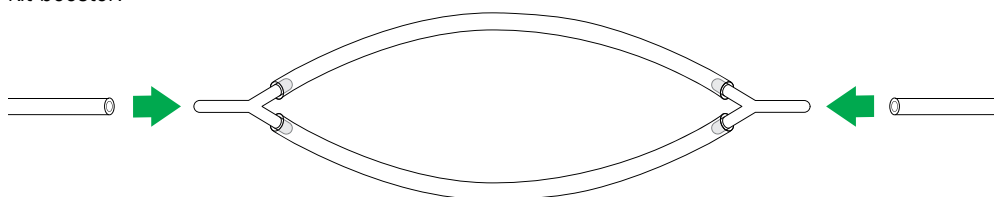
Figure 17

### 3.3.4 Distribution assembly

#### GL45 bottle:

Figure 18

- Place the tube weight at the bottom of the bottle and screw on the GL45 cap.
- Place the tubings paying attention to the direction of the flow. This direction is indicated by the arrows on the housing.
- If you are using tubings with an internal diameter smaller than 6.4 mm (delivered as standard), please refer to chapter 3.3.5 "Adjusting the pump and insertion of the tubings".
- Place the tubing in the tubing holder.
- Remove the protection cap and insert the nozzle into the nozzle holder (see Figure 18).

#### Broth bag:


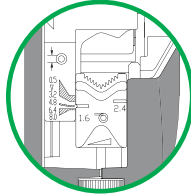
Figure 19

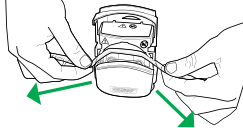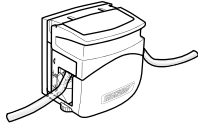
- Screw in the integral stand for broth bags (delivered with 4 screws and a screwdriver).
- Replace the GL45 cap, the small tubing and its weight with a trocar.
- Insert the trocar into the pouch and install the pouch on its stand.
- Place the tubings paying attention to the direction of the flow, which is indicated by the arrows on the housing of the Diluflux Pro.
- If you are using tubings with an internal diameter smaller than 6.4 mm (delivered as standard), please refer to chapter 3.3.5 “Adjusting the pump and insertion of the tubings”.
- Place the tubing in the tubing holder.
- Remove the cap and insert the nozzle into the nozzle holder (see Figure 19).

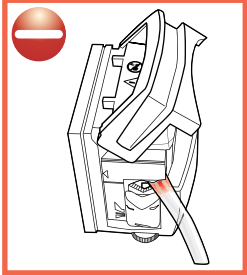
### 3.3.5 Adjusting the pump and inserting of the tubings

- 1 When the pump is opened, the factory setting is 6.4 (or depending on the tubing diameter).  
Turn the button on the side of the pump to obtain the right setting.

- 2 Insert the tubing into the pump.

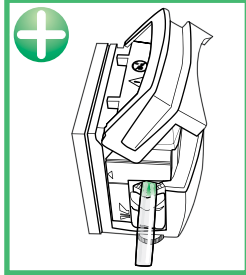
- 3 Keep the tubing stretched while you close the pump.

- 4

**WRONG**

The tubing is pinched.  
It can be pierced.

**RIGHT**

The tubing is not pinched.  
There is no risk.

# 4

## 4. How to use

### 4.1 Safety precautions

- The bag + sample + diluent must not exceed the authorized weight on the weighing platform. For the DiluFlux Pro-3: 3000 g, for the DiluFlux Pro-5 dilutor: 5000 g.
- Avoid placing any excess pressure on the weighing cell.
- Leave a free space of at least 5 cm behind the back of the DiluFlux Pro dilutor to allow for ventilation.
- Only use your finger to press the keyboard (no pens, sharp nails...)
- Use only accessories provided by Thermo Fisher Scientific (foot pedal, USB cable...)

**Before use, check that the drip tray is correctly positioned. If it is not so, it may cause malfunction of the weighing platform.**

- Be sure that there are no visible defaults (e.g.: unstable zero, zero drift...)
- Check that there is no obstacle which may interfere with the weighing cell.
- For the DiluFlux Pro-5 dilutor, check that there are no obstacles that could interfere with the robotic arm.

**Do not activate the robotic arm manually. The arm must not undergo any kind of shock or pressure.**

### 4.2 Keyboard

The interface has:

- 10 keys indicating an action
- 14 control LEDs indicating the current status of the process and the result of the dilution/distribution process.
- Two control screens showing the weight and dilution rate (dilution mode), the adjustable weight targeted by the user (distribution mode)


Figure 20




**Note: The unit automatically goes into sleep mode after 30 minutes of inactivity. Press the  key on the keyboard to resume activity.**
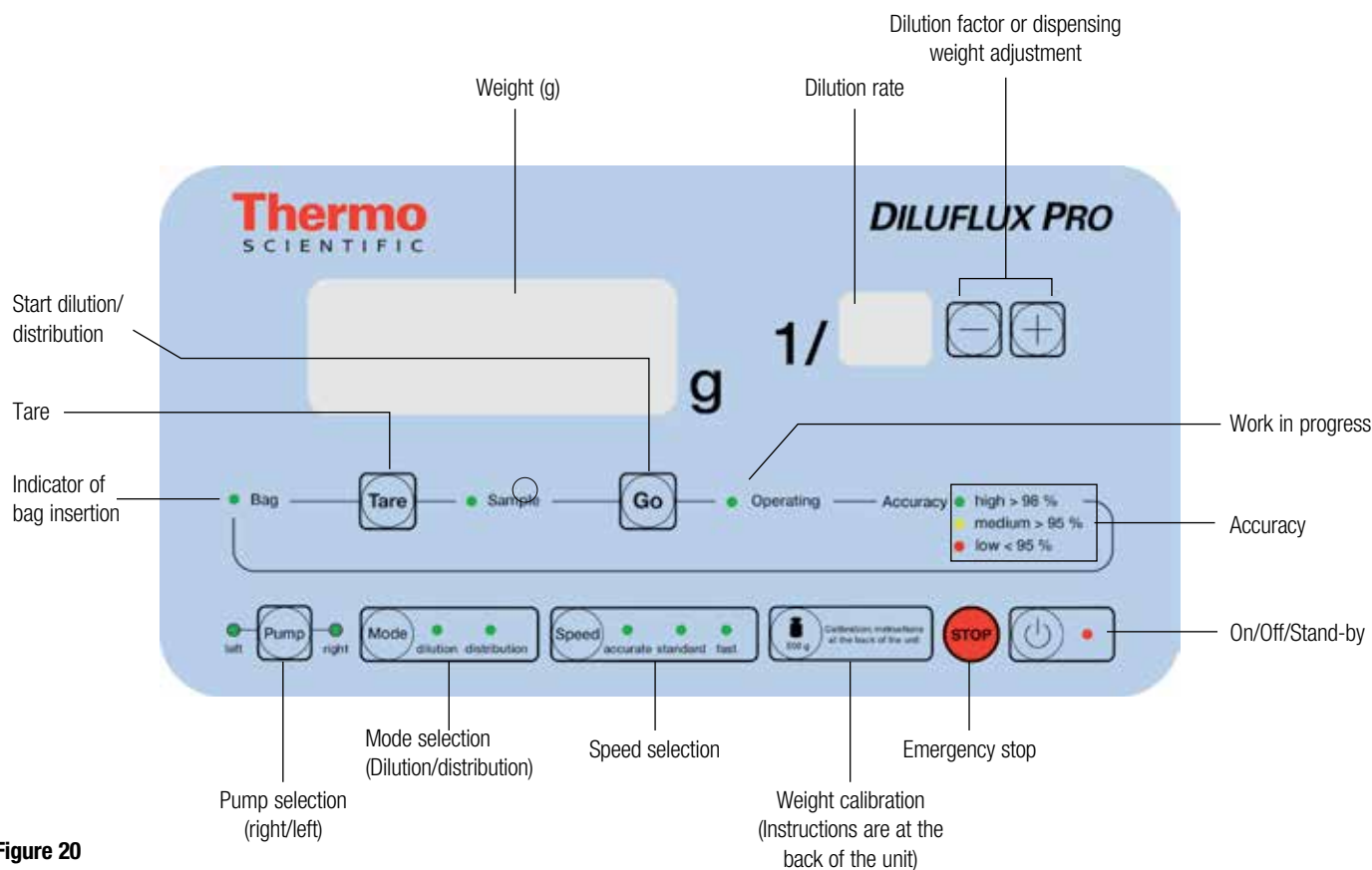
## 4.3 Dilution

Plug the power cord into an electrical outlet and flip the switch at the back of the unit in "1" position to turn it on.

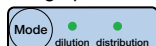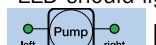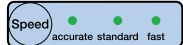A green light glows under the device, indicating that it is ready to function. DiluFlux Pro-5 dilutor: the arm is initialized.

**To ensure greater accuracy of the weighing platform, it is recommended to wait 10 minutes before the first use so that the weighing cell reaches room temperature. Then, go to dilution mode by pressing the  key, tare by pressing , put a certified standard weight of 500 g on the BagOpen™ and check that the displayed weight corresponds to  $500 \pm 0.5$  g. If not, proceed to a calibration (see chapter 4.5 "Calibration").**

- Place the BagOpen™ on the unit. Insert a sterile bag.
- With the  key, select the "Dilution" mode. Two beeps will sound. The "Dilution" LED should light up.
- With the  key, select the pump (right or left).
- With the  key select one of the 3 modes:
  - FAST: Fast flow with lower accuracy
  - STANDARD: Standard flow – standard accuracy
  - ACCURATE: Slow flow for optimal accuracy

**The ACCURATE mode is recommended to prime the DiluFlux Pro dilutor to guarantee an accuracy superior to 98% for the first dilution, after opening the pumps, changing the tubings or changing the container.**

- Select your dilution rate with the 2 keys  .

**NOTE: The Dilution/Distribution mode, the selected pump, the dilution rate and the chosen speed are memorized even when the unit is switched off.**

- Place the bag in the BagOpen™. Press the  key. A beep will sound.
- Once the LED “Place sample” is lit, place the sample in the bag. The weight of the sample shows up on the screen. Check the color of the light under the unit:
  - Blue indicates that the weighing cell is not stable. Wait for the unit to stabilize before proceeding.
  - Green indicates that the weighing platform is stable. You may proceed with your operation.
- Press the  key or press on the pedal to start the dilution. The LED “in progress” lights up, indicating that the dilution is proceeding. The lights under the DiluFlux Pro dilutor become orange and will flash slowly. For the DiluFlux Pro 5 dilutor the robotic arm moves simultaneously forward.

The dilution can be stopped anytime by pressing the  key. A beep will sound.

To resume the dilution, press  key.

It is possible to repeat this operation as many times as necessary, during the dilution. You may, during these interruptions, change the diluent, the tubings, the pump, (right/left) and speed before ending the current dilution.

To abort the run and start again, press a second time on the  key. Three beeps will sound.

- Once the dilution is complete, a beep will sound. The final weight is displayed and one of the 3 LEDs “high” (green), “medium” (orange) or “low” (red) indicates the obtained accuracy of the dilution. The lights under the unit also flash quickly showing the accuracy with a color code. For the DiluFlux Pro 5 dilutor, the robotic arm goes backward.
- To repeat the operation again with the same type of bag, just press the  key again once the bag is in the BagOpen™.

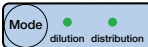
This functionality enables automatic tare of the bag by memorizing the weight of the empty bag. It is very useful for successive dilutions with bags of the same weight: it enables you to prepare a set of samples in their bags and dilute them sequentially. Please check that the weight of the bags is always the same.

## 4.4 Distribution

Plug the power cord in to an electrical outlet and put the switch at the back of the unit to the “I” position to turn it on.

A green light illuminates under the unit, indicating it is ready for use. DiluFlux Pro-5 dilutor: the arm initializes itself.

**To ensure greater accuracy of the weighing platform, it is best to wait 10 minutes before the first use, so that the weighing cell adjusts to room temperature. Next, switch to dilution mode by pressing the  key, tare by pressing the  key, place a 500 g calibration weight on the BagOpen™ and check that the displayed weight corresponds to 500 ± 0.5 g. If not, proceed to a calibration (see chapter 4.5 Calibration).**

- Place the BagOpen™ on the unit. Insert a sterile bag. If you are using another container, remove the BagOpen™ and place the container on the suitable support.
- With the  key select the “Distribution”. Two beeps will sound. The “Distribution” LED should be on.

In the “Distribution” mode, the small display, at the top right, is off.

- The main display shows the target weight. The target weight value can be changed by pressing the   keys.
- Select the pump (right or left) by pressing the  key.
- With the  key, select one of the 3 modes: accurate, standard or fast.
  - FAST: Fast flow with lower accuracy
  - STANDARD: Standard flow – standard accuracy
  - ACCURATE: Slow flow for optimal accuracy

**The ACCURATE mode is recommended to prime the DiluFlux Pro dilutor to guarantee an accuracy superior to 98% for the first dilution, after opening the pumps, changing the tubings or container.**

**NOTE : The chosen Dilution/Distribution mode, the selected pump, the amount of liquid to distribute and the chosen speed are memorized even after turning off the unit.**

- Press the  key or press the foot pedal to start the dispensing. The “in progress” LED lights up, indicating that the dispensing is in progress. The lights under the DiluFlux Pro- become orange and will flash slowly. For the DiluFlux Pro-5 dilutor the robotic arm will move simultaneously forward.

The distribution can be interrupted anytime by pressing  key. A beep will sound.

To resume the distribution, press the  key.

It is possible to repeat the operation as many times as necessary during the dispensing. You may, during these interruptions, change the diluent, the tubings, the pump (right/left) and speed before ending the current dispensing.

To abort the run and start again press again on . Three beeps will sound.

- Once the dispensing is over, a beep will sound. The final weight is displayed and one of the three LEDs “high” (green), “medium” (orange) or “low” (red) indicates the accuracy of the dispensing. The lighting under the unit flashes also quickly with the color of the accuracy. For the DiluFlux Pro-5 dilutor the robotic arm goes backwards automatically.
- To do the operation again with the same kind of container, just place the container and press the  key again.

## 4.5 Calibration

### 4.5.1 How to do a calibration

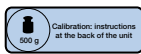
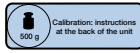
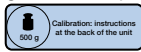
A calibration may be necessary if, after taring, the displayed weight does not show the following level of accuracy:

- $\pm 0.05$  g for weights inferior to 100 g
- $\pm 0.1$  % for weights superior to 100 g

Please find the calibration instructions at the back of the unit.

- Put the Diluflux Pro dilutor into “Dilution” mode

**ATTENTION: To calibrate the unit, use a certified weight of 500 g. The unit must NOT be calibrated with another weight. Thermo Fisher can provide you with a certified COFRAC weight.**

- Press the  key for two seconds.
- “0.00 g” flashes on the screen and three beeps will sound.
- Check that the BagOpen™ or the platform is clean and empty and press the  again.
- One beep sounds and then “500.0 g” flashes on the screen. Carefully place your standard 500 g weight on the platform.
- Press the  key again. The DiluFlux Pro dilutor beeps 3 times.
- Remove the standard weight, the DiluFlux Pro dilutor is calibrated and ready for use.

### 4.5.2 Metrological data

- Minimal weight of the sample: 1 g
- Maximum authorized weight: 3000 g for DiluFlux Pro-3 dilutor  
5000 g for DiluFlux Pro-5 dilutor
- Ideal functioning temperature for optimal accuracy of the weighing platform: 10°C / + 40°C
- Accuracy of the unit from 1 to 100 g:  $\pm 0.05$  g

- Accuracy of the unit from 100 to 5000 g:  $\pm 0.1 \%$

## 4.6 Traceability

### 4.6.1 Printer

Zebra™ printer is optional (Ref: DB4102A) and can be connected directly to DiluFlux Pro dilutor via the RS232 port on the back of the unit. It is delivered with a RS232 cable and a power cord.

Before using the printer, you need to load paper or labels.

#### 4.6.1.1 Loading roll paper / label roll

- Open the printer by pulling the two green buttons (on the sides of the printer) toward the front of the printer (see Figure 21).
- Open the paper roll holders. Pull the paper guides open with your free hand and place the paper roll on the roll holders and release the guides. Orient the paper roll so that its printing surface will face up as it passes over the platen (drive) roller.
- Pull the paper so that it extends out of the front of the printer. Make sure that the roll turns freely. The roll must not sit in the bottom of the paper compartment. Verify that the printing surface is facing up (see Figure 22).
- Push the paper under both of the media guides. (Figure 23)
- Close the printer. Press down until the cover snaps closed.

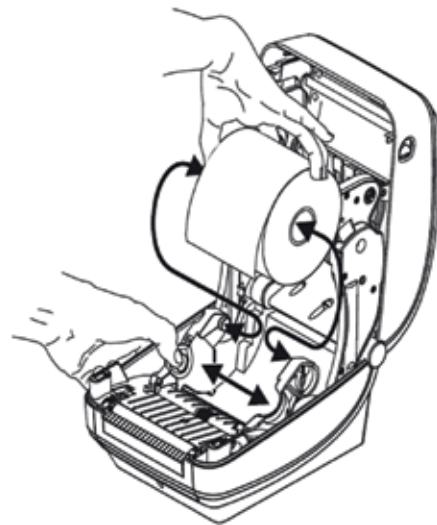
Figure 21

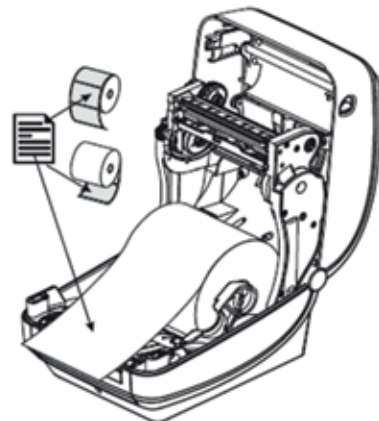
Figure 22


Figure 23

### 4.6.1.2 How to use

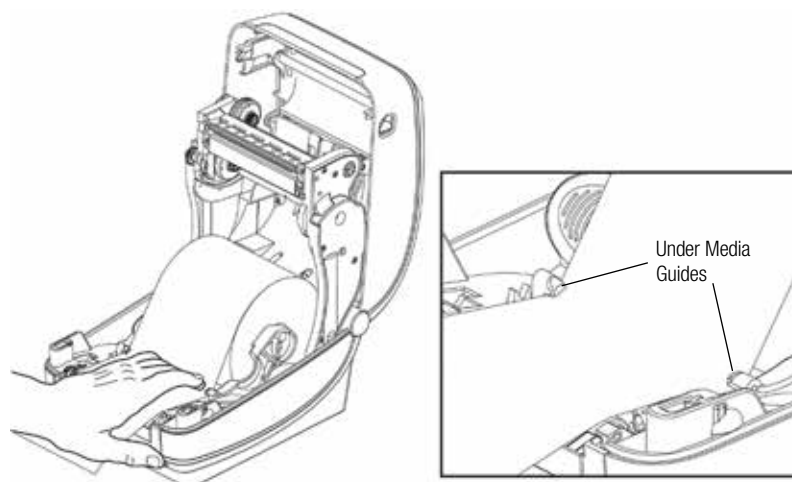Once it is connected to the Diluflux Pro, plugged to the mains and loaded with paper or labels, the printer issues a report on paper or label after each dilution or distribution.

#### WARNING:

**The DiluFlux Pro dilutor is set by default for paper rolls. For label rolls, a reconfiguration of the DiluFlux Pro dilutor is necessary.**

**Put the DiluFlux Pro dilutor in standby mode by pressing the standby key . Then press successively the ,  and  keys.**

**Turn the DiluFlux Pro dilutor back on by pressing the standby key .**

**The configuration is saved on the machine. If you want to switch back to rolls of continuous paper, put the machine in standby mode with the  key. Then press successively the ,  and  keys.**

**Turn the DiluFlow back on by pressing the standby key .**  
**The configuration is saved on the machine.**

Dilution information printed on the report:

- Right or left pump
- Dilution factor
- Sample weight
- Total weight
- Dilution accuracy

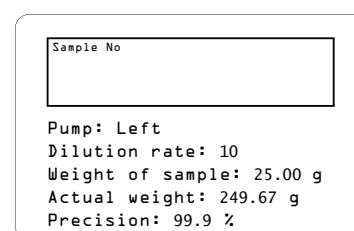
Figure 24

Dispensing information printed on the report:

- Right or left pump
- Target weight
- Dispensed weight
- Dispensing accuracy


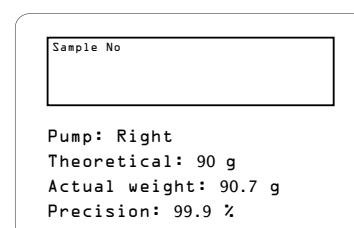
Figure 25

**NOTE: More information may be printed if using the monitoring software (See 4.6.2.3 Stickers).**

### 4.6.1.3 Loading transfer film

**NOTE: The printer is delivered with a transfer film in it. This transfer film can be ordered separately for future needs (Ref. DB4101A).**

Prepare the ribbon by removing its wrapping and pulling its adhesive strip free.

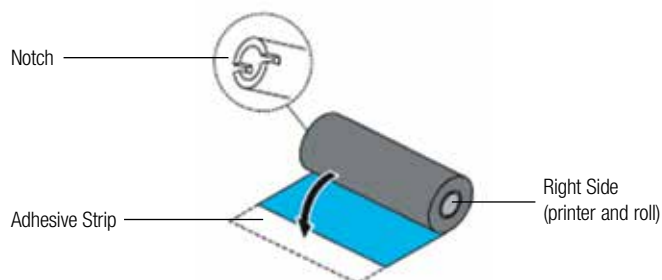
**Figure 26**

- With the printer open, place a new ribbon roll on printer's lower ribbon supply spindles. Rotate the roll until the notches align and lock into the left side of the supply hub. (See Figure 27)
- Place an empty ribbon core on printer's take-up spindles. Rotate the ribbon core until the notches align and lock into the left side of the take-up hub.
- You can find your first ribbon take-up core in the packing box. Subsequently, use the empty supply core to take up the next roll of ribbon. (See Figure 28)
- Pull the transfer ribbon's leader off the roll and attach it with the adhesive strip on the leader to the empty ribbon core on the supply spindle. Center the ribbon on the ribbon core. (See Figure 29)
- Rotate the thumb wheel on the left side of the supply spindle until the ribbon is pulled tight across the print head. (See Figure 30)

**Figure 27**

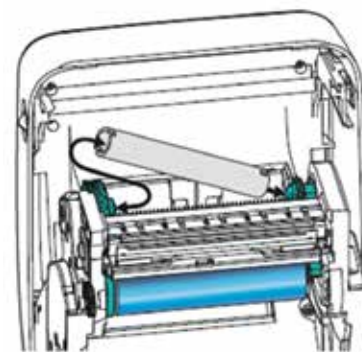**Figure 28**

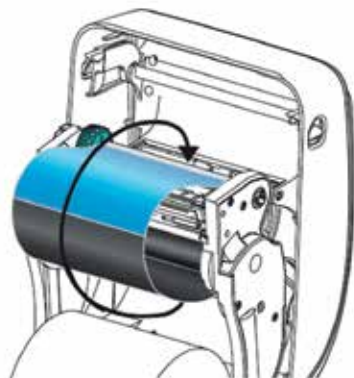
Figure 29

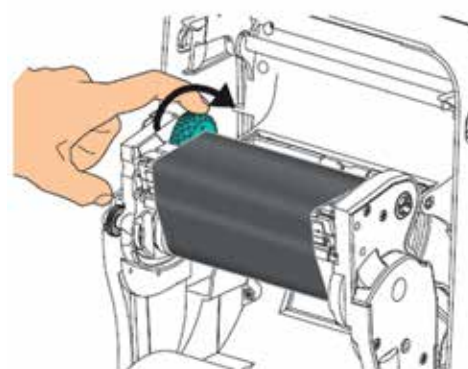
Figure 30

## 4.6.2 Monitoring software

The “Diluflux Pro” monitoring software enables the traceability of dilution and distribution via a spreadsheet (Microsoft® Excel® or OpenOffice™ or LibreOffice Calc™) and the printing of results.

### 4.6.2.1 Download

- Connect your computer to the DiluFlux Pro dilutor with the supplied USB cable.
- Insert the CD supplied with the unit and launch the “DiluFlow.exe” software.

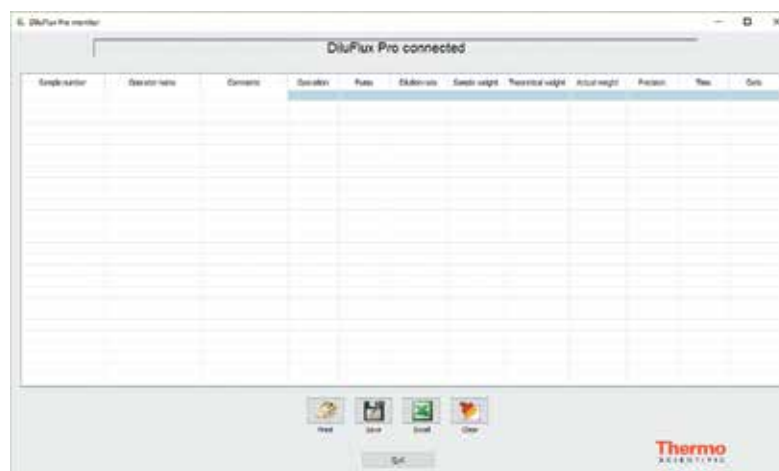
Figure 31

- When this screen (Figure 31) is displayed, the DiluFlux Pro dilutor is connected.

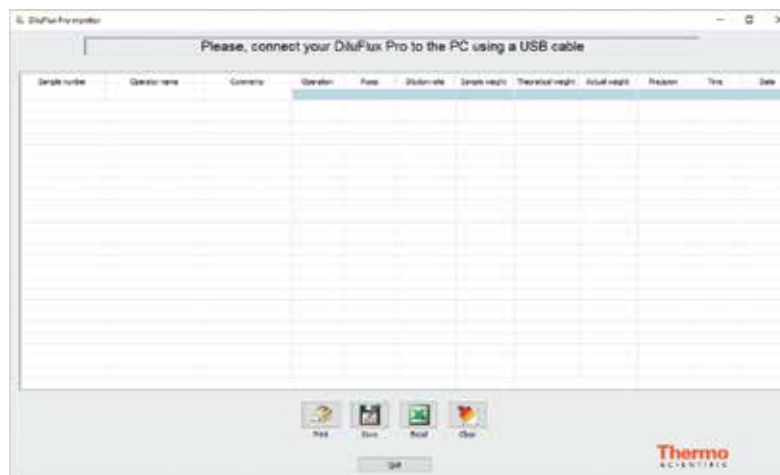
Figure 32

- If a screen such as Figure 32 is displayed, check that:
  - The DiluFlux Pro dilutor is turned on
  - The cable is connected
  - Relaunch the “DiluFlow.exe” software

## 4.6.2.2 Use of software

### Dilution/distribution

This function enables you to have records of your dilutions and distributions. Launch the process (dilution or distribution).

For each operation, a new line will be displayed, showing:

- Sample Number: as entered in the computer keyboard or the barcode reader
- Operator name: as entered in the computer keyboard or the barcode reader
- Comments: as entered on computer keyboard
- Operation: can take two values, i.e. “Dispensing” or “Dilution”
- Pump: can take two values, i.e. “Left pump” or “Right pump”
- Dilution rate: displayed when the operation performed is a dilution. This is a full number : the number N when the dilution ratio is 1/N
- Sample weight: in grams to 2 decimal places
- Theoretical weight: when the operation performed is a distribution, in grams, without decimal points
- Actual weight: in grams to two decimal places
- Precision: percentage, with one decimal point
- Time: computer time
- Date: computer date

#### NOTE:

- the titles of the columns are in English
- the separator fields of the CSV file are commas or semicolons depending on the country and computer
- the decimal separator is a full stop or comma depending on the country and the computer language
- dates and times are formatted according to the country and the computer language

The first 3 parameters (first 3 columns) should be filled in by the user. The other columns on green background are filled in automatically after each DiluFlux Pro dilutor operation.

Sample number	Operator name	Comments	Operation	Rate	Dilution	Sample weight	Theoretical weight	Actual weight	Precision	Time	Date
2345-12345	John Doe	OK	Dilution	1000	10	100.00	100.00	99.98	0.02	15:30:15	2023-10-27
10-1000-012	Marie Curie	N.A.S.	Dilution	1000	10	100.00	100.00	99.95	0.05	15:30:15	2023-10-27
10-1000-012	Marie Curie	N.A.S.	Dilution	1000	10	100.00	100.00	99.95	0.05	15:30:15	2023-10-27

Figure 33

To complete missing data (Number of sample, name of operator, comments), click on the corresponding cell, fill in the data and press Enter on your computer's keyboard. The cursor then moves to the next input field. To stop the editing process, press the "Esc" on your computer's keyboard or simply click elsewhere on the screen.

### Examples

The data in the first columns can be entered after operations, as shown in Figure 33.

As each new operation will automatically be recorded below the previous ones, it is recommended to enter the data in the first columns beforehand, as shown in Figure 34.

Sample number	Operator name	Comments	Operation	Rate	Dilution	Sample weight	Theoretical weight	Actual weight	Precision	Time	Date
1111-11111	John Doe										
1111-11111	John Doe										
10-1000-012	Marie Curie										

Figure 34

### Buttons at the bottom of the screen

To save your operations in a file, click on  and save in the directory and the desired file. The format will be CSV: values separated by commas or semicolons, depending on your country.

It is possible to export data under Excel® or to OpenOffice™ or LibreOffice Calc™ by clicking on  provided that one of this software is already installed in your computer.

Click on  to erase all the content of the sheet.

### 4.6.2.3 Stickers

#### Introduction

The DiluFlux Pro dilutor monitoring software enables you to print a label per operation (dilution or distribution) on a Zebra™ printer (optional), to apply to your sample. The key parameters of your dilution or distribution are printed on this label.

The operator's name must be written in the 2nd column of the sheet and may contain 30 characters. The name will be printed on the labels.

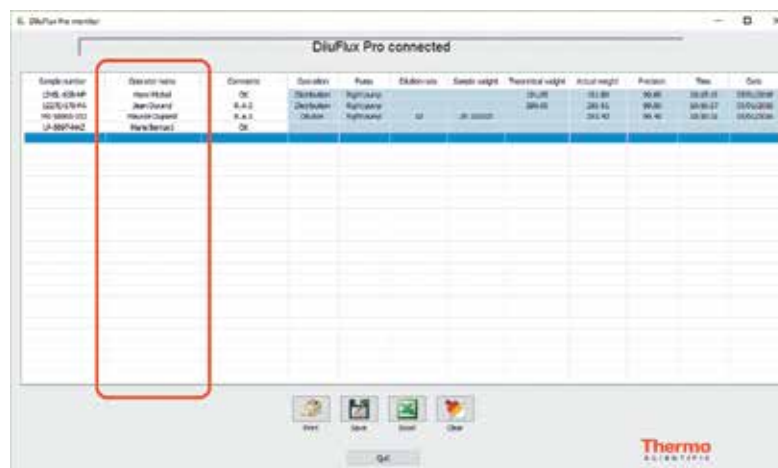
Figure 35

#### How to use

When a printer is connected to the Diluflux Pro dilutor, the software automatically prints a label after each operation (dilution or distribution).

You must enter the sample number and the operator name before the operation. The sticker is printed as below:

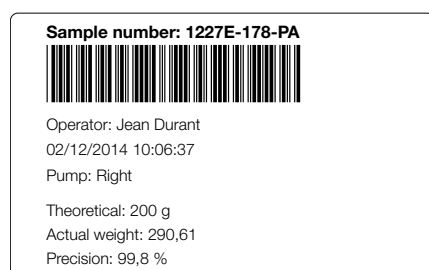
Figure 36

If the sample number is not specified, the DiluFlux Pro dilutor prints a label with an empty frame allowing you to write in the missing information.

<p><b>Sample No:</b></p> <p><b>Operator:</b></p>
<p>Pump: Right  Dilution rate: 10  Sample weight: 29,16 g  Actual weight: 293,43 g  Précision: 99,4 %</p>

**Figure 37**

To print a new label again, as long as the sample number is entered, you must select the corresponding line and click . To select a line, click on its right, green, part. The selected line then becomes dark blue.

## 4.7 Error messages

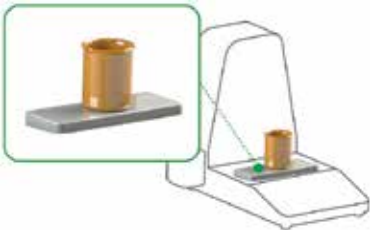
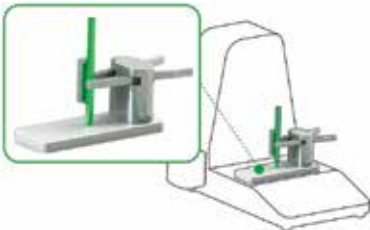
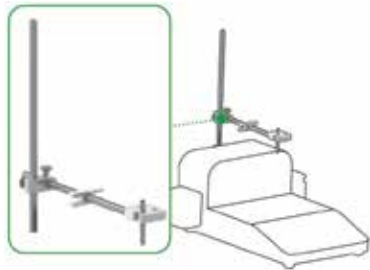
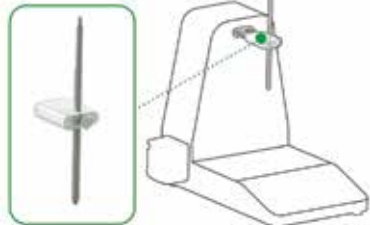
If operation fails, the device displays “Err” followed by a number:

- Err1:** (DiluFlux Pro-5 dilutor only) The robotic arm was unable to operate. Check that nothing is in the way and repeat the operation.
- Err2:** For a dilution, the weight of the sample is too low (lower than 0.5 gram) or too high (the total weight after dilution should not exceed the weight capacity of the Diluflux Pro dilutor). Replace the sample and start over.
- Err3:** The liquid does not flow. Check that the tubing is correctly placed in the pump and nothing is preventing the liquid from flowing in the tubing and then press again the .
- Err4:** At the end of a calibration, one of the two measures, either the empty weight or the 500 g calibrated weight is wrong. Check that nothing is resting on the weighing platform that could disturb the measurement.

# 5

## 5. Accessories (Optional)

Ref	Description	Figures	N°	Autoclavable	DiluFlux Pro-3 dilutor	DiluFlux Pro-5 dilutor
DB4103A	<b>BagOpen™ 400 for Diluflux Pro</b>		1		✓	✓
DB4104A	<b>Weighing platform</b>		2		✓	✓
DB4105A	<b>Tongs platform (Ø 10-60 mm)</b>		3		✓	✓
DB4106A	<b>Adjustable distribution system for DiluFlux Pro-3 dilutor</b>		4		✓	
DB4106A	<b>Adjustable distribution system for DiluFlux Pro-5 dilutor</b> (Incl. nozzle holder + 4 mm inner Ø (L: 23.5 cm) + 5 mm inner Ø (L: 12 cm) + Allen wrench + screw)		5			✓

Ref	Description	Figures	N°	Autoclavable	DiluFlux Pro-3 dilutor	DiluFlux Pro-5 dilutor
DB4107A	<b>Jumbo kit</b> (Incl. BagOpen™ 3500 + 3500 mL drip tray + 2 sets of adhesive gums + 1 protection cover)		6			✓
DB4108A	<b>Set of 10 adhesives for BagOpen™</b>		7		✓	✓
DB4109A	<b>Integral stand for broth bags</b> (Incl. 4 screws + 1 Torx screwdriver).		8	✓	✓	✓
DB4110A	<b>Booster kit</b> (Incl. 2 pump tubings (6.4 mm inner Ø L: 20 cm) + Y)		9		✓	✓
DB4113A	<b>Independent stand for broth bags</b>			✓	✓	✓

Ref	Description	Figures	N°	Autoclavable	DiluFlux Pro-3 dilutor	DiluFlux Pro-5 dilutor
DB4002A	<b>316 L stainless steel trocar for broth bags</b>			✓	✓	✓
DB4114A	<b>Set of 5 autoclavable PTFE filters</b> 0.2 µm Ø 25 mm			✓	✓	✓
DB4009A	<b>Set of 5 stainless steel nozzles for 6.4 mm tubings</b> (inner Ø 4.8 mm to 8 mm)			✓	✓	✓
DB4010A	<b>Set of 5 tube weights</b> (Ø 6 mm)			✓	✓	✓
DB4007A	<b>Foot pedal</b>				✓	✓
DB4102A	<b>Printer</b> (Incl. 1 USB cable + RS232- RS232 cable + 1 power cord + 1 printer)				✓	✓
DB4101A	<b>Film for printer</b>				✓	✓
DB4016A	<b>Printer labels</b> (x 2)				✓	✓

Ref	Description	Figures	N°	Autoclavable	DiluFlux Pro-3 dilutor	DiluFlux Pro-5 dilutor
DB4017A	<b>Printer paper</b> (x 2)				✓	✓
DB4118A	<b>Dispensing assembly</b> (inner Ø 6.4 mm) <b>for GL45 bottle</b>			✓	✓	✓
DB4116A	<b>Booster kit dispensing assembly</b> (inner Ø 6.4 mm) <b>for GL45 bottle</b>			✓	✓	✓
DB4119A	<b>Set of 5 silicone tubings</b> (inner Ø 6.4 mm)			✓	✓	✓
DB4117A	<b>Set of 5 silicone tubings for booster kit</b> (inner Ø 6.4 mm)			✓	✓	✓
DB4123A	<b>Set of 6 silicone tubings for pump</b> (inner Ø 6.4 mm)			✓	✓	✓

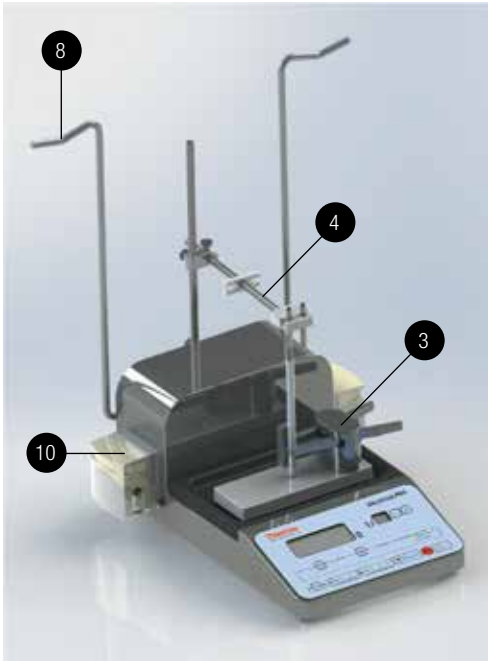
Figure 38

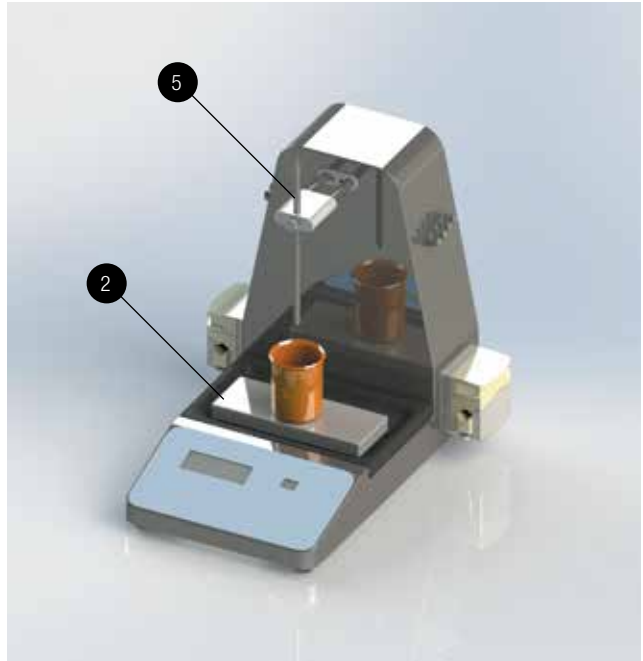
Figure 39

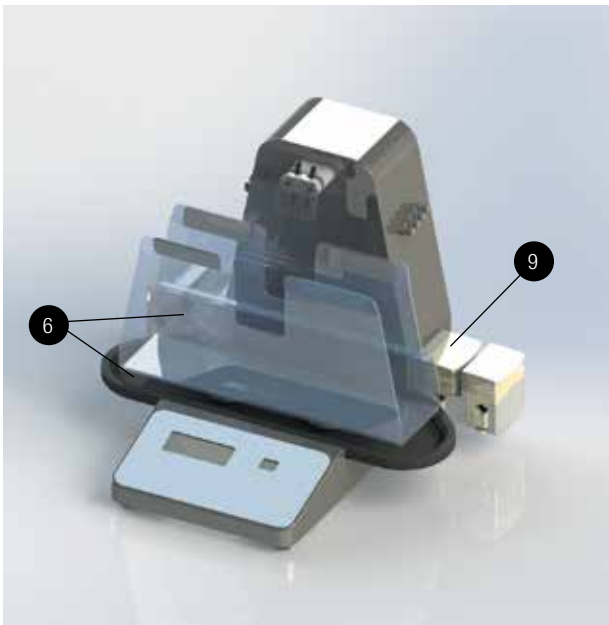
Figure 40

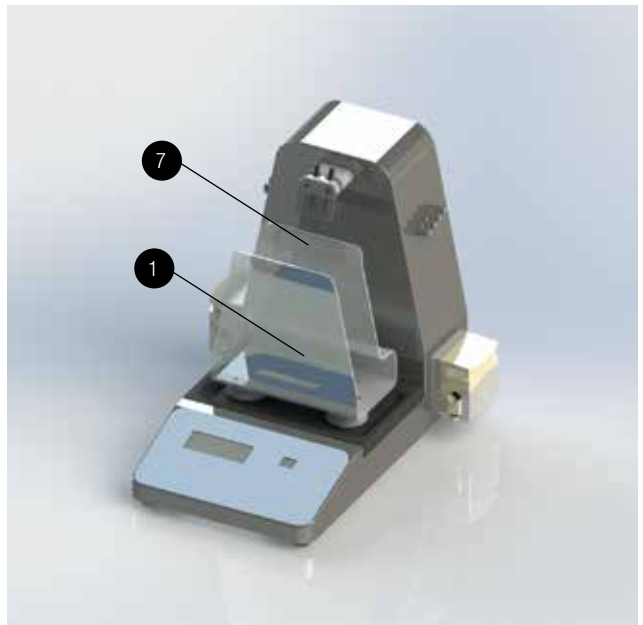
Figure 41

# 6

## 6. Maintenance

DiluFlux Pro dilutor requires little maintenance as long as leakages are cleaned quickly and thoroughly.

**Autoclavable items: dispensing assemblies, filter, bottle.**

**Non-autoclavable items: the unit itself, pumps, BagOpen™, DripTray, fixed arm, the nozzle holder, protecting plate for nozzles.**

**Accessories: refer to the chart 5 “Accessories (optional)”**


**ATTENTION: Do not use bleach to clean the housing of the DiluFlux Pro dilutor as it could corrode the stainless steel.**

- During cleaning, do not let any liquid drip into the unit.
- Autoclave all the dispensing assemblies every day when working in sterile conditions.
- Replace the silicone tubings whenever the tubing is not air-tight or becomes sticky or stiff.
- Every 20 autoclavings: replace the air filter (Ref DB4114A).
- To rinse the tubing using the DiluFlux Pro dilutor:
  - Take a bottle filled with water or another rinsing fluid.
  - Place the tube weight and its tubing in the bottle.
  - Pull the nozzle out of the nozzle holder by leaving the tubing in the pump and place it in the rinsing bottle to close the circuit.
  - Select the distribution mode by pressing , choose 100 g to dispense and press .
  - The tubing is rinsed.
  - Do this operation again as many times as necessary.

## 6.1 Pump heads maintenance

The pump head is not autoclavable. To clean it, unplug the unit, wipe out the dust and dirt with a soft cloth and then all surfaces with a disinfecting wipe.

Once a week, grease the pump head with a universal multi-purpose grease to the friction areas (see diagram below) to prevent premature wearing.

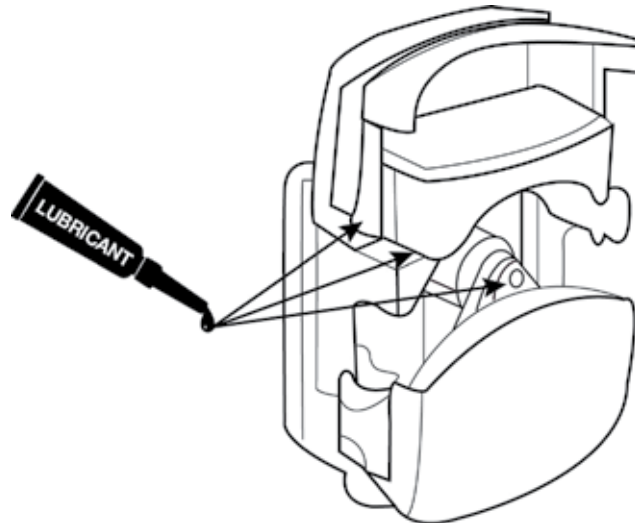**Figure 42**

Signs of wearing of pump head are:

- Difficult to prime
- Excessive to distribute
- Unsuccessful trial with a new tubing
- Gap between the cover and the body of the pump

If you notice these symptoms, please replace the pump head. The frequency of replacement depends on use and handling conditions. For information, peristaltic heads have been tested to undergo approximately 10,000 openings.

## 6.2 DiluFlux Pro-5 dilutor: cleaning of robotic arm

Figure 43

A daily cleaning of the two metallic rods of the arm is mandatory to ensure that the unit functions correctly. (see Figure 43).

- Switch the unit to standby mode by pressing the  key.
- Press the  button twice. The arm initializes and stops.
- Clean the 2 rods with ethanol.
- Turn the unit on again by pressing the  key. The arm initializes and retracts.

## 6.3 DiluFlux Pro-5 dilutor: nozzle holder cleaning

- Put the machine in standby mode by pressing the  key.
- Press the  button twice. The arm initializes and stops.
- Unscrew the nozzle holder using the Allen key supplied with the instrument. The screw is located under the holder. Take the holder out.
- Clean the holder (Caution: roller is not autoclavable).
- Put the holder back into place and screw closed.
- Turn the unit on by pressing the  key. The arm initializes and retracts.

# 7

## 7. Troubleshooting

Before contacting Thermo Fisher Technical Support, please check that your problem is not described here:

PROBLEM	CHECK-UP
<b>This display shows Err n°</b>	See chapter 4.7 Error messages
<b>DiluFlux Pro dilutor does not start</b>	<ul style="list-style-type: none"> <li>• Check that the switch is on "I"</li> <li>• Check that the power cord is correctly plugged in to the power supply and to the Diluflux Pro.</li> <li>• Check that the 2 fuses at the back of the unit are OK (value: T3.15A).</li> </ul>
<b>There are air bubbles in the diluent</b>	Check that you have correctly inserted the tubing in the peristaltic pump according to the direction of the flow indicated on the side of the unit. If it is not the case, open the pump and insert the tubing the right way.
<b>The indicated weight does not remain stable</b>	<ul style="list-style-type: none"> <li>• Check that the unit has been turned on for 10 minutes.</li> <li>• Check that the DripTray is well positioned.</li> <li>• Check that the weighing cell holes are free.</li> <li>• Check that there are no vibrations on the bench where you have installed the Diluflux Pro dilutor.</li> <li>• Check that the unit is horizontal.</li> <li>• Check that there is no airflow affecting the weighing.</li> <li>• Switch the unit off and on to re-initialize it.</li> <li>• Repeat a calibration.</li> </ul>
<b>The motor makes a noise but the pump does not start</b>	<ul style="list-style-type: none"> <li>• Check that the tubing is correctly positioned in the pump (see 3.3.5 "Adjusting the pump and insertion of the tubings").</li> <li>• Check that the tubing along its whole length is not pinched.</li> <li>• Check that the rollers of the pump are not blocked.</li> <li>• Check that the pumps are set correctly (see 3.3.5 "Adjusting the pump and insertion of the tubings").</li> <li>• If all these elements have been checked, stretch the tubings on both sides without opening the pump.</li> </ul>
<b>The pump perforates the tubings</b>	<ul style="list-style-type: none"> <li>• Check that the tubings have been correctly inserted and center them in the pump (see 3.3.5 "Adjusting the pump and insertion of the tubings")</li> <li>• Check that the pumps are set correctly (see 3.3.5 "Adjusting the pump and insertion of the tubings").</li> </ul>

PROBLEM	CHECK-UP
<b>The pump does not take up the liquid</b>	<ul style="list-style-type: none"> <li>• Check that the pumps are correctly set (see 3.3.5 "Adjusting the pump and insertion of the tubings").</li> <li>• Check that you have inserted the tubing in the peristaltic pump in the correct direction, according to the direction flow indicated at the side of the unit. If it is not the case, open the pump and insert the tubing the proper way.</li> </ul>
<b>The liquid flows back to the bottle or the pouch</b>	<ul style="list-style-type: none"> <li>• Check that the pumps are correctly set (see 3.3.5 "Adjusting the pump and insertion of the tubings").</li> <li>• Check that the tubings are not pinched.</li> </ul>
<b>STANDARD or ACCURATE Mode: The pump slows down before the end of the dilution or dispensing</b>	This could take place when mechanical vibrations disturb the device. Place the DiluFlux Pro dilutor on a table with a rigid and heavy platform.
<b>FAST Mode: The DiluFlux Pro dilutor dispenses too much or not enough liquid and the accuracy is not satisfactory</b>	An error occurred during auto-calibration of the pump flow. Proceed to a new dispensing/dilution operation.
<b>The printer does not switch on</b>	<ul style="list-style-type: none"> <li>• Check that the ON/OFF switch is well positioned.</li> <li>• Check that the power cord is connected to the power source and to the printer.</li> </ul>
<b>The printer does not work and an orange circle flashes around the green key of the printer</b>	<ul style="list-style-type: none"> <li>• There is no more paper. Please insert some more.</li> <li>• The cover of the printer is not shut.</li> <li>• Check the connection printer/Diluflux Pro dilutor.</li> <li>• Press on the green key until the flashing stops and replace the housing.</li> </ul>
<b>The computer does not display dilution or dispensing reports</b>	Check that the Diluflux Pro dilutor is connected to the USB cable. To do this, disconnect and then reconnect the cable and check that the computer makes a sound every time (the sound must be active). If the computer does not have a sound card, after the connection, check in the list displayed by the device manager that a new device has been added.
<b>The bag adhesive is no longer effective</b>	<ul style="list-style-type: none"> <li>• Clean it with water and soap</li> <li>• After several cleanings, replace it with a new cut (a set of 10 is provided with the unit). You may also order them (Ref: DB4018A).</li> <li>• If the sticker is not adapted to your application, use the adhesive gum provided with the unit.</li> </ul>

# 8

## 8. Appendix

### 8.1 Technical specifications

	DiluFlux Pro-3 Double pump	DiluFlux Pro-5 dilutor Double pump
Reference	DB4100A	DB4150A
<b>Weighing range</b>	0.1 g to 3 000 g	0.1 g to 5 000 g
<b>Weighing accuracy</b>	0 to 100 g: ± 0.05 g 100 g to 3 000 g: ± 0,1 %	0 to 100 g: ± 0.05 g 100g to 5 000 g: ± 0,1 %
<b>Minimum weight for dilution</b>	5 g	
<b>Minimum weight for distribution</b>	5 g	
<b>Maximum weight without stand</b>	4 500 g	7 000 g
<b>Dilution factor</b>	1/2 to 1/99	
<b>Dilution accuracy</b>	> 98 %	> 99 %*
<b>Dilution time for a 25 g sample diluted at 1/10</b>	< 9 s**	
<b>Dilution time for a 375 g sample diluted at 1/10</b>	-	< 80 s**
<b>Dispensing time for 90 mL</b>	< 7 s**	
<b>Dispensing time for 225 mL</b>	< 10 s**	
<b>Max number of pumps</b>	2 integrated	
<b>Dispensing/distribution mode</b>	Fast / standard / accurate	
<b>Traceability</b>	Printer / Monitoring software (data export to Excel™, OpenOffice™, LibreOffice Calc™, printer)	
<b>Exported data</b>	Sample weight / Total weight / Dilution factor / Dilution accuracy / Right or left pump + with monitoring software: Sample number / Name of operator / Date / Time	
<b>Connection</b>	RS232 output: To connect a printer USB output: To connect a computer Jack output: To connect a foot pedal	
<b>Weight</b>	13 kg	15 kg
<b>Dimensions (w x d x h)</b>	37 x 44 x 38 cm	37 x 42 x 38 cm
<b>In compliance with</b>	ISO 7218 / ISO 6887-1 / FDA-BAM / EI 61010-1: 2010	

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\* en mode STANDARD ou ACCURATE

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[thermofisher.com/microbiology](http://thermofisher.com/microbiology)

© 2016 Thermo Fisher Scientific Inc. All trademarks are the property of Thermo Fisher Scientific unless otherwise specified. BagFilter, BagPage, BagLight, BagOpen, RollBag are trademarks of INTERSCIENCE. Microsoft Excel is a trademark of Microsoft Corporation. OpenOffice is a trademark of Apache Software Foundation. Zebra is a trademark of ZIH Corp.

**Contact Information**

microbiology@thermofisher.com  
USA +800 255 6730  
International +44 (0) 1256 841144

May 2016  
LT2279A  
993-235

**Thermo**  
S C I E N T I F I C  
A Thermo Fisher Scientific Brand