

Alexa Fluor[®] Antibody Labeling Kits

Revision Date 27 July 2012

MAN0006869

Detailed protocol is available online at www.lifetechnologies.com/manuals.

Quick Protocol

- 1. Prepare a 1 M sodium bicarbonate solution by resuspending Component B in 1 mL of deionized water.
- Adjust the antibody concentration to 1.0 mg/mL, then add 1/10th volume of the 1 M sodium bicarbonate solution (from step 1).
- Add 100 μL of the antibody (from step 2) to the vial of Alexa Fluor[®] dye. Invert to mix and incubate for 1 hour at room temperature.
- Assemble the Purification Column. Prepare 1.5 mL of resin bed and centrifuge the column at 1100 x g for 3 minutes.
- 5. Add the reaction mix (from step 3) and centrifuge the column at $1100 \times g$ for 5 minutes to collect the labeled antibody.

For research use only. Not for use in diagnostic procedures.



Alexa Fluor® Antibody Labeling Kit	Ex/Em (nm)	Cat. no.
Alexa Fluor [®] 350 Antibody Labeling Kit	346/442	A20180
Alexa Fluor [®] 488 Antibody Labeling Kit	494/519	A20181
Alexa Fluor® 532 Antibody Labeling Kit	530/554	A20182
Alexa Fluor® 546 Antibody Labeling Kit	554/570	A20183
Alexa Fluor® 555 Antibody Labeling Kit	555/565	A20187
Alexa Fluor® 568 Antibody Labeling Kit	577/603	A20184
Alexa Fluor® 594 Antibody Labeling Kit	590/617	A20185
Alexa Fluor [®] 647 Antibody Labeling Kit	650/668	A20186
Alexa Fluor [®] 680 Antibody Labeling Kit	679/702	A20188
Alexa Fluor® 790 Antibody Labeling Kit	785/810	A20189

Limited Product Warranty

Life Technologies Corporation and/or its affiliate(s) warrant their products as set forth in the Life Technologies' General Terms and Conditions of Sale found on Life Technologies' website at www.lifetechnologies.com/termsandconditions. If you have any questions, please contact Life Technologies at www.lifetechnologies.com/support.

© 2012 Life Technologies Corporation. All rights reserved. The trademarks mentioned herein are the property of Life Technologies Corporation or their respective owners.

LIFE TECHNOLOBIES CORPORATION AND/OR ITS AFFLIATESI DISCLAM ALL WARRANTES WITH RESPECT TO THIS DOCUMENT, EXPRESSIO OR MOVELO, INCLUDING BUT NOT UNITED TO THOSE OF MERCHANTABILITY, FITHASES FOR A APATICULAR PURPOSE, OR NON-INFRRIGEMENT. TO THE EVET ALLOWED BY LAW, IN NO EVENT SHALL LIFE TECHNOLOBIES AND/OR ITS AFFLIATESI BE LABLE, WHETHET IN CONTRACT, TORT, WARRANT, OR UNDER ANY STATUTE OR ON ANY OTHER BASIS FOR SPECIAL, INDERNIAL, INDIRECT, PUNITIVE, MULTIPLE OR CONSEQUENTIAL DAMAGES IN CONNECTION WITH OR ARISING FORM THIS DOCUMENT, INCLUMING BUT NOT UNITED TO THE USE THEREOF.

For support visit www.lifetechnologies.com/support

www.lifetechnologies.com