LIVE/DEAD[™] Fixable Dead Cell Stain Kits

Pub. No. MAN0002416 (MP34955) Rev. A.0

Table 1. Contents and storage

Material	Amount	Storage	Stability	
Individual Kits: Blue, violet, aqua, yellow-, green, red, far red, or near-IR fluorescent reactive dye (Component A)	2 vials, each (80 assays); 5 vials, each (200 assays); 10 vials, each (400 assays)			
Sampler Kit: Blue, violet, aqua, yellow, green, red, far red, and near-IR fluorescent reactive dye (Components A–H)	1 vial of each dye (8 total)	 ≤-20°C Desiccate Protect from light 	When stored as directed, kit components are stable for at least 6 months.	
Dimethylsulfoxide (DMSO), anhydrous (Component B in Individual Kits, Component I in Sampler Kit)	500 µL	igne		
Number of assays: Each vial of the reactive dye provides sufficient material for approximately 40 flow cytometry assays, for a total of 80, 200, or 400 assays depending on the kit configuration (see Table 2). The Sampler Kit contains one vial of each fluorescent reactive dye with 40 assays per vial of the reactive dye, for a total of 320 assays per kit.				

Approximate fluorescence excitation/emission maxima: See Table 2.

Introduction

The LIVE/DEAD[™] Fixable Dead Cell Stain Kits use a novel method to evaluate the viability of mammalian cells by flow cytometry. These assays are based on the reaction of a fluorescent reactive dye with cellular amines. The reactive dye can permeate the compromised membranes of necrotic cells and react with free amines both in the interior and on the cell surface, resulting in intense fluorescent staining. In contrast, only the cell-surface amines of viable cells are available to react with the dye, resulting in relatively dim staining (Figure 1, page 3). The difference in intensity between the live and dead cell populations is typically greater than 50-fold (Figure 2, page 3). The discrimination is completely preserved following fixation of the sample by formaldehyde, under conditions that inactivate pathogens. Moreover, these single-color assays use only one channel of a flow cytometer, leaving the other channels available for multicolor experiments. The assays can also be used to detect dead cells by microscopy; however, the difference in fluorescence intensity of the live and dead cells can be appreciable, making it relatively difficult to simultaneously photograph the two populations.



The single-color LIVE/DEAD[™] Fixable Dead Cell Stain Kits are identical except for the fluorescent color of the included dye—blue, violet, aqua, yellow, green, red, far red, or near-IR (infra red).

- Cells labeled by green fluorescent or red fluorescent reactive dye are excited by the 488 nm line of an argon-ion laser; green fluorescence is typically detected in the green channel of the flow cytometer (530/30 nm) and red fluorescence is detected in the red channel (630/30 nm). The blue fluorescent reactive dye requires UV (350–360 nm) excitation with fluorescence emission read at ~450 nm.
- The violet fluorescent reactive dye requires violet (~405 nm) excitation with fluorescence emission read at ~440 nm.
- The aqua fluorescent reactive dye is efficiently excited with ~405 nm light (or UV light) and has fluorescence emission monitored at ~525 nm.
- The yellow fluorescent reactive dye requires violet (~405 nm) excitation with fluorescence emission read at ~575 nm (appropriate channels for the violet-excitable reactive dyes may vary depending on the instrument).
- The far red and near-IR fluorescent reactive dyes are excited at 633/635 nm with fluorescence emission monitored at 665 nm and 775 nm, respectively.
- The LIVE/DEAD[™] Fixable Dead Cell Stain Sampler Kit (Cat. no. L34960) contains one vial of each of the eight different fluorescent reactive dyes.

We also offer a LIVE/DEAD[™] Reduced Biohazard Cell Viability Kit (Cat. no. L7013) for determining cell viability. It is based on a different staining principle using two dyes of different colors, and is described separately. For additional information, visit **www.thermofisher.com**.

Reactive dye	Number of assays (Cat. no.)	Excitation source	Ex*	Em*
	80 (L34961)			
Blue fluorescent reactive dye	200 (L23105)	UV	350 nm	450 nm
	400 (L34962)			
	80 (L34963)			
Violet fluorescent reactive dye	200 (L34955)	405 nm	416 nm	451 nm
	400(L34964)			
	80 (L34965)			
Aqua fluorescent reactive dye	200 (L34957)	405 nm	367 nm	526 nm
	400 (L34966)			
Yellow fluorescent reactive dye	80 (L34967)			
	200 (L34959)	405 nm	400 nm	575 nm
	400 (L34968)			
Green fluorescent reactive dye	80 (L34969)			
	200 (L23101)	488 nm	495 nm	520 nm
	400 (L34970)			
Red fluorescent reactive dye	80 (L34971)			
	200 (L23102)	488 nm	595 nm	615 nm
	400 (L34972)			
Far red fluorescent reactive dye	80 (L34973)			
	200 (L10120)	633/635 nm	650 nm	665 nm
	400 (L34974)			
	80 (L34975)			
Near-IR fluorescent reactive dye	200 (L10119)	633/635 nm	750 nm	775 nm
	400 (L34976)			
*Approximate fluorescence excitation	(Ex) and emission (Em) m	axima, in nm.		

Table 2. Approximate fluorescence excitation and emission maxima and kit sizes for the LIVE/DEAD[™] Fixable Dead Cell Stain single-color dyes.



Figure 1. Principle of the LIVE/DEAD[™] Fixable Dead Cell Stain Kits. Live cells (left) react with the kit's fluorescent reactive dye only on their surface to yield weakly fluorescent cells. Cells with compromised membranes (right) react with the dye throughout their volume, yielding brightly stained cells. In both cases, the excess reactive dye is subsequently washed away.



Figure 2. Live and dead cells distinguished by flow cytometry. Each of the LIVE/DEAD[™] Fixable Dead Cell Stain Kits was used to differentially stain a mixture of live (left peak) and heat-treated Jurkat cells (right peak) according to the protocol provided in this document. (A) LIVE/DEAD[™] Fixable Blue Stain Kit with UV excitation and ~440 nm emission; (B) LIVE/DEAD[™] Fixable Violet Stain Kit with violet excitation and ~440 nm emission; (C) LIVE/DEAD[™] Fixable Aqua Dead Cell Stain Kit using 405 nm excitation and ~525 nm emission; (D) LIVE/DEAD[™] Fixable Yellow Stain Kit with 405 nm excitation and ~575 nm emission; (E) LIVE/DEAD[™] Fixable Green Stain Kit with 488 nm excitation and ~525 nm emission; (F) LIVE/DEAD[™] Fixable Red Stain Kit with 488 nm excitation and ~585 nm emission; (G) LIVE/DEAD[™] Fixable Far Red Stain Kit with 633 nm excitation and ~660 nm emission; and (H) LIVE/DEAD[™] Fixable Near-IR Stain Kit with 633 nm excitation and ~780 nm emission. Following the staining reaction, the cells were fixed in 3.7% formaldehyde and analyzed by flow cytometry. The live cell population is easily distinguished from the killed population, and nearly identical results were obtained using unfixed cells (data not shown).

Allow the reagents to warm to room temperature before opening the vials.

Materials required but	
not provided	Phosphate buffered saline (PBS)
	• PBS with 1% bovine serum albumin (BSA)
	• Formaldehyde
	 AbC[™] Total Antibody Compensation Bead Kit (if calculating compensation in multicolor immunophenotyping experiments using mouse, rat, rabbit, or hamster antibodies)
	 ArC[™] Amine Reactive Compensation Bead Kit (if calculating compensation in multicolor experiments using a LIVE/DEAD[™] Fixable Dead Cell Stain)
Caution	DMSO is hazardous; avoid contact with skin and eyes and do not swallow. Handle reagents containing DMSO using equipment and practices appropriate for the hazards posed by such materials.

Experimental Protocols

The following protocol has been used successfully in our laboratories to differentially stain live and dead cells and then to fix the cells in formaldehyde for subsequent analysis by flow cytometry. Excellent results have been obtained with a variety of cell types, including: Jurkat, MDCK, COLO205, CHO-K1, K-562, HeLa, P3X, 3T3, B35, BPAC, CHO-M1, A549, MMM, MRC5, U205, U266, 293 MSR, and RAW tissue-culture cell lines and human peripheral blood lymphocytes. If another staining reaction is to be performed on the same sample, determine the optimal staining sequence for the two procedures and whether or not the additional staining reaction will tolerate fixation by formaldehyde.

Dye preparation For convenience, the reactive dye in each individual kit is supplied in five separate vials, while the sampler kit contains one vial of each reactive dye. Each vial provides sufficient material for staining at least 40 cell samples. However, once reconstituted, the DMSO solution of reactive dye is somewhat unstable, especially if exposed to moisture. Unused portions may be used for up to 2 weeks if stored at ≤–20°C, protected from light and moisture.

- **1.1** Bring one vial of the fluorescent reactive dye (Component A) and the vial of anhydrous DMSO (Component B) to room temperature before removing the caps.
- **1.2** Add 50 μ L of DMSO to the vial of reactive dye. Mix well and visually confirm that all of the dye has dissolved.
- **1.3** Use the solution of reactive dye as soon as possible (see below), ideally within a few hours of reconstitution.

- **Cell staining** Buffers appropriate for cell staining include phosphate-buffered saline (PBS), Hanks' Balanced Salt Solution (HBSS), and Dulbecco's PBS (D-PBS), without extraneous proteins such as bovine serum albumin or serum. When using an amino-reactive dye, for example the violet fluorescent reactive dye, Tris buffers and solutions containing sodium azide or extraneous protein should not be used for cell resuspension and washing.
 - **2.1** Centrifuge a sample of cells in suspension containing at least 1×10^{6} cells. Discard the supernatant.
 - 2.2 Wash the cells once with 1 mL of PBS.
 - **2.3** Resuspend the cells in 1 mL of PBS.
 - **2.4** Count the cells and adjust the density with PBS to 1×10^6 cells in a 1 mL volume.
 - **2.5** Add 1 μ L of the reconstituted fluorescent reactive dye (from step 1.3) to 1 mL of the cell suspension and mix well.
 - 2.6 Incubate at room temperature or on ice for 30 minutes, protected from light.

Note: If fixation is not required, then you can skip the steps 2.7–2.10 below. Instead, wash the cells twice with 1 mL of PBS with 1% bovine serum albumin, and resupend in 1 mL of PBS with 1% bovine serum albumin.

- 2.7 Wash the cells once with 1 mL of PBS, and resuspend the cells in 900 µL of PBS.
- **2.8** Add 100 µL of 37% formaldehyde.
- 2.9 Incubate at room temperature for 15 minutes.
- **2.10** Wash once with 1 mL of PBS with 1% bovine serum albumin, and resuspend the cells in 1 mL of PBS with 1% bovine serum albumin.
- **2.11** Analyze the fixed cell suspension by flow cytometry using the appropriate excitation and detection channel (see Table 3, below).

Note: The appropriate excitation and detection channels may vary depending on instrument used.

Table 3. Appropriate excitation and detection channels to use with the amine-reactive LIVE/DEAD	fixable
dead cell stains	

Reactive dye	Excitation	Detection*	Recommended filters
Blue fluorescent reactive dye	UV	~450 nm	450/50 nm or similar
Violet fluorescent reactive dye	405 nm	~450 nm	450/50 nm or similar
Aqua fluorescent reactive dye	405 nm	~525 nm	525/50 nm or similar
Yellow fluorescent reactive dye	405 nm	~575 nm	575/26 nm, 585/42 nm, or similar
Green fluorescent reactive dye	488 nm	~530 nm	530/30 nm or similar
Red fluorescent reactive dye	488 nm	~585 nm	585/42 nm, 610/20 nm, or similar
Far red fluorescent reactive dye	633/635 nm	~660 nm	
Near-IR fluorescent reactive dye	633/635 nm	~780 nm	
*For the approximate fluorescence excitation and emission maxima of the amine-reactive LIVE/DEAD [™] dyes, see Table 2 (page 2) and Figure 2 (page 3).			

- **Intracellular staining** This procedure is completely compatible with common fixation and permeabilization methods for performing intracellular staining for flow cytometry. The procedure for staining with a LIVE/DEAD[™] Fixable Dead Cell Stain combined with immunophenotyping is as follows:
 - **3.1** Centrifuge a sample of cells in suspension containing at least 1×10^6 cells. Discard the supernatant.
 - 3.2 Wash the cells once with 1 mL of PBS.
 - 3.3 Resuspend the cells in 1 mL of PBS.
 - **3.4** Count the cells and adjust the density with PBS to 1×10^6 cells in a 1 mL volume.
 - **3.5** Add 1 µL of the reconstituted fluorescent reactive dye (from step 1.3) to 1 mL of the cell suspension and mix well.
 - 3.6 Incubate at room temperature or on ice for 30 minutes, protected from light.
 - 3.7 Wash the cells with 1 mL of PBS and resuspend in 100 µl of PBS.
 - 3.8 Stain as usual for surface markers, and incubate for desired time for antibody staining.
 - 3.9 Wash the cells with PBS and resuspend in 900 µl of PBS
 - **3.10** Add 100 µL of 37% formaldehyde.
 - 3.11 Incubate at room temperature for 15 minutes.
 - 3.12 Wash once with 1 mL of PBS with 1% bovine serum albumin, and resuspend the cells in 100 μ L of PBS with 1% bovine serum albumin.
 - **3.11** Add permeabilization reagent and stain as usual for intracellular markers. Incubate for the desired time for antibody staining.
 - **3.12** Wash once with 1 mL of PBS with 1% bovine serum albumin, and resuspend the cells in 1 mL of PBS with 1% bovine serum albumin.
 - **3.13** Analyze the fixed cell suspension by flow cytometry using the appropriate excitation and detection channel (from step 2.11)

Compensation Using ArC[™] Amine Reactive Beads

The ArC^{$^{\text{M}}$} Amine Reactive Compensation Bead Kit is designed to facilitate compensation when using any of the LIVE/DEAD^{$^{\text{M}}$} fixable dead cell stains, providing a consistent, accurate and simple-to-use technique for the setting of flow cytometry compensation. The ArC^{$^{\text{M}}$} Amine Reactive Compensation Bead Kit includes two types of specially modified polystyrene microspheres to allow easy compensation of the LIVE/DEAD^{$^{\text{M}}$} fixable stains: the ArC^{$^{\text{M}}$} reactive beads (Component A), which bind any of the amine-reactive dyes, and the ArC^{$^{\text{M}}$} negative beads (Component B), which have no reactivity. After incubation with any amine-reactive dye, the two kit components will provide distinct positive and negative populations of beads that can be used to set compensation. We recommend the following protocol for using ArC^{$^{\text{M}}</sup> Amine Reactive Compensation:</sup>$

- **4.1.** Gently vortex ArC[™] Amine Reactive Compensation Bead Kit components for 30 seconds to completely resuspend before use.
- **4.2.** Add 1 drop of ArC^{TM} reactive beads (Component A) to a labeled sample tube.
- **4.3.** Allow ArC[™] reactive beads to sit in the tube for 5 minutes to warm to room temperature.
- **4.4.** Prepare fluorescent amine-reactive dye according to instructions included in the LIVE/DEAD[™] Fixable Dead Cell Kit. For optimal performance of the ArC[™] reactive beads, use freshly prepared amine-reactive dye. Do not use previously frozen dye solution.
- **4.5.** Add the amount of LIVE/DEAD[™] fixable dead cell stain listed in Table 4, below, to the bead suspension and mix well. Make sure to deposit the amine-reactive dye directly to the bead suspension.

Table 4. Amount of amine-reactive LIVE/DEAD[™] fixable dead cell stain for use with ArC[™] reactive beads

Amine-reactive dye for use with ArC [™] Reactive Beads	Amount
LIVE/DEAD [™] Fixable Blue stain	3 µL
LIVE/DEAD [™] Fixable Violet stain	1 µL
LIVE/DEAD [™] Fixable Aqua stain	3 µL
LIVE/DEAD [™] Fixable Yellow stain	3 µL
LIVE/DEAD [™] Fixable Green stain	3 µL
LIVE/DEAD [™] Fixable Red stain	1 µL
LIVE/DEAD [™] Fixable Far Red stain	3 µL
LIVE/DEAD [™] Fixable Near-IR stain	1 µL

- **4.6.** Incubate for 30 minutes at room temperature, **protected from light**.
- **4.7.** Add 3 mL of PBS or other buffer to sample tube. Centrifuge for 5 minutes at $300 \times g$.
- **4.8.** Carefully remove all the supernatant from tube.

Note: If using the red fluorescent reactive dye, repeat step 4.7.

- **4.9.** Resuspend bead pellet by adding 0.5 mL of buffer to sample tube
- **4.10.** Add one drop of ArC^{TM} negative beads (Component B) to sample tube. Mix thoroughly.
- **4.11.** Vortex tubes before analyzing using flow cytometry.
- **4.12.** Perform manual or automatic compensation according to the preferred procedure for the flow cytometer in use. Gate on the bead singlet population based on FSC and SSC characteristics.

Combining ArC[™] and AbC[™] Kits The AbC[™] Total Antibody Compensation Bead Kit (Cat. nos. A10497, A105213) provides a consistent, accurate, and simple-to-use technique for the setting of flow cytometry compensation when using fluorochrome-conjugated mouse antibodies. The kit contains two types of specially modified polystyrene microspheres: the AbC[™] capture beads (Component A) that bind all isotypes of mouse, rat, rabbit, and hamster immunoglobulin, and the negative beads (Component B) that have no antibody binding capacity. After incubating with a fluorochrome-conjugated antibody, the two components provide distinct positive and negative populations of beads that you can use to set compensation. You can use the AbC[™] Total Antibody Compensation Bead Kit and the ArC[™] Amine Reactive Compensation Bead Kit together to calculate compensation in multicolor immunophenotyping experiments that incorporate a LIVE/DEAD[™] fixable dye by following the protocol outlined below:

- **5.1.** Gently vortex the ArC[™] Amine Reactive Compensation Bead Kit and the AbC[™] Total Antibody Compensation Bead Kit components for 30 seconds to completely resuspend before use.
- **5.2.** Label a sample tube for the amine-reactive dye you are using and add 1 drop of ArC[™] reactive beads (Component A in the ArC[™] Amine Reactive Compensation Bead Kit) to the labeled sample tube. Allow ArC[™] reactive beads to sit in the tube for 5 minutes to warm to room temperature.
- **5.3.** Prepare fluorescent reactive dye according to kit instructions included in the LIVE/ DEAD[™] Fixable Dead Cell Stain Kit. For optimal performance of ArC[™] reactive beads, always use freshly prepared amine-reactive dye. Do **not** use previously frozen dye solution.
- **5.4.** Add the amount of LIVE/DEAD[™] fixable dead cell stain listed in Table 4 (page 7) to the bead suspension and mix well. Make sure to deposit the amine-reactive dye directly to the bead suspension.
- **5.5.** Label another sample tube for each fluorochrome-conjugated antibody you are using, and add 1 drop of AbC[™] capture beads (Component A in the AbC[™] Total Antibody Compensation Bead Kit) to each labeled tube.
- **5.6.** Add a pre-titrated amount of antibody conjugate to the appropriate tube and mix well. Make sure to deposit the antibody directly to the bead suspension.
- 5.7. Incubate for 30 minutes at room temperature, protected from light.
- **5.8.** Add 3 mL of PBS or other buffer to each sample tube. Centrifuge at $300 \times g$ for 5 minutes to collect beads.
- 5.9. Carefully remove all supernatant from each tube.

Note: If using the red fluorescent reactive dye, repeat step 5.8 for that tube.

- 5.10. Resuspend bead pellet by adding 0.5 mL of staining buffer to each sample tube.
- **5.11.** Add one drop of negative beads (Component B in the AbC[™] Total Antibody Compensation Bead Kit) to sample tube(s) containing the AbC[™] capture beads.
- **5.12.** Add one drop of ArC[™] negative beads (Component B in the ArC[™] Amine Reactive Compensation Bead Kit) to sample tube(s) containing the ArC[™] reactive beads.
- 5.13. Vortex tubes before analyzing using flow cytometry.
- **5.14.** Perform manual or automatic compensation according to the preferred procedure for the flow cytometer in use. Gate on the bead singlet population based on FSC and SSC characteristics.

1. J Immunol Methods 313, 199 (2006). 2. Blood 111, 1344 (2008). 3. J Immunol 180, 774 (2008).

Product List Current prices may be obtained from our website or from our Customer Service Department.

Cat. no.	Product Name
L34961	LIVE/DEAD [™] Fixable Blue Dead Cell Stain Kit *for UV excitation* *80 assays*
L23105	LIVE/DEAD [™] Fixable Blue Dead Cell Stain Kit *for UV excitation* *200 assays*
L34962	LIVE/DEAD [™] Fixable Blue Dead Cell Stain Kit *for UV excitation* *400 assays*
L34963	LIVE/DEAD [™] Fixable Violet Dead Cell Stain Kit *for 405 nm excitation* *80 assays*
L34955	LIVE/DEAD [™] Fixable Violet Dead Cell Stain Kit *for 405 nm excitation* *200 assays*
L34964	LIVE/DEAD [™] Fixable Violet Dead Cell Stain Kit *for 405 nm excitation* *400 assays*
L34965	LIVE/DEAD [™] Fixable Aqua Dead Cell Stain Kit *for 405 nm excitation* *80 assays*
L34957	LIVE/DEAD [™] Fixable Aqua Dead Cell Stain Kit *for 405 nm excitation* *200 assays*
L34966	LIVE/DEAD [™] Fixable Aqua Dead Cell Stain Kit *for 405 nm excitation* *400 assays*
L34967	LIVE/DEAD [™] Fixable Yellow Dead Cell Stain Kit *for 405 nm excitation* *80 assays*
L34959	LIVE/DEAD [™] Fixable Yellow Dead Cell Stain Kit *for 405 nm excitation* *200 assays*
L34968	LIVE/DEAD [™] Fixable Yellow Dead Cell Stain Kit *for 405 nm excitation* *400 assays*
L34969	LIVE/DEAD [™] Fixable Green Dead Cell Stain Kit *for 488 nm excitation* *80 assays [*]
L23101	LIVE/DEAD [™] Fixable Green Dead Cell Stain Kit *for 488 nm excitation* *200 assays*
L34970	LIVE/DEAD [™] Fixable Green Dead Cell Stain Kit *for 488 nm excitation* *400 assays*
L34791	LIVE/DEAD [™] Fixable Red Dead Cell Stain Kit *for 488 nm excitation* *80 assays [*]
L23102	LIVE/DEAD [™] Fixable Red Dead Cell Stain Kit *for 488 nm excitation* *200 assays*
L34972	LIVE/DEAD [™] Fixable Red Dead Cell Stain Kit *for 488 nm excitation* *400 assays*
L34973	LIVE/DEAD [™] Fixable Far Red Dead Cell Stain Kit *for 633/635 nm excitation* *80 assays*
L10120	LIVE/DEAD [™] Fixable Far Red Dead Cell Stain Kit *for 633/635 nm excitation* *200 assays*
L34974	LIVE/DEAD [™] Fixable Far Red Dead Cell Stain Kit *for 633/635 nm excitation* *400 assays*
L34975	LIVE/DEAD [™] Fixable Near-IR Dead Cell Stain Kit *for 633/635 nm excitation* *80 assays*
L10119	LIVE/DEAD [™] Fixable Near-IR Dead Cell Stain Kit *for 633/635 nm excitation* *200 assays*
L34976	LIVE/DEAD [™] Fixable Near-IR Dead Cell Stain Kit *for 633/635 nm excitation* *400 assays*
L34960	LIVE/DEAD [™] Fixable Dead Cell Stain Sampler Kit *for flow cytometry* *320 assays*
Related Produ	cts
A10497	AbC [™] Total Antibody Compensation Bead Kit *for flow cytometry* *100 tests*
A10513	AbC [™] Total Antibody Compensation Bead Kit *for flow cytometry* *25 tests*
A10344	AbC [™] Anti-Mouse Bead Kit *for mouse antibody capture* *for flow cytometry* *100 tests*
A10346	ArC [™] Amine Reactive Compensation Bead Kit *for use with amine reactive dyes * *for flow cytometry compensation*
	100 tests
L7013	LIVE/DEAD [™] Reduced Biohazard Cell Viability Kit #1 *green and red fluorescence* *100 assays*
GAS-003	Fixation and Permeabilization, 1 × 5 mL *for 50 tests*
GAS-004	Fixation and Permeabilization, 4 × 5 mL *for 200 tests*
GAS001S-100	Fixation Medium - Bulk, (MEDIUM A),1 × 100 mL Fixation Medium
GAS002S-100	Permeabilization Medium - Bulk, (MEDIUM B),1 × 100 mL Permeabilization Medium
FB001	IC Fixation Buffer
PB001	IC Permeabilization Buffer

Purchaser notification

These high-quality reagents and materials must be used by, or directly under the supervision of, a technically qualified individual experienced in handling potentially hazardous chemicals. Read the Safety Data Sheet provided for each product; other regulatory considerations may apply.

Obtaining support

For the latest services and support information for all locations, go to thermofisher.com/support.

- At the website, you can:
- · Access worldwide telephone and fax numbers to contact Technical Support and Sales facilities
- Search through frequently asked questions (FAQs)
- Submit a question directly to Technical Support (thermofisher.com/support)
- Search for user documents, SDSs, vector maps and sequences, application notes, formulations, handbooks, certificates of analysis, citations, and other product support
- documents

 Obtain information about customer training
- Obtain information about customer training
 Download software updates and patches

SDS

Safety Data Sheets (SDSs) are available at thermofisher.com/support.

Certificate of Analysis

The Certificate of Analysis provides detailed quality control and product qualification information for each product. Certificates of Analysis are available on our website. Go to **thermofisher.com/support** and search for the Certificate of Analysis by product lot number, which is printed on the product packaging (tube, pouch, or box).

Limited Product Warranty

Life Technologies Corporation and/or its affiliate(s) warrant their products as set forth in the Life Technologies' General Terms and Conditions of Sale found on Life Technologies' website at www.lifetechnologies.com/termsandconditions. If you have any questions, please contact Life Technologies at www.lifetechnologies.com/support.

For Research Use Only. Not for use in diagnostic procedures.

Disclaimer

TO THE EXTENT ALLOWED BY LAW, LIFE TECHNOLOGIES AND/OR ITS AFFILIATE(S) WILL NOT BE LIABLE FOR SPECIAL, INCIDENTAL, INDIRECT, PUNITIVE, MULTIPLE OR CONSEQUENTIAL DAMAGES IN CONNECTION WITH OR ARISING FROM THIS DOCUMENT, INCLUDING YOUR USE OF IT.

Important Licensing Information

These products may be covered by one or more Limited Use Label Licenses. By use of these products, you accept the terms and conditions of all applicable Limited Use Label Licenses.

Corporate entity

Life Technologies I Carlsbad, CA 92008 USA I Toll free in USA 1.800.955.6288

All trademarks are the property of Thermo Fisher Scientific and its subsidiaries, unless otherwise specified.

©2016 Thermo Fisher Scientific Inc. All rights reserved.

