

# invitrogen™

by *life* technologies™

## One Shot® *ccdB* Survival™ 2 T1<sup>R</sup> Chemically Competent Cells

Cat. no. A10460

Size 10 reactions

Store at **-85°C to -68°C**

Doc. Part no. 100003536

Pub. no. MAN0000761

Rev. 2.0

### Description

The One Shot® *ccdB* Survival™ 2 T1<sup>R</sup> strain is resistant to the toxic effects of the *ccdB* gene (Bernard and Couturier, 1992; Bernard et al., 1993) and it may be used to propagate and maintain vectors containing the *ccdB* gene (e.g. Gateway® Technology vectors). The transformation efficiency of One Shot® *ccdB* Survival™ 2 T1<sup>R</sup> chemically competent cells is greater than  $1 \times 10^9$  cfu/ $\mu$ g pUC19 Control DNA.

Component	Amount
<i>ccdB</i> Survival™ 2 T1 <sup>R</sup> Competent Cells	11 × 50 $\mu$ L
pUC19 Control DNA (10 pg/ $\mu$ L)	50 $\mu$ L
S.O.C. Medium	6 mL

### Genotype

F *mcrA*  $\Delta$ (*mrr-hsdRMS-mcrBC*)  $\Phi$ 80*lacZ* $\Delta$ M15  $\Delta$ *lacX74 recA1 ara* $\Delta$ 139  
*Δ(ara-leu)*7697 *galU galK rpsL* (Str<sup>R</sup>) *endA1 nupG fhuA::IS2*

### Certificate of Analysis

The Certificate of Analysis provides detailed quality control and product qualification information for each product. Certificates of Analysis are available on our website. Go to [www.lifetechnologies.com/support](http://www.lifetechnologies.com/support) and search for the Certificate of Analysis by product lot number, which is printed on the box.

For research use only. Not for use in diagnostic procedures.

## General Guidelines

- Handle competent cells gently; they are highly sensitive to changes in temperature and mechanical lysis caused by pipetting.
- Thaw the competent cells on ice, and transform them immediately. Add DNA to tube and mix contents by gentle swirling or tapping. *Do not mix cells by pipetting.*

## Cloning Unstable Inserts

When cloning unstable inserts such as lentiviral DNA containing direct repeats (e.g. ViraPower™ Lentiviral Expression Kits), use the following modifications to reduce the chance of recombination between direct repeats:

- Select and culture transformants at 25°C–30°C.
- **Do not use** “rich” bacterial medias, because they tend to give rise to a greater number of unwanted recombinants.
- If your plasmid confers chloramphenicol resistance, select and culture transformants using LB medium containing 15–30 µg/mL chloramphenicol *in addition to* the antibiotic appropriate for selection of your plasmid.

## Transforming Competent Cells

Before starting the transformation procedure:

- Equilibrate a water bath to 42°C.
- Warm the vial of S.O.C. Medium (supplied with the kit) and LB Medium (if needed) to room temperature.
- Warm the selective plates in a 37°C incubator for 30 minutes (use 1 or 2 plates for each transformation). If you are including the pUC19 control, warm one LB agar plate containing 100 µg/mL ampicillin.

## Transformation Procedure

Use this procedure to transform One Shot<sup>®</sup> *ccdB* Survival<sup>™</sup> 2 T1<sup>R</sup> chemically competent cells. Include the pUC19 control plasmid DNA to verify the transformation efficiency. *Do not* use these cells for electroporation.

1. Thaw one vial of One Shot<sup>®</sup> cells on ice for each transformation.
2. Add 1–5  $\mu\text{L}$  of DNA (10 pg to 100 ng) into a vial of One Shot<sup>®</sup> cells and mix gently. *Do not mix by pipetting up and down.* For the pUC19 control, add 1  $\mu\text{L}$  (10 pg) of DNA into a separate vial of One Shot<sup>®</sup> cells and mix gently.
3. Incubate the vials on ice for 30 minutes.
4. Heat-shock the cells for 30 seconds at 42°C without shaking.
5. Remove the vials from the 42°C bath and place them on ice for 2 minutes.
6. Add 250  $\mu\text{L}$  of pre-warmed S.O.C. Medium to each vial.
7. Cap the vials tightly and shake horizontally at 37°C for 1 hour at 225 rpm in a shaking incubator.\*
8. Spread 25–100  $\mu\text{L}$  from each transformation on a pre-warmed selective plate. Plate two different volumes to ensure that at least one plate will have well-spaced colonies. For the pUC19 control, dilute the transformation mix 1:10 into LB Medium and plate 25–100  $\mu\text{L}$ .
9. Store the remaining transformation mix at 4°C. Additional cells may be plated out the next day, if desired.
10. Incubate plates overnight at 37°C. \*

\* **Note:** When cloning unstable inserts, incubate the cells at 25°C–30°C.

## Calculating Transformation Efficiency

Use the following formula to calculate the transformation efficiency as the number of transformants (in cfu) per  $\mu\text{g}$  of plasmid DNA.

$$\frac{\text{\# of colonies}}{\text{pg pUC19 DNA}} \times \frac{10^6 \text{ pg}}{\mu\text{g}} \times \frac{\text{volume of transformants}}{X \mu\text{L plated}} \times \text{dilution factor}$$

For example, if transformation of 10 pg of pUC19 DNA yields 50 colonies when 25  $\mu\text{L}$  of a 1:10 dilution is plated, then the transformation efficiency is:

$$\frac{50 \text{ colonies}}{10 \text{ pg DNA}} \times \frac{10^6 \text{ pg}}{\mu\text{g}} \times \frac{300 \mu\text{L}}{25 \mu\text{L plated}} \times 10 = 6 \times 10^8$$

## References

Bernard, P., Couturier, M., "Cell killing by the F plasmid CcdB protein involves poisoning of DNA-topoisomerase II complexes". *J. Mol. Biol.* (1992) Aug 5; 226(3):735-45.

Bernard, P., Kezdy, K.E., Van Melderen, L., Steyaert, J., Wyns, L., Pato, M.L., *et al.* (1993) The F plasmid CcdB protein induces efficient ATP-dependent DNA cleavage by gyrase. *J Mol Biol* 234: 534-541.

## Limited Product Warranty

Life Technologies Corporation and/or its affiliate(s) warrant their products as set forth in the Life Technologies' General Terms and Conditions of Sale found on Life Technologies' website at [www.lifetechnologies.com/termsandconditions](http://www.lifetechnologies.com/termsandconditions). If you have any questions, please contact Life Technologies at [www.lifetechnologies.com/support](http://www.lifetechnologies.com/support).

©2013 Life Technologies Corporation. All rights reserved. The trademarks mentioned herein are the property of Life Technologies Corporation or their respective owners.

**DISCLAIMER** - LIFE TECHNOLOGIES CORPORATION AND/OR ITS AFFILIATE(S) DISCLAIM ALL WARRANTIES WITH RESPECT TO THIS DOCUMENT, EXPRESSED OR IMPLIED, INCLUDING BUT NOT LIMITED TO THOSE OF MERCHANTABILITY, FITNESS FOR A PARTICULAR PURPOSE, OR NON-INFRINGEMENT. TO THE EXTENT ALLOWED BY LAW, IN NO EVENT SHALL LIFE TECHNOLOGIES AND/OR ITS AFFILIATE(S) BE LIABLE, WHETHER IN CONTRACT, TORT, WARRANTY, OR UNDER ANY STATUTE OR ON ANY OTHER BASIS FOR SPECIAL, INCIDENTAL, INDIRECT, PUNITIVE, MULTIPLE OR CONSEQUENTIAL DAMAGES IN CONNECTION WITH OR ARISING FROM THIS DOCUMENT, INCLUDING BUT NOT LIMITED TO THE USE THEREOF.

For support visit [lifetechnologies.com/support](http://lifetechnologies.com/support) or email [techsupport@lifetech.com](mailto:techsupport@lifetech.com)

[lifetechnologies.com](http://lifetechnologies.com)

