

Stealth™ RNAi Positive Control Duplexes (Human)

Cat. No. 12935-140

Cat. No. 12935-141

Cat. No. 12935-142

Stealth™ RNAi GAPDH Positive Control

Stealth™ RNAi Actin Positive Control

Stealth™ RNAi Cyclophilin B Positive Control

Description

Stealth™ RNAi Positive Control Duplexes (Human) are ideal for use in RNA interference (RNAi) experiments as a control for quantitatively assessing transfection efficiency and optimizing RNAi experiments following Stealth™ RNAi delivery in any human vertebrate cell line. Each Stealth™ RNAi Positive Control Duplex is functionally tested and guaranteed to knockdown their intended targets. The Stealth™ RNAi Positive Control Duplexes are supplied in a ready-to-use format. 1X RNA Annealing/Dilution Buffer is included for dilution of the Stealth™ RNAi stock solution, if desired.

Components

Item	Composition	Amount
Stealth™ RNAi Positive Control Duplex	20 μM in 1X RNA Annealing/Dilution Buffer	250 μl
1X RNA Annealing/Dilution Buffer	10 mM Tris-HCl, pH 8.0 20 mM NaCl 1 mM EDTA, pH 8.0	1 ml

Storage

Store the Stealth™ RNAi Positive Control Duplexes and the 1X RNA Annealing/Dilution Buffer at -20°C.

Stealth™ RNAi

Stealth™ RNAi is chemically modified dsRNA developed to overcome the limitations of traditional siRNA. Using Stealth™ RNAi for RNAi analysis offers the following advantages:

- Obtain effective target gene knockdown
- Eliminates sense strand off-target effects for higher specificity
- Exhibits enhanced stability for greater flexibility in RNAi analysis
- Avoids induction of cellular stress response pathways

For more information about Stealth™ RNAi, visit www.invitrogen.com/rnai.

Amount of Stealth™ RNAi Positive Control Duplex to Transfect

The amount of Stealth™ RNAi Positive Control Duplex required to achieve quantitative knockdown should be determined experimentally for each human cell line. **As a starting point, we recommend using 40 nM Stealth™ RNAi Positive Control Duplex for transfection.** To optimize transfection conditions, vary transfection reagent concentrations and the final concentrations of Stealth™ RNAi from 10 to 200 nM, as necessary for your cell line. Use the 1X RNA Annealing/Dilution Buffer supplied with the kit to dilute the Stealth™ RNAi Positive Control Duplex stock solution, if necessary.

Handling the Stealth™ RNAi Positive Control Duplexes

- Thaw Stealth™ RNAi Positive Control Duplex stock solutions on ice or at room temperature. After use, return to -20°C storage.
- Multiple freeze/thaw cycles are permitted without loss of activity if stock solutions are handled properly.
- Ensure that the stock solutions do not become contaminated with RNase.

General Guidelines for Transfection

- Use a transfection reagent suitable for delivery of Stealth™ RNAi to mammalian cells. **For optimal results, we recommend using Lipofectamine™ 2000 Reagent (Cat. no. 11668-027) or Lipofectamine™ RNAiMax (Cat. no. 13778-075) available from Invitrogen.** For a protocol to transfect Stealth™ RNAi into cells using Lipofectamine™ 2000 or Lipofectamine™ RNAiMax, see the RNAi resource page at www.invitrogen.com/rnai; click on Protocols. **Note:** Do not use the DNA transfection protocol included with the product.
- Use low-passage cells, and make sure that cells are healthy and greater than 90% viable before transfection.
- Transfect cells at the density recommended by the manufacturer of your transfection reagent.
- Assay for target gene knockdown at a suitable time period (typically 24 to 72 hours) after transfection.
- For a convenient tool to assess transfection efficiency and Stealth™ RNAi uptake, we recommend using the BLOCK-iT™ Fluorescent Oligo (Cat. no. 2013) available from Invitrogen.

Limited Use Label License No. 196: Stealth™ RNAi

The purchase of this product conveys to the buyer the non-transferable right to use the purchased amount of the product and components of the product in research conducted by the buyer (whether the buyer is an academic or for-profit entity). The buyer cannot sell or otherwise transfer (a) this product (b) its components or (c) materials made using this product or its components to a third party or otherwise use this product or its components or materials made using this product or its components for Commercial Purposes. The buyer may transfer information or materials made through the use of this product to a scientific collaborator, provided that such transfer is not for any Commercial Purpose, and that such collaborator agrees in writing (a) not to transfer such materials to any third party, and (b) to use such transferred materials and/or information solely for research and not for Commercial Purposes. Commercial Purposes means any activity by a party for consideration and may include, but is not limited to: (1) use of the product or its components in manufacturing; (2) use of the product or its components to provide a service, information, or data; (3) use of the product or its components for therapeutic, diagnostic or prophylactic purposes; or (4) resale of the product or its components, whether or not such product or its components are resold for use in research. Invitrogen Corporation will not assert a claim against the buyer of infringement of patents owned by Invitrogen and claiming this product based upon the manufacture, use or sale of a therapeutic, clinical diagnostic, vaccine or prophylactic product developed in research by the buyer in which this product or its components was employed, provided that neither this product nor any of its components was used in the manufacture of such product. If the purchaser is not willing to accept the limitations of this limited use statement, Invitrogen is willing to accept return of the product with a full refund. For information on purchasing a license to this product for purposes other than research, contact Licensing Department, Invitrogen Corporation, 1600 Faraday Avenue, Carlsbad, California 92008. Phone (760) 603-7200. Fax (760) 602-6500.

©2007 Invitrogen Corporation. All rights reserved.

For research use only. Not intended for any animal or human therapeutic or diagnostic use.