

ProcartaPlex™ Human SARS-CoV2 Spike Omicron (B.1.1.529) Neutralizing Ab Simplex USER GUIDE

Using Magnetic Beads for Serum and Plasma

Catalog Number EPX010-16019N-901

Publication Number MAN0026539

Revision A.0 (30)



Bender MedSystems GmbH | Campus Vienna Biocenter 2 | 1030 Vienna, Austria

For descriptions of symbols on product labels or product documents, go to [thermofisher.com/symbols-definition](https://www.thermofisher.com/symbols-definition).

The information in this guide is subject to change without notice.

DISCLAIMER: TO THE EXTENT ALLOWED BY LAW, THERMO FISHER SCIENTIFIC INC. AND/OR ITS AFFILIATE(S) WILL NOT BE LIABLE FOR SPECIAL, INCIDENTAL, INDIRECT, PUNITIVE, MULTIPLE, OR CONSEQUENTIAL DAMAGES IN CONNECTION WITH OR ARISING FROM THIS DOCUMENT, INCLUDING YOUR USE OF IT.

Revision history: Pub. No. MAN0026539

Revision	Date	Description
A.0	7 March 2022	New document for ProcartaPlex™ Human SARS-CoV2 Spike Omicron (B.1.1.529) Neutralizing Ab Simplex.

End User License Agreement for Luminex Assay: Notice to Purchaser: By opening the packaging containing this Assay Product (which contains fluorescently labeled microsphere beads authorized by Luminex Corporation) or using this Assay Product in any manner, you are consenting and agreeing to be bound by the End User Terms and Conditions and the End User License Agreement available at <https://www.luminexcorp.com/enduser-terms-and-conditions/#sec3>. If you do not agree to all of the terms and conditions, you must promptly return this Assay Product for a full refund prior to using it in any manner.

TRADEMARKS: All trademarks are the property of Thermo Fisher Scientific and its subsidiaries unless otherwise specified. A subset of reagents are sourced from ACROBiosystems Inc. xMAP, Luminex 100/200, MAGPIX, INTELLIFLEX and FLEXMAP 3D are trademarks of Luminex Corp. BioPlex is a trademark of Bio-Rad Laboratories Inc. Bio-Rad is a registered trademark of Bio-Rad Laboratories Inc.

©2022 Thermo Fisher Scientific Inc. All rights reserved.

Contents

■	CHAPTER 1	Product information	5
		Product description	5
		Technology overview	5
		Contents and storage	6
		Required materials not supplied	6
■	CHAPTER 2	Methods	7
		Workflow	7
		Procedural guidelines	9
		Prepare samples	9
		Prepare plasma samples	9
		Prepare serum samples	9
		Dilute samples	10
		Prepare the reagents	10
		Prepare 1X wash buffer	10
		Prepare the positive control	10
		Prepare 1X Biotinylated Detection Reagent	11
		Perform the assay	11
		Define the plate map	11
		Add magnetic beads to the 96-well Flat Bottom plate	11
		Wash the magnetic beads	11
		Add Assay Diluent, controls, and samples, then incubate	12
		Add the 1X Biotinylated Detection Reagent, then incubate	12
		Add Streptavidin-PE (SA-PE), then incubate	12
		Add Reading Buffer, then run the plate on a Luminex™ instrument	13
		Set up the instrument	13
		Re-read the plate after a failed run	14
		Analyze the results	14
■	APPENDIX A	Troubleshooting	15
■	APPENDIX B	Performance characteristics	17
		Performance characteristics	17

- **APPENDIX C Recommended plate layouts** 18
 - Recommended plate layouts 18
- **APPENDIX D Documentation and support** 19
 - Customer and technical support 19
 - Limited product warranty 19



Product information



WARNING! Read the Safety Data Sheets (SDSs) and follow the handling instructions. Wear appropriate protective eyewear, clothing, and gloves. Safety Data Sheets (SDSs) are available from [thermofisher.com/support](https://www.thermofisher.com/support).

Product description

This user manual is for a ProcartaPlex™ Immunoassay Kit to perform multiplexed protein measurements from serum and plasma samples using Luminex™ magnetic bead technology. Other biological samples may be suitable for use in the assay.

IMPORTANT! ProcartaPlex™ Human SARS-CoV2 Spike Omicron (B.1.1.529) Neutralizing Ab Simplex (Cat. No. [EPX010-16019N-901](#)) **cannot** be combined with other ProcartaPlex™ Panels or Simplex Kits.

For the most current version of user documentation, visit our website www.thermofisher.com.

Technology overview

ProcartaPlex™ Immunoassays incorporate magnetic microsphere technology licensed from Luminex™ to enable the simultaneous detection and quantitation of multiple protein targets in diverse matrices. The platform allows the simultaneous detection from a single sample of up to 80 protein targets on the Luminex™ 200™ and FLEXMAP 3D™ platforms, and 50 protein targets on the MAGPIX™ platform.

This assay is designed to detect the level of SARS-CoV2 neutralizing antibodies of the Omicron variant (B.1.1.529) in serum and plasma. It allows direct comparison of neutralizing potential of antibodies towards five described variants of SARS-CoV2, namely wild-type and variants B.1.1.7 (α, originating in the United Kingdom), B.1.351 (β, originating in South Africa), P1 (γ, originating in Brazil), and B.1.617.2 (δ, originating in India). The capture beads are coupled with spike S1 protein from the SARS-CoV2 Omicron variant (B.1.1.529), containing the known relevant mutations in the spike protein. Samples with neutralizing antibodies compete with an excess amount of biotinylated ACE2 (Biotinylated Detection Reagent). Any biotinylated ACE2 that binds to the proteins on the beads will produce a median fluorescence intensity (MFI) signal. Signals are inversely proportional to the level of neutralizing antibodies since this is a competitive immunoassay. The negative control (Assay Diluent without neutralizing antibodies) will give the highest MFI values. Neutralizing antibodies in the sample or in the positive control will reduce the signal. The neutralization (%) for samples are calculated from the MFI of the samples and the MFI of the negative control, see “Analyze the results” on page 14.

Contents and storage

ProcartaPlex™ Immunoassay Kits contain the components listed below. For quantities and details of components supplied, see the Certificate of Analysis and Table 1. Store kit at 2°C–8°C. See the expiration date on the label. Do not use expired product.

Contents	Amount
Bead Mix Neutr Ab (magnetic beads, 1X) ^[1]	1 × 5 mL
Biotinylated Detection Reagent (lyophilized)	2 vials
Detection AB Diluent	1 × 3 mL
Control Neutr Ab (lyophilized)	2 vials
Streptavidin-PE (SA-PE, 1X)	1 × 5 mL
Wash Buffer (10X) ^[1]	1 × 25 mL
Assay Diluent ^[1]	1 × 60 mL
Reading Buffer ^[1]	1 × 40 mL
96-well Flat Bottom Plate	1 plate
Black Microplate Lid	1 lid
Plate Seals	8 seals

^[1] Contains sodium azide. Read the Safety Data Sheets (SDSs) and follow the handling instructions. Safety Data Sheets (SDSs) are available from [thermofisher.com/support](https://www.thermofisher.com/support).

Table 1 Bead mix neutralizing antibodies

Analyte	Beadsets
SARS-CoV2 o B.1.1.529 neutralizing antibodies	57

Required materials not supplied

- MAGPIX™, Luminex™ 200™, FLEXMAP 3D™, or equivalent Luminex™-based instrument
- Glass-distilled or deionized water
- Adjustable single and multichannel pipettes with disposable tips
- Multichannel pipette reservoir
- Beakers, flasks, and cylinders for preparation of reagents
- Hand-held magnetic plate washer
- Vortex mixer
- Microtiter™ plate shaker

Workflow

Assay protocol

Prepare samples

Prepare plasma samples and serum samples.

Prepare the reagents

Prepare 1X wash buffer and positive control.

Define the plate map

Mark sample and control wells (see page 18).

Add magnetic beads to the 96-well Flat Bottom plate

1. Vortex the capture beads for 30 seconds, then add 50 μL of the capture beads to each well.
2. Remove the liquid.

Note: Wash the plate after adding the beads.

Add Assay Diluent, controls, and samples, then incubate

1. Add the following reagents according to the sample type.
 - a. **Positive control:** 25 μL of Assay Diluent, followed by 25 μL of the reconstituted positive control.
 - b. **Negative control:** 50 μL of 1X Assay Diluent.
 - c. **Serum and plasma samples:** 25 μL of 1X Assay Diluent, followed by 25 μL of pre-diluted samples.
2. Seal the plate, then incubate with shaking at room temperature for 120 minutes.
3. Wash the plate twice.

Add the 1X Biotinylated Detection Reagent, then incubate

1. Add 25 μL of prediluted 1X Detection Antibody mixture to each well.
2. Seal the plate, then incubate with shaking at room temperature for 30 minutes.
3. Wash the plate twice.

Assay protocol

Add Streptavidin-PE (SA-PE), then incubate

1. Add 50 μL of SA-PE solution to each well.
2. Seal the plate, then incubate with shaking at room temperature for 30 minutes.
3. Wash the plate twice.

Add Reading Buffer, then run the plate on a Luminex™ instrument

1. Add 120 μL of Reading Buffer to each well.
2. Seal the plate, then incubate with shaking at room temperature for 5 minutes.
3. Remove the Plate Seal, then run the plate on a Luminex™ instrument.

Analyze the results

Analyze the results using the neutralization (%) equation.

Procedural guidelines

- Thoroughly read this user manual and the product insert that is included with the assay kit. The product insert contains specific instructions for proper use of your kit.
- For Luminex™ 200™ and FLEXMAP 3D™ instruments, initiate the startup protocol to warm up the lasers for at least 30 minutes.
 - Ensure that the Luminex™ machine is calibrated according to the manufacturer's instructions.
 - The MAGPIX™ instrument does not require additional warm up.
- When working with samples and standards, change the pipette tips after every transfer and avoid creating bubbles when pipetting.
- During the incubation steps, cover the 96-well Flat Bottom Plate with the Black Microplate Lid provided in the kit to minimize exposure of the beads to light.
- Do not invert the 96-well Flat Bottom Plate during the assay or allow contents from one well to mix with another well.
- Use a multichannel pipette and reagent reservoirs whenever possible to achieve optimal assay precision.

Prepare samples

- For frozen samples: thaw samples on ice and mix well by vortexing, followed by centrifugation at $10,000 \times g$ for 5–10 minutes to remove particulates.

IMPORTANT! Avoid multiple freeze/thaw cycles.

- For samples with high lipid content: centrifuge at $10,000 \times g$ for 10 minutes and transfer contents to a new tube.

Prepare plasma samples

1. Collect samples in sodium citrate or EDTA tubes.

When using heparin as an anticoagulant, no more than 10 IU of heparin per mL of blood collected should be used because an excess of heparin may give falsely high values of some of the analytes.
2. Centrifuge samples at $1,000 \times g$ at 4°C for 10 minutes within 30 minutes of collection.
3. Collect the plasma fraction. Use immediately or aliquot and store at –80°C.

Prepare serum samples

For best results, centrifuge serum samples immediately before running the assay.

1. Allow blood to clot for 20–30 minutes at 20°C–25°C.
2. Centrifuge at $1,000 \times g$ for 10 minutes at 20°C–25°C.
3. Use immediately or aliquot and store at –80°C.

Dilute samples

Levels may vary between different samples. You may need to further dilute your samples if the analyte concentration exceeds the assay upper limit of quantitation (ULOQ). When preparing dilution of serum and plasma samples, use Assay Diluent. The suggested dilution for serum/plasma is 1:100.

1. Prepare a 1:100 external dilution of serum or plasma samples in Assay Diluent (e.g., add 5 μ L of sample to 495 μ L of Assay Diluent).
2. Mix diluted samples gently.

Prepare the reagents

Prepare 1X wash buffer

Bring the Wash Buffer Concentrate (10X) to room temperature and vortex for 15 seconds. Mix 20 mL of the Wash Buffer Concentrate (10X) with 180 mL ddH₂O. Store 1X wash buffer at 2°C to 8°C for up to 6 months if needed.

Note: Wash Buffer Concentrate (10X) volume might not be sufficient if using an automated plate washer. For bulk orders use Cat. No. [EPX-66666-001](#).

Prepare the positive control

Lyophilized neutralizing antibodies are used as the positive control. Assay Diluent is used as the negative control.

Note: After use, discard any remaining positive controls.

1. Centrifuge the control vial at 2,000 $\times g$ for 10 seconds.
2. Add Assay Diluent to the control vial (see the label for the appropriate reconstitution volume).
3. Gently vortex the vial for 10 seconds and centrifuge at 2,000 $\times g$ for 10 seconds to collect contents at the bottom of the vials.
4. Incubate at room temperature for 10 minutes to ensure complete reconstitution.
The reconstituted positive control is ready to use, no pre-dilution is needed.

Prepare 1X Biotinylated Detection Reagent

Reconstitute the Biotinylated Detection Reagent (lyophilized) 15 minutes before use.

Note: The reconstituted Biotinylated Detection Reagent should be kept on ice.

1. Centrifuge the Biotinylated Detection Reagent (lyophilized) vial at $2,000 \times g$ for 10 seconds.
2. Add Assay Diluent into the vial (see the label for the appropriate reconstitution volume).
3. Gently vortex the vial for 10 seconds and centrifuge at $2,000 \times g$ for 10 seconds to collect contents at the bottom of the vials.
4. Incubate for 10 minutes to ensure complete reconstitution.
5. Add 60 μL of reconstituted Biotinylated Detection Reagent concentrate (50X) to the mixing bottle, then add the Detection AB Diluent to a final volume of 3 mL if assaying a whole plate (otherwise scale the volume accordingly).

Perform the assay

Define the plate map

Mark sample and control wells using the plate map at the end of this manual (see page 18).

Add magnetic beads to the 96-well Flat Bottom plate

The panel is provided with magnetic beads at a 1X working concentration.

1. Vortex the bead vials for 30 seconds.
2. Pour the bead mix into a disposable reservoir, then add 50 μL of bead solution to each well of the plate. Use a multichannel pipette for this step as well as for the steps that follow.

Wash the magnetic beads

1. Securely insert the 96-well Flat Bottom Plate into the hand-held magnetic plate washer, ensure that the plate is held in place by the tabs, and wait 2 minutes to allow the beads to accumulate on the bottom of each well.
2. Remove the liquid in the wells by quickly inverting the hand-held magnetic plate washer and 96-well Flat Bottom Plate assembly over a sink or waste container. Do not remove the 96-well Flat Bottom Plate from the hand-held magnetic plate washer.
3. Blot the inverted assembly onto several layers of paper towels or absorbent surface to remove any residual solution.
4. Add 150 μL of 1X wash buffer into each well and wait 30 seconds to allow the beads to accumulate on the bottom of each well.

5. Remove the wash buffer in the wells by quickly inverting the hand-held magnetic plate washer and 96-well Flat Bottom Plate assembly over a sink or waste container. Do not remove the 96-well Flat Bottom Plate from the hand-held magnetic plate washer.
6. Blot the inverted assembly onto several layers of paper towels or absorbent surface to remove any residual solution.
7. Remove the 96-well Flat Bottom Plate from the hand-held magnetic plate washer and proceed to the next step.

Add Assay Diluent, controls, and samples, then incubate

1. For the positive control, add 25 μ L of Assay Diluent to the dedicated wells of the plate, followed by 25 μ L of the reconstituted positive control.

Note: The reconstituted positive control is ready to use, no pre-dilution is needed.

2. For the negative control, add 50 μ L of 1X Assay Diluent to the dedicated wells.
3. For serum and plasma samples, add 25 μ L of 1X Assay Diluent to the appropriate wells, followed by 25 μ L of pre-diluted samples.
4. Seal the plate with the provided Plate Seal. Cover the plate with the Black Microplate Lid, then incubate at room temperature for 120 minutes on a plate shaker set at 500 rpm.
5. Wash the 96-well plate twice. Follow “Wash the magnetic beads” on page 11.

Add the 1X Biotinylated Detection Reagent, then incubate

1. Add 25 μ L of prediluted 1X Biotinylated Detection Reagent mixture to each well.
2. Seal the plate with a new Plate Seal, cover the plate with the Black Microplate Lid, then incubate at room temperature for 30 minutes on a plate shaker set at 500 rpm.
3. Wash the 96-well plate twice. Follow “Wash the magnetic beads” on page 11.

Add Streptavidin-PE (SA-PE), then incubate

1. Add 50 μ L of SA-PE solution to each well.
2. Seal the plate with a new Plate Seal, cover the plate with the Black Microplate Lid, then incubate at room temperature for 30 minutes on a plate shaker set at 500 rpm.
3. Wash the 96-well plate twice. Follow “Wash the magnetic beads” on page 11.

Add Reading Buffer, then run the plate on a Luminex™ instrument

1. Add 120 µL of Reading Buffer into each well.
2. Seal the plate with a new Plate Seal, cover the plate with the Black Microplate Lid, then incubate at room temperature for 5 minutes on a plate shaker set at 500 rpm.
3. Remove the Plate Seal, then run the plate on a Luminex™ instrument.

Note: For details about instrument settings, see the next section.

Set up the instrument

Instrument	Acquisition volume	Timeout (optional)	Bead type	DD gate	Reporter gain	Minimum bead count
MAGPIX™	50 µL	N/A	N/A	N/A	Standard PMT	50
INTELLIFLEX™	30 µL	40 seconds	MagPlex™	4,000–13,000	Standard PMT	50
FLEXMAP 3D™ Luminex™ 200™	50 µL	60 seconds	MagPlex™	7,500–25,000	Standard PMT	50
Bio-Rad™ BioPlex™	50 µL	60 seconds	MagPlex™	5,000–25,000	Standard PMT	50

Before running the assay, ensure that the probe height has been calibrated with the 96-well Flat Bottom Plate that is supplied with the kit. Failure to adjust the probe height can cause damage to the instrument or low bead count. The Luminex™ system allows for calibration of low and high RP1 target values. We recommend RP1 low target value settings for ProcartaPlex™ immunoassays. When entering the information into the Luminex™ Acquisition Software, refer to the Certificate of Analysis provided with the kit for bead region.

For Luminex™ 200™ and FLEXMAP 3D™ instruments, initiate the startup protocol to warm up the lasers for at least 30 minutes.

- Ensure that the Luminex™ instrument is calibrated according to the manufacturer's instructions.
- The MAGPIX™ instrument does not require additional warm-up.

Re-read the plate after a failed run

If there is a malfunction of the Luminex™ instrument or software during the run, the 96-well Flat Bottom Plate can be re-read using the following instructions:

1. Remove the 96-well Flat Bottom Plate from the instrument, then insert the 96-well Flat Bottom Plate into the hand-held magnetic plate washer.
2. Wait 2 minutes.
3. Remove the buffer in the wells by quickly inverting the 96-well Flat Bottom Plate over a sink or waste container. Do not remove the 96-well Flat Bottom Plate from the hand-held magnetic plate washer.
4. Blot the assembly onto several layers of paper towels to remove any residual solution.
5. Resuspend the beads in 120 µL of Reading Buffer.
6. Remove from the hand-held magnetic plate washer.
7. Seal the 96-well Flat Bottom Plate with a new Plate Seal and lid.
8. Shake at 500 rpm for 5 minutes at room temperature.

Note: The assayed samples may take longer to read because there will be less beads in the wells.

Analyze the results

Since this is a competitive assay, the negative control will give the highest MFI values without any reduction of the signal due to neutralizing antibodies. See “Performance characteristics” on page 17 for expected MFI levels for the negative and positive controls.

Calculate the ratio of values using the following equation:

$$\text{Neutralization (\%)} = (1 - (\text{MFI of samples} / \text{MFI of negative control})) \times 100$$

- >25% = Positive
- <25% = Negative

The cut-off level was determined at 25% after screening 160 healthy samples.



Troubleshooting

Observation	Possible cause	Recommended action
Low flow rate	Samples/beads were stuck in flow cell.	Remove the 96-well plate, then perform a wash and rinse cycle.
High CVs	Contamination occurred from re-using the Plate Seal.	Use a new Plate Seal for each incubation step.
	Washing was incomplete.	After adding the controls and samples, it is very important that any excess controls are removed during the wash step.
	Cross-well contamination occurred.	Avoid splashing the wash buffer into adjacent wells during wash steps.
	Poor pipetting techniques were used.	Use a multichannel pipettor and careful pipette techniques. Avoid touching pipette tips to the sides of the wells when adding wash buffer.
Limited dynamic range for BioPlex™ software users	The instrument was calibrated at high PMT settings.	Calibrate the instrument using the CAL2 Low RP1 target value.
Low bead count	The volume of bead solution used was too low.	Add 120 µL Reading Buffer to each well, then shake at 500 rpm for 5 minutes at room temperature to resuspend the beads before reading on the Luminex™ instrument.
	The beads aggregated.	Vortex the bead suspension well before using in the assay and ensure that the beads are properly mixed during the incubation steps.
	The dyes contained in the beads were photo-bleached from overexposure to light.	Store the bead solution and the 96-well plate in the dark.
	Samples caused the instrument to clog.	Remove the 96-well Flat Bottom Plate, then perform a wash and rinse of the instrument. Rerun the assay with further diluted samples.
	The probe height was incorrect.	See the Luminex™ manual for instructions to adjust the needle to the proper height.
	The instrument needle was partially clogged.	Replace or clean the needle according to the recommendations of the manufacturer.
	Beads adhered to the bottom of the plate.	Confirm that the plate shaker is set to 500 rpm. Shake the plate for at least 5 minutes before reading.
	An air bubble formed in the sample loop.	See the Luminex™ manual for instructions to remove the air bubble.



Observation	Possible cause	Recommended action
Low signal or sensitivity of the negative control	Shaking was insufficient, or the Biotinylated Detection Reagent was not reconstituted and diluted properly.	Prepare fresh Biotinylated Detection Reagent following the instructions provided (see “Prepare 1X Biotinylated Detection Reagent” on page 11).
High signal or sensitivity of the positive control, or no difference compared to the negative control	The positive control was not reconstituted or diluted properly.	Prepare fresh controls following the instructions provided (see “Prepare the positive control” on page 10).



Performance characteristics

Performance characteristics

The data presented in this section are examples of typical data, and levels can vary depending upon the instrument used for measurement, and according to the conditions under which the assay is performed (e.g., operator, pipetting technique, washing technique, or temperature effects).

The positive control (Neutralizing Antibodies (lyophilized) reconstituted with Assay Diluent) and the negative control (Assay Diluent) provided in the kit were run in duplicate according to the protocol. The results were measured with a Luminex™ 200™ instrument.

Table 2 Performance of positive control and negative control (assay diluent)

Analyte	Negative control MFI	Positive control MFI	Neutralization
SARS-CoV2 o B.1.1.529 neutralizing antibodies	2321	784	66%

Neutralization of the positive control is expected to be at least 30%.



Recommended plate layouts

Recommended plate layouts

Table 3 Layout for using the negative control (no competition, high MFI signal) and positive control

	1	2	3	4	5	6	7	8	9	10	11	12
A	Neg. Control	Neg. Control										
B	Pos. Control	Pos. Control										
C												
D												
E												
F												
G												
H												



Documentation and support

Customer and technical support

Visit [thermofisher.com/support](https://www.thermofisher.com/support) for the latest service and support information.

- Worldwide contact telephone numbers
- Product support information
 - Product FAQs
 - Software, patches, and updates
 - Training for many applications and instruments
- Order and web support
- Product documentation
 - User guides, manuals, and protocols
 - Certificates of Analysis
 - Safety Data Sheets (SDSs; also known as MSDSs)

Note: For SDSs for reagents and chemicals from other manufacturers, contact the manufacturer.

Limited product warranty

Life Technologies Corporation and/or its affiliate(s) warrant their products as set forth in the Life Technologies' General Terms and Conditions of Sale at www.thermofisher.com/us/en/home/global/terms-and-conditions.html. If you have any questions, please contact Life Technologies at www.thermofisher.com/support.

