

Disposable Column Trial Pack

29925

0031.3

Number

29925

Description**Disposable Column Trial Pack**, 2 each of 3 different column sizes, plus accessories**Trial Pack Contents:****Disposable Polystyrene Columns**, 2 columns, appropriate for 0.5-2.0mL gel beds**Disposable Polypropylene Columns**, 2 columns, appropriate for 1-5mL gel beds**Disposable Polypropylene Columns**, 2 columns, appropriate for 2-10mL gel beds**Porous Polyethylene Discs**, 3 × 5 discs (5 discs each of 7, 10 and 15mm diameter for small, medium and large column sizes, respectively), 30µm pore size, for use above and below the packed gel bed**Top Caps**, 3 × 2 push-on, water-tight polypropylene caps for the column tops**Bottom Caps**, 3 × 2 push-on, water-tight polypropylene caps for the column tips**Serum Separators**, 3 units, each 10cm long but different diameters for insertion of porous discs into small, medium and large columns, respectively**Disposable Polypropylene Funnel**, 1 unit, for use as a column extender to extend volume of column reservoir by an additional 25mL**Storage:** Upon receipt store at room temperature. Product shipped at ambient temperature.

Introduction

The Thermo Scientific Disposable Column Trial Pack is an economical and convenient set of three different columns sizes and accessories for packing gel supports for gravity-flow affinity and size exclusion chromatography procedures. Components in the trial pack are used in many of our affinity purification products. The durable columns have leak-proof top and bottom caps that enable end-over-end mixing of reaction solutions and long-term storage of packed columns without risk of dehydration.

The porous polyethylene discs, when positioned above as well as below the packed gel bed (Figure 1, see last page), confer a unique stop-flow action to the column. Solutions applied to a column will automatically stop dripping from the column tip when the surface of solution drains down to the top disc, thereby keeping air bubbles from being drawn into the gel bed and preventing the gel from drying out if the column is left unattended for brief periods of time. Also, use of a top porous disc prevents suspension of the gel when pouring or pipetting solutions into the column and ensures that the quantity of solution applied to the column always equals the volume of eluate emerging from the tip.

The reservoir (wide-mouth) portion of each column holds 3-4mL of solution. The volume of this reservoir can be increased to nearly 30mL using the included Disposable Polypropylene Funnel, which fits onto each column size. Additional funnels are available in packages of 50 units, and each of the column sizes and their matched accessories are available in packages of 100 units (see Related Thermo Scientific Products).

Procedure for Packing Gel into a Column

1. Equilibrate column, degassed 50% gel slurry and degassed buffer solution (or water) to room temperature.
2. Secure a bottom cap on the column tip and place the column in a 16 × 125mm glass test tube (in a test tube rack) or clamp the column upright in a laboratory stand.
3. Add a sufficient volume of degassed buffer/water to the column to fill it up to the reservoir (wide-mouth) portion, then gently tap the end and side of the column to dislodge any air bubbles.
Note: Adding detergent such as Thermo Scientific Triton[®] X-100 Surfact-Amps Solution (Product No. 28314) to the buffer solution to a concentration of 0.03-0.05% will make it easier to remove air bubbles, especially from the conical tip of the column.
4. Float the appropriate diameter porous disc on top of the liquid within the column.
5. Using the open end of the appropriate diameter serum separator, push the disc evenly to the bottom of the column.
6. Decant most of the liquid from the empty column, being sure to avoid getting air bubbles in the tip region of the column below the inserted disc. Place the column back in its stand with bottom cap still in place.
7. Add sufficient volume of degassed gel slurry to obtain the desired settled gel volume.
8. Allow gel to settle in the column for at least 30 minutes.
9. Position a second porous disc on top of the settled gel bed by floating it on the liquid within the column and pushing it down to just above the settled gel. Leave 1-2mm of space between the top of the gel bed and the top disc; do not compress the gel bed.
10. Wash the inside top part of the column with buffer/water to remove residual gel that may have remained along the sides during packing.
11. The packed column is now ready for storage or use.
 - Store the packed column upright at 4°C with the gel bed submerged under 1-2mL of buffer/water and a top column cap securely in place. Sodium azide added to the storage buffer/water to a concentration of 0.02% will help to prevent microbial growth in carbohydrate-containing gel supports.
 - Always remove the top cap before the bottom cap to avoid drawing air bubbles down into the gel bed.
 - Prevent air bubbles from forming in the gel bed by using only degassed buffer and sample solutions. Degassing involves subjecting a solution to vacuum to “boil” off excess dissolved air (see Additional Information section).

Additional Information

A. Please visit the web site for additional information about this products, including the following items:

- Tech Tip #29: Degas buffers for use in affinity and gel filtration columns
- Tech Tip #7: Remove air bubbles from columns to restore flow rate

B. Chemical Compatibility and Physical Properties of Plastic Columns

Polystyrene and Polypropylene Columns are compatible with dilute or weak acids, aliphatic alcohols and bases. Strong or concentrated acids can be used for brief periods. The columns are not compatible with aldehydes, esters, hydrocarbons, ketones or strong oxidizing agents. Polypropylene Columns (Product No. 29922, 29924) generally are more resistant to organic solvents, at least for brief periods of exposure, than Polystyrene Columns (Product No. 29920). Disposable Plastic Columns are intended for use between 4°C and 50°C, and they will deform above 90°C. The polyethylene porous discs used in the columns are more resistant to chemicals of all kinds than the column plastic.

Related Thermo Scientific Products

89896	Pierce® Centrifuge Columns, 2mL
89897	Pierce Centrifuge Columns, 5mL
29920	Disposable Polystyrene Columns, 100 columns plus accessories, appropriate for 0.5-2.0mL gel beds
29922	Disposable Polypropylene Columns, 100 columns plus accessories, appropriate for 1-5mL gel beds
29924	Disposable Polypropylene Columns, 100 columns plus accessories, appropriate for 2-10mL gel beds
69700	Spin Cup Columns, microcentrifuge cup columns for 50-200µL chromatography gel
69705	Mini-Spin Columns Plus Accessories, microcentrifuge columns for 20-400µL chromatography gel
69710	Resin Separators

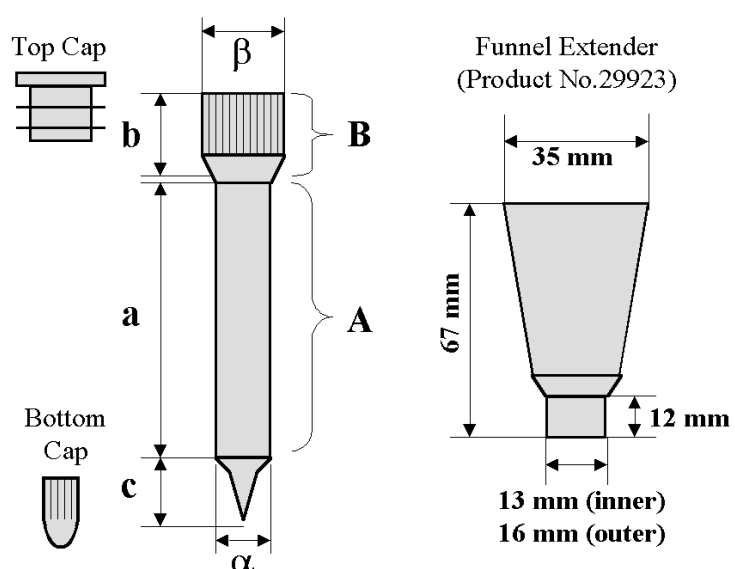
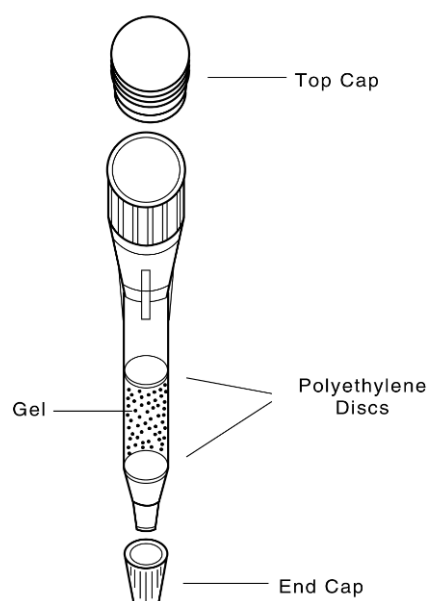


Figure 1. Assembly of Disposable Columns.

Figure 2. Column parameters described in Table 1.

Table 1. Dimensions (in millimeters) and Volumes (in milliliters) of Pierce Disposable Columns.

Column Product No.	Column Bed			Column Reservoir			Tip	Total Column			
	a Height	α Diameter Inner Outer		A Volume	b Height	β Diameter Inner Outer		B Volume	c Height	Height (w/o cap)	Height (with cap)
29920	60	7	9	2.3	20	16	19	4.0	17	102	107
29922	70	10	13	5.5	14	16	19	2.8	20	107	110
29924	70	15	18	12.3	14	16	19	2.8	20	107	110

This product ("Product") is warranted to operate or perform substantially in conformance with published Product specifications in effect at the time of sale, as set forth in the Product documentation, specifications and/or accompanying package inserts ("Documentation") and to be free from defects in material and workmanship. Unless otherwise expressly authorized in writing, Products are supplied for research use only. No claim of suitability for use in applications regulated by FDA is made. The warranty provided herein is valid only when used by properly trained individuals. Unless otherwise stated in the Documentation, this warranty is limited to one year from date of shipment when the Product is subjected to normal, proper and intended usage. This warranty does not extend to anyone other than the original purchaser of the Product ("Buyer").

No other warranties, express or implied, are granted, including without limitation, implied warranties of merchantability, fitness for any particular purpose, or non infringement. Buyer's exclusive remedy for non-conforming Products during the warranty period is limited to replacement of or refund for the non-conforming Product(s).

There is no obligation to replace Products as the result of (i) accident, disaster or event of force majeure, (ii) misuse, fault or negligence of or by Buyer, (iii) use of the Products in a manner for which they were not designed, or (iv) improper storage and handling of the Products.

Current product instructions are available at www.thermoscientific.com/pierce. For a faxed copy, call 800-874-3723 or contact your local distributor.

© 2011 Thermo Fisher Scientific Inc. All rights reserved. Unless otherwise indicated, all trademarks are property of Thermo Fisher Scientific Inc. and its subsidiaries. Printed in the USA.