

Stealth™ RNAi Reporter Control Duplexes

Cat. No. 12935-145

Cat. No. 12935-146

Cat. No. 12935-147

Cat. No. 12935-148

GFP Reporter Control

Luciferase Reporter Control

LacZ Reporter Control

β-lactamase Reporter Control

Description

Stealth™ RNAi Reporter Control Duplexes are ideal for use in RNA interference (RNAi) experiments to help you optimize your transfection conditions in any vertebrate cell line. Each Stealth™ RNAi Reporter Control Duplex is designed to efficiently knock-down its intended target and minimize sequence homology to any other known vertebrate transcript. The Stealth™ RNAi Reporter Control Duplexes are supplied in a ready-to-use format and 1X RNA Annealing/Dilution Buffer is included for dilution of the Stealth™ RNAi stock solution, if desired.

Components

Each Stealth™ RNAi Negative Control Duplex includes the appropriate Stealth™ RNAi Reporter Control Duplex and 1 ml of 1X RNA Annealing/Dilution Buffer, with the following compositions:

Item	Composition	Amount
Stealth™ RNAi Reporter Control Duplex	20 μM in 1X RNA Annealing/Dilution Buffer	250 μl
1X RNA Annealing/Dilution Buffer	10 mM Tris-HCl, pH 8.0 20 mM NaCl 1 mM EDTA, pH 8.0	1 ml

Storage

Store the Stealth™ RNAi Reporter Control Duplexes and the 1X RNA Annealing/Dilution Buffer at -20°C.

Part No.: 12935.rep.pps

Rev. Date: 19 April 2007

Stealth™ RNAi

Stealth™ RNAi is chemically modified dsRNA developed to overcome the limitations of traditional siRNA. Using Stealth™ RNAi for RNAi analysis offers the following advantages:

- Obtain effective target gene knockdown
- Eliminates sense strand off-target effects for higher specificity
- Exhibits enhanced stability for greater flexibility in RNAi analysis
- Avoids induction of cellular stress response pathways

For more information about Stealth™ RNAi, see www.invitrogen.com/rnai.

Stealth™ RNAi Reporter Control Duplex Target Genes

The Stealth™ RNAi Reporter Control Duplex target the following genes:

- **Stealth™ RNAi GFP Reporter Control** targets the following fluorescent proteins: Emerald Green Fluorescent Protein (EmGFP), Enhanced GFP, Enhanced Cyan Fluorescent Protein (ECFP), Yellow Fluorescent Protein (YFP), and Topaz fluorescent protein (Tpz).
- **Stealth™ RNAi Luciferase Reporter Control** targets the firefly (*Photinus pyralis*) luciferase gene, commonly used in reporter vectors such as pGL2 and pGL3. The Stealth™ RNAi Luciferase Reporter Control does not knock-down luciferase genes from other species, such as *Renilla* luciferase.
- **Stealth™ RNAi LacZ Reporter Control** targets the *E. coli* *LacZ* gene, encoding β -galactosidase, which is a frequently used reporter gene and fusion protein.
- **Stealth™ RNAi β -lactamase Reporter Control** targets the *bla* gene encoding the β -lactamase enzyme, a regularly used reporter gene, for example in the GeneBLAzer® Technology from Invitrogen.

Handling the Stealth™ RNAi Reporter Control Duplexes

- Thaw Stealth™ RNAi Reporter Control Duplex stock solutions on ice or at room temperature. After use, return to -20°C storage.
- Multiple freeze/thaw cycles are permitted without loss of activity if stock solutions are handled properly.
- Ensure that the stock solutions do not become contaminated with RNase.

Amount of Stealth™ RNAi Reporter Control Duplex to Transfect

The amount of Stealth™ RNAi duplex or the Stealth™ RNAi Reporter Control Duplex required to achieve optimal target gene knockdown or minimal knockdown, as appropriate, should be determined experimentally for each human cell line. **As a starting point, we recommend using 10 nM Stealth™ RNAi duplex or Stealth™ RNAi Reporter Control Duplex for transfection.**

To optimize transfection conditions, vary transfection reagent concentrations and the final concentrations of Stealth™ RNAi from 1 to 200 nM, as necessary for your cell line. Use the 1X RNA Annealing/Dilution Buffer supplied with the kit to dilute the Stealth™ RNAi Reporter Control Duplex stock solution, if necessary.

General Guidelines for Transfection

- Use a transfection reagent suitable for delivery of Stealth™ RNAi to mammalian cells. **For optimal results, we recommend using Lipofectamine™ RNAiMAX (Cat. No. 13778-075) available from Invitrogen.** For a protocol to transfect Stealth™ RNAi into cells using Lipofectamine™ RNAiMAX, see the RNAi resource page at www.invitrogen.com/rnai; click on Protocols.
- Use cells that express the appropriate target gene for the Stealth™ RNAi Reporter Control Duplex you are using (see **Stealth™ RNAi Reporter Control Duplex Target Genes**), for example through stable integration or viral transduction of the reporter construct. Alternatively, co-transfect the Stealth™ RNAi duplex with the appropriate reporter expression construct.
Note: For co-transfection experiments with a plasmid, we recommend using Lipofectamine™ 2000, since Lipofectamine™ RNAiMAX was specifically designed for use with Stealth™ RNAi or siRNA duplexes
- Use low-passage cells, and make sure that cells are healthy and greater than 90% viable before transfection.
- Transfect cells at the density recommended by the manufacturer of your transfection reagent (*e.g.* 30-50% confluence at time of transfection if using Lipofectamine™ RNAiMAX).
- Assay for target gene knockdown at a suitable time period (typically 24 to 72 hours) after transfection.
- For a convenient tool to assess transfection efficiency and Stealth™ RNAi uptake, we recommend using the BLOCK-iT™ Alexa Fluor® Red Fluorescent Control (Cat. No. 14750-100) available from Invitrogen.

Quality Control

- The identity and concentration of each corresponding single-stranded RNA oligo is verified by mass spectrometry and optical density reading, respectively.
- After annealing, the Stealth™ RNAi Negative Control Duplex is analyzed by polyacrylamide gel electrophoresis to verify its integrity and to confirm the absence of RNA degradation.

Limited Use Label License No. 196: Stealth™ RNAi

The purchase of this product conveys to the buyer the non-transferable right to use the purchased amount of the product and components of the product in research conducted by the buyer (whether the buyer is an academic or for-profit entity). The buyer cannot sell or otherwise transfer (a) this product (b) its components or (c) materials made using this product or its components to a third party or otherwise use this product or its components or materials made using this product or its components for Commercial Purposes. The buyer may transfer information or materials made through the use of this product to a scientific collaborator, provided that such transfer is not for any Commercial Purpose, and that such collaborator agrees in writing (a) not to transfer such materials to any third party, and (b) to use such transferred materials and/or information solely for research and not for Commercial Purposes. Commercial Purposes means any activity by a party for consideration and may include, but is not limited to: (1) use of the product or its components in manufacturing; (2) use of the product or its components to provide a service, information, or data; (3) use of the product or its components for therapeutic, diagnostic or prophylactic purposes; or (4) resale of the product or its components, whether or not such product or its components are resold for use in research. Invitrogen Corporation will not assert a claim against the buyer of infringement of patents owned by Invitrogen and claiming this product based upon the manufacture, use or sale of a therapeutic, clinical diagnostic, vaccine or prophylactic product developed in research by the buyer in which this product or its components was employed, provided that neither this product nor any of its components was used in the manufacture of such product. If the purchaser is not willing to accept the limitations of this limited use statement, Invitrogen is willing to accept return of the product with a full refund. For information on purchasing a license to this product for purposes other than research, contact Licensing Department, Invitrogen Corporation, 1600 Faraday Avenue, Carlsbad, California 92008. Phone (760) 603-7200. Fax (760) 602-6500.

©2007 Invitrogen Corporation. All rights reserved. Stealth™ is a registered trademark of Invitrogen corporation.

For research use only. Not intended for any animal or human therapeutic or diagnostic use.