

PRODUCT INFORMATION

SP6 RNA

Polymerase

Pub. No. MAN0012020

Rev. Date 16 September 2016 (Rev. B.00)

#_ **Lot:** _ **Expiry Date:** _

Store at -20 °C

Components	#EP0131	#EP0133
SP6 RNA Polymerase	20 U/ μ L 2000 U	200 U/ μ L HC, 5000 U
5X Transcription Buffer	0.8 mL	1.25 mL

BSA included

www.thermofisher.com

For Research Use Only. Not for use in diagnostic procedures.

Description

Bacteriophage SP6 RNA Polymerase is a DNA-dependent RNA polymerase with strict specificity for its respective double-stranded promoter. The enzyme catalyzes the 5' \rightarrow 3' synthesis of RNA on either single-stranded DNA or double-stranded DNA downstream from the promoter.

SP6 RNA Polymerase accepts modified nucleotides (e.g., biotin-, digoxigenin-, fluorescein-labeled nucleotides) as substrates for RNA synthesis.

Applications

Synthesis of unlabeled and labeled RNA that can be used (see protocols on back page):

- for hybridization (1), *in vitro* RNA translation (2);
- as aRNA (3), siRNA (4), substrate in RNase protection assays (5), template for genomic DNA sequencing (6);
- in studies of RNA secondary structure and RNA-protein interactions (7), RNA splicing (8).

Source

E.coli cells with a cloned gene encoding the SP6 RNA Polymerase.

Molecular Weight

99 kDa monomer.

Rev.8

Definition of Activity Unit

One unit of the enzyme incorporates 1 nmol of AMP into a polynucleotide fraction in 60 minutes at 37 °C.

Storage Buffer

The enzyme is supplied in: 50 mM Tris-HCl (pH 8.0), 150 mM NaCl, 5 mM DTT, 0.1 mg/mL BSA, 0.03% (v/v) ELUGENT Detergent and 50% (v/v) glycerol.

5X Transcription Buffer

200 mM Tris-HCl (pH 7.9 at 25 °C), 30 mM MgCl₂, 50 mM DTT, 50 mM NaCl and 10 mM spermidine.

Inhibition and Inactivation

- Inhibitors: metal chelators, enzyme activity is reduced by 50% at NaCl or KCl concentration above 150 mM.
- Inactivated by heating at 70 °C for 10 min or by addition of EDTA.

CERTIFICATE OF ANALYSIS

Endodeoxyribonuclease Assay

No detectable degradation was observed after incubation of supercoiled plasmid DNA with SP6 RNA Polymerase.

Ribonuclease Assay

No detectable degradation was observed after incubation of [3H]-RNA with SP6 RNA Polymerase.

Endo-, exodeoxyribonuclease and phosphatase Assay

No detectable degradation after incubation of single stranded and double stranded radiolabeled oligonucleotides with SP6 RNA Polymerase.

Functional Assay

SP6 RNA Polymerase was tested in *in vitro* transcription reaction.

Quality authorized by:

 Jurgita Zilinskiene

(continued on back page)

Protocol for *in vitro* transcription

1. Linearize template DNA with a restriction enzyme. Extract DNA with phenol/chloroform, then with chloroform/isoamyl alcohol, and precipitate with ethanol. Dissolve DNA in DEPC-treated Water (#R0601).
2. Prepare the following reaction mixture:

5X Transcription buffer	10 μ L
ATP/GTP/CTP/UTP Mix, 10 mM each	10 μ L (2 mM final concentration)
Linear template DNA	1 μ g
Thermo Scientific RiboLock RNase Inhibitor (#EO0381)	1.25 μ L (50 U)
SP6 RNA Polymerase	30 U
DEPC-treated water (#R0601)	to 50 μ L

3. Incubate at 37 °C for 2 hours.
4. Optional: To remove template DNA add 2 μ L (2 U) of DNase I, RNase-free (#EN0521), mix and incubate at 37 °C for 15 min.
5. Inactivate DNase I by phenol/chloroform extraction.

Note

- The transcription reaction should be performed under conditions that exclude contamination with RNases. The tips, tubes and water should be nuclease free. All the solutions should be made up in nuclease free water. Wearing gloves is advisable.
- The reaction mixture should be prepared at room temperature, since DNA may precipitate in the presence of spermidine at 4°C.
- Under the conditions described above, more than 10 μ g RNA per 1 μ g template DNA is obtained.
- The yield of proper length transcripts decreases if the template DNA is incompletely linearized due to a read-through reaction and accumulation of longer transcripts of a variable length.
- The reaction mixture can be scaled up or down.

Protocol for synthesis of radiolabeled RNA probes of high specific activity

1. Linearize template DNA with a restriction enzyme. Extract DNA with phenol/chloroform, then with chloroform/isoamyl alcohol, and precipitate with ethanol. Dissolve DNA in DEPC-treated Water (#R0601).
2. Prepare the following reaction mixture:

5X Transcription buffer	4 μ L
3 NTP mix, 10 mM each	1 μ L (0.5 mM final concentration)
100 μM CTP (#R0451)	2.4 μ L (12 μ M final concentration)
[α-³²P]-CTP, ~30 TBq/mmol (800 Ci/mmol)	1.85 MBq (50 μ Ci)
Linear template DNA	0.2-1.0 μ g
RiboLock™ RNase Inhibitor (#EO0381)	0.4 μ L (20 U)
SP6 RNA Polymerase	20 U
DEPC-treated water (#R0601)	to 20 μ L

3. Incubate at 37 °C for 2 hours.
4. Stop the reaction by cooling at -20 °C.
5. Determine the percentage of label incorporated into RNA.

Note

- RNA synthesized under the conditions described above usually has a specific activity of 3-5 x10⁸dpm/ μ g.
- RNA can be radiolabeled with [³²P], [³⁵S] or [³H]-ribonucleotides. The use of 1.85 MBq (50 μ Ci) of 5'-[α -³²P]-CTP, ~30 TBq/mmol (800 Ci/mmol), 11.1 MBq (300 μ Ci) of 5'-[α -³⁵S]-UTP, >37 TBq/mmol (>1000 Ci/mmol), 0.925 MBq (25 μ Ci) of 5,6-[³H]-UTP, 1.1-2.2 TBq/mmol (30-60 Ci/mmol) for 20 μ L reaction mixture is recommended.
- The yield of full-length transcripts is reduced when the final concentration of labeled NTP is below 12 μ M.

References

1. Melton, D.A., et al., Efficient *in vitro* synthesis of biologically active RNA and RNA hybridization probes from plasmids containing a bacteriophage SP6 promoter, *Nucleic Acids Res.*, 12, 7035-7056, 1984.
2. Krieg, P.A., Melton, D.A., Functional messenger RNAs are produced by SP6 *in vitro* transcription of cloned cDNAs, *Nucleic Acids Res.*, 12, 7057-7070, 1984.
3. Melton, D.A., Injected antisense RNAs specifically block messenger RNA translation *in vivo*, *Proc. Natl. Acad. Sci. USA*, 82, 144-148, 1985.
4. Bernstein, E., et al., Role for bidentate ribonuclease in the initiation step of RNA interference, *Nature*, 409, 363-366, 2001.
5. Peebles, C.L., et al., A self-splicing RNA excises an intron lariat, *Cell*, 44, 213-223, 1986.
6. Church, G.M., Gilbert, W., Genomic sequencing, *Proc. Natl. Acad. Sci. USA*, 81, 1991-1995, 1984.
7. Witherell, G.W., et al., Cooperative binding of R17 coat protein to RNA, *Biochemistry*, 29, 11051-11057, 1990.
8. Krainer, A.R., et al., Normal and mutant human β -globin pre-mRNAs are faithfully and efficiently spliced *in vitro*, *Cell*, 36, 993-1005, 1984.
9. Jorgensen, E.D., et al., Specific contacts between the bacteriophage T3, T7, and SP6 RNA polymerases and their promoters, *J. Biol. Chem.*, 266, 645-651, 1991.

LIMITED USE LABEL LICENSE: Internal Research and Development Use Only.

The purchase of this product conveys to the buyer the limited, non-exclusive, non-transferable right (without the right to resell, repackage, or further sublicense) to use this product for internal research and development purposes. No other license is granted to the buyer whether expressly, by implication, by estoppel or otherwise. In particular, the purchase of the product does not include or carry any right or license to use, develop, or otherwise exploit this product commercially and no rights are conveyed to the buyer to use the product or components of the product for purposes including but not limited to provision of services to a third party, generation of commercial databases or clinical diagnostics. This product is sold pursuant to authorization from Thermo Fisher Scientific and Thermo Fisher Scientific reserves all other rights. For information on purchasing a license for uses other than internal research and development purposes, please contact outlicensing@lifetech.com or Out Licensing, Life Technologies Inc., 5791 Van Allen Way, Carlsbad, California 92008.

PRODUCT USE LIMITATION

This product is developed, designed and sold exclusively for research purposes and *in vitro* use only. The product was not tested for use in diagnostics or for drug development, nor is it suitable for administration to humans or animals.

Please refer to www.thermofisher.com for Material Safety Data Sheet of the product.

© 2016 Thermo Fisher Scientific Inc. All rights reserved. All trademarks are the property of Thermo Fisher Scientific Inc. and its subsidiaries.