

Human Adiponectin ELISA Kit

Enzyme-linked Immunosorbent Assay for quantitative detection of human Adiponectin

Catalog Numbers BMS2032-2 or BMS2032-2TEN

Pub. No. MAN0016492 **Rev.** A.0 (30)

WARNING! Read the Safety Data Sheets (SDSs) and follow the handling instructions. Wear appropriate protective eyewear, clothing, and gloves. Safety Data Sheets (SDSs) are available from thermofisher.com/support.

Product description

The Human Adiponectin ELISA Kit is an enzyme-linked immunosorbent assay for the quantitative detection of human Adiponectin.

Summary

Adiponectin (also referred to as Acrp30, apM1) is a protein hormone that modulates a number of metabolic processes, including glucose regulation and fatty acid catabolism. Adiponectin is exclusively secreted from adipose tissue into the bloodstream and is very abundant in plasma relative to many hormones. Levels of the hormone are inversely correlated with body mass index (BMI). The hormone plays a role in the suppression of the metabolic derangements that may result in type 2 diabetes, obesity, atherosclerosis and non-alcoholic fatty liver disease (NAFLD).

Adiponectin was first characterized in mice as a transcript overexpressed in preadipocytes (precursors of fat cells) differentiating into adipocytes. The human homologue was identified as the most abundant transcript in adipose tissue. Contrary to expectations, despite being produced in adipose tissue, adiponectin was found to be decreased in obesity. This downregulation has not been fully explained. The gene was localized to chromosome 3p27, a region highlighted as affecting genetic susceptibility to type 2 diabetes and obesity. Adiponectin is a 244-amino-acid-long polypeptide. There are four distinct regions of adiponectin. The first is a short signal sequence that targets the hormone for secretion outside the cell; next is a short region that varies between species; the third is a 65-amino acid region with similarity to collagenous proteins; the last is a globular domain.

Adiponectin is secreted into the bloodstream where it accounts for approximately 0.01% of all plasma protein. Plasma concentrations reveal a sexual dimorphism, with females having higher levels than males. Levels of adiponectin are reduced in diabetics compared to non-diabetics. Weight reduction significantly increases circulating levels. Adiponectin automatically self-associates into larger structures.

Adiponectin binds to a number of receptors. So far, two receptors have been identified, with homology to G protein-coupled receptors. These have distinct tissue specificities within the body and have different affinities to the various forms of adiponectin. The receptors affect the downstream target AMP kinase, an important cellular metabolic rate control point. Expression of the receptors are correlated with insulin levels, as well as reduced in mouse models of diabetes, particularly in skeletal muscle and adipose tissue.

Adiponectin exerts some of its weight reduction effects via the brain. This is similar to the action of leptin, but the two hormones perform complementary actions, and can have additive effects.

Adiponectin affects glucose flux (gluconeogenesis, glucose uptake), lipid catabolism (β -oxidation, triglyceride clearance), endothelial dysfunction, insulin sensitivity, and weight loss.

For literature update refer to our website.

For Research Use Only. Not for use in diagnostic procedures.

Principles of the test

An anti-human Adiponectin coating antibody is adsorbed onto microwells.

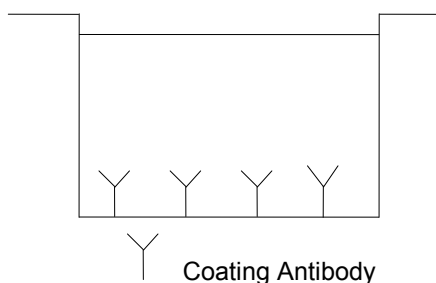


Fig. 1 Coated microwell.

Human Adiponectin present in the sample or standard binds to antibodies adsorbed to the microwells. A biotin-conjugated anti-human Adiponectin antibody is added and binds to human Adiponectin captured by the first antibody.

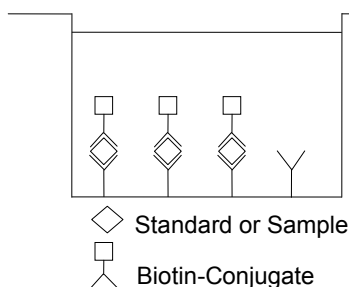


Fig. 2 First incubation.

Following incubation unbound biotin-conjugated anti-human Adiponectin antibody is removed during a wash step. Streptavidin-HRP is added and binds to the biotin-conjugated anti-human Adiponectin antibody.

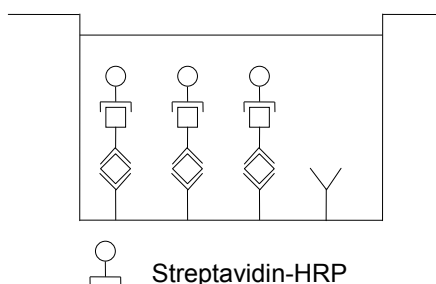


Fig. 3 Second incubation.

Following incubation unbound Streptavidin-HRP is removed during a wash step, and substrate solution reactive with HRP is added to the wells.

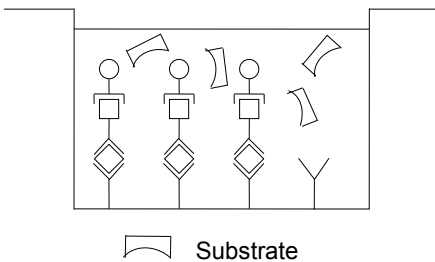


Fig. 4 Third incubation.

A colored product is formed in proportion to the amount of human Adiponectin present in the sample or standard. The reaction is terminated by addition of acid and absorbance is measured at 450 nm. A standard curve is prepared from 7 human Adiponectin standard dilutions and human Adiponectin sample concentration determined.

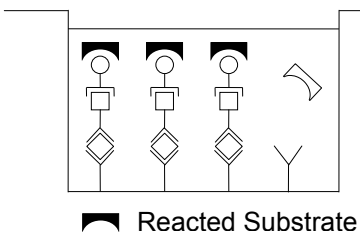


Fig. 5 Stop reaction.

Reagents provided

Reagents for human Adiponectin ELISA BMS2032-2 (96 tests)

- 1 aluminum pouch with a Microwell Plate (12 strips of 8 wells each) coated with monoclonal antibody to human Adiponectin
- 1 vial (70 μ L) Biotin-Conjugate anti-human Adiponectin monoclonal antibody
- 1 vial (150 μ L) Streptavidin-HRP
- 2 vials human Adiponectin Standard lyophilized, 30 ng/mL upon reconstitution
- 2 vial (5 mL) Assay Buffer Concentrate 20x (PBS with 1% Tween™ 20, 10% BSA)
- 1 bottle (50 mL) Wash Buffer Concentrate 20x (PBS with 1% Tween™ 20)
- 1 vial (15 mL) Substrate Solution (tetramethyl-benzidine)
- 1 vial (15 mL) Stop Solution (1M Phosphoric acid)
- 4 Adhesive Films

Reagents for human Adiponectin ELISA BMS2032-2TEN (10x96 tests)

- 10 aluminum pouches with a Microwell Plate (12 strips of 8 wells each) coated with monoclonal antibody to human Adiponectin
- 10 vials (70 μ L) Biotin-Conjugate anti-human Adiponectin monoclonal antibody
- 10 vials (150 μ L) Streptavidin-HRP
- 10 vials human Adiponectin Standard lyophilized, 30 ng/mL upon reconstitution
- 2 vials (50 mL) Assay Buffer Concentrate 20x (PBS with 1% Tween™ 20, 10% BSA)
- 6 bottles (50 mL) Wash Buffer Concentrate 20x (PBS with 1% Tween™ 20)
- 10 vials (15 mL) Substrate Solution (tetramethyl-benzidine)
- 1 vial (100 mL) Stop Solution (1M Phosphoric acid)
- 20 Adhesive Films

Storage instructions – ELISA kit

Store kit reagents between 2°C and 8°C.

Immediately after use remaining reagents should be returned to cold storage (2°C to 8°C). Expiry of the kit and reagents is stated on labels.

Expiry of the kit components can only be guaranteed if the components are stored properly, and if, in case of repeated use of one component, this reagent is not contaminated by the first handling.

Sample collection and storage instructions

Cell culture supernatant, serum and plasma (citrate, heparin) were tested with this assay. Other biological samples might be suitable for use in the assay. Remove serum or plasma from the clot or cells as soon as possible after clotting and separation.

Pay attention to a possible *Hook Effect* due to high sample concentrations (see “Calculation of results” on page 5).

Samples containing a visible precipitate must be clarified prior to use in the assay. Do not use grossly hemolyzed or lipemic samples.

Samples should be aliquoted and must be stored frozen at –20°C to avoid loss of bioactive human Adiponectin. If samples are to be run within 24 hours, they may be stored at 2°C to 8°C (for sample stability refer to “Sample stability” on page 6).

Avoid repeated freeze-thaw cycles. Prior to assay, the frozen sample should be brought to room temperature slowly and mixed gently.

Materials required but not provided

- 5 mL and 10 mL graduated pipettes
- 5 μ L to 1000 μ L adjustable single channel micropipettes with disposable tips
- 50 μ L to 300 μ L adjustable multichannel micropipette with disposable tips
- Multichannel micropipette reservoir
- Beakers, flasks, cylinders necessary for preparation of reagents
- Device for delivery of wash solution (multichannel wash bottle or automatic wash system)
- Microwell strip reader capable of reading at 450 nm (620 nm as optional reference wave length)
- Glass-distilled or deionized water
- Statistical calculator with program to perform regression analysis

Precautions for use

- All reagents should be considered as potentially hazardous. We therefore recommend that this product is handled only by those persons who have been trained in laboratory techniques and that it is used in accordance with the principles of good laboratory practice. Wear suitable protective clothing such as laboratory overalls, safety glasses and gloves. Care should be taken to avoid contact with skin or eyes. In the case of contact with skin or eyes wash immediately with water. See material safety data sheet(s) and/or safety statement(s) for specific advice.
- Reagents are intended for research use only and are not for use in diagnostic or therapeutic procedures.
- Do not mix or substitute reagents with those from other lots or other sources.
- Do not use kit reagents beyond expiration date on label.
- Do not expose kit reagents to strong light during storage or incubation.
- Do not pipet by mouth.
- Do not eat or smoke in areas where kit reagents or samples are handled.
- Avoid contact of skin or mucous membranes with kit reagents or samples.
- Rubber or disposable latex gloves should be worn while handling kit reagents or samples.
- Avoid contact of substrate solution with oxidizing agents and metal.
- Avoid splashing or generation of aerosols.

- To avoid microbial contamination or cross-contamination of reagents or samples that may invalidate the test, use disposable pipette tips and/or pipettes.
- Use clean, dedicated reagent trays for dispensing the conjugate and substrate reagent.
- Exposure to acid inactivates the conjugate.
- Glass-distilled water or deionized water must be used for reagent preparation.
- Substrate solution must be at room temperature prior to use.
- Decontaminate and dispose samples and all potentially contaminated materials as if they could contain infectious agents. The preferred method of decontamination is autoclaving for a minimum of 1 hour at 121.5°C.
- Liquid wastes not containing acid and neutralized waste may be mixed with sodium hypochlorite in volumes such that the final mixture contains 1.0% sodium hypochlorite. Allow 30 minutes for effective decontamination. Liquid waste containing acid must be neutralized prior to the addition of sodium hypochlorite.

Preparation of reagents

1. Buffer Concentrates should be brought to room temperature and should be diluted before starting the test procedure.
2. If crystals have formed in the Buffer Concentrates, warm them gently until they have completely dissolved.

Wash buffer (1x)

1. Pour entire contents (50 mL) of the Wash Buffer Concentrate (20x) into a clean 1000 mL graduated cylinder. Bring to final volume of 1000 mL with glass-distilled or deionized water. Mix gently to avoid foaming.
2. Transfer to a clean wash bottle and store at 2°C to 25°C. The Wash Buffer (1x) is stable for 30 days.
3. Wash Buffer (1x) may also be prepared as needed according to the following table:

Number of Strips	Wash Buffer Concentrate (20x) (mL)	Distilled Water (mL)
1-6	25	475
1-12	50	950

Assay buffer (1x)

1. For BMS2032-2:
 - a. Pour the entire contents (5 mL) of the Assay Buffer Concentrate (20x) into a clean 100 mL graduated cylinder. Bring to final volume of 100 mL with distilled water. Mix gently to avoid foaming.
 - b. Store at 2°C to 8°C. The Assay Buffer (1x) is stable for 30 days.
2. For BMS2032-2TEN:
 - a. Pour the entire contents (50 mL) of the Assay Buffer Concentrate (20x) into a clean 1000 mL graduated cylinder. Bring to final volume of 1000 mL with distilled water. Mix gently to avoid foaming.
 - b. Store at 2°C to 8°C. The Assay Buffer (1x) is stable for 30 days.
3. Assay Buffer (1x) may also be prepared as needed according to the following table:

Number of Strips	Assay Buffer Concentrate (20x) (mL)	Distilled Water (mL)
1-6	2.5	47.5
1-12	5.0	95.0

Biotin-Conjugate

Note: The Biotin-Conjugate should be used within 30 minutes after dilution.

Make a 1:100 dilution of the concentrated Biotin-Conjugate solution with Assay Buffer (1x) in a clean plastic tube as needed according to the following table:

Number of Strips	Biotin-Conjugate (mL)	Assay Buffer (1x) (mL)
1-6	0.03	2.97
1-12	0.06	5.94

Streptavidin-HRP

Note: The Streptavidin-HRP should be used within 30 minutes after dilution.

Make a 1:400 dilution of the concentrated Streptavidin-HRP solution with Assay Buffer (1x) in a clean plastic tube as needed according to the following table:

Number of Strips	Streptavidin-HRP (mL)	Assay Buffer (1x) (mL)
1-6	0.015	5.985
1-12	0.030	11.97

Human Adiponectin standard

1. Reconstitute human Adiponectin standard by addition of distilled water.
2. Reconstitution volume is stated on the label of the standard vial. Swirl or mix gently to insure complete and homogeneous solubilization (concentration of reconstituted standard = 30 ng/mL).
3. Allow the standard to reconstitute for 10-30 minutes. Mix well prior to making dilutions.

The standard has to be used immediately after reconstitution and cannot be stored.

External standard dilution

1. Label 7 tubes, one for each standard point: S1, S2, S3, S4, S5, S6, S7.
2. Prepare 1:2 serial dilutions for the standard curve as follows: Pipet 225 µL of Assay Buffer (1x) into each tube.
3. Pipet 225 µL of reconstituted standard (concentration of standard = 30 ng/mL) into the first tube, labeled S1, and mix (concentration of standard 1 = 15 ng/mL).
4. Pipet 225 µL of this dilution into the second tube, labeled S2, and mix thoroughly before the next transfer.
5. Repeat serial dilutions five more times thus creating the points of the standard curve (see Figure 6).
Assay Buffer (1x) serves as blank.

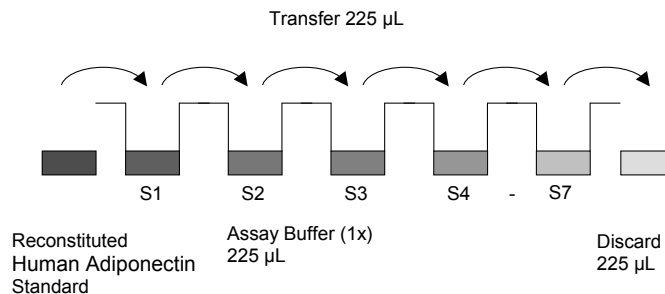


Fig. 6 Dilute standards - tubes.

Test protocol

Note: In case of incubation without shaking the obtained O.D. values may be lower than indicated below. Nevertheless the results are still valid.

1. Predilute your samples before starting with the test procedure. Dilute serum, plasma, and cell culture samples 1:2500 with Assay Buffer (1x) according to the following scheme:
 - Dilution 1: 10 μ L sample + 490 μ L Assay Buffer (1x)
 - Dilution 2: 10 μ L of dilution 1 + 490 μ L Assay Buffer (1x)
2. Determine the number of microwell strips required to test the desired number of samples plus appropriate number of wells needed for running blanks and standards. Each sample, standard, blank and optional control sample should be assayed in duplicate. Remove extra microwell strips from holder and store in foil bag with the desiccant provided at 2°C to 8°C sealed tightly.
3. Wash the microwell strips twice with approximately 400 μ L Wash Buffer per well with thorough aspiration of microwell contents between washes. Allow the Wash Buffer to sit in the wells for about 10–15 seconds before aspiration. Take care not to scratch the surface of the microwells.
4. After the last wash step, empty wells and tap microwell strips on absorbent pad or paper towel to remove excess Wash Buffer. Use the microwell strips immediately after washing. Alternatively microwell strips can be placed upside down on a wet absorbent paper for not longer than 15 minutes. Do not allow wells to dry.
5. Standard dilution on the microwell plate (alternatively the standard dilution can be prepared in tubes, see “External standard dilution” on page 3):

Add 100 μ L of Assay Buffer (1x) in duplicate to all standard wells. Pipet 100 μ L of prepared standard (see Preparation of Standard 0, concentration = 30.0 ng/mL) in duplicate into well A1 and A2 (see Table 1). Mix the contents of wells A1 and A2 by repeated aspiration and ejection (concentration of standard 1, S1 = 15.0 ng/mL), and transfer 100 μ L to wells B1 and B2, respectively (see Figure 7). Take care not to scratch the inner surface of the microwells. Continue this procedure 5 times, creating two rows of human Adiponectin standard dilutions ranging from 15.0–0.23 ng/mL. Discard 100 μ L of the contents from the last microwells (G1, G2) used.

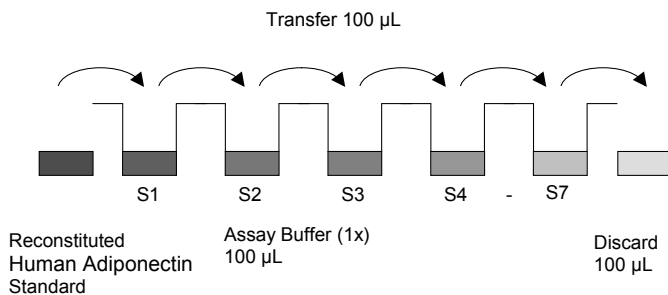


Fig. 7 Dilute standards - microwell plate.

In case of an external standard dilution (see “External standard dilution” on page 3), pipet 100 μ L of these standard dilutions (S1–S7) in the standard wells according to Table 1.

Table 1 Example of the arrangement of blanks, standards, and samples in the microwell strips.

	1	2	3	4
A	Standard 1 15.0 ng/mL	Standard 1 15.0 ng/mL	Sample 1	Sample 1
B	Standard 2 7.5 ng/mL	Standard 2 7.5 ng/mL	Sample 2	Sample 2
C	Standard 3 3.75 ng/mL	Standard 3 3.75 ng/mL	Sample 3	Sample 3
D	Standard 4 1.88 ng/mL	Standard 4 1.88 ng/mL	Sample 4	Sample 4
E	Standard 5 0.94 ng/mL	Standard 5 0.94 ng/mL	Sample 5	Sample 5
F	Standard 6 0.47 ng/mL	Standard 6 0.47 ng/mL	Sample 6	Sample 6
G	Standard 7 0.23 ng/mL	Standard 7 0.23 ng/mL	Sample 7	Sample 7
H	Blank	Blank	Sample 8	Sample 8

6. Add 100 μ L of Assay Buffer (1x) in duplicate to the blank wells.
7. Add 50 μ L of Assay Buffer (1x) to the sample wells.
8. Add 50 μ L of each prediluted sample in duplicate to the sample wells.
9. Prepare Biotin-Conjugate (see “Biotin-Conjugate” on page 3).
10. Add 50 μ L of Biotin-Conjugate to all wells.
11. Cover with an adhesive film and incubate at room temperature (18°C to 25°C) for 2 hours on a microplate shaker.
12. Prepare Streptavidin-HRP (see “Streptavidin-HRP” on page 3).
13. Remove adhesive film and empty wells. Wash microwell strips 6 times according to step 3. Proceed immediately to the next step.
14. Add 100 μ L of diluted Streptavidin-HRP to all wells, including the blank wells.
15. Cover with an adhesive film and incubate at room temperature (18°C to 25°C) for 1 hour on a microplate shaker.
16. Remove adhesive film and empty wells. Wash microwell strips 6 times according to step 3. Proceed immediately to the next step.
17. Pipet 100 μ L of TMB Substrate Solution to all wells.
18. Incubate the microwell strips at room temperature (18°C to 25°C) for about 30 minutes. Avoid direct exposure to intense light.

The color development on the plate should be monitored and the substrate reaction stopped (see next step) before positive wells are no longer properly recordable. Determination of the ideal time period for color development has to be done individually for each assay.

It is recommended to add the stop solution when the highest standard has developed a dark blue color. Alternatively the color development can be monitored by the ELISA reader at 620 nm. The substrate reaction should be stopped as soon as Standard 1 has reached an OD of 0.9–0.95.

19. Stop the enzyme reaction by quickly pipetting 100 μ L of Stop Solution into each well. It is important that the Stop Solution is spread quickly and uniformly throughout the microwells to completely inactivate the enzyme. Results must be read immediately after the Stop Solution is added or within 1 hour if the microwell strips are stored at 2°C to 8°C in the dark.
20. Read absorbance of each microwell on a spectrophotometer using 450 nm as the primary wave length (optionally 620 nm as the reference wave length; 610–650 nm is acceptable). Blank the plate reader according to the manufacturer's instructions by using the blank wells. Determine the absorbance of both the samples and the standards.

Calculation of results

- Calculate the average absorbance values for each set of duplicate standards and samples. Duplicates should be within 20% of the mean value.
- Create a standard curve by plotting the mean absorbance for each standard concentration on the ordinate against the human Adiponectin concentration on the abscissa. Draw a best fit curve through the points of the graph (a 5-parameter curve fit is recommended).
- To determine the concentration of circulating human Adiponectin for each sample, first find the mean absorbance value on the ordinate and extend a horizontal line to the standard curve. At the point of intersection, extend a vertical line to the abscissa and read the corresponding human Adiponectin concentration.
- If instructions in this protocol have been followed, samples have been diluted 1:5000 (1:2500 external predilution, 1:2 dilution on the plate: 50 μ L sample + 50 μ L Assay Buffer (1x)) and the concentration read from the standard curve must be multiplied by the dilution factor (\times 5000).
- Calculation of 1:5000 prediluted samples with a concentration exceeding standard 1 will result in incorrect, low human Adiponectin levels (Hook Effect). Such samples require further external predilution according to expected human Adiponectin values with Assay Buffer (1x) in order to precisely quantitate the actual human Adiponectin level.
- It is suggested that each testing facility establishes a control sample of known human Adiponectin concentration and runs this additional control with each assay. If the values obtained are not within the expected range of the control, the assay results may be invalid.
- A representative standard curve is shown in Figure 8.

Note: Do not use this standard curve to derive test results. Each laboratory must prepare a standard curve for each group of microwell strips assayed.

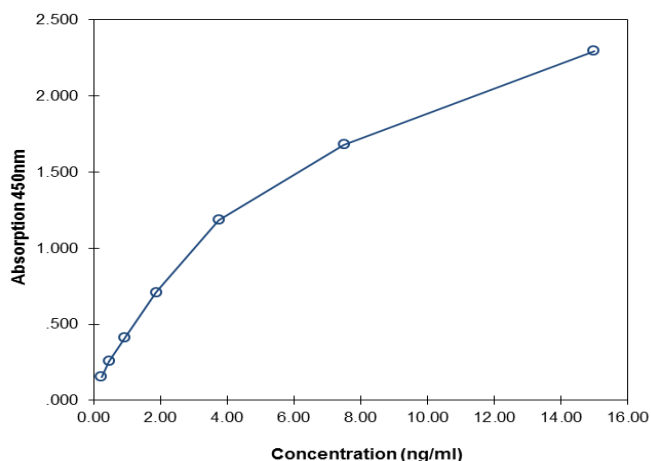


Fig. 8 Representative standard curve for Human Adiponectin ELISA Kit. Human Adiponectin was diluted in serial 2-fold steps in Assay Buffer (1x).

Table 2 Typical data using the Human Adiponectin ELISA Kit.
Measuring wavelength: 450 nm
Reference wavelength: 620 nm

Standard	Human Adiponectin concentration (ng/mL)	O.D. at 450 nm	Mean O.D. at 450 nm	C.V. (%)
1	15.0	2.314 2.275	2.295	0.9
2	7.50	1.698 1.663	1.680	1.0
3	3.75	1.185 1.183	1.184	0.1
4	1.88	0.712 0.711	0.711	0.1
5	0.94	0.418 0.410	0.414	1.0
6	0.47	0.251 0.262	0.256	2.2
7	0.23	0.160 0.147	0.154	4.4
Blank	0.00	0.026 0.024	0.025	2.8

The OD values of the standard curve may vary according to the conditions of assay performance (e.g., operator, pipetting technique, washing technique, or temperature effects). Furthermore, shelf life of the kit may affect enzymatic activity and thus color intensity. Values measured are still valid.

Limitations

- Because exact conditions may vary from assay to assay, a standard curve must be established for every run.
- Bacterial or fungal contamination of either screen samples or reagents or cross-contamination between reagents may cause erroneous results.
- Disposable pipette tips, flasks, or glassware are preferred, reusable glassware must be washed and thoroughly rinsed of all detergents before use.
- Improper or insufficient washing at any stage of the procedure will result in either false positive or false negative results. Empty wells completely before dispensing fresh wash solution, fill with Wash Buffer as indicated for each wash cycle, and do not allow wells to sit uncovered or dry for extended periods.
- The use of radioimmunotherapy has significantly increased the number of patients with human anti-mouse IgG antibodies (HAMA). HAMA may interfere with assays utilizing murine monoclonal antibodies leading to both false positive and false negative results. Serum samples containing antibodies to murine immunoglobulins can still be analyzed in such assays when murine immunoglobulins (serum, ascitic fluid, or monoclonal antibodies of irrelevant specificity) are added to the sample.

Performance characteristics

Sensitivity

The limit of detection of human Adiponectin defined as the analyte concentration resulting in an absorbance significantly higher than that of the dilution medium (mean plus 2 standard deviations) was determined to be 0.012 ng/mL (mean of 6 independent assays).

Reproducibility

Intra-assay

Reproducibility within the assay was evaluated in 3 independent experiments. Each assay was carried out with 6 replicates of 8 serum samples containing different concentrations of human Adiponectin. Two standard curves were run on each plate. Data below show the mean human Adiponectin concentration and the coefficient of

variation for each sample. The calculated overall intra-assay coefficient of variation was 4.2%.

Table 3 The mean human Adiponectin concentration and the coefficient of variation for each sample.

Sample	Experiment	Mean human Adiponectin concentration (ng/mL)	Coefficient of variation (%)
1	1	3341.47	4.5
	2	3183.93	2.0
	3	3256.38	2.4
2	1	3725.36	7.1
	2	3543.81	5.3
	3	3713.56	5.2
3	1	7394.21	5.0
	2	7442.24	5.4
	3	7633.32	1.9
4	1	9973.84	4.1
	2	10,238.92	3.9
	3	10,298.49	2.2
5	1	8618.62	5.7
	2	8794.77	6.3
	3	9123.28	2.1
6	1	5571.56	7.0
	2	5693.79	3.5
	3	5907.67	1.9
7	1	1263.71	5.0
	2	1334.75	2.4
	3	1337.23	2.0
8	1	637.49	6.0
	2	681.93	8.2
	3	736.65	2.4

Inter-assay

Assay-to-assay reproducibility within one laboratory was evaluated in 3 independent experiments. Each assay was carried out with 6 replicates of 8 serum containing different concentrations of human Adiponectin. Two standard curves were run on each plate. Data below show the mean human Adiponectin concentration and the coefficient of variation calculated on 18 determinations of each sample. The calculated overall inter-assay coefficient of variation was 3.1%.

Table 4 The mean human Adiponectin concentration and the coefficient of variation of each sample.

Sample	Mean human Adiponectin concentration (ng/mL)	Coefficient of variation (%)
1	3260.59	2.4
2	3660.91	2.8
3	7489.92	1.7
4	10,170.42	1.7
5	8845.56	2.9
6	5724.34	3.0
7	1311.90	3.2
8	685.36	7.2

Spike recovery

The spike recovery was evaluated by spiking 2 levels of human Adiponectin into serum, plasma and cell culture supernatant. Recoveries were determined with 4 replicates each. The amount of endogenous human Adiponectin in unspiked samples was subtracted from the spike values. Recoveries were shown to depend on the samples used.

Sample matrix ^[1]	Spike medium (%)	Spike low (%)
Serum	110	121
Plasma (citrate)	92	98
Plasma (heparin)	119	130
Cell culture supernatant	111	122

^[1] Due to high sample dilution, data for high spikes could not be evaluated.

Dilution parallelism

Serum, plasma, and cell culture supernatant samples with different levels of human Adiponectin were analyzed at serial 2-fold dilutions with 4 replicates each.

Sample matrix	% recovery of exp. val.	
	Range	Mean
Serum	76–117	90
Plasma (citrate)	74–105	90
Plasma (heparin)	80–113	90
Cell culture supernatant	92–108	98

Sample stability

Freeze-thaw stability

Aliquots of serum samples (spiked or unspiked) were stored at –20°C and thawed five times, and the human Adiponectin levels determined. There was no significant loss of human Adiponectin immunoreactivity detected by freezing and thawing.

Storage stability

Aliquots of serum samples (spiked or unspiked) were stored at –20°C, 2°C to 8°C, room temperature, and at 37°C, and the human Adiponectin level determined after 24 hours. There was no significant loss of human Adiponectin immunoreactivity detected during storage under above conditions.

Specificity

The cross-reactivity and interference of circulating factors of the immune system was evaluated by spiking these proteins at physiologically relevant concentrations into a human Adiponectin positive sample. No cross-reactivity or interference was detected.

Expected values

Panels of 40 serum as well as citrate and heparin plasma samples from randomly selected apparently healthy donors (males and females) were tested for human Adiponectin. The levels measured may vary with the sample collection used.

Sample matrix	Number of samples evaluated	Range (ng/mL)	Mean (ng/mL)
Serum	40	1309–13,929	5377
Plasma (citrate)	40	1401–8,377	4301
Plasma (heparin)	40	764–16,893	4238

Reagent preparation summary

Wash buffer (1x)

Add Wash Buffer Concentrate 20x (50 mL) to 950 mL distilled water.

Number of Strips	Wash Buffer Concentrate (mL)	Distilled Water (mL)
1–6	25	475
1–12	50	950

Assay buffer (1x)

- For BMS2032-2: Add Assay Buffer Concentrate 20x (5 mL) to 95 mL distilled water.
- For BMS2032-2TEN: Add Assay Buffer Concentrate 20x (50 mL) to 950 mL distilled water.

Number of Strips	Assay Buffer Concentrate (mL)	Distilled Water (mL)
1–6	2.5	47.5
1–12	5.0	95.0

Biotin-Conjugate

Make a 1:100 dilution of Biotin-Conjugate in Assay Buffer (1x).

Number of Strips	Biotin-Conjugate (mL)	Assay Buffer (1x) (mL)
1–6	0.03	2.97
1–12	0.06	5.94

Streptavidin-HRP

Make a 1:400 dilution of Streptavidin-HRP in Assay Buffer (1x):

Number of Strips	Streptavidin-HRP (mL)	Assay Buffer (1x) (mL)
1–6	0.015	5.985
1–12	0.030	11.97

Human Adiponectin standard

Reconstitute lyophilized human Adiponectin standard with distilled water. (Reconstitution volume is stated on the label of the standard vial.)

Test protocol summary

Note: If instructions in this protocol have been followed, samples have been diluted 1:5000 (50 µL 1:2500 prediluted sample + 50 µL Assay Buffer (1x)) and the concentration read from the standard curve must be multiplied by the dilution factor (x 5000).

- Predilute sample with Assay Buffer (1x) 1:2500.
- Determine the number of microwell strips required.
- Wash microwell strips twice with Wash Buffer.
- Standard dilution on the microwell plate: Add 100 µL Assay Buffer (1x), in duplicate, to all standard wells. Pipette 100 µL prepared standard into the first wells and create standard dilutions by transferring 100 µL from well to well. Discard 100 µL from the last wells.

Alternatively, external standard dilution in tubes (see “External standard dilution” on page 3): Pipette 100 µL of these standard dilutions in the microwell strips.

- Add 100 µL Assay Buffer (1x), in duplicate, to the blank wells.
- Add 50 µL Assay Buffer (1x) to sample wells.
- Add 50 µL prediluted sample in duplicate, to designated sample wells.
- Prepare Biotin-Conjugate.
- Add 50 µL Biotin-Conjugate Conjugate to all wells.
- Cover microwell strips and incubate 2 hours at room temperature (18°C to 25°C).
- Prepare Streptavidin-HRP.
- Empty and wash microwell strips 6 times with Wash Buffer.
- Add 100 µL diluted Streptavidin-HRP to all wells.
- Cover microwell strips and incubate 1 hour at room temperature (18°C to 25°C).
- Empty and wash microwell strips 6 times with Wash Buffer.
- Add 100 µL of TMB Substrate Solution to all wells.
- Incubate the microwell strips for about 30 minutes at room temperature (18°C to 25°C).
- Add 100 µL Stop Solution to all wells.
- Blank microwell reader and measure color intensity at 450 nm.

Customer and technical support

Visit [thermofisher.com/support](http://www.thermofisher.com/support) for the latest service and support information.

- Worldwide contact telephone numbers
- Product support information
 - Product FAQs
 - Software, patches, and updates
 - Training for many applications and instruments
- Order and web support
- Product documentation
 - User guides, manuals, and protocols
 - Certificates of Analysis
 - Safety Data Sheets (SDSs; also known as MSDSs)

Note: For SDSs for reagents and chemicals from other manufacturers, contact the manufacturer.

Limited product warranty

Life Technologies Corporation and/or its affiliate(s) warrant their products as set forth in the Life Technologies' General Terms and Conditions of Sale at www.thermofisher.com/us/en/home/global/terms-and-conditions.html. If you have any questions, please contact Life Technologies at www.thermofisher.com/support.



Bender MedSystems GmbH | Campus Vienna Biocenter 2 | 1030 Vienna, Austria

For descriptions of symbols on product labels or product documents, go to thermofisher.com/symbols-definition.

The information in this guide is subject to change without notice.

DISCLAIMER: TO THE EXTENT ALLOWED BY LAW, THERMO FISHER SCIENTIFIC INC. AND/OR ITS AFFILIATE(S) WILL NOT BE LIABLE FOR SPECIAL, INCIDENTAL, INDIRECT, PUNITIVE, MULTIPLE, OR CONSEQUENTIAL DAMAGES IN CONNECTION WITH OR ARISING FROM THIS DOCUMENT, INCLUDING YOUR USE OF IT.

Important Licensing Information: These products may be covered by one or more Limited Use Label Licenses. By use of these products, you accept the terms and conditions of all applicable Limited Use Label Licenses.

©2019 Thermo Fisher Scientific Inc. All rights reserved. All trademarks are the property of Thermo Fisher Scientific and its subsidiaries unless otherwise specified. All other trademarks are the property of their respective owners.

