

Mouse TNF alpha High Sensitivity ELISA Kit

Enzyme-linked Immunosorbent Assay for quantitative detection of mouse TNF- α

Catalog Number BMS607HS

Pub. No. MAN0016886 Rev. B.0 (31)

WARNING! Read the Safety Data Sheets (SDSs) and follow the handling instructions. Wear appropriate protective eyewear, clothing, and gloves. Safety Data Sheets (SDSs) are available from thermofisher.com/support.

Product description

The Mouse TNF alpha High Sensitivity ELISA Kit is an enzyme-linked immunosorbent assay for the quantitative detection of mouse TNF- α .

Summary

TNF- α is a multifunctional cytokine involved in many different pathways, in homeostasis and pathophysiology of mammals. It can show opposing biological effects suggesting complex regulatory mechanisms.

TNF- α , also known as cachectin, was first detected as a cytotoxic factor inducing lysis of certain tumor cells. The TNF- α gene is member 2 of the TNF-superfamily (consisting of at least 20 distinct members).

TNF- α release is mainly triggered by viral infections and endotoxins, lipopolysaccharides or other bacterial components, by tissue injury, DNA-damage and by IL-1, PDGF and TNF- α itself. It is primarily expressed in macrophages, but also in monocytes, neutrophils, NK-cells, mast-cells, endothelial cells and activated lymphocytes. TNF- α expression in endothelial cells and fibroblasts can be induced by IL-17.

The expression of other cytokines, chemokines, reactive oxygen intermediates, nitric oxide and prostaglandins is stimulated by TNF- α . The initially membrane bound TNF- α is enzymatically cleaved by TACE (= ADAM17). The soluble monomers aggregate to homotrimers and are secreted into blood and other biological fluids. The membrane bound and the soluble form are biologically active and bind to the TNF-receptors TNFR1 (= TNFRSF1A, p55-60) and TNFR2 (= TNFRSF1B, TNFR2, p75-80).

Upon ligand binding, receptors form trimers leading to conformational changes, protein dissociation (SODD = silencer of death domains, BAG4, Bcl2-associated athanogene 4) and association (TRADD = TNF-R1 associated death domain protein) and yielding the following biological activities:

- transcription of anti-apoptotic factors and proteins involved in cell proliferation and inflammation via binding of TRAF2 (TNF-R associated factor 2) and RIPK1 (TNF-R interacting serine-threonine kinase 1) and activation of the transcription factor NF- κ B.
- cell proliferation, differentiation but also apoptosis via TRAF2 binding, kinases activation, activation of c-Jun and ATF2 (JNK-MAPK-pathway).
- apoptosis via the binding of FADD (Fas associating protein with death domain) to TRADD and activation of caspases (including caspase 8 = FLICE).
- necrosis, a caspase independent cell death, mediated by NADPH oxidases, which form a complex with TRADD and RIPK1, leading to the generation of oxygen species.

TNF-R2 contains no DD (death domain), but exhibits its function via direct TRAF binding.

Thus the multiple biological functions of TNF- α comprise cellular proliferation and differentiation, tumorigenesis, apoptotic or necrotic

cell death (including certain tumor cell lines), immunoregulatory activities, lipid metabolism, coagulation and endothelial function. It promotes local or systemic inflammation (TNF- α is a potent pyrogen) and stimulates the acute phase response. Very high expressions of TNF- α after infection can lead to septic shock (TNF- α is highly cytotoxic.), whereas sustained low levels induce cachexia and inflammation.

Dysregulation of TNF- α is involved in many diseases.

For literature update refer to our website.

Principles of the test

An anti-mouse TNF- α coating antibody is adsorbed onto microwells.

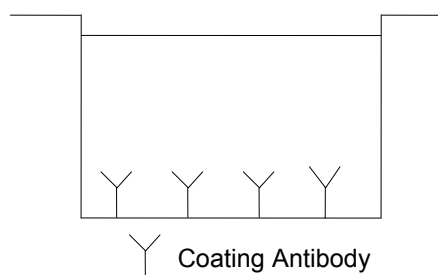


Fig. 1 Coated microwell

Mouse TNF- α present in the sample or standard binds to antibodies adsorbed to the microwells. A biotin-conjugated anti-mouse TNF- α antibody is added and binds to mouse TNF- α captured by the first antibody.

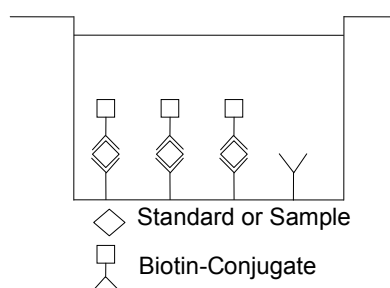


Fig. 2 First incubation

Following incubation unbound biotin-conjugated anti-mouse TNF- α antibody is removed during a wash step. Streptavidin-HRP is added and binds to the biotin-conjugated anti-mouse TNF- α antibody.

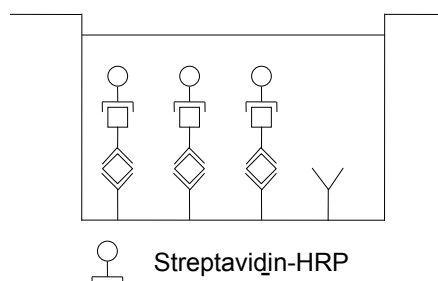


Fig. 3 Second incubation

Following incubation unbound Streptavidin-HRP is removed during a wash step, and amplification reagent I (Biotinyl-Tyramide) is added to the wells.

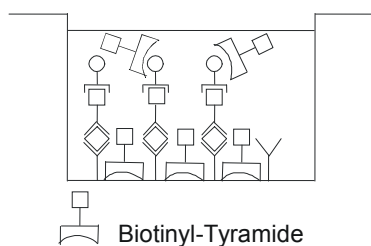


Fig. 4 Third incubation

Following incubation unbound amplification reagent I is removed during a wash step and amplification reagent II (Streptavidin-HRP) is added.

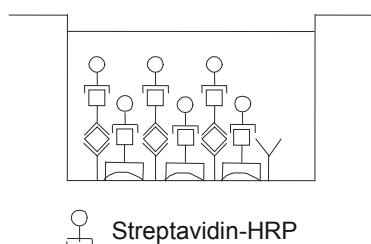


Fig. 5 Fourth incubation

Following incubation unbound amplification reagent II is removed during a wash step and substrate solution reactive with HRP is added.

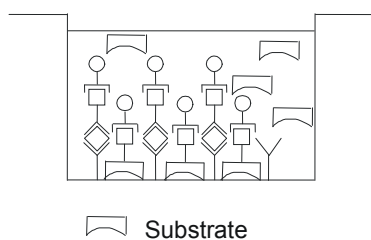


Fig. 6 Fifth incubation

A colored product is formed in proportion to the amount of mouse TNF- α present in the sample or standard. The reaction is terminated by addition of acid and absorbance is measured at 450 nm. A standard curve is prepared from 7 mouse TNF- α standard dilutions and mouse TNF- α sample concentration determined.

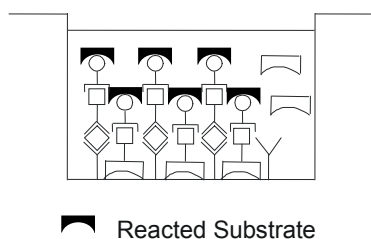


Fig. 7 Stop reaction

Principle of amplification reaction

The amplification reaction is based upon PerkinElmer Life Sciences TSA™ technology.

Amplification reagent I contains biotinyl-tyramide. HRP converts multiple biotinyl-tyramide molecules into highly reactive derivatives (free radicals). These free radicals bind covalently to any protein in the well.

Thus, the amount of reacted biotinyl-tyramide is proportional to the amount of HRP in the well.

Following incubation unbound biotinyl-tyramide is removed during a wash step. Amplification reagent II contains Streptavidin-HRP, which binds to the biotin sites created during the biotinyl-tyramide reaction,

thus multiplying the HRP molecules available at the surface for the substrate reaction.

Reagents provided

1 aluminum pouch with a Microwell Plate (12 strips of 8 wells each) coated with monoclonal antibody to mouse TNF- α

1 vial (70 μ L) Biotin-Conjugate anti-mouse TNF- α monoclonal antibody

1 vial (150 μ L) Streptavidin-HRP

2 vials mouse TNF- α Standard lyophilized, 200 pg/mL upon reconstitution

1 bottle (5 mL) Calibrator Diluent

1 bottle (12 mL) Sample Diluent

1 vial (5 mL) Assay Buffer Concentrate 20x (PBS with 1% Tween 20 and 10% BSA)

1 vial (7 mL) Amplification Diluent Concentrate (2x)

1 vial (75 μ L) Amplification Reagent I

Note: reagent contains ethyl alcohol

2 vials (15 μ L) Amplification Reagent II

2 bottles (50 mL) Wash Buffer Concentrate 20x (PBS with 1% Tween 20)

1 vial (15 mL) Substrate Solution (tetramethyl-benzidine)

1 vial (15 mL) Stop Solution (1M Phosphoric acid)

8 Adhesive Films

Storage instructions – ELISA kit

Store kit reagents between 2° and 8°C.

Immediately after use remaining reagents should be returned to cold storage (2° to 8°C). Expiry of the kit and reagents is stated on labels.

Expiry of the kit components can only be guaranteed if the components are stored properly, and if, in case of repeated use of one component, this reagent is not contaminated by the first handling.

Sample collection and storage instructions

Cell culture supernatant, serum and EDTA plasma were tested with this assay. Other biological samples might be suitable for use in the assay.

Samples containing a visible precipitate must be clarified prior to use in the assay. Do not use grossly hemolyzed or lipemic specimens.

Samples should be aliquoted and must be stored frozen at -20°C to avoid loss of bioactive mouse TNF- α . If samples are to be run within 24 hours, they may be stored at 2° to 8°C (refer to “Sample stability” on page 6).

Avoid repeated freeze-thaw cycles. Prior to assay, the frozen sample should be brought to room temperature slowly and mixed gently.

Materials required but not provided

- 5 mL and 10 mL graduated pipettes
- 5 μ L to 1000 μ L adjustable single channel micropipettes with disposable tips
- 50 μ L to 300 μ L adjustable multichannel micropipette with disposable tips
- Multichannel micropipette reservoir
- Beakers, flasks, cylinders necessary for preparation of reagents
- Device for delivery of wash solution (multichannel wash bottle or automatic wash system)
- Microplate shaker
- Microwell strip reader capable of reading at 450 nm (620 nm as optional reference wave length)
- Glass-distilled or deionized water
- Statistical calculator with program to perform regression analysis

Precautions for use

- All chemicals should be considered as potentially hazardous. We therefore recommend that this product is handled only by those persons who have been trained in laboratory techniques and that it is used in accordance with the principles of good laboratory practice. Wear suitable protective clothing such as laboratory overalls, safety glasses, and gloves. Care should be taken to avoid contact with skin or eyes. In the case of contact with skin or eyes wash immediately with water. See material safety data sheet(s) and/or safety statement(s) for specific advice.
- Reagents are intended for research use only and are not for use in diagnostic or therapeutic procedures.
- Do not mix or substitute reagents with those from other lots or other sources.
- Do not use kit reagents beyond expiration date on label.
- Do not expose kit reagents to strong light during storage or incubation.
- Do not pipet by mouth.
- Do not eat or smoke in areas where kit reagents or samples are handled.
- Avoid contact of skin or mucous membranes with kit reagents or samples.
- Rubber or disposable latex gloves should be worn while handling kit reagents or samples.
- Avoid contact of substrate solution with oxidizing agents and metal.
- Avoid splashing or generation of aerosols.
- To avoid microbial contamination or cross-contamination of reagents or samples that may invalidate the test, use disposable pipette tips and/or pipettes.
- Use clean, dedicated reagent trays for dispensing the conjugate and substrate reagent.
- Exposure to acid inactivates the conjugate.
- Glass-distilled water or deionized water must be used for reagent preparation.
- Substrate solution must be at room temperature prior to use.
- Decontaminate and dispose samples and all potentially contaminated materials as if they could contain infectious agents. The preferred method of decontamination is autoclaving for a minimum of 1 hour at 121.5°C.
- Liquid wastes not containing acid and neutralized waste may be mixed with sodium hypochlorite in volumes such that the final mixture contains 1.0% sodium hypochlorite. Allow 30 minutes for effective decontamination. Liquid waste containing acid must be neutralized prior to the addition of sodium hypochlorite.

Preparation of reagents

- Buffer Concentrates should be brought to room temperature and should be diluted before starting the test procedure.
- If crystals have formed in the Buffer Concentrates, warm them gently until they have completely dissolved.

Wash buffer (1x)

- Pour entire contents (50 mL) of the Wash Buffer Concentrate (20x) into a clean 1000 mL graduated cylinder. Bring to final volume of 1000 mL with glass-distilled or deionized water. Mix gently to avoid foaming.
- Transfer to a clean wash bottle and store at 2° to 25°C. Please note that Wash Buffer (1x) is stable for 30 days.
- Wash Buffer (1x) may also be prepared as needed according to the following table:

Number of Strips	Wash Buffer Concentrate (20x) (mL)	Distilled Water (mL)
1 - 6	25	475
1 - 12	50	950

Assay buffer (1x)

- Pour the entire contents (5 mL) of the Assay Buffer Concentrate (20x) into a clean 100 mL graduated cylinder. Bring to final volume of 100 mL with distilled water. Mix gently to avoid foaming.
- Store at 2° to 8°C. Please note that the Assay Buffer (1x) is stable for 30 days.
- Assay Buffer (1x) may also be prepared as needed according to the following table:

Number of Strips	Assay Buffer Concentrate (20x) (mL)	Distilled Water (mL)
1 - 6	2.5	47.5
1 - 12	5.0	95.0

Biotin-Conjugate

Note: The Biotin-Conjugate should be used within 30 minutes after dilution.

Make a 1:100 dilution of the concentrated Biotin-Conjugate solution with Assay Buffer (1x) in a clean plastic tube as needed according to the following table:

Number of Strips	Biotin-Conjugate (mL)	Assay Buffer (1x) (mL)
1 - 6	0.03	2.97
1 - 12	0.06	5.94

Streptavidin-HRP

Note: The Streptavidin-HRP should be used within 30 minutes after dilution.

Make a 1:200 dilution of the concentrated Streptavidin-HRP solution with Assay Buffer (1x) in a clean plastic tube as needed according to the following table:

Number of Strips	Streptavidin-HRP (mL)	Assay Buffer (1x) (mL)
1 - 6	0.03	5.97
1 - 12	0.06	11.94

Mouse TNF- α standard

- Reconstitute mouse TNF- α standard by addition of Calibrator Diluent (for subsequent measurement of serum or plasma samples) or Sample Diluent (for subsequent measurement of cell culture supernatant samples).
- Reconstitution volume is stated on the label of the standard vial. Swirl or mix gently to insure complete and homogeneous solubilisation (concentration of reconstituted standard = 200 pg/mL). Allow the standard to reconstitute for 10-30 minutes. Mix well prior to making dilutions.
- After usage remaining standard cannot be stored and has to be discarded.

External standard dilution

- Label 6 tubes, one for each standard point: S2, S3, S4, S5, S6, S7. The reconstituted standard serves as S1.
- Prepare 1:2 serial dilutions for the standard curve as follows: Pipette 150 μ L of Calibrator Diluent (for subsequent measurement of serum or plasma samples) or Sample Diluent (for subsequent measurement of cell culture supernatant samples) into each tube.
- Pipette 150 μ L of reconstituted standard (concentration of the standard S1 = 200 pg/mL) into the tube labeled with S2 and mix (concentration of the standard S2 = 100 pg/mL).
- Pipette 150 μ L of this dilution into the second tube labeled S3 and mix thoroughly before the next transfer.

- Repeat serial dilutions 4 more times thus creating the points of the standard curve (see Figure 8).

Calibrator Diluent (serum or plasma samples) or Sample Diluent (cell culture supernatant samples) serves as blank.

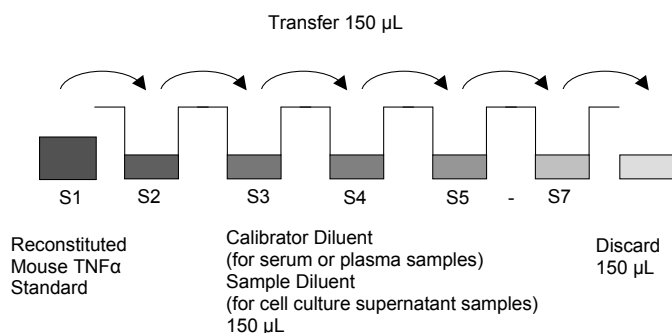


Fig. 8 External standard dilution

Amplification diluent (1x)

Preparation of Amplification Diluent (1x) has to be done immediately prior to use. Make a 1:2 dilution of the concentrated Amplification Diluent (2x) as needed according to the following table:

Number of Strips	Amplification Diluent (2x) (mL)	Distilled Water (mL)
1 - 6	3	3
1 - 12	6	6

Amplification solution I

- Prepare Amplification Solution I immediately prior to application on the plate.
- Dilute Amplification Reagent I in Amplification Diluent (1x) as indicated in the Certificate of Analysis.
- Discard immediately any prediluted Amplification Solution I after usage.

Amplification solution II

- Prepare Amplification Solution II immediately prior to application on the plate.
- Centrifuge vial for a few seconds in a microcentrifuge before opening to collect liquid trapped in the lid.
- Dilute Amplification Reagent II in Assay Buffer (1x) as indicated in the Certificate of Analysis.
- Discard immediately any prediluted Amplification Solution II after usage.

Test protocol

IMPORTANT!

- Because this ELISA is a high sensitivity system, it is extremely important to stick exactly to the manual (washing procedure, chronology of and preparation of solutions, incubation time) to obtain optimal test performance.
- Amplification Solutions must be prepared immediately prior to application on the plate! It is extremely important to wash the wells properly to obtain a good test performance.
- Shaking is absolutely necessary for an optimal test performance. Protect microwell plate from light during incubation steps

- Determine the number of microwell strips required to test the desired number of samples plus appropriate number of wells needed for running blanks and standards. Each sample, standard, blank and optional control sample should be assayed in duplicate. Remove extra microwell strips from holder and store in foil bag with the desiccant provided at 2°-8°C sealed tightly.
- Prepare Biotin-Conjugate (see "Biotin-Conjugate" on page 3).

- Wash the microwell strips twice with exactly 400 μ L Wash Buffer per well with thorough aspiration of microwell contents between washes. Allow the Wash Buffer to sit in the wells for about 10 – 15 seconds before aspiration. Soaking is highly recommended between the washes to obtain a good test performance! Take care not to scratch the surface of the microwells.

After the last wash step, empty wells and tap microwell strips on absorbent pad or paper towel to remove excess Wash Buffer. Use the microwell strips immediately after washing. Do not allow wells to dry.

- External standard dilution (see "External standard dilution" on page 3):
- Add 50 μ L of Sample Diluent in duplicate to all wells.
- Add 50 μ L of extern diluted Standards (refer to "External standard dilution" on page 3) in duplicate to the corresponding standard well (see Table 1).

Table 1 Example of the arrangement of blanks, standards and samples in the microwell strips.

	1	2	3	4
A	Standard 1 200 pg/mL	Standard 1 200 pg/mL	Sample 1	Sample 1
B	Standard 2 100 pg/mL	Standard 2 100 pg/mL	Sample 2	Sample 2
C	Standard 3 50 pg/mL	Standard 3 50 pg/mL	Sample 3	Sample 3
D	Standard 4 25 pg/mL	Standard 4 25 pg/mL	Sample 4	Sample 4
E	Standard 5 12.5 pg/mL	Standard 5 12.5 pg/mL	Sample 5	Sample 5
F	Standard 6 6.25 pg/mL	Standard 6 6.25 pg/mL	Sample 6	Sample 6
G	Standard 7 3.13 pg/mL	Standard 7 3.13 pg/mL	Sample 7	Sample 7
H	Blank	Blank	Sample 8	Sample 8

- Add 50 μ L of each sample in duplicate to the sample wells.
- Add 50 μ L of Calibrator Diluent (for serum or plasma samples) or Sample Diluent (for cell culture supernatant samples) in duplicate to the blank wells.
- Add 50 μ L of Biotin Conjugate to all wells.
- Cover with an adhesive film and incubate at room temperature (18° to 25°C) for 2 hours on a microplate shaker. (Shaking is absolutely necessary for an optimal test performance.)
- Prepare Streptavidin-HRP (refer to "Streptavidin-HRP" on page 3).
- Remove adhesive film and empty wells. Wash microwell strips 6 times according to point 3 of the test protocol. Proceed immediately to the next step.
- Add 100 μ L of diluted Streptavidin-HRP to all wells, including the blank wells.
- Cover with an adhesive film and incubate at room temperature (18° to 25°C) for 1 hour on a microplate shaker. (Shaking is absolutely necessary for an optimal test performance.)
- Prepare Amplification Solution I diluted in Amplification Diluent (1x) (see "Amplification solution I" on page 4) immediately prior to use.
- Remove adhesive film and empty wells. Wash microwell strips 6 times according to point 3 of the test protocol. Proceed immediately to the next step.
- Add 100 μ L of Amplification Solution I to all wells, including the blank wells.
- Cover with an adhesive film and incubate at room temperature (18° to 25°C) for exactly 15 minutes. (Shaking is absolutely necessary for an optimal test performance.)
- Prepare Amplification Solution II diluted in Assay Buffer Concentrate(1X) (see "Amplification solution II" on page 4) immediately prior to use.

20. Remove adhesive film and empty wells. Wash microwell strips 6 times according to point 3 of the test protocol. Proceed immediately to the next step.
21. Add 100 μ L of Amplification Solution II to all wells, including the blank wells.
22. Cover with an adhesive film and incubate at room temperature (18° to 25°C) for exactly 30 minutes on a microplate shaker. (Shaking is absolutely necessary for an optimal test performance.)
23. Remove adhesive film and empty wells. Wash microwell strips 6 times according to point 3 of the test protocol. Proceed immediately to the next step.
24. Pipette 100 μ L of TMB Substrate Solution to all wells.
25. Incubate the microwell strips at room temperature (18° to 25°C) for about 10-20 minutes. Avoid direct exposure to intense light.

The color development on the plate should be monitored and the substrate reaction stopped (see next point of this protocol) before positive wells are no longer properly recordable. Determination of the ideal time period for color development has to be done individually for each assay.

It is recommended to add the stop solution when the highest standard has developed a dark blue color. Alternatively the color development can be monitored by the ELISA reader at 620 nm. The substrate reaction should be stopped as soon as Standard 1 has reached an OD of 0.9 – 0.95.

26. Stop the enzyme reaction by quickly pipetting 100 μ L of Stop Solution into each well. It is important that the Stop Solution is spread quickly and uniformly throughout the microwells to completely inactivate the enzyme. Results must be read immediately after the Stop Solution is added or within one hour if the microwell strips are stored at 2 - 8°C in the dark.
27. Read absorbance of each microwell on a spectro-photometer using 450 nm as the primary wave length (optionally 620 nm as the reference wave length; 610 nm to 650 nm is acceptable). Blank the plate reader according to the manufacturer's instructions by using the blank wells. Determine the absorbance of both the samples and the standards.

Note: If instructions in this protocol have been followed samples have not been diluted, the concentration read from the standard curve must not be multiplied by a dilution factor.

Calculation of results

- Calculate the average absorbance values for each set of duplicate standards and samples. Duplicates should be within 20 percent of the mean value.
- Create a standard curve by plotting the mean absorbance for each standard concentration on the ordinate against the mouse TNF- α concentration on the abscissa. Draw a best fit curve through the points of the graph (a 5-parameter curve fit is recommended).
- To determine the concentration of circulating mouse TNF- α for each sample, first find the mean absorbance value on the ordinate and extend a horizontal line to the standard curve. At the point of intersection, extend a vertical line to the abscissa and read the corresponding mouse TNF- α concentration.
- Calculation of samples with a concentration exceeding standard 1 may result in incorrect, low mouse TNF- α levels. Such samples require further external predilution according to expected mouse TNF- α values with Sample Diluent in order to precisely quantitate the actual mouse TNF- α level.
- It is suggested that each testing facility establishes a control sample of known mouse TNF- α concentration and runs this additional control with each assay. If the values obtained are not within the expected range of the control, the assay results may be invalid.

- A representative standard curve is shown in Figure 9. This curve cannot be used to derive test results. Each laboratory must prepare a standard curve for each group of microwell strips assayed.

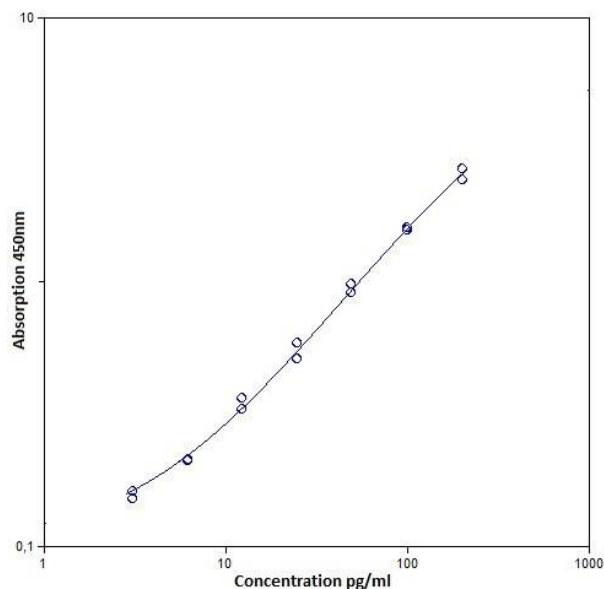


Fig. 9 Representative standard curve for mouse TNF- α ELISA. Mouse TNF- α was diluted in serial 2-fold steps in Sample Diluent. Do not use this standard curve to derive test results. A standard curve must be run for each group of microwell strips assayed.

Table 2 Typical data using the mouse TNF- α ELISA

Measuring wavelength: 450 nm

Reference wavelength: 620 nm

Standard	Mouse TNF- α Concentration (pg/mL)	O.D. at 450 nm	Mean O.D. at 450 nm	C.V. (%)
1	200.00	2.669	2.414	5.0
2	100.00	1.600	1.557	1.3
3	50.00	0.975	0.902	3.9
4	25.00	0.585	0.510	6.9
5	12.50	0.361	0.326	5.0
6	6.25	0.211	0.208	0.7
7	3.13	0.150	0.160	3.1
Blank	0.00	0.100	0.120	9.0

The OD values of the standard curve may vary according to the conditions of assay performance (e.g. operator, pipetting technique, washing technique or temperature effects). Furthermore shelf life of the kit may affect enzymatic activity and thus color intensity. Values measured are still valid.

Limitations

- Since exact conditions may vary from assay to assay, a standard curve must be established for every run.
- Bacterial or fungal contamination of either screen samples or reagents or cross-contamination between reagents may cause erroneous results.
- Disposable pipette tips, flasks or glassware are preferred, reusable glassware must be washed and thoroughly rinsed of all detergents before use.

- Improper or insufficient washing at any stage of the procedure will result in either false positive or false negative results. Empty wells completely before dispensing fresh wash solution, fill with Wash Buffer as indicated for each wash cycle and do not allow wells to sit uncovered or dry for extended periods.

Performance characteristics

Sensitivity

The limit of detection of mouse TNF- α defined as the analyte concentration resulting in an absorbance significantly higher than that of the dilution medium (mean plus 2 standard deviations) was determined to be 0.75 pg/mL (mean of 4 independent assays).

Reproducibility

Intra-assay

Reproducibility within the assay was evaluated in 3 independent experiments. Each assay was carried out with 6 replicates of 7 serum, samples containing different concentrations of mouse TNF- α . 2 standard curves were run on each plate. Data below show the mean mouse TNF- α concentration and the coefficient of variation for each sample (see Table 3). The calculated overall intra-assay coefficient of variation was 9.2%.

Table 3 The mean mouse TNF- α concentration and the coefficient of variation for each sample

Sample	Experiment	Mean Mouse TNF- α Concentration (pg/mL)	Coefficient of Variation (%)
1	1	114.8	6.6%
	2	105.9	14.9%
	3	114.4	11.0%
2	1	94.5	5.4%
	2	91.7	11.3%
	3	87.1	6.8%
3	1	83.2	7.5%
	2	72.8	10.6%
	3	73.9	8.6%
4	1	67.7	8.4%
	2	61.9	9.9%
	3	70.5	8.2%
5	1	44.9	8.5%
	2	43.1	7.9%
	3	48.7	6.3%
6	1	45.2	9.8%
	2	43.6	12.8%
	3	43.9	8.3%
7	1	43.4	4.4%
	2	41.6	13.9%
	3	43.2	12.4%

Inter-assay

Assay to assay reproducibility within one laboratory was evaluated in 3 independent experiments. Each assay was carried out with 6 replicates of 7 serum samples containing different concentrations of mouse TNF- α . 2 standard curves were run on each plate. Data below show the mean mouse TNF- α concentration and the coefficient of

variation calculated on 18 determinations of each sample (see Table 4). The calculated overall inter-assay coefficient of variation was 4.7%.

Table 4 The mean mouse TNF- α concentration and the coefficient of variation of each sample

Sample	Mean Mouse TNF- α Concentration (pg/mL)	Coefficient of Variation (%)
1	112	4.5%
2	91	4.1%
3	77	7.5%
4	67	6.6%
5	46	6.3%
6	44	1.9%
7	43	2.4%

Spike recovery

The spike recovery was evaluated by spiking 3 levels of mouse TNF- α into serum, EDTA plasma and cell culture supernatant. Recoveries were determined with 4 replicates each. The amount of endogenous mouse TNF- α in unspiked sample was subtracted from the spike values.

Sample matrix	Spike high (%)	Spike medium (%)	Spike low (%)
Serum	105	90	94
Plasma (EDTA)	120	99	103
Cell culture supernatant	114	104	104

Dilution parallelism

Serum, plasma (EDTA) and cell culture supernatant samples with different levels of mouse TNF- α were analyzed at serial 2 fold dilutions with 4 replicates each.

Sample matrix	Recovery of Exp. Val.		
	Dilution	Mean (%)	Range (%)
Serum	1:4	102	96-109
	1:8	112	103-126
	1:16	111	89-129
Plasma (EDTA)	1:4	96	-
	1:8	106	-
	1:16	103	-
Cell culture supernatant	1:4.5	100	92-107
	1:13.5	95	86-104
	1:40.5	91	56-126

Sample stability

Freeze-Thaw stability

Aliquots of spiked serum samples were stored at -20°C and thawed 5 times, and the mouse TNF- α levels determined. There was no significant loss of mouse TNF- α immunoreactivity detected by freezing and thawing.

Storage stability

Aliquots of spiked serum samples were stored at -20°C, 2-8°C, room temperature, and at 37°C, and the mouse TNF- α level determined after 24 hours. There was no significant loss of mouse TNF- α immunoreactivity detected during storage at -20°C and 2-C.

Specificity

The interference of circulating factors of the immune system was evaluated by spiking these proteins at physiologically relevant concentrations into a mouse TNF- α positive serum.

There was no cross reactivity observed with mouse IL-2, GM-CSF, IL-5, IL-6, IFN-g, IL-4, IL-10, IL-17, IL-1a, TNFR II and CD40L.

Calibration

The immunoassay is calibrated with highly purified recombinant mouse TNF- α which has been evaluated against the international Reference Standard NIBSC 88/532.

NIBSC/WHO 88/532 approximate value is quantified in arbitrarily units. 1 Arbitrarily Unit corresponds to 2.5 pg.

Reagent preparation summary

Wash buffer (1x)

Add Wash Buffer Concentrate 20x (50 mL) to 950 mL distilled water.

Number of Strips	Wash Buffer Concentrate (20x) (mL)	Distilled Water (mL)
1 - 6	25	475
1 - 12	50	950

Assay buffer (1x)

Add Assay Buffer Concentrate 20x (5 mL) to 95 mL distilled water.

Number of Strips	Assay Buffer Concentrate (20x) (mL)	Distilled Water (mL)
1 - 6	2.5	47.5
1 - 12	5.0	95.0

Biotin-Conjugate

Make a 1:100 dilution of Biotin-Conjugate in Assay Buffer (1x):

Number of Strips	Biotin-Conjugate (mL)	Assay Buffer (1x) (mL)
1 - 6	0.03	2.97
1 - 12	0.06	5.94

Streptavidin-HRP

Make a 1:200 dilution of Streptavidin-HRP in Assay Buffer (1x):

Number of Strips	Streptavidin-HRP (mL)	Assay Buffer (1x) (mL)
1 - 6	0.03	5.97
1 - 12	0.06	11.94

Mouse TNF- α standard

1. Reconstitute lyophilized mouse TNF- α standard by addition of Calibrator Diluent (for subsequent measurement of serum or plasma samples) or Sample Diluent (for subsequent measurement of cell culture supernatant samples). Reconstitution volume is stated on the label of the standard vial.
2. The reconstituted mouse TNF- α standard serves as S1.

Amplification diluent (1x)

Prepare Amplification Diluent (1x) immediately prior to use.

Number of Strips	Amplification Diluent (2x) (mL)	Distilled Water (mL)
1 - 6	3	3
1 - 12	6	6

Amplification solution I

Dilute Amplification Reagent I in Amplification Diluent (1x) immediately prior to application on the plate as indicated in the Certificate of Analysis.

Amplification solution II

Centrifuge vial for a few seconds in a micro-centrifuge before opening to collect liquid trapped in the lid. Dilute Amplification Solution II in Assay Buffer (1x) immediately prior to application on the plate as indicated in the Certificate of Analysis.

Test protocol summary

Note: Prepare Amplification Solutions immediately prior to application on the plate. It is extremely important to wash the wells properly to obtain a good test performance.

1. Determine the number of microwell strips required.
2. Prepare Biotin-Conjugate.
3. Wash microwell strips twice with Wash Buffer.
4. Add 50 μ L Sample Diluent in duplicate, to the all wells.
5. Add 50 μ L of prepared Standard dilutions (see "External standard dilution" on page 3) in duplicate to standard wells.
6. Add 50 μ L samples in duplicate, to designated sample wells.
7. Add 50 μ L of Calibrator Diluent (for serum or plasma samples) or Sample Diluent (for cell culture supernatant samples) in duplicate to the blank wells.
8. Add 50 μ L Biotin-Conjugate to all wells.
9. Cover microwell strips and incubate 2 hours at room temperature (18° to 25°C). (Shaking is absolutely necessary for an optimal test performance.)
10. Prepare Streptavidin-HRP.
11. Empty and wash microwell strips 6 times with Wash Buffer.
12. Add 100 μ L diluted Streptavidin-HRP to all wells.
13. Cover microwell strips and incubate 1 hour at room temperature (18° to 25°C). (Shaking is absolutely necessary for an optimal test performance.)
14. Prepare Amplification Solution I diluted in Amplification Diluent (1x) immediately prior to application on the plate.
15. Empty and wash microwell strips 6 times with Wash Buffer.
16. Add 100 μ L Amplification Solution I to all wells.
17. Cover microwell strips and incubate for exactly 15 minutes at room temperature (18° to 25°C). (Shaking is absolutely necessary for an optimal test performance.)
18. Prepare Amplification Solution II diluted in Assay Buffer immediately prior to application on the plate.
19. Empty and wash microwell strips 6 times with Wash Buffer.
20. Add 100 μ L Amplification Solution II to all wells.
21. Cover microwell strips and incubate for exactly 30 minutes at room temperature (18° to 25°C).
22. Empty and wash microwell strips 6 times with Wash Buffer.
23. Add 100 μ L of TMB Substrate Solution to all wells.
24. Incubate the microwell strips for about 10-20 minutes at room temperature (18° to 25°C).
25. Add 100 μ L Stop Solution to all wells.
26. Blank microwell reader and measure color intensity at 450 nm.

If instructions in this protocol have been followed samples have not been diluted, the concentration read from the standard curve must not be multiplied by a dilution factor.

Customer and technical support

Visit thermofisher.com/support for the latest service and support information.

- Worldwide contact telephone numbers
- Product support information
 - Product FAQs
 - Software, patches, and updates
 - Training for many applications and instruments
- Order and web support
- Product documentation
 - User guides, manuals, and protocols
 - Certificates of Analysis
 - Safety Data Sheets (SDSs; also known as MSDSs)

Note: For SDSs for reagents and chemicals from other manufacturers, contact the manufacturer.

Limited product warranty

Life Technologies Corporation and/or its affiliate(s) warrant their products as set forth in the Life Technologies' General Terms and

Conditions of Sale at www.thermofisher.com/us/en/home/global/terms-and-conditions.html. If you have any questions, please contact Life Technologies at www.thermofisher.com/support.



Bender MedSystems GmbH | Campus Vienna Biocenter 2 | 1030 Vienna, Austria

For descriptions of symbols on product labels or product documents, go to thermofisher.com/symbols-definition.

The information in this guide is subject to change without notice.

DISCLAIMER: TO THE EXTENT ALLOWED BY LAW, THERMO FISHER SCIENTIFIC INC. AND/OR ITS AFFILIATE(S) WILL NOT BE LIABLE FOR SPECIAL, INCIDENTAL, INDIRECT, PUNITIVE, MULTIPLE, OR CONSEQUENTIAL DAMAGES IN CONNECTION WITH OR ARISING FROM THIS DOCUMENT, INCLUDING YOUR USE OF IT.

Important Licensing Information: These products may be covered by one or more Limited Use Label Licenses. By use of these products, you accept the terms and conditions of all applicable Limited Use Label Licenses.

©2019 Thermo Fisher Scientific Inc. All rights reserved. All trademarks are the property of Thermo Fisher Scientific and its subsidiaries unless otherwise specified. TSA is a trademark of PerkinElmer Life Sciences.