Pierce[™] Dye and Biotin Removal Plate

96-well plates for processing 40–100 µL sample volumes

Catalog Numbers A44304 and A44305

Doc. Part No. 2162738 Pub. No. MAN0018805 Rev. A.0

WARNING! Read the Safety Data Sheets (SDSs) and follow the handling instructions. Wear appropriate protective eyewear, clothing, and gloves. Safety Data Sheets (SDSs) are available from **thermofisher.com/support**.

Product description

The Thermo Scientific[™] Pierce[™] Dye and Biotin Removal Plate contains a ready-to-use resin that is uniquely designed for rapid removal of non-conjugated fluorescent dyes, biotin, reducing agents, and crosslinkers with exceptional protein recovery. One plate can process 96 samples in five minutes, enabling high-throughput sample purification for downstream analysis.

Removal of free dye after a labeling reaction is essential for the accurate determination of dye-to-protein ratios. The Pierce^T Dye and Biotin Removal Plate can be used to remove most fluorescent dyes. For optimal protein recovery and dye removal, ensure that the appropriate amount of sample and buffer conditions are used.

This product is recommended for the removal of unreacted biotin, fluorescent dyes, crosslinkers, and reducing agents from proteins >7 kDa.

Contents and storage

Product	Cat. No.	Amount ^[1]	Storage
Pierce [™] Dye and Biotin Removal Plate, 96-well ^[2]	A44304	2 filter plates	4°C
	A44305	4 filter plates	4 U

^[1] Two wash plates and two collection plates are included in each package.

 $^{[2]}$ Each well contains approximately 420 μL of settled resin in a 0.1 N NaCl/0.05% sodium azide solution.

Required materials not supplied

Unless otherwise indicated, all materials are available through **thermofisher.com**. MLS: Fisher Scientific (**fisherscientific.com**) or other major laboratory supplier.

Item	Source
Equipment	
Variable-speed centrifuge with rotor and adaptor capable of handling stacked plates [4.4 cm height] at 1,000 $\times g$	MLS
Multichannel pipette	MLS

ltem	Source	
Consumables		
Pipette tips (as recommended by pipette manufacturer)	thermofisher.com/ pipettetips	

Procedural guidelines

- Do not reuse the purification resin.
- Limit DMF and other organic solvents to $\leq 10\%$.
- Do not use this product for desalting or buffer exchange.

Prepare the plate

- 1. Equilibrate the filter plate to room temperature.
- 2. Use the tab to remove the seal from the bottom of the plate, then place the filter plate on top of the wash plate.
- 3. Remove the seal from the top of the filter plate.
- 4. Place the plate assembly into a 96-well plate-carrier rotor, then centrifuge at $1,000 \times g$ for 2 minutes to remove the storage buffer. Discard the flow-through.
- 5. (*Optional*) Add 300–400 μ L of equilibration buffer to each well of the filter plate, then centrifuge at 1,000 × *g* for 2 minutes. Discard the flow-through.

Rinse the wash plate three times with deionized water, allow the plate to dry, then store for future use.

Process the samples

- 1. Place the prepared filter plate on top of a blue collection plate, aligning the alphanumeric grids of the plates.
- Apply 40–100 μL of sample to the center of the settled resin. Touch the pipette tip to the resin to expel the entire sample.
- **3.** Centrifuge the plate assembly at $1,000 \times g$ for 2 minutes.

The sample is in the blue collection plate.



Troubleshooting

Observation	Possible cause	Recommended action
Sample or buffer does not flow A problem occurred during		Ensure that the centrifuge is in proper working condition.
through the resin	centrifugation.	Ensure that the bottom plug of the column is removed and the top cap is loosened.
Insufficient removal of free dye	Low resin-to-free dye ratio.	Repeat the procedure using a higher resin-to-free dye ratio.
Sample contains contaminants The sample was not loaded on the column properly.		Add the sample directly to the center of the settled resin, then carefully touch the pipette tip to the resin to expel the entire sample.
		Avoid contact with the sides of the column.
	A problem occurred during centrifugation.	Do not exceed the recommended centrifugation speed or time.

Related products

Product	Cat. No.
Pierce [™] Biotin Quantitation Kit	28005
Pierce [™] Fluorescence Biotin Quantitation Kit	46610
Pierce™ Rapid Gold BCA Protein Assay Kit	A53225
Slide-A-Lyzer™ G2 Dialysis Cassettes, 20K MWCO, 70 mL, 6 units	87738
EZ-Link™ Sulfo-NHS-SS-Biotin, No-Weigh™ Format	A39258
EZ-Link™ Hydrazide-PEG4-Biotin	21360
Bond-Breaker™ TCEP Solution, Neutral pH	77720
Pierce™ TCEP-HCl, No-Weigh™ Format	A35349
BS3 (bis(sulfosuccinimidyl)suberate), No-Weigh™ Format	A39266
Alexa Fluor™ 660 NHS Ester (Succinimidyl Ester)	A20007
DyLight™ 488 NHS Ester	46402
Pierce™ DTT (Dithiothreitol)	A39255
DTNB (Ellman's Reagent) (5,5-dithio-bis-(2- nitrobenzoic acid)	22582
Pierce™ BCA Protein Assay Kit	23225

Limited product warranty

Life Technologies Corporation and/or its affiliate(s) warrant their products as set forth in the Life Technologies' General Terms and Conditions of Sale at www.thermofisher.com/us/en/home/global/terms-and-conditions.html. If you have any questions, please contact Life Technologies at www.thermofisher.com/support.

Pierce Biotechnology, Inc. | Thermo Fisher Scientific | 3747 N. Meridian Road | Rockford, Illinois 61101 USA

For descriptions of symbols on product labels or product documents, go to **thermofisher.com/symbols-definition**.

The information in this guide is subject to change without notice.

DISCLAIMER: TO THE EXTENT ALLOWED BY LAW, THERMO FISHER SCIENTIFIC INC. AND/OR ITS AFFILIATE(S) WILL NOT BE LIABLE FOR SPECIAL, INCIDENTAL, INDIRECT, PUNITIVE, MULTIPLE, OR CONSEQUENTIAL DAMAGES IN CONNECTION WITH OR ARISING FROM THIS DOCUMENT, INCLUDING YOUR USE OF IT.

Revision history: Pub. No. MAN0018805

Revision	Date	Description
A.0	6 August 2019	New document for new product.

Important Licensing Information: These products may be covered by one or more Limited Use Label Licenses. By use of these products, you accept the terms and conditions of all applicable Limited Use Label Licenses.

©2019 Thermo Fisher Scientific Inc. All rights reserved. All trademarks are the property of Thermo Fisher Scientific and its subsidiaries unless otherwise specified.

