



High-performance quantitative RT-PCR in one step


To achieve success in quantitative RT-PCR, you need the right RT and PCR technologies. With the Platinum® Quantitative RT-PCR ThermoScript™ One-Step System, you get both with accurate real-time quantification of RNA. Obtain the high specificity, yield and sensitivity essential to your experiments, with one-step convenience.

One-step qRT-PCR protocol

The Platinum® Quantitative RT-PCR ThermoScript™ One-Step System is optimized to deliver accuracy and reliability. It combines two high-performance enzymes, ThermoScript™ reverse transcriptase and Platinum® *Taq* DNA polymerase, in a pre-mixed formulation. You use just one tube, where cDNA synthesis and amplification occur sequentially. Simply add total or poly(A)⁺ RNA, gene-specific reverse transcription and amplification primers, and fluorogenic

probes to your enzyme mixture and you're ready to go. This one-step protocol reduces reaction variability, lowers risk of contamination, and provides highly sensitive and efficient qRT-PCR. The convenient format makes it well suited for high throughput applications. As Figure 1 shows, the system provides exceptional sensitivity—you can achieve precise quantification from as low as 10 molecules of RNA template or 1 pg of total RNA.

Figure 1 - One-Step Quantitative RT-PCR

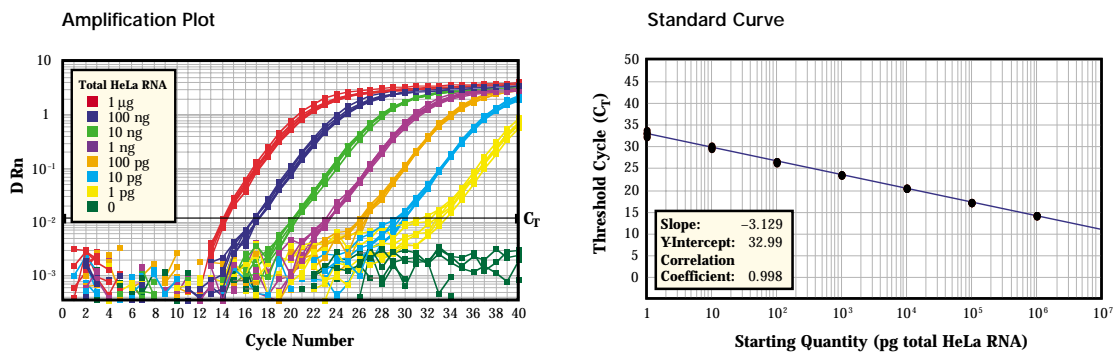


Quantitation of β -actin mRNA. Amplification of six replicates of various amounts of HeLa total RNA was performed with the Platinum® Quantitative RT-PCR ThermoScript™ One-Step System and detected using a 6-FAM-labeled *TaqMan*® probe on an ABI PRISM® 7700. TAMRA was used as a passive reference for normalization of reporter signal (ΔRn).

Thermostability for enhanced cDNA synthesis

ThermoScript™ RT, a cloned avian RNase H⁻ reverse transcriptase, is engineered to provide highly specific cDNA synthesis at temperatures up to 70°C. It offers the same performance advantages of SuperScript™ II with added thermostability. With its reduced RNase H activity,

ThermoScript™ improves the yields of full-length cDNA. With its high thermostability, you can overcome the challenges of high GC content and secondary structure to achieve highly efficient cDNA synthesis at elevated temperatures (Figure 2).

Figure 2 - High yield RT-PCR of GC-rich RNA with extensive secondary structure

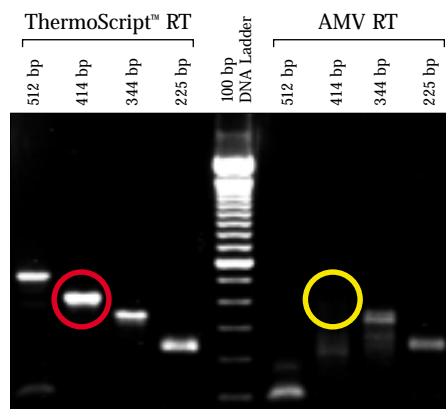


cDNA was synthesized from 10 ng total RNA using 15 units of ThermoScript™ RT or 15 units AMV RT according to manufacturers' recommendations. One tenth of each reaction was amplified with Platinum® *Taq* DNA Polymerase High Fidelity.

Platinum® automatic hot-start for high specificity

Along with ThermoScript™ Plus RT, you get the power of Platinum® technology, which confers a high level of specificity in PCR. Platinum® *Taq* DNA polymerase provides antibody-mediated automatic hot-start capability. With significantly reduced mispriming and non-specific amplification, you get higher yields, vastly improved specificity, and greater sensitivity than with other DNA polymerases (Figure 3).

Figure 3 - Complete inhibition of DNA polymerase activity at room temperature

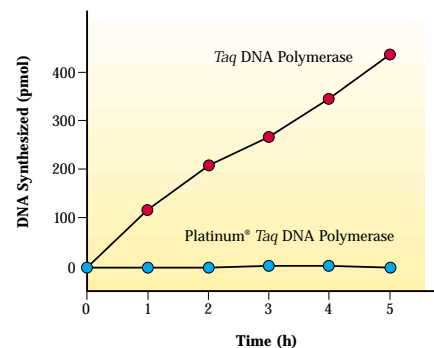


Activity assays were run at 37°C for 5 hours with *Taq* DNA polymerase and Platinum® *Taq* DNA Polymerase.

The right combination to surpass the competition

Together, the combination of RT and PCR technologies gives you the elements you need—yield, specificity, and sensitivity—for high performance qRT-PCR. Figure 4 compares the results achieved with the Platinum® Quantitative RT-PCR ThermoScript™

One-Step System to the leading competitor's. The Platinum® kit demonstrates lower Ct (threshold cycle) and gives higher yield, making it more sensitive to detect low copies of mRNA transcripts.

Figure 4 - High yield with Platinum® Quantitative RT-PCR ThermoScript™ One-Step System

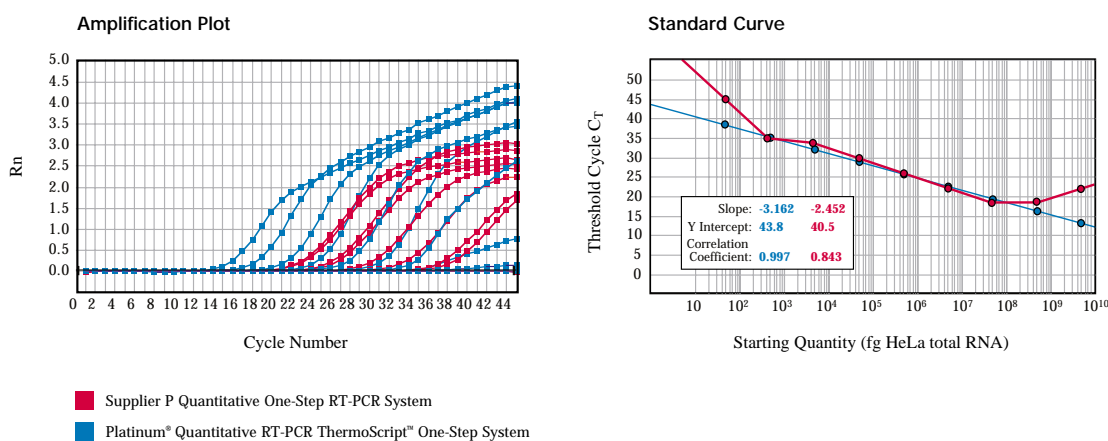

qRT-PCR of a 145 bp β -actin fragment was performed with the Platinum® Quantitative RT-PCR ThermoScript™ One-Step System (blue plot) or Supplier P Quantitative One-Step RT-PCR System (red plot) according to manufacturers' recommendations. Starting RNA varied from 5 μ g to 50 fg in 10-fold serial dilutions. Normalized relative fluorescence (Δ Rn) was collected using TAMRA as the passive reference. **Panel A:** Linear scale amplification plots. **Panel B:** Standard curve plots.

More options for detection and analysis

Choose from TaqMan® probes, Molecular Beacons, or whatever fluorescent detection system you desire. Complete your analysis on any number of instruments, such as the ABI PRISM® 7700/7000/7900, Bio-Rad's iCycler, Stratagene MX4000™, Cepheid

SmartCycler®, etc. With the Platinum® Quantitative RT-PCR ThermoScript™ One-Step System, you don't need to limit yourself to one particular detection method or one real-time PCR machine.

One step to accurate qRT-PCR

The Platinum® Quantitative RT-PCR ThermoScript™ One-Step System provides the leading RT and PCR technologies in real-time quantification. Order today, and you'll get accurate quantitative RT-PCR, one-step convenience, and the flexibility to use the fluorescent detection method and instrument of your choice.

Ordering information

Product	Quantity	Cat. no.
Platinum® Quantitative RT-PCR ThermoScript™ One-Step System	100 rxns	11731-015
	500 rxns	11731-023
Rox Reference Dye*	500 rxns	12223-012

* There is no additional charge for Rox reference dye when ordered in conjunction with the Platinum® Quantitative RT-PCR ThermoScript™ One-Step System.

ABI Prism® is a registered trademarks of Applied BioSystems. TaqMan® is a registered trademark of Roche Molecular Systems. iCycler™ is a trademark of Bio-Rad Laboratories, Inc. Mx4000™ is a trademark of Stratagene. SmartCycler® is a registered trademark of Cepheid. Platinum® Quantitative RT-PCR ThermoScript™ One-Step System is covered by Limited Label License Nos. 2, 13, 14, and 24. Please refer to the Invitrogen web site or catalog for the corresponding Limited Label License statements.



Printed in the U.S.A. ©2002 Invitrogen Corporation. Reproduction forbidden without permission.

Corporate headquarters:

1600 Faraday Avenue • Carlsbad, CA 92008 USA • Tel: 760 603 7200 • Fax: 760 602 6500 • Toll Free Tel: 800 955 6288 • E-mail: tech_service@invitrogen.com • www.invitrogen.com

European headquarters:

Invitrogen Ltd, 3 Fountain Drive • Inchinnan Business Park • Paisley PA4 9RF, UK • Tel: +44 (0) 41 814 6100 • Fax: +44 (0) 141 814 6260 • E-mail: eurotech@invitrogen.com