INSTRUCTIONS



Pierce[™] Chromatography Cartridges Protein A

89924 89925

1969.2

Number	Description
89924	Pierce Chromatography Cartridge Protein A, 2 × 1mL
89925	Pierce Chromatography Cartridge Protein A, 1 × 5mL
	Binding Capacity: \geq 35mg human IgG/ml of resin bed
	Note: The Protein A is immobilized on crosslinked 6% beaded agarose supplied in 0.05% sodium azide in water. Each product contains an accessory pack (1 female Luer-Lok TM Adapter, 1 connector fitting, 1 column plug and 1 or 2 bottom caps).
	Storage: Upon receipt store at 4-8°C. Product is shipped at ambient temperature. Do not freeze.

Introduction

The Thermo ScientificTM PierceTM Chromatography Cartridges Protein A are convenient, ready-to-use pre-packed devices for the isolation and purification of polyclonal and monoclonal antibodies. Protein A binds specifically and with high affinity to the Fc region of immunoglobulins, particularly IgG. These cartridges are effective for routine affinity purification of polyclonal and monoclonal IgG from serum, ascites and cell culture supernatant.

Pierce Cartridges are compatible with the major automated liquid-chromatography systems or for manual syringe processing (see Table 1 for general properties of the cartridges). The cartridges attach directly to ÄKTATM or FPLC Systems without additional connectors. An accessory pack, included with each product, readily adapts cartridges for use with Luer-Lok Syringe Fittings or 1/16" tubing. The Protein A cartridges provide fast, easy and reproducible chromatographic separations and can be regenerated for multiple uses.

Support	Crosslinked 6% beaded agarose
Ligand	Immobilized Protein A, MW 46,700
Binding Capacity	\geq 35mg human IgG/mL of resin bed
Cartridge Dimensions	0.7×2.7 cm (1mL column); 1.3×3.8 cm (5mL column)
Particle Size	45-165µm
Void Volume	0.32mL (1mL column); 1.5mL (5mL column)

Table 1. Properties of the Thermo Scientific Pierce Protein A Cartridges.

Recommended Flow Rate	1mL/min (1mL column); 1-5mL/min (5mL column)	
Maximum Recommended Flow Rate	4mL/min (1mL column); 5mL/min (5mL column)	
pH Limits	3-10	
Maximum Operating Pressure	0.3MPa, 43.5psi or 3bar	
Cartridge Material	Polypropylene	
Frit	Polyethylene, 10μm	
Storage Solution	0.02-0.05% Sodium azide	
Accessory Pack	Luer-Lok Adapter to 10-32 male Finger-tight 10-32 connector fitting for 1/16" OD tubing Plug for 10-32 coned port Cap 1/16 male	



Important Product Information

- Use high-purity buffers and water. For best results, degas or filter buffers through a 0.45µm filter.
- Serum samples, ascites fluid, plasma or tissue culture supernatant may be used with this product. The interaction between antibodies and Protein A is not equivalent for all species or for all subgroups of IgG. Refer to Tech Tip #34: Binding characteristics of Protein A, Protein G, Protein A/G and Protein L for further information on species and subgroup specificities of Protein A.
- For optimal recovery, use a sample size such that the expected IgG load on the column is less than 80% of the maximum binding capacity. The total IgG content of serum is approximately 10-15mg/mL. The concentration of antibody in tissue culture supernatant varies considerably among hybridoma clones. Be aware that antibodies from fetal bovine serum (FBS) culture media supplement will be purified along with the antibody of interest.
- Do not freeze the resin as this will cause irreversible damage to the bead structure.
- Pierce Cartridges may be used singly or connected in series (2-3 columns) to increase capacity. Back pressure will be greater when columns are used in a series than when used as single columns.
- To monitor protein as it emerges from the column, measure the UV absorbance at 280nm.

Additional Materials Required

- Suitable liquid chromatography system (LC procedure only) with 1/16" tubing or syringes
- Binding Buffer: 100mM sodium phosphate, 150mM sodium chloride, pH 7.2. Alternatively, prepare a Thermo Scientific[™] BupH[™] Phosphate Buffered Saline Pack (Product No. 28372) with high-purity water according to instructions. Sodium azide (0.02%) may be added to the buffer solution as a preservative. For best results, process the buffer through a 0.45µm filter before use in LC applications.
- Elution Buffer: 0.1M glycine, pH 2-3, or Pierce IgG Elution Buffer (Product No. 21004). Sodium azide (0.02%) may be added to the buffer solution as a preservative. For best results, process the buffer through a 0.45µm filter before use in LC applications.
- Neutralization Buffer: High ionic strength alkaline buffer such as 1M phosphate or 1M Tris at pH 7.5-9
- Pierce Chromatography Cartridge Desalting (Product No. 89934 or 89935), Thermo Scientific[™] Zeba[™] Spin Desalting Columns (Product No. 89891, 89892, 89893 or 89894) or Thermo Scientific[™] Slide-A-Lyzer[™] Dialysis Cassettes for optional buffer exchange.
- Additional connectors and fittings are required to attach to the Bio-Rad BioLogic[™] System.

Procedure for Antibody Purification using a Liquid Chromatography System

- 1. Equilibrate the cartridge and all buffers to room temperature. Ensure all solutions are degassed.
- 2. Prepare the LC system by filling tubing with buffer. Remove top plug from cartridge and carefully snap off the end-tab (do not twist). To avoid introducing air into the system, let a few drops of buffer flow from tubing into column top then connect cartridge top to the tubing; allow a few drops to emerge from the column before connecting to the LC inlet port.
- 3. Equilibrate the resin with 5-10 column volumes of the Binding Buffer at a flow rate of 1mL/minute for the 1mL column or 5mL/minute for the 5mL column.
- 4. For maximum binding, adjust the sample to the ionic strength and pH of the Binding Buffer by diluting it at least 1:1 before applying to the column. Alternatively, buffer exchange the sample against the Binding Buffer. If the sample contains insoluble matter, centrifuge or filter (0.45µm filter) it before use.

Note: Plasma may become hazy upon dilution with the Binding Buffer because of lipoprotein precipitation. Centrifuge the diluted sample at $10,000 \times g$ for 20 minutes and apply the supernatant to the equilibrated column.

5. Apply the diluted sample to the column. For maximum binding, apply at a flow rate of 1mL/minute for the 1mL column and 1-2mL/minute for the 5mL column. Apply any volume of sample provided the total amount of antibody is < 80% of Protein A resin capacity.

Note: If the sample exceeds the binding capacity of the resin (or contains an antibody type that does not bind), non-bound antibody can be recovered in the flow-through and analyzed by antibody-specific assays.



- 6. Wash the resin with 5-10 column volumes of Binding Buffer or until the absorbance approaches baseline.
- Elute with approximately 2-5 column volumes of Elution Buffer and collect 0.5-1mL fractions. Immediately adjust eluted fractions to physiological pH by adding 100µL of the Neutralization Buffer per 1mL of eluate. Alternatively, add the Neutralization Buffer to the collection tubes before eluting.
- 8. Analyze the purified fractions directly by SDS-PAGE, or dialyze or desalt the sample into a buffer that is compatible with the specific downstream application.
- 9. The resin may be regenerated and reused multiple times without significant loss of binding capacity. To prevent crosscontamination, reuse columns with identical antibodies. Regenerate the resin with 3-5 column volumes of Elution Buffer and re-equilibrate with Binding Buffer until the pH returns to the buffer value.
- 10. For storage, wash the resin with five column volumes of water containing 0.02-0.05% sodium azide. Attach supplied bottom cap followed by the top plug. Store the column at 4°C.

Procedure for Antibody Purification using a Syringe

Note: The void volumes are 0.320mL for the 1mL columns and 1.5mL for the 5mL columns.

- 1. Equilibrate the cartridge and all solutions to room temperature. Ensure all solutions are degassed.
- 2. Fill a syringe with 5-10 column volumes of buffer.
- 3. Attach the syringe to the Luer-Lok Adapter included in the accessory pack. Remove top plug from cartridge and carefully snap off the end-tab. To avoid introducing air into the system, allow a few drops to emerge from the Luer-Lok Adapter and then connect to the cartridge top. Securely tighten the connection.
- 4. Equilibrate the Protein A resin with 5-10 column volumes of buffer at a flow rate of ~1mL/minute for the 1mL column or ~5 ml/minute for 5 ml column. Remove syringe from the Luer-Lok Adapter.
- 5. For maximum binding, adjust the sample to the ionic strength and pH of the Binding Buffer by diluting it at least 1:1 before applying to the column. Alternatively, buffer-exchange the sample against the Binding Buffer. If the sample contains insoluble matter, centrifuge or filter (0.45µm filter) it before use.

Note: Plasma may become hazy upon dilution with the Binding Buffer because of lipoprotein precipitation. Centrifuge the diluted sample at $10,000 \times g$ for 20 minutes and apply the supernatant to the equilibrated column.

6. Fill an appropriately sized syringe with the diluted sample and connect it to the Luer-Lok Adapter. Depress the syringe plunger to pass the sample through the column. For maximum binding, apply at a flow rate of 1mL/minute for the 1mL column and 1-2mL/minute for the 5mL column. Any volume may be applied provided the total amount of antibody is less than 80% of the column capacity.

Note: If the sample exceeds the binding capacity of the resin (or contains an antibody type that does not bind), non-bound antibody can be recovered in the flow-through and analyzed by antibody-specific assays.

- 7. Change the syringe and wash the resin with 5-10 column volumes of Binding Buffer.
- Change syringe and elute with approximately 2-5 column volumes of Elution Buffer and collect 0.5-1mL fractions. Immediately adjust eluted fractions to physiological pH by adding 100µL of the Neutralization Buffer per 1mL of eluate. Alternatively, add the Neutralization Buffer to the collection tube before eluting.
- 9. Analyze the purified fractions directly by SDS-PAGE, or dialyze or desalt the sample into a buffer that is compatible with the specific downstream application.
- 10. The resin may be regenerated and reused multiple times without significant loss of binding capacity. To prevent crosscontamination, reuse columns with identical antibodies. Regenerate the resin in 3-5 column volumes of Elution Buffer and re-equilibrate in Binding Buffer until the pH returns to the buffer value.
- 11. For storage, wash resin with five column volumes of water containing 0.02-0.05% sodium azide. Attach supplied bottom cap followed by the top plug. Store the cartridge at 4°C.



Troubleshooting

Problem	Possible Cause	Solution	
Sample in flow-through fractions	Exceeded resin's binding capacity	Make sure the total amount of protein loaded is < 80% of the resin's binding capacity	
	Sample or resin not at proper pH or correct ionic strength for optimum binding	Dilute sample 1:1 with Binding Buffer or perform a buffer exchange so the ionic strength and pH is optimal for binding and make sure to equilibrate the resin with the Binding Buffer	
	Sample being applied too fast for proper binding	Slow the flow rate to 0.1-1mL/minute during sample loading	
No antibody detected in any elution fraction	Sample devoid of antibody species or subclass that binds to Protein A	Refer to the Pierce catalog or website to select the best antibody-binding protein for a specific species	
Considerable antibody purified,	Antibody of interest is at low	Use serum-free medium for cell supernatant samples	
but no specific antibody of interest detected	concentration	Use the specific antigen coupled to an affinity support such as Thermo Scientific TM AminoLink TM Plus Immobilization Kit (Product No. 44894) to purify the antibody	
Antibody of interest purified, but it is degraded (determined	Antibody is sensitive to low-pH Elution Buffer	Try Gentle Ag/Ab Elution Buffer (see Related Thermo Scientific Products)	
by lack of function in downstream assay)	Downstream application is sensitive to neutralized Elution Buffer	Desalt or dialyze eluted sample into suitable buffer	

Related Thermo Scientific Products

89971	Accessory Pack (1 female Luer-Lok Adapter, 1 connector fitting, 1 column plug and 1 bottom cap)
28372	BupH TM Phosphate Buffered Saline Pack, 40 packs
21004	IgG Elution Buffer, 1L
21027	Gentle Ag/Ab Elution Buffer, 500mL
89935	Pierce Chromatography Cartridge Desalting, 5 × 5mL
89806	Protein Stabilizing Cocktail (4X), 10mL

General References

Ey, P.L., *et al.* (1978). Isolation of pure IgG₁, IgG_{2a}, and IgG_{2b} immunoglobulins from mouse serum using protein A-Sepharose. *Immunochem* **15**:429-36. Goding, J.W. (1978). Use of Staphylococcal protein A as an immunological reagent. *J Immunol Methods* **20**:241-53

Hjelm, H., *et al.* (1975). Immunologically active and structurally similar fragments of protein A from *Staphylococcus aureus*. *Eur J Biochem* 57:395-403.
Kronvall, G. (1970). Phylogenetic insight into evolution of mammalian Fc fragment of gamma G globulin using staphylococcal protein A. *J Immunol* 104:140-7.

Kronvall, G., et al. (1970). Protein A reactivity with mouse immunoglobulins. J Immunol 105:1116-23.

Kronvall, G. and Williams, R.C. (1969). Differences in anti-protein A activity among IgG subgroups. J Immunol 103(4):828-33.

Lindmark, R., *et al.* (1983). Binding of immunoglobulins to protein A and immunoglobulin levels in mammalian sera. *J Immunol Methods* **62:**1-13. Surolia, A., *et al.* (1982). Protein A: Nature's universal anti-body. *TIBS* **7:**74-6.

Products are warranted to operate or perform substantially in conformance with published Product specifications in effect at the time of sale, as set forth in the Product documentation, specifications and/or accompanying package inserts ("Documentation"). No claim of suitability for use in applications regulated by FDA is made. The warranty provided herein is valid only when used by properly trained individuals. Unless otherwise stated in the Documentation, this warranty is limited to one year from date of shipment when the Product is subjected to normal, proper and intended usage. This warranty does not extend to anyone other than Buyer. Any model or sample furnished to Buyer is merely illustrative of the general type and quality of goods and does not represent that any Product will conform to such model or sample.

NO OTHER WARRANTIES, EXPRESS OR IMPLIED, ARE GRANTED, INCLUDING WITHOUT LIMITATION, IMPLIED WARRANTIES OF MERCHANTABILITY, FITNESS FOR ANY PARTICULAR PURPOSE, OR NON INFRINGEMENT. BUYER'S EXCLUSIVE REMEDY FOR NON-CONFORMING PRODUCTS DURING THE WARRANTY PERIOD IS LIMITED TO REPAIR, REPLACEMENT OF OR REFUND FOR THE NON-CONFORMING PRODUCT(S) AT SELLER'S SOLE OPTION. THERE IS NO OBLIGATION TO REPAIR, REPLACE OR REFUND FOR PRODUCTS AS THE RESULT OF (I) ACCIDENT, DISASTER OR EVENT OF FORCE MAJEURE, (II) MISUSE, FAULT OR NEGLIGENCE OF OR BY BUYER, (III) USE OF THE PRODUCTS IN A MANNER FOR WHICH THEY WERE NOT DESIGNED, OR (IV) IMPROPER STORAGE AND HANDLING OF THE PRODUCTS.

Unless otherwise expressly stated on the Product or in the documentation accompanying the Product, the Product is intended for research only and is not to be used for any other purpose, including without limitation, unauthorized commercial uses, in vitro diagnostic uses, ex vivo or in vivo therapeutic uses, or any type of consumption by or application to humans or animals.

Pierce Biotechnology	PO Box 117	(815) 968-0747	www.thermoscientific.com/pierce
3747 N. Meridian Road	Rockford, IL 61105 USA	(815) 968-7316 fax	



Current product instructions are available at <u>www.thermoscientific.com/pierce</u>. For a faxed copy, call 800-874-3723 or contact your local distributor. © 2013 Thermo Fisher Scientific Inc. All rights reserved. ÄKTA is a trademark of GE Healthcare. BioLogic is a trademark of Bio-Rad Laboratories, Inc. All (other) trademarks are the property of Thermo Fisher Scientific Inc. and its subsidiaries. Printed in the USA.