



00-2019 500 ml
TMB Solution
(Ready-to-use)
Lot No.

TMB SOLUTION (READY-TO-USE) for IMMUNOBLOT

SUBSTRATE FOR PEROXIDASE

FORM:

500 ml. TMB solution. (Ready-to-use)

Tetramethyl Benzidine (TMB) is an insoluble substrate for peroxidase which is especially good for staining blots.

BLOTTING PROCEDURE:

Add 10-20 ml of TMB solution or enough to cover blots completely.
Shake for 10-30 minutes or until background begins to develop.

STORAGE:

Store at 2-8°C and protect from light.

REMARKS:

For research use only. Not for human or drug use. No preservative added.

www.invitrogen.com

Invitrogen Corporation • 542 Flynn Rd • Camarillo • CA 93012 • Tel: 800.955.6288

E-mail: techsupport@invitrogen.com

PI002019

(Rev 10/08) DCC-08-1089

Important Licensing Information - These products may be covered by one or more Limited Use Label Licenses (see the Invitrogen Catalog or our website, www.invitrogen.com). By use of these products you accept the terms and conditions of all applicable Limited Use Label Licenses. Unless otherwise indicated, these products are for research use only and are not intended for human or animal diagnostic, therapeutic or commercial use.