

Corning® Rocker Cell Culture Bag


CORNING

Corning® single-use rocker cell culture bags feature a multi-layered film with an inert PE (Polyethylene) fluid contact layer. The bag is designed for use with commercial rocking-motion bioreactor chambers. Corning offers this single-use cell culture bag in 2L to 50L configurations. The bag dimensions, port locations, and stability bars located on each side, allow commercially available bioreactors to accommodate the cell culture bag.

The rocking-motion of the bioreactor platform creates effective, non-invasive mixing of the cell culture/media with efficient gas transfer. Rocking-motion prevents cells from settling at the bottom of the rocker cell culture bag while generating large amounts of recombinant protein in both mammalian and insect cell lines.

Features

- ▶ Provides non-invasive mixing of the cell culture/media with efficient gas transfer
- ▶ Eliminates the risk of cross-contamination
- ▶ Used for basic research to large-scale cell culture applications
- ▶ Available in the following sizes: 2L, 10L, 20L, 22L, and 50L

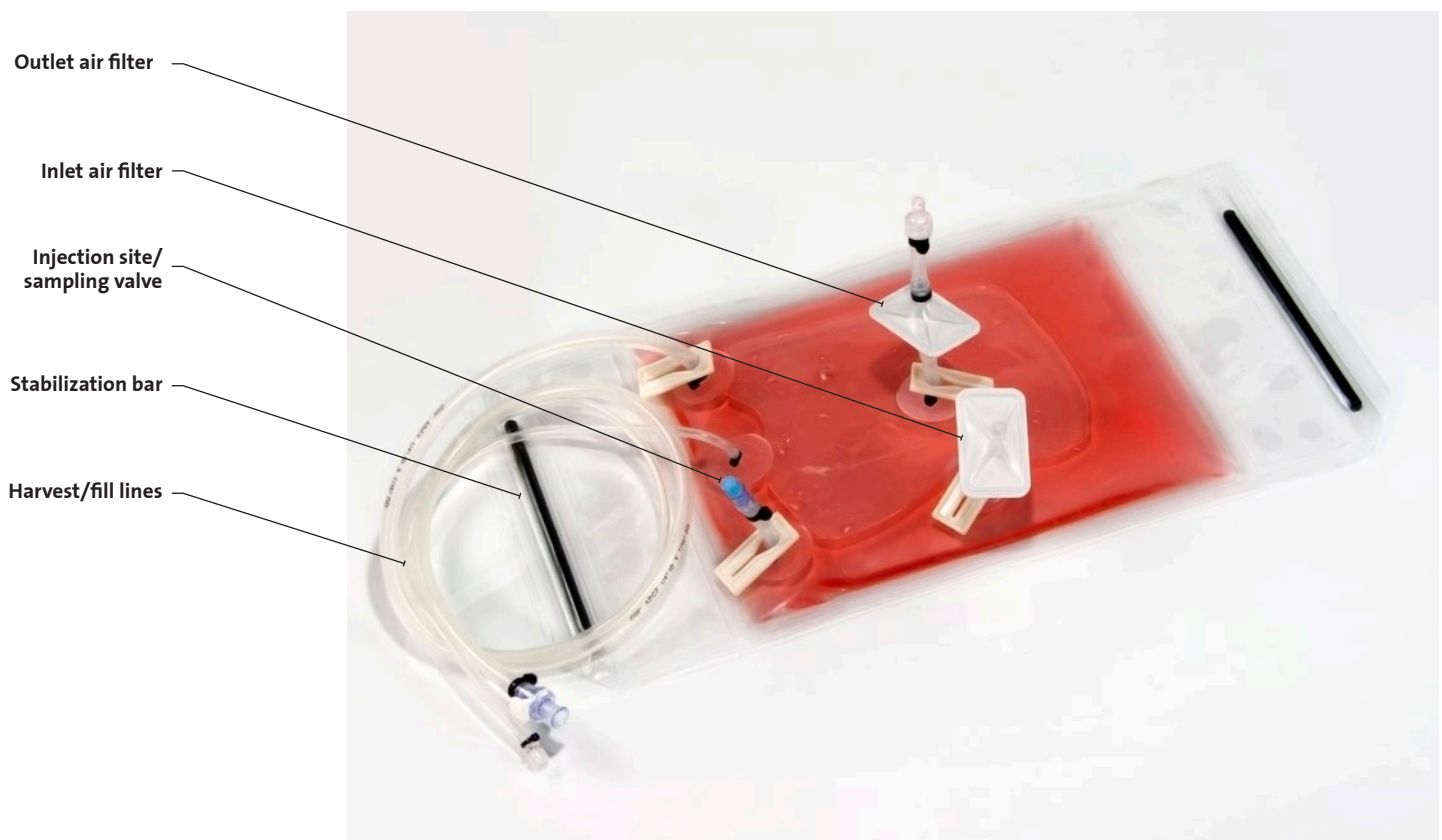


Corning® Rocker Bags

2L Rocker Bag

Cat. No.	Description	Dimensions	Film Type	Unit Size	Qty
91-200-80	Rocker Bag	22.25" x 10.75"	EVOH/ULDPE	2L	1

Port Configuration

	Check valve with 1/4" barb
P1	Pall 8004022 filter 2" length 3/16" I.D. x 3/8" O.D. PCMGS tube
P2	Pall 8004022 filter 2" length 3/16" I.D. x 3/8" O.D. PCMGS tube
P3	Female MPC connector with 1/4" barb Male MPC plug 40" length 1/4" I.D. x 7/16" O.D. chemically resistant, heat sealable flexible tube
P4	Female Luer with 1/4" barb Male Luer plug 40" length 1/8" I.D. x 1/4" O.D. chemically resistant, heat sealable flexible tube
P5	Female Luer with 1/4" barb Swabbable valve Male Luer plug without stem

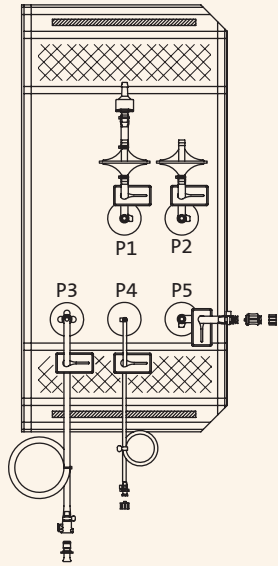


10L Rocker Bag

Cat. No.	Description	Dimensions	Film Type	Unit Size	Qty
91-200-79	Rocker Bag	22" x 13"	EVOH/ULDPE	10L	1

Port Configuration

	Check valve with 1/4" barb
P1	Pall 8004022 filter 2" length 3/16" I.D. x 3/8" O.D. PCMGS tube
P2	Pall 8004022 filter 2" length 3/16" I.D. x 3/8" O.D. PCMGS tube
P3	Female MPC connector with 1/4" barb Male MPC plug 40" length 1/4" I.D. x 7/16" O.D. chemically resistant, heat sealable flexible tube
P4	Female Luer with 1/4" barb Male Luer plug 40" length 1/8" I.D. x 1/4" O.D. chemically resistant, heat sealable flexible tube
P5	Female Luer with 1/4" barb Swabbable valve Male Luer plug without stem

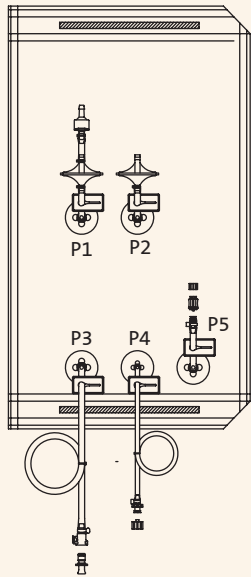



20L Rocker Bag

Cat. No.	Description	Dimensions	Film Type	Unit Size	Qty
91-200-78	Rocker Bag	26" x 22"	EVOH/ULDPE	20L	1

Port Configuration

	Check valve with 1/4" barb
P1	Pall 8004022 filter 2" length 3/16" I.D. x 3/8" O.D. PCMGS tube
P2	Pall 8004022 filter 2" length 3/16" I.D. x 3/8" O.D. PCMGS tube
P3	Female MPC connector with 1/4" barb Male MPC plug 40" length 1/4" I.D. x 7/16" O.D. chemically resistant, heat sealable flexible tube
P4	Female Luer with 1/4" barb Male Luer plug 40" length 1/8" I.D. x 1/4" O.D. chemically resistant, heat sealable flexible tube
P5	Female Luer with 1/4" barb Swabbable valve Male Luer plug without stem

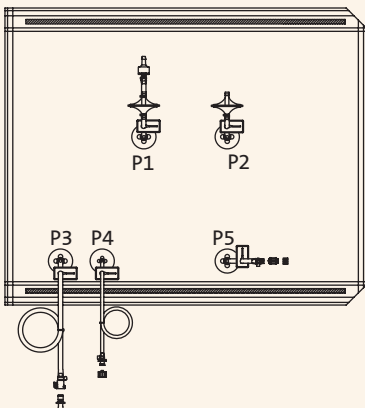

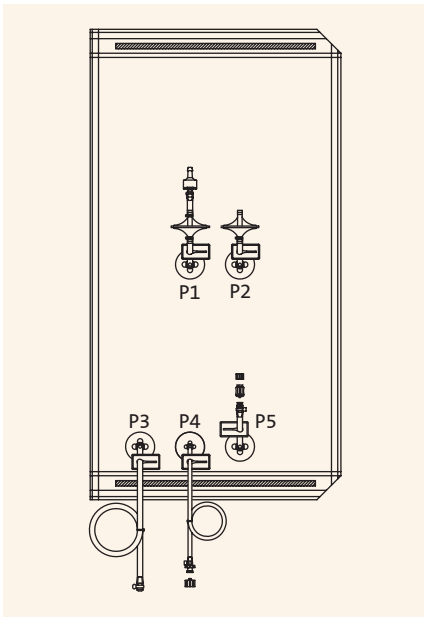


22L Rocker Bag

Cat. No.	Description	Dimensions	Film Type	Unit Size	Qty
91-200-92	Rocker Bag	28.25" x 15.063"	EVOH/ULDPE	22L	1

Port Configuration

	Check valve with 1/4" barb
P1	Pall 8004022 filter 2" length 3/16" I.D. x 3/8" O.D. PCMGs tube
P2	Pall 8004022 filter 2" length 3/16" I.D. x 3/8" O.D. PCMGs tube
P3	Plug with 1/4" barb Male MPC plug 48" length 1/4" I.D. x 7/16" O.D. chemically resistant, heat sealable flexible tube
P4	Female Luer with 1/4" barb Male Luer plug 48" length 1/8" I.D. x 1/4" O.D. chemically resistant, heat sealable flexible tube
P5	Female Luer with 1/4" barb Swabbable valve Male Luer plug without stem

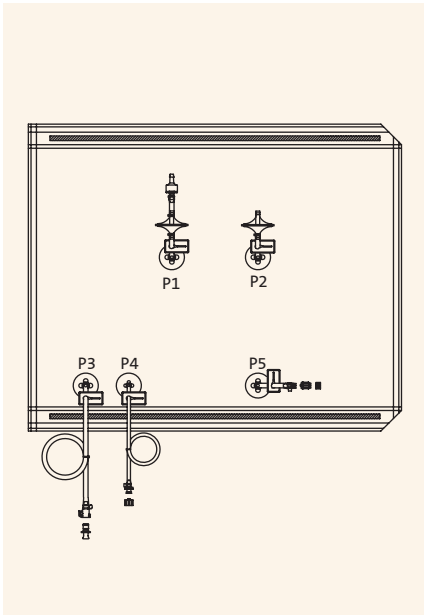

50L Rocker Bag


Cat. No.	Description	Dimensions	Film Type	Unit Size	Qty
91-200-81	Rocker Bag	29.25" x 28.25"	EVOH/ULDPE	50L	1

Port Configuration

	Check valve with 1/4" barb
P1	Pall 8004022 filter 2" length 3/16" I.D. x 3/8" O.D. PCMGs tube
P2	Pall 8004022 filter 2" length 3/16" I.D. x 3/8" O.D. PCMGs tube
P3	Female MPC connector with 1/4" barb Male MPC plug 40" length 1/4" I.D. x 7/16" O.D. chemically resistant, heat sealable flexible tube
P4	Female Luer with 1/4" barb Male Luer plug 40" length 1/8" I.D. x 1/4" O.D. chemically resistant, heat sealable flexible tube
P5	Female Luer with 1/4" barb Swabbable valve Male Luer plug without stem

For more specific information on claims, visit the Certificates page at www.corning.com/lifesciences.

Warranty/Disclaimer: Unless otherwise specified, all products are for research use only. Not intended for use in diagnostic or therapeutic procedures. Not for use in humans. Corning single-use containers are not intended for use in parenteral applications. Corning Life Sciences makes no claims regarding the performance of these products for clinical or diagnostic applications.

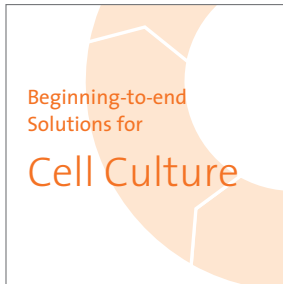

www.corning.com/lifesciences/solutions

At Corning, cells are in our culture. In our continuous efforts to improve efficiencies and develop new tools and technologies for life science researchers, we have scientists working in Corning R&D labs across the globe, doing what you do every day. From seeding starter cultures to expanding cells for assays, our technical experts understand your challenges and your increased need for more reliable cells and cellular material.

It is this expertise, plus a 160-year history of Corning innovation and manufacturing excellence, that puts us in a unique position to offer a beginning-to-end portfolio of high-quality, reliable cell culture consumables.

For additional product or technical information, visit www.corning.com/lifesciences or call 800.492.1110. Outside the United States, call +1.978.442.2200 or contact your local Corning sales office.

Corning Incorporated
Life Sciences

836 North St.
Building 300, Suite 3401
Tewksbury, MA 01876
t 800.492.1110
t 978.442.2200
f 978.442.2476

www.corning.com/lifesciences

**Worldwide
Support Offices**

ASIA/PACIFIC

Australia/New Zealand
t 61 427286832

China
t 86 21 3338 4338
f 86 21 3338 4300

India
t 91 124 4604000
f 91 124 4604099

Japan
t 81 3-3586 1996
f 81 3-3586 1291

Korea
t 82 2-796-9500
f 82 2-796-9300

Singapore
t 65 6572-9740
f 65 6861-2913

Taiwan
t 886 2-2716-0338
f 886 2-2516-7500

EUROPE
France
t 0800 916 882
f 0800 918 636

Germany
t 0800 101 1153
f 0800 101 2427

The Netherlands
t 020 655 79 28
f 020 659 76 73

United Kingdom
t 0800 376 8660
f 0800 279 1117

**All Other European
Countries**
t 31 (0) 20 659 60 51
f 31 (0) 20 659 76 73

LATIN AMERICA
grupoLA@corning.com
Brasil
t (55-11) 3089-7400
Mexico
t (52-81) 8158-8400

CORNING | **FALCON** | **AXYGEN** | **GOSELIN** | **PYREX**

For a listing of trademarks, visit www.corning.com/clstrademarks.
All other trademarks are the property of their respective owners.