

# StemPro®-34 SFM Complete Medium

## Description

StemPro®-34 SFM is a serum-free medium specifically formulated to support the growth of human hematopoietic progenitor cells (for example, CD34+) in culture. StemPro®-34 SFM can be used to study hematopoietic cells from bone marrow, peripheral blood, or neonatal cord blood. StemPro®-34 SFM consists of a basal liquid medium and frozen 40X nutrient supplement (StemPro®-34 Nutrient Supplement) which is added at the time of use. StemPro®-34 SFM is manufactured without cytokines and hematopoietic growth factors, giving the investigator the freedom to use any factor or combination of factors required in ones' studies. StemPro®-34 SFM contains components of human, recombinant or synthetic origin and is performance-tested on normal human bone marrow.

Product	Catalog no.	Amount	Storage	Shelf life*
StemPro®-34 SFM Kit	10639-011	1 kit		
Contains:				
StemPro®-34 SFM (1X), liquid	10640-019	500 mL	2°C to 8°C; Protect from light	18 months
StemPro®-34 Nutrient Supplement (40X)**	10641-025*	13 mL	–20°C to –5°C; Protect from light	30 months

<sup>\*</sup> Shelf life duration is determined from Date of Manufacture.

#### Product use

For Research Use Only. Not for use in diagnostic procedures.

## Important information

- Do not use this product if the product packaging appears compromised, if StemPro®-34 Nutrient Supplement is received thawed, or if StemPro®-34 SFM appears cloudy or a visible precipitate is observed.
- Complete StemPro®-34 SFM has a shelf life of 30 days when stored in the dark at 2° to 8°C.
- Thawed StemPro®-34 SFM Nutrient Supplement is stable for 14 days when stored in the dark at 2° to 8°C.

#### Safety information

Read the Safety Data Sheets (SDSs) and follow the handling instructions. Wear appropriate protective eyewear, clothing, and gloves.

# StemPro®-34 Nutrient Supplement (40X) only

Caution: Human origin materials are non-reactive (donor level) for anti-HIV 1 & 2, anti-HCV, and  $HB_sAg$ . Handle in accordance with established bio-safety practices.

## Prepare medium

Supplement StemPro®-34 SFM Basal Medium with StemPro®-34 Nutrient Supplement and L-glutamine:

- 1. Thaw StemPro®-34 Nutrient Supplement overnight at 4°C. Thoroughly mix by gently inverting the vial several times.
- 2. Aseptically add the entire contents (13 mL) of StemPro®-34 Nutrient Supplement directly to a 500 mL bottle of StemPro®-34 SFM Basal Medium before use. Swirl to mix.

**Note:** StemPro®-34 Nutrient Supplement may be thawed, aliquoted and refrozen once. Avoid repeated freeze/thaw cycles and use immediately once thawed.

- 3. Aseptically add 5 mL 200 mM L-glutamine (2 mM final) to 500 mL of StemPro $^{\circ}$ -34 SFM Basal Medium before use.
- Supplement Complete StemPro®-34 SFM with cytokines or growth factors for extended support of progenitor cells.
  Note: For hematopoietic stem cell (HSC) culture, we recommend adding the cytokines SCF (100 ng/mL), IL-3 (50 ng/mL), and GM-CSF (25 ng/mL) to Complete

StemPro®-34 SFM immediately before use.

#### Culture conditions

Media: Complete StemPro®-34 SFM

Cells: Human hematopoietic progenitor cells

Temperature range: 36°C to 38°C

**Incubator atmosphere:** Humidified atmosphere of 4–6% CO<sub>2</sub> in air. Ensure proper gas exchange and minimize exposure of

cultures to light.

Culture conditions: Stationary suspension Recommended culture vessels: T-Flasks

Our research data has shown that normal bone marrow cells can be held at 2°C to 8°C, in complete StemPro $^{\odot}$ -34 SFM, without added growth factors, for 5–7 days without significant loss of cell viability as determined by trypan blue dye exclusion.

#### Use

For procedures detailing StemPro®-34 SFM use, refer to publication number MAN0005808 included with StemPro®-34 Medium and CD34+ Cell Kit (Cat. no. A14509) available at www.lifetechnologies.com/manuals.

#### Related products

Product	Catalog no.	
L-Glutamine, 200 mM (100X), liquid	25030	
GlutaMAX <sup>™</sup> -I, 200 mM (100X), liquid	35050	
SCF (c-kit Ligand) Recombinant Human	PHC2111	
IL-3 Recombinant Human	PHC0033	
GM-CSF Recombinant Human	PHC2013	
StemPro®-34 Medium and CD34⁺ Cells Kit	A14509	
Dynal® CD34 Progenitor Cell Selection System	113-01D	
Trypan Blue Solution, 0.4%	15250	
Countess® Automated Cell Counter	C10227	

Publication Number MAN0007286 Rev. 1.00

<sup>\*\*</sup>StemPro 34® SFM Nutrient Supplement is not sold separately.

## Explanation of symbols and warnings

The symbols present on the product label are explained below:

<u></u>	i	*	$\downarrow$			
Caution, consult accompanying documents	Consult instructions for use	Keep away from light	Temperature Limitation			
MM-TOTY	LOT	REF	STERILE A			
Use By:	Batch code	Catalog number	Sterilized using aseptic processing techniques			

## Limited product warranty

Life Technologies Corporation and/or its affiliate(s) warrant their products as set forth in the Life Technologies' General Terms and Conditions of Sale found on Life Technologies' website at www.lifetechnologies.com/termsandconditions. If you have any questions, please contact Life Technologies at www.lifetechnologies.com/support.

## Important licensing information

These products may be covered by one or more Limited Use Label Licenses. By use of these products, you accept the terms and conditions of all applicable Limited Use Label Licenses.

For additional technical information such as Safety Data Sheets (SDS), Certificates of Analysis, visit www.lifetechnologies.com/support For further assistance, email **techsupport@lifetech.com** 

All trademarks are the property of Thermo Fisher Scientific and its subsidiaries unless otherwise specified. ©2014 Thermo Fisher Scientific Inc. All rights reserved.

DISCLAIMER - LIFE TECHNOLOGIES CORPORATION AND/OR ITS AFFILIATE(S) DISCLAIM ALL WARRANTIES WITH RESPECT TO THIS DOCUMENT, EXPRESSED OR IMPLIED, INCLUDING BUT NOT LIMITED TO THOSE OF MERCHANTABILITY, FITNESS FOR A PARTICULAR PURPOSE, OR NON-INFRINGEMENT. TO THE EXTENT ALLOWED BY LAW, IN NO EVENT SHALL LIFE TECHNOLOGIES AND/OR ITS AFFILIATE(S) BE LIABLE, WHETHER IN CONTRACT, TORT, WARRANTY, OR UNDER ANY STATUTE OR ON ANY OTHER BASIS FOR SPECIAL, INCIDENTAL INDIRECT, PUNITIVE, MULTIPLE OR CONSEQUENTIAL DAMAGES IN CONNECTION WITH OR ARISING FROM THIS DOCUMENT, INCLUDING BUT NOT LIMITED TO THE USE THEREOF.

