



Netwell™ Carrier and Handle Instructions

CORNING

To Load the Netwell Carrier

An easy method to load the Netwell inserts into the reusable carrier is to invert the plate, leaving the Netwell inserts in the plate cover, then place the carrier over the Netwells, and reinvert the assembly.

To Use the Handle

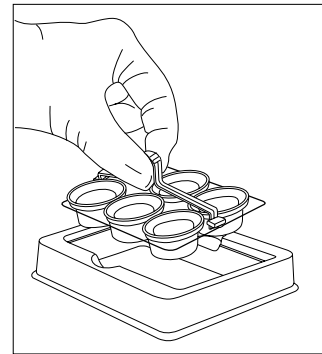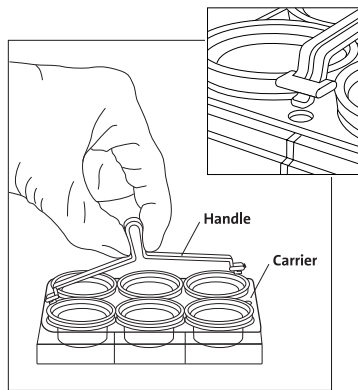

1. Remove the plate lid.
2. Grasp the handle along its axis between the thumb and forefinger and squeeze.
3. Insert both ends of the handle into the carrier (see insert drawing), and release the handle.
4. Change the grip on the handle 90° (this will prevent you from accidentally squeezing the handle and releasing the Netwell inserts).
5. Pick up the Netwell inserts and transport to the next workstation.

For more specific information on claims, visit www.corning.com/certificates.

Warranty/Disclaimer: Unless otherwise specified, all products are for research use only. Not intended for use in diagnostic or therapeutic procedures. Not for use in humans. Corning Life Sciences makes no claims regarding the performance of these products for clinical or diagnostic applications.

For additional product or technical information, visit www.corning.com/lifesciences or call 800.492.1110. Outside the United States, call +1.978.442.2200 or contact your local Corning sales office.

CORNING

Corning Incorporated
Life Sciences

836 North St.
Building 300, Suite 3401
Tewksbury, MA 01876
t 800.492.1110
t 978.442.2200
f 978.442.2476

www.corning.com/lifesciences

ASIA/PACIFIC
Australia/New Zealand
t 61 427286832

Chinese Mainland
t 86 21 3338 4338
f 86 21 3338 4300

India
t 91 124 4604000
f 91 124 4604099

Japan
t 81 3-3586 1996
f 81 3-3586 1291

Korea
t 82 2-796-9500
f 82 2-796-9300

Singapore
t 65 6572-9740
f 65 6735-2913

Taiwan
t 886 2-2716-0338
f 886 2-2516-7500

EUROPE
CSEurope@corning.com

France
t 0800 916 882
f 0800 918 636

Germany
t 0800 101 1153
f 0800 101 2427

The Netherlands
t 020 655 79 28
f 020 659 76 73

United Kingdom
t 0800 376 8660
f 0800 279 1117

All Other European Countries
t +31 (0) 206 59 60 51
f +31 (0) 206 59 76 73

LATIN AMERICA
grupoLA@corning.com

Brazil
t 55 (11) 3089-7400
Mexico
t (52-81) 8158-8400