

Pierce™ Cypridina-Firefly Luciferase Dual Assay Kit

16183 16184

2370.2

Number	Description
16183	<p>Pierce Cypridina-Firefly Luciferase Dual Assay Kit, sufficient reagents to perform 100 assays for <i>Cypridina</i>-firefly luciferase activity in cultured cell lysate</p> <p>Kit Contents:</p> <p>Cypridina-Firefly Dual Assay Buffer, 5mL, store at 4°C</p> <p>Vargulin (100X), 50µL, store at -20°C</p> <p>D-Luciferin, Lyophilized, 3mg, store at 4°C</p> <p>2X Cell Lysis Buffer, 6mL, store at room temperature</p>
16184	<p>Pierce Cypridina-Firefly Luciferase Dual Assay Kit, sufficient reagents to perform 1000 assays for <i>Cypridina</i>-firefly luciferase activity in cultured cell lysate</p> <p>Kit Contents:</p> <p>Cypridina-Firefly Dual Assay Buffer, 50mL, store at 4°C</p> <p>Vargulin (100X), 0.5mL, store at -20°C</p> <p>D-Luciferin, Lyophilized, 30mg, store at 4°C</p> <p>2X Cell Lysis Buffer, 60mL, store at room temperature</p> <p>Storage: Upon receipt store kit at -20°C or store individual components as indicated above. Kit is shipped on dry ice.</p>

Table of Contents

Introduction	2
Procedure Summary.....	3
Important Product Information	4
Additional Materials Required.....	4
Material Preparation	4
Assay Protocol.....	5
A. Cell Transfection	5
B. Cell Lysis.....	5
C. Luciferase Dual Assay.....	5
Troubleshooting.....	6
Related Thermo Scientific Products	6
General Reference	7

Introduction

The Thermo Scientific™ Pierce™ Cypridina-Firefly Luciferase Dual Assay Kit provides a highly sensitive system for detecting intracellular luciferase activity from promoter or pathway activation in mammalian cell culture experiments. The activity of *Cypridina* luciferase expressed from *Cypridina* luc reporter plasmid is normalized by the activity of red firefly luciferase from the red firefly luc control plasmid. Additionally, red firefly luciferase works as a second experimental reporter for monitoring secondary regulatory element activity.

The *Cypridina* luciferase bioluminescence signal has greater stability and brightness than firefly and native *Renilla* luciferase. The bioluminescent signal ($\lambda_{\text{max}} = 463\text{nm}$) produced by *Cypridina* luciferase results from the oxidation of Vargulin (Figure 1). This reaction does not require adenosine triphosphate (ATP) or other cofactors. The light output correlates with the amount of *Cypridina* protein expressed, which is used to determine promoter activity for *Cypridina* expression.

The second assay uses red firefly luciferase, which is a mutant form of the Italian firefly luciferase from *Luciola cruciata*. This luciferase produces a red-shifted emission spectrum ($\lambda_{\text{max}} = 613\text{nm}$) that results from the oxidation of D-luciferin in the presence of ATP (Figure 2). Red firefly and *Cypridina* luciferases are used in the dual assay because the light output is spectrally resolvable (Figure 3). *Cypridina* luciferase plasmid acts as an experimental reporter coupled with red firefly expression plasmid (CMV-Red Firefly) as a normalization control. This reporter and control combination enables simultaneous monitoring of experimental reporter and control luciferase activity in a single-read assay without the need for two-step addition of substrate reagents or quenching (Figure 3). Specificity analysis of the two luciferase emissions has less than 1% of *Cypridina* signal crossover to the 640nm LP channel (Figure 4).

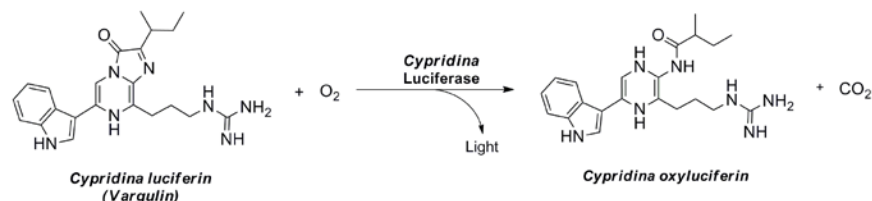


Figure 1. Chemical reaction of Vargulin and *Cypridina* luciferase. Light, with an emission maximum of 463nm, is produced from the oxidation of Vargulin by *Cypridina* luciferase.



Figure 2. Chemical reaction of luciferin and red firefly luciferase. Light, with an emission maximum of 613nm, is produced from the oxidation of D-luciferin by red firefly luciferase in an ATP-dependent reaction.

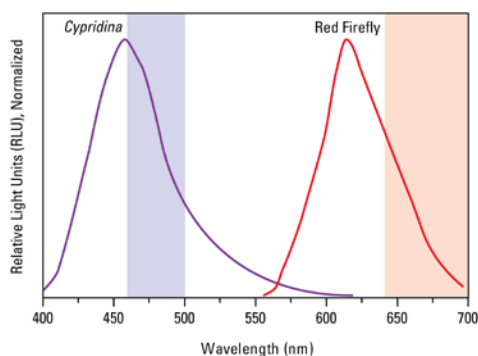


Figure 3. Thermo Scientific Pierce Cypridina-Firefly Dual Luciferase Assay emission spectra profile. The shaded boxes represent the filter ranges. See Table 1 for the filter details.

Table 1. Filter requirement for the Thermo Scientific Pierce Cypridina-Firefly Luciferase Dual Assay Kit.

Luciferase	Emission Range (nm)	Emission Maximum (nm)	Recommended Filter* (nm)
<i>Cypridina</i>	400-600	463	480±20 BP
Red Firefly	560-700	613	640 LP

*The 480±20nm bandpass (BP) is designed to capture light wavelengths ranging from 460 to 500nm (Figure 3). Similarly, the 640nm longpass (LP) filter collects wavelengths above 640nm. Each luminometer requires specific filters; filter availability and dimensions are available from Omega Optical, Inc. or Chroma Technology Corp.


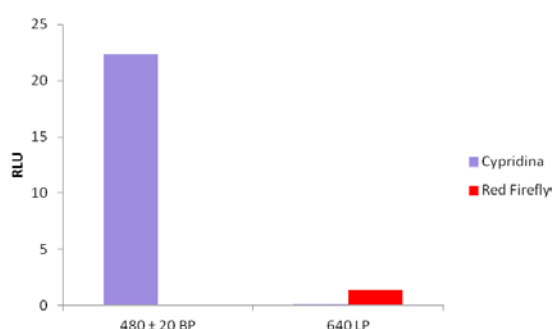


Figure 4. Specificity analysis of the two luciferase emissions. *Cypridina* luciferase was measured using a 480±20nm BP filter and red firefly luciferase was measured using a 640nm LP filter on a Thermo Scientific™ Luminoskan™ Ascent Luminometer.

Procedure Summary

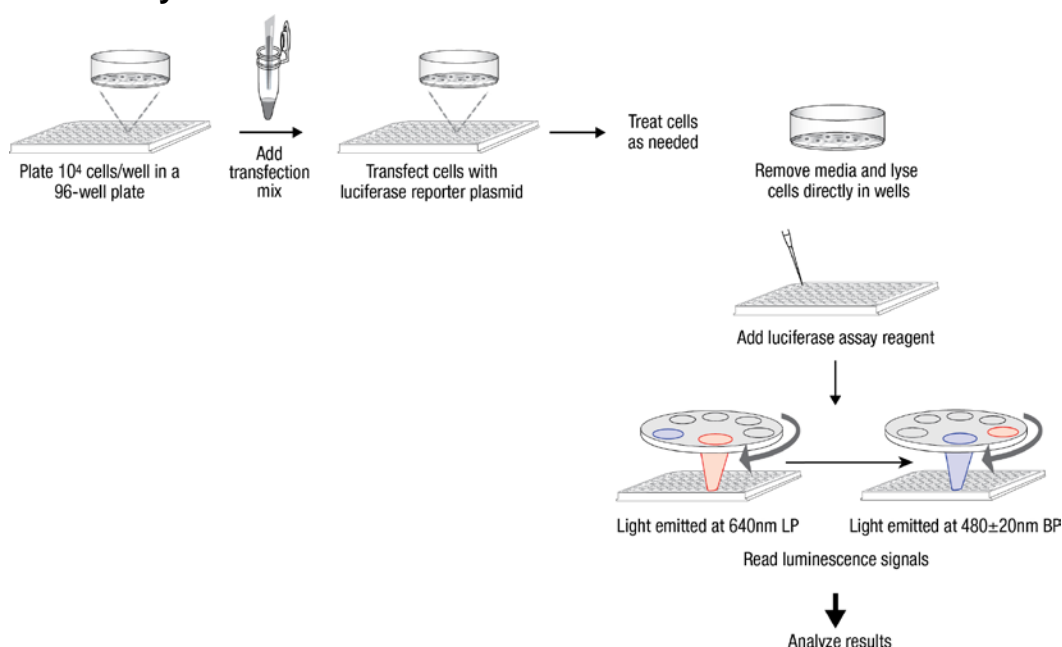

Important Product Information

- For long-term use, store Vargulin (100X) at -20°C and D-Luciferin (lyophilized) at 4°C.
- Briefly centrifuge tubes of Vargulin (100X) before use. Vargulin (100X) is volatile; seal tube tightly after use.
- Cypridina-Firefly Luciferase Dual Assay Working Solution (Working Solution), Vargulin (100X) and D-Luciferin must always be stored protected from light.
- *Cypridina* luciferase protein is significantly more stable than red firefly luciferase protein in cell culture lysate. Red firefly luciferase in cell lysate is subject to degradation. When working with red firefly luciferase in cell lysates, store the container on ice and perform luciferase assays immediately after cell lysis. Adding protease inhibitors (see Related Thermo Scientific Products) to the cell lysis buffer increases the stability of red firefly luciferase.
- To avoid cross-contamination, use a new disposable pipette tip for each transfer. If using a multichannel pipette, always use a new disposable reagent reservoir.
- Avoid exposing reagents to excessive heat or light during storage and incubation.
- Do not mix reagents from different lots. Do not discard unused working solutions after assay completion. Do not combine leftover reagents with those reserved for additional plates.
- Individual components might contain corrosives and/or preservatives. Wear gloves while performing the assay to avoid contact with samples and reagents. Please follow proper disposal procedures.
- Dispense and equilibrate to room temperature only the reagent volumes needed for the number of plates being used.

Additional Materials Required

- Reagents and equipment for propagating mammalian cells in culture
- Reagents (e.g., Thermo Scientific™ TurboFect™ Transfection Reagent, Product No. R0533) and equipment for transfecting plasmid DNA into mammalian cells
- Laboratory platform shaker
- Pipettes and/or liquid handling equipment
- Luminometer or other luminescence-monitoring instrument equipped with reagent injectors and filter wheel
- 480±20nm BP and 640nm LP filters (see Figure 3 and Table 1)
- White or black opaque, 96- or 384-well microplates

Material Preparation

100X D-Luciferin Stock Solution	Reconstitute lyophilized D-Luciferin pellet in 50µL (100 reaction kit) or 0.5mL (1000 reaction kit) of Cypridina-Firefly Luciferase Dual Assay Buffer. Store at -20°C for up to two months.
Working Solution	For 100 reactions, add 50µL of 100X Vargulin and 50µL of 100X reconstituted D-Luciferin to 5mL of Cypridina-Firefly Luciferase Dual Assay Buffer. Use 50µL of the Working Solution per reaction. Use within 2 hours or store in single-use volumes at -20°C for up to two months.
1X Cell Lysis Buffer	Dilute 2X Cell Lysis Buffer with an equal volume of ultrapure water.

Assay Protocol

A. Cell Transfection

1. Plate ~10,000 cells/well in a 96-well plate. Incubate cells overnight at 37°C in 5% CO₂. If using a different plate size, adjust the cell number accordingly. Use only cells growing in log phase at a passage number ≤ 15.

Note: Plate sufficient wells to perform the experiment in triplicate; include appropriate controls, such as nonspecific signal (i.e., non-transfected cells).

2. Use a standard protocol to transfect mammalian cells with two plasmids; *Cypridina* luciferase driven by an experimental promoter and red firefly luciferase driven by a strong constitutive promoter (CMV).

Note: Empirically determine the optimum ratio for the two plasmids. When using TurboFect Transfection Reagent (Product No. R0533), use an equal molar ratio of two plasmids up to 100ng total per well of a 96-well plate (e.g., 50ng of *Cypridina* luc reporter plasmid: 50ng of red firefly luc control plasmid).

Note: For optimal results, use a visual transfection control. For example, transfect cells in a separate well with a constitutively expressed GFP plasmid and observe GFP expression using a fluorescence microscope.

3. Incubate cells for 16-72 hours at 37°C in 5% CO₂ in a cell culture incubator.
4. Proceed with the individual experimental protocol for cell treatment.

B. Cell Lysis

1. Aspirate media from the cells, rinse once with 50-100µL/well of 1X DPBS buffer (Thermo Scientific™ BupH™ Modified Dulbecco's PBS, Product No. 28374), aspirate DPBS and add 100µL/well of 1X Cell Lysis Buffer. Do not disturb the cell monolayer during the DPBS rinse.
2. Shake the plate on a platform shaker at moderate speed for 15 minutes. Check for complete cell lysis using a light microscope. If lysis is incomplete, continue shaking the plate for 15 additional minutes.

C. Luciferase Dual Assay

1. Program the luminometer and injector and prime the system with Working Solution.
2. Add 10-20µL of cell lysate to a black flat-bottomed 96-well plate.
3. Using the luminometer's injectors, inject 50µL of Working Solution into each well containing cell lysate.
Note: For best results, use injectors for >24 wells.
4. Immediately after adding the reagent, program the luminometer to detect the light output using a 640nm LP filter to capture the red firefly luciferase signal; then immediately read the sample again using a 480±20nm BP filter to capture the *Cypridina* luciferase signal.
5. Repeat Steps 3 and 4 for all samples in the plate. Use the red firefly luciferase signal as a normalization control for the *Cypridina* luciferase signal.

Note: Follow the manufacturer's recommendations for using injector velocity to obtain a uniform coating of liquid in the well. Adjust the detector's integration time to achieve a signal within the linear range of the instrument.

Troubleshooting

Problem	Possible Cause	Solution
No signal	Low transfection efficiency	Optimize transfection conditions using a visual transfection control (e.g., a plasmid over-expressing a fluorescent protein)
		Verify plasmid DNA quality using restriction digestion and agarose gel electrophoresis Note: Most high-quality plasmid DNA should be supercoiled.
		Use actively dividing, low-passage cells
		Use a different cell type
	No promoter induction	Incubate cells using promoter-specific inducing conditions
		Incubate the cells for a longer time after treatment
		Change growth conditions to improve expression
		Use a different promoter
	Vargulin or D-Luciferin auto-oxidized	Protect substrate from light and air
		Maintain 100X Vargulin and D-Luciferin at -20°C
Low signal in lysate	Low luciferase expression	Lyse cells in a smaller volume of 1X Cell Lysis Buffer
		Use a different promoter or growth conditions to improve expression
		Increase the integration time on the instrument
		Scale-up the sample volume and reagent per well
High signal	High luciferase expression	Reduce incubation time before collecting samples
		Decrease the integration time on the instrument
		Dilute the sample Note: A low sample volume can increase assay variability. Dilute the sample and use the recommended volume of 10-20µL per assay
High background signal	Nonspecific oxidation of Vargulin and D-Luciferin	Avoid repeated freezing and thawing of the sample
	Control sample was contaminated	Change pipette tips after each well
		Reduce shaker speed during the cell lysis step to avoid contaminating the wells

Related Thermo Scientific Products

See our website for a complete list of related luciferase products.

16149	pMCS-Cypridina Luc
16150	PCMV-Cypridina Luc
16155	pMCS-Red Firefly Luc
16156	pCMV-Red Firefly Luc
16151	pTK-Cypridina Luc
16157	pTK-Red Firefly Luc
16189	Pierce Luciferase Cell Lysis Buffer (2X)
28374	BupH Modified Dulbecco's PBS Packs, 40 packs

78425 **Halt™ Protease Inhibitor Single-Use Cocktail, EDTA-free (100X), 24 × 100µL**
R0533 **TurboFect Transfection Reagent, 200µL**

General Reference

Osamu, S. (2006). Bioluminescence: Chemical Principles and Methods. World Scientific Publishing Co. Pte. Ltd, Hakensack NJ.

Products are warranted to operate or perform substantially in conformance with published Product specifications in effect at the time of sale, as set forth in the Product documentation, specifications and/or accompanying package inserts ("Documentation"). No claim of suitability for use in applications regulated by FDA is made. The warranty provided herein is valid only when used by properly trained individuals. Unless otherwise stated in the Documentation, this warranty is limited to one year from date of shipment when the Product is subjected to normal, proper and intended usage. This warranty does not extend to anyone other than Buyer. Any model or sample furnished to Buyer is merely illustrative of the general type and quality of goods and does not represent that any Product will conform to such model or sample.

NO OTHER WARRANTIES, EXPRESS OR IMPLIED, ARE GRANTED, INCLUDING WITHOUT LIMITATION, IMPLIED WARRANTIES OF MERCHANTABILITY, FITNESS FOR ANY PARTICULAR PURPOSE, OR NON INFRINGEMENT. BUYER'S EXCLUSIVE REMEDY FOR NON-CONFORMING PRODUCTS DURING THE WARRANTY PERIOD IS LIMITED TO REPAIR, REPLACEMENT OF OR REFUND FOR THE NON-CONFORMING PRODUCT(S) AT SELLER'S SOLE OPTION. THERE IS NO OBLIGATION TO REPAIR, REPLACE OR REFUND FOR PRODUCTS AS THE RESULT OF (I) ACCIDENT, DISASTER OR EVENT OF FORCE MAJEURE, (II) MISUSE, FAULT OR NEGLIGENCE OF OR BY BUYER, (III) USE OF THE PRODUCTS IN A MANNER FOR WHICH THEY WERE NOT DESIGNED, OR (IV) IMPROPER STORAGE AND HANDLING OF THE PRODUCTS.

Unless otherwise expressly stated on the Product or in the documentation accompanying the Product, the Product is intended for research only and is not to be used for any other purpose, including without limitation, unauthorized commercial uses, in vitro diagnostic uses, ex vivo or in vivo therapeutic uses, or any type of consumption by or application to humans or animals.

Current product instructions are available at www.thermoscientific.com/pierce. For a faxed copy, call 800-874-3723 or contact your local distributor.

© 2013 Thermo Fisher Scientific Inc. All rights reserved. All (other) trademarks are the property of Thermo Fisher Scientific Inc. and its subsidiaries. Printed in the USA.