INSTRUCTIONS

Pierce Premium Grade TCEP•HCl

PG82080	PG82089	PG82090	2544.0
Number	Description		
PG82080	Premium Grade TCEP•HCl, Tris(2-carboxyethyl)phosphine hydrochloride, 1g		
PG82089	Premium Grade TCEP•HCl, Tris(2-carboxyethyl)phosphine hydrochloride, 10g		
PG82090	Premium Grade TCEP•HO	Cl, Tris(2-carboxyethyl)phosphine hydrochloride, 100g	
	Molecular Weight: 286.65		
	CAS #51805-45-9		
	Storage: Upon receipt store Product is shipped at ambien	product at room temperature in sealed container to prevent oxic at temperature.	lation.

Introduction

Thermo Scientific[™] Pierce[™] Premium Grade Reagents are high-quality formulations of selected chemical modification reagents, specially characterized for applications where product integrity and risk minimization are critical. Compared to standard grade equivalents, Pierce Premium Grade Reagents provide more clearly defined quality and product support by including: (a) increased analytical testing and product characterization, (b) greater batch-specific information and quality assurance review, (c) extensive lot sample retention, and (d) change control notification.

The Thermo ScientificTM PierceTM Premium Grade TCEP is a potent, versatile, odorless, thiol-free reducing agent with broad application to protein and other research involving reduction of disulfide bonds (Figure 1). The unique compound is easily soluble and very stable in many aqueous solutions. TCEP reduces disulfide bonds as effectively as dithiothreitol (DTT), but unlike DTT and other thiol-containing reducing agents, TCEP does not have to be removed before certain sulfhydryl-reactive crosslinking reactions.

The ability and virtues of trialkylphosphine compounds to reduce protein disulfide bonds have been known for many years.^{1,2} Phosphines are stable in aqueous solution, selectively reduce disulfide bonds and are essentially non-reactive toward other functional groups commonly found in proteins.² However, widespread adoption of trialkylphosphines as reductants for protein research was hindered by their disagreeable odor and poor water solubility. These obstacles were overcome by discovery of Tris(2-carboxyethyl)phosphine (TCEP).³

TCEP selectively and completely reduces even the most stable water-soluble alkyl disulfides over a wide pH range.⁴ Reductions frequently require less than five minutes at room temperature. TCEP is non-volatile, odorless, and unlike most other reducing agents, is resistant to air oxidation. Compared to DTT, TCEP is more stable, more effective and able to reduce disulfide bonds at lower pHs.⁵

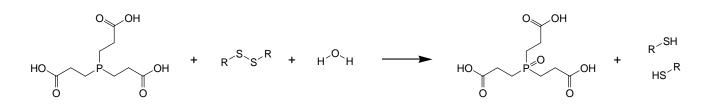


Figure 1. Reduction of organic disulfide bonds with TCEP.





Product Information

Solubility

• The hydrochloride salt (TCEP•HCl, MW 286.65) has a solubility in water of 310g/L (1.08M). Being hydrophilic, TCEP is generally very soluble in aqueous buffers at nearly any pH. Therefore, working concentrations and 10X stock solutions may be readily prepared in most aqueous buffers (see subsequent points about stability). TCEP has only minimal solubility in organic solvents, including methanol and ethanol.

Stability in Solution

- TCEP is stable in aqueous, acidic and basic solutions. When TCEP is dissolved directly in water, the resulting pH is approximately 2.5. Studies indicate that no change in concentration of TCEP occurs after 24-hour incubation at room temperature in 100mM HCl, 100mM NaOH, or in any of the following 50mM buffers: Tris•HCl (pH 7.5, 8.5 and 9.5), HEPES (pH 6.8 and 8.2), borate (pH 8.2 and 10.2) and CAPS (pH 9.7 and 11.1).⁵ Even after three weeks in these buffers, less than 20% of the TCEP was oxidized.
- TCEP is not particularly stable in phosphate buffers, especially at neutral pH. Experiments indicate that TCEP completely oxidizes within 72 hours in 0.35M phosphate-buffered saline (PBS), pH 7.0. Approximately 50% oxidation occurs in the same amount of time in 0.15M PBS, pH 8.0. Only minimal oxidation occurs in PBS at pH > 10.5 or pH < 6.0. Therefore, if using TCEP in PBS buffers, prepare the working solution immediately before use.

Effective Reducing pH

• TCEP effectively reduces disulfide bonds over a broad pH range.⁵ In one experiment, TCEP completely reduced 2,2'dithiodipyridine (2,2'-DTDP) within 30 seconds at 1.5 < pH < 9.0. Above pH 9.0, only 50% reduction occurred. TCEP is a more effective than DTT at pH < 8.0; TCEP will even reduce oxidized DTT.⁵

Working Concentration

For most applications, 5-50mM TCEP provides sufficient molar excess to effectively reduce peptide or protein disulfide bonds within a few minutes at room temperature. Han demonstrated complete reduction of 2,2'-DTDP (20μM) by TCEP (30μM) at pH 1.5 to 8.5, within 40 seconds.⁵ When molar equivalents of TCEP are used, the reaction time is much longer. Mery *et al.* reported that a 1:1 ratio of TCEP to disulfides required nearly one hour to complete the reduction.⁶

Compatible Applications

- Because TCEP does not contain thiols, it does not have to be removed from solutions before performing reactions involving maleimide labeling or crosslinking reagents. In most situations, TCEP concentrations < 10-20mM are compatible with maleimide reaction chemistry.
- TCEP may be used as a substitute for DTT or 2-mercaptoethanol (2-ME) in reducing sample loading buffer for SDS-PAGE; use a final concentration of 50mM TCEP (see also Related Thermo Scientific Products). However, because TCEP is charged in solution, it is not compatible for use in isoelectric focusing (IEF), which is the first dimension of 2D electrophoresis.
- One strategy for labeling antibodies involves reducing disulfide bonds in the hinge region of IgG molecules without also reducing those connecting heavy and light chains. This "partial" reduction is often accomplished with 2-mercapto-ethylamine (2-MEA, see Related Thermo Scientific Products). Bläuenstein *et al.* demonstrated that TCEP can be used for this purpose⁷: Add TCEP at a final concentration of 3.8-4.0mM to 10mg/mL IgG in 0.1M phosphate buffer (pH 4.6-7.5); incubate for 20-30 minutes at room temperature, then use a desalting column to remove the TCEP products.

Related Thermo Scientific Products

24090	Pierce TCEP•HCl, Tris(2-carboxyethyl)phosphine hydrochloride, 1g
24091	Pierce TCEP•HCl, Tris(2-carboxyethyl)phosphine hydrochloride, 10g
77720	Bond-Breaker TM TCEP Solution, Neutral pH, 5mL
77712	Immobilized TCEP Disulfide Reducing Gel, 5mL
20408	2-Mercaptoethylamine•HCl (2-MEA), 6 × 6mg
20290	Dithiothreitol (DTT), 5g
20291	Dithiothreitol (DTT), No-Weigh TM Format, 48 tubes × 7.7mg



References

- 1. Ruegg, U.T. and Rudinger, J. (1977). Reductive cleavage of cystine disulfides with tributylphosphine. Methods Enzymol 47:111-26.
- 2. Kirley, T.L. (1989). Reduction and fluorescent labeling of cyst(e)ine-containing proteins for subsequent structural analysis. Anal Biochem 180:231.
- Burns, J.A., et al. (1991). Selective reduction of disulfides by tris-(2-carboxyethyl)-phosphine. J Org Chem 56:2648-50. 3.
- Levison, M.E., et al. (1969). Reduction of biological substances by water-soluble phosphines: Gamma-globulin. Experentia 25:126-7. 4
- Han, J.C., et al. (1994). A procedure for quantitative determination of tris(2-carboxyethyl)phosphine, an odorless reducing agent more stable and 5. effective than dithiothreitol. Anal Biochem 220:5-10.
- Mery, J., et al. (1993). Disulfide linkage to polyacrylic resin for automated Fmoc peptide synthesis. Immunochemical applications of peptide resins and 6. mercaptoamide peptides. Int J Peptide Protein Res 42:44-52.
- Blauenstein, P., et al. (1995). Experience with the iodine-123 and technetium-99m labeled anti-granulocyte antibody Mab47: a comparison of labeling 7 methods. Eur J Nuclear Med 22:690-8.

Products are warranted to operate or perform substantially in conformance with published Product specifications in effect at the time of sale, as set forth in the Product documentation, specifications and/or accompanying package inserts ("Documentation"). No claim of suitability for use in applications regulated by FDA is made. The warranty provided herein is valid only when used by properly trained individuals. Unless otherwise stated in the Documentation, this warranty is limited to one year from date of shipment when the Product is subjected to normal, proper and intended usage. This warranty does not extend to anyone other than Buyer. Any model or sample furnished to Buyer is merely illustrative of the general type and quality of goods and does not represent that any Product will conform to such model or sample.

NO OTHER WARRANTIES, EXPRESS OR IMPLIED, ARE GRANTED, INCLUDING WITHOUT LIMITATION, IMPLIED WARRANTIES OF MERCHANTABILITY, FITNESS FOR ANY PARTICULAR PURPOSE, OR NON INFRINGEMENT, BUYER'S EXCLUSIVE REMEDY FOR NON-CONFORMING PRODUCTS DURING THE WARRANTY PERIOD IS LIMITED TO REPAIR, REPLACEMENT OF OR REFUND FOR THE NON-CONFORMING PRODUCT(S) AT SELLER'S SOLE OPTION. THERE IS NO OBLIGATION TO REPAIR. REPLACE OR REFUND FOR PRODUCTS AS THE RESULT OF (I) ACCIDENT, DISASTER OR EVENT OF FORCE MAJEURE, (II) MISUSE, FAULT OR NEGLIGENCE OF OR BY BUYER, (III) USE OF THE PRODUCTS IN A MANNER FOR WHICH THEY WERE NOT DESIGNED, OR (IV) IMPROPER STORAGE AND HANDLING OF THE PRODUCTS.

Unless otherwise expressly stated on the Product or in the documentation accompanying the Product, the Product is intended for research only and is not to be used for any other purpose, including without limitation, unauthorized commercial uses, in vitro diagnostic uses, ex vivo or in vivo therapeutic uses, or any type of consumption by or application to humans or animals.

Current product instructions are available at www.thermoscientific.com/pierce. For a faxed copy, call 800-874-3723 or contact your local distributor. © 2013 Thermo Fisher Scientific Inc. All rights reserved. Unless otherwise indicated, all trademarks are property of Thermo Fisher Scientific Inc. and its subsidiaries. Printed in the USA.