

FAK [pY397] ELISA Kit

Catalog Number KHO0441 (96 tests)

Pub. No. MAN0015690 Rev. 2.0 (30)

CAUTION! This kit contains materials with small quantities of sodium azide. Sodium azide reacts with lead and copper plumbing to form explosive metal azides. Upon disposal, flush drains with a large volume of water to prevent azide accumulation. Avoid ingestion and contact with eyes, skin and mucous membranes. In case of contact, rinse affected area with plenty of water. Observe all federal, state, and local regulations for disposal.

Note: For safety and biohazard guidelines, see the “Safety” appendix in the *ELISA Technical Guide* (Pub. no. MAN0006706). Read the Safety Data Sheets (SDSs) and follow the handling instructions. Wear appropriate protective eyewear, clothing, and gloves.

Product description

The Invitrogen™ FAK [pY397] ELISA Kit is a solid-phase sandwich Enzyme-Linked Immunosorbent Assay (ELISA). This assay is designed to detect and quantify level of human and mouse FAK phosphorylated at tyrosine residue 397 (FAK [pY397] in cell lysates. The assay recognizes both natural and recombinant FAK [pY397].

Focal Adhesion Kinase (FAK), also known as pp125FAK and FADK 1 (EC,2.7.1.112) is a non-receptor protein-tyrosine kinase that localizes to focal adhesions. FAK appears to be ubiquitously expressed among all mammalian tissues, with highest expression levels observed in brain tissue.

FAK is activated by phosphorylation events in response to several stimuli. In response to integrin engagement, FAK is autophosphorylated at tyrosine residue 397. Phosphorylation also occurs at tyrosine residues 407, 576, 577, 861, and 925.

A FAK (Total) ELISA kit (Cat. No. KHO0431), which is independent of phosphorylation status, is available for normalizing FAK in samples.

Contents and storage

Upon receipt, store the kit at 2°C to 8°C.

Contents	Cat. No. KHO0441 (96 tests)
FAK [pY397] Standard, lyophilized; contains 0.1% sodium azide	2 vials
Standard Diluent Buffer; contains 0.1% sodium azide	25 mL
Antibody Coated Plate, 96-well plate	1 plate
FAK [pY397] Detection Antibody; contains 0.1% sodium azide and blue dye ^[1]	6.0 mL
Anti-Rabbit IgG HRP (100X); contains 3.3 mM thymol	0.125 mL
HRP Diluent; contains 3.3 mM thymol and yellow dye ^[1]	25 mL
Wash Buffer Concentrate (25X)	100 mL
Stabilized Chromogen, Tetramethylbenzidine (TMB)	25 mL
Stop Solution	25 mL
Plate Covers, adhesive strips	3

^[1] In order to help our customers avoid any mistakes in pipetting the ELISAs, we provide colored *Standard Diluent Buffer*, *Detection Antibody*, and *HRP Diluent* to help monitor the addition of solution to the reaction well. This does not in any way interfere with the test results.

Required materials not supplied

- Distilled or deionized water
- Microtiter plate reader with software capable of measurement at or near 450 nm
- Plate washer—automated or manual (squirt bottle, manifold dispenser, or equivalent)
- Calibrated adjustable precision pipettes and glass or plastic tubes for diluting solutions; beakers, flask and cylinders for preparation of reagents
- Cell Extraction Buffer (Cat. No. FNN0011, or see “Prepare Cell Extraction Buffer”)

Before you begin

IMPORTANT! Reagents are lot-specific. Do not mix or interchange different reagent lots from various kit lots.

- Review the **Procedural guidelines** and **Plate washing directions** in the *ELISA Technical Guide* available at thermofisher.com.
- Allow reagents to reach room temperature before use. Mix to redissolve any precipitated salts.

Prepare 1X Wash Buffer

1. Dilute 16 mL of Wash Buffer Concentrate (25X) with 384 mL of deionized or distilled water. Label as 1X Wash Buffer.
2. Store the concentrate and 1X Wash Buffer in the refrigerator. Use the diluted buffer within 14 days.

Prepare Cell Extraction Buffer

Note: See the *ELISA Technical Guide* for detailed information on preparing Cell Extraction Buffer.

1. Prepare Cell Extraction Buffer.

Cell Extraction Buffer consists of 10 mM Tris (pH 7.4), 100 mM NaCl, 1 mM EDTA, 1 mM EGTA, 1 mM NaF, 20 mM Na₄P₂O₇, 2 mM Na₃VO₄, 1% Triton™ X-100, 10% glycerol, 0.1% SDS, and 0.5% deoxycholate.

2. Immediately before use, add PMSF (0.3 M stock in DMSO) to 1 mM and 50 µL protease inhibitor cocktail (e.g., Sigma Cat. No. P-2714) for each 1 mL of Cell Extraction Buffer.

Prepare cell lysate

1. Collect cells by centrifugation (non-adherent cells) or scraping from culture flasks (adherent cells), then wash cells twice with cold PBS.
2. Remove and discard the supernatant and collect the cell pellet. The pellet can be stored at -80°C and lysed at a later date if desired.
3. Lyse the cell pellet in Cell Extraction Buffer for 30 minutes, on ice. Vortex at 10-minute intervals.

Note: The volume of Cell Extraction Buffer used depends on the number of cells in the cell pellet, and expression levels of FAK [pY397]. Researchers must optimize the extraction procedures for their own applications.

4. Transfer the lysate into microcentrifuge tubes and centrifuge at 13,000 rpm for 10 minutes at 4°C.
5. Transfer the supernatant into clean microcentrifuge tubes. Samples can be stored at -80°C (avoid multiple freeze-thaw cycles).

Pre-dilute samples

Sample concentrations should be within the range of the standard curve. Because conditions may vary, each investigator should determine the optimal dilution for each application.

- Perform sample dilutions with Standard Diluent Buffer (serum/plasma) or with the corresponding cell culture medium (cell culture supernatant).
- Dilute use of 1-10 µL of the clarified cell extract to a volume of 50 µL/well in Standard Diluent Buffer (See Assay Method) is sufficient for the detection of FAK [pY397].
- Dilute samples prepared in Cell Extraction Buffer 1:5 or greater in Standard Diluent Buffer (e.g., 10 µL sample into 40 µL buffer).


This dilution is necessary to reduce the matrix effect of the Cell Extraction Buffer. SDS concentration should be less than 0.01% before adding to the plate. While a 1:5 sample dilution has been found to be satisfactory, higher dilutions such as 1:10 or 1:20 may be optimal.

Dilute standards

Note: Use glass or plastic tubes for diluting standards.

Note: The FAK [pY397] Standard is prepared using purified, full length, recombinant, phosphorylated FAK protein. One Unit of standard is equivalent to the amount of FAK [pY397] autophosphorylated from 300 pg of total FAK protein.

1. Reconstitute FAK [pY397] Standard to 100 Units/mL with Standard Diluent Buffer. Refer to the standard vial label for instructions. Swirl or mix gently and allow the contents to sit for 10 minutes to ensure complete reconstitution. Label as 100 Units/mL FAK [pY397]. **Use the standard within 1 hour of reconstitution.**
2. Add 150 µL Standard Diluent Buffer to each of 7 tubes labeled as follows: 50, 25, 12.5, 6.25, 3.12, 1.6, and 0 Units/mL FAK [pY397].
3. Make serial dilutions of the standard as shown in the following dilution diagram. Mix thoroughly between steps.
4. Remaining reconstituted standard should be discarded or frozen in aliquots at -80°C for further use. Standard can be frozen and thawed one time only without loss of immunoreactivity.

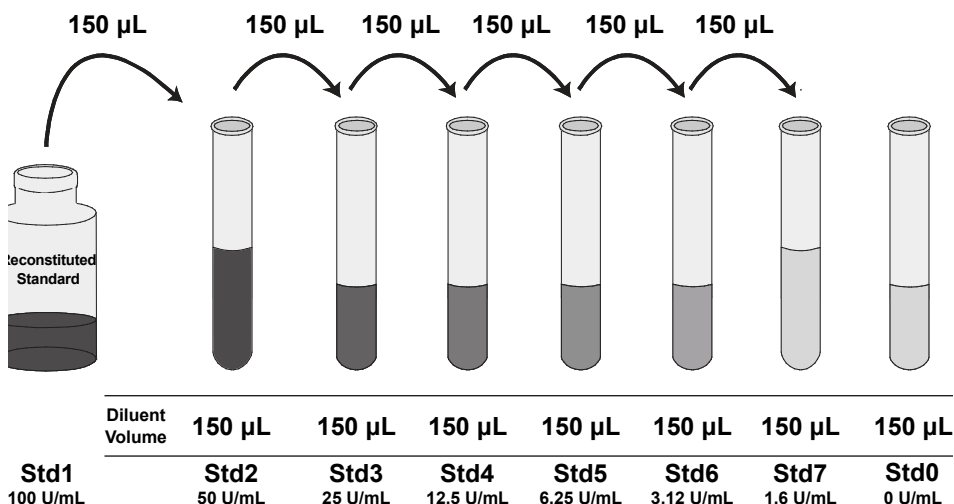

Prepare 1X Anti-Rabbit IgG HRP solution

Note: Prepare 1X Anti-Rabbit IgG HRP solution within 15 minutes of usage.

The Anti-Rabbit IgG HRP (100X) is in 50% glycerol, which is viscous. To ensure accurate dilution:

1. For each 8-well strip used in the assay, pipet 10 µL Anti-Rabbit IgG HRP (100X) solution, wipe the pipette tip with clean absorbent paper to remove any excess solution, and dispense the solution into a tube containing 1 mL of HRP Diluent. Mix thoroughly.
2. Return the unused Anti-Rabbit IgG HRP (100X) solution to the refrigerator.

Perform ELISA (Total assay time: 4 hours)

IMPORTANT! Perform a standard curve with each assay.

- Allow all components to reach room temperature before use. Mix all liquid reagents prior to use.
- Determine the number of 8-well strips required for the assay. Insert the strips in the frames for use. Re-bag any unused strips and frames, and store at 2°C to 8°C for future use.



1	Bind antigen and add detector	<ol style="list-style-type: none"> Add 50 μL of standards, controls, or samples (see “Pre-dilute samples” on page 2) to the appropriate wells. Leave the wells for chromogen blanks empty. Add 50 μL of FAK [pY397] Detection Antibody solution into each well except the chromogen blanks. Tap the side of the plate to mix. Cover the plate with a plate cover and incubate 3 hours at room temperature. Thoroughly aspirate the solution and wash wells 4 times with 1X Wash Buffer.
2	Add IgG HRP	<ol style="list-style-type: none"> Add 100 μL 1X Anti-Rabbit IgG HRP solution into each well except the chromogen blanks. Cover the plate with plate cover and incubate for 30 minutes at room temperature. Thoroughly aspirate the solution and wash wells 4 times with 1X Wash Buffer.
3	Add Stabilized Chromogen	<ol style="list-style-type: none"> Add 100 μL Stabilized Chromogen to each well. The substrate solution begins to turn blue. Incubate for 30 minutes at room temperature in the dark. <p>Note: TMB should not touch aluminum foil or other metals.</p>
4	Add Stop Solution	Add 100 μ L Stop Solution to each well. Tap the side of the plate to mix. The solution in the wells changes from blue to yellow.

Read the plate and generate the standard curve

1. Read the absorbance at 450 nm. Read the plate within 2 hours after adding the Stop Solution.
2. Use curve-fitting software to generate the standard curve. A 4 parameter algorithm provides the best standard curve fit. Optimally, the background absorbance may be subtracted from all data points, including standards, unknowns and controls, prior to plotting.
3. Read the concentrations for unknown samples and controls from the standard curve. Multiply value(s) obtained for sample(s) by the appropriate factor to correct for the sample dilution.

Note: Dilute samples producing signals greater than the upper limit of the standard curve in Standard Diluent Buffer (serum/plasma) or with the corresponding cell culture medium (cell culture supernatant) and reanalyze. Multiply the concentration by the appropriate dilution factor.

Performance characteristics

Standard curve example

The following data were obtained for the various standards over the range of 0 to 100 Units/mL FAK [pY397].

Std. FAK [pY397] (U/mL)	Optical Density (450 nm)	Std. FAK [pY397] (U/mL)	Optical Density (450 nm)
100	2.82	6.25	0.36
50	1.60	3.12	0.29
25	0.94	1.6	0.24
12.5	0.61	0	0.19

Intra-assay precision

Samples of known FAK [pY397] concentration were assayed in replicates of 16 to determine precision within an assay.

Parameters	Sample 1	Sample 2	Sample 3
Mean (U/mL)	71.6	21.5	6.8
Standard Deviation	2.9	0.86	0.25
% Coefficient of Variation	4.1	4.1	3.7

Inter-assay precision

Samples were assayed 48 times in multiple assays to determine precision between assays.

Parameters	Sample 1	Sample 2	Sample 3
Mean (U/mL)	71.5	23.0	7.1
Standard Deviation	5.8	1.7	0.6
% Coefficient of Variation	8.1	7.4	8.5


Linearity of dilution

3T3L1 cells grown in cell culture medium containing 10% fetal calf serum were treated with 1 mM sodium orthovanadate for 5 hours and lysed with cell extraction buffer. The lysate was diluted in Standard Diluent Buffer over the range of the assay and measured for FAK [pY397]. Linear regression analysis of sample values versus the expected concentration yielded a correlation coefficient of 0.99.


Dilution	Cell Lysate		
	Measured (U/mL)	Expected (U/mL)	
			%
Neat	67.1	67.1	100
1/2	36.6	33.6	108.9
1/4	16.7	16.8	99.4
1/8	8.1	8.4	96.4
1/16	5.2	4.2	124

Specificity


The FAK [pY397] ELISA Kit recognizes **human, rat, and mouse** FAK. Specificity for FAK [pY397] was confirmed by peptide competition. The data shows that only the tyrosine 397 phosphopeptide blocks the ELISA signal. The non-phosphorylated peptide sequence or other FAK phosphopeptides do not block the signal.

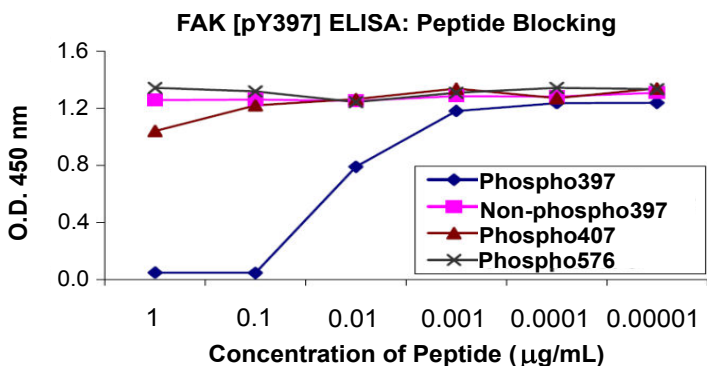

A comparison of FAK [pY397] ELISA Kit and FAK (Total) ELISA Kits was also used to confirm specificity. Lysates from 3T3L1 cells treated with sodium orthovanadate (1 mM for 5 hours) and untreated controls, and phosphorylated and non-phosphorylated versions of recombinant FAK protein were assayed. The FAK [pY397] ELISA Kit detected phosphorylated FAK protein and phosphorylated FAK in treated cells, but not non-phosphorylated FAK protein or non-phosphorylated FAK in untreated cells.

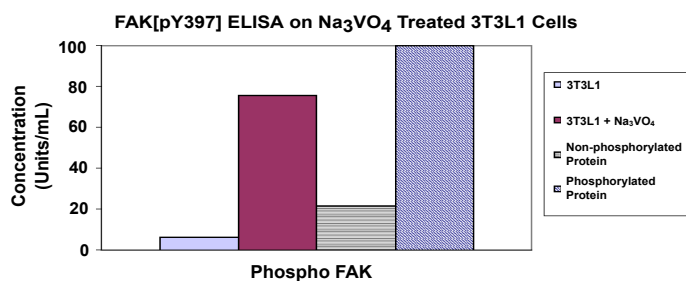

In contrast, the FAK [pY397] ELISA Kit detected both phosphorylated and non-phosphorylated FAK proteins, and FAK in both treated and untreated cells.

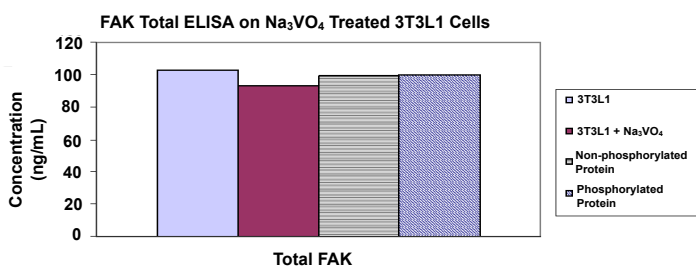


Sensitivity

The analytical sensitivity of the assay is <0.9 Units/mL FAK [pY397]. This was determined by adding two standard deviations to the mean O.D. obtained from 30 assays of the zero standard. This level of sensitivity is equivalent to the detection of FAK [pY397] in 5,000 sodium orthovanadate treated 3T3L1 cells.


The sensitivity of this ELISA is ~1:4 greater than that of western blotting when tested against known quantities of FAK [pY397]. The western blot was developed using rabbit anti-FAK [pY397], and an alkaline phosphatase conjugated anti-rabbit IgG followed by chemiluminescent substrate and autoradiography.

Detection of FAK [pY397] by ELISA vs Western Blot: 3T3L1 Cells Stimulated by Na₃VO₄

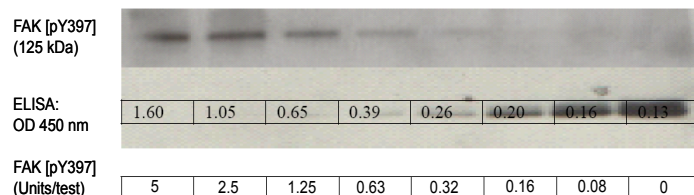

Parallelism

Natural FAK [pY397] from lysates of sodium orthovanadate-treated 3T3L1 cells were serially diluted in Standard Diluent Buffer. The optical density of each dilution was plotted against the FAK [pY397] standard curve. The standard accurately reflects FAK [pY397] content in samples.

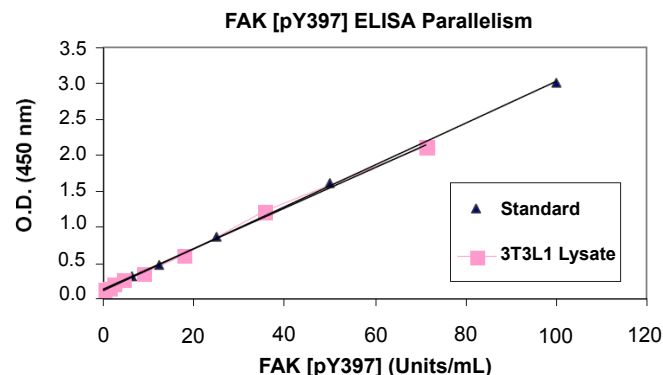

Recovery

FAK [pY397] ELISA Kit Standard was spiked into 200 µg/mL 3T3L1 cell lysate at 3 different concentrations to evaluate recovery. The average recovery was 97%.








High-dose hook effect

Samples spiked with FAK [pY397] Standard up to 400 Units/mL give responses higher than that obtained from the last standard point.

Limited product warranty

Life Technologies Corporation and/or its affiliate(s) warrant their products as set forth in the Life Technologies' General Terms and Conditions of Sale found on Life Technologies' website at www.thermofisher.com/us/en/home/global/terms-and-conditions.html. If you have any questions, please contact Life Technologies at www.thermofisher.com/support.

Product label explanation of symbols and warnings

 REF	Catalog Number	 LOT	Batch code		Temperature limitation		Use by		Manufacturer		Consult instructions for use		Caution, consult accompanying documents
--	----------------	---	------------	---	------------------------	---	--------	---	--------------	---	------------------------------	---	---

Manufacturer's address: Bender MedSystems GmbH | Campus Vienna Biocenter 2 | 1030 Vienna, Austria

The information in this guide is subject to change without notice.

DISCLAIMER: TO THE EXTENT ALLOWED BY LAW, LIFE TECHNOLOGIES AND/OR ITS AFFILIATE(S) WILL NOT BE LIABLE FOR SPECIAL, INCIDENTAL, INDIRECT, PUNITIVE, MULTIPLE, OR CONSEQUENTIAL DAMAGES IN CONNECTION WITH OR ARISING FROM THIS DOCUMENT, INCLUDING YOUR USE OF IT.

Important Licensing Information: These products may be covered by one or more Limited Use Label Licenses. By use of these products, you accept the terms and conditions of all applicable Limited Use Label Licenses.

©2017 Thermo Fisher Scientific Inc. All rights reserved. All trademarks are the property of Thermo Fisher Scientific and its subsidiaries unless otherwise specified.