

Jasplakinolide (J7473)

Quick Facts

Storage upon receipt:

• ≤-20°C

Desiccate

· Protect from light

Molecular weight: 709.68

Solvent for stock: DMSO or methanol

Absorbance in methanol: 278 ± 3 nm

Introduction

Jasplakinolide (Figure 1) was the first cyclodepsipeptide to be isolated from a marine sponge, specifically the Indo-Pacific sponge *Jaspis johnstoni*. It is a nonfluorescent, cell-permeant F-actin probe composed of three amino acids, including the rare D- β -tyrosine and D-2-bromoabrine. A potent inducer of actin polymerization and stabilization, jasplakinolide ($K_d = 15 \text{ nM}$) competes with phallotoxins for actin binding and exhibits fungicidal, insecticidal, and anticancer properties. In studies with *Candida albicans*, the MIC and MLC for jasplakinolide were established at 25 μ g/mL. The compound shows no effect upon a broad range of bacteria. ^{2,3}

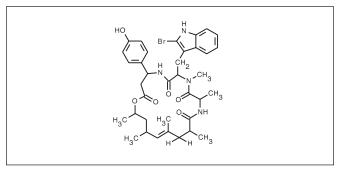


Figure 1. The structure of jasplakinolide.

Materials

Jasplakinolide is supplied in a unit size of 100 μ g. Upon receipt, it should be stored frozen at \leq -20°C, desiccated and protected from light. Properly stored, it is stable for up to 2 years.

To prepare a 1 mM stock solution of jasplakinolide, reconstitute in 140 μ L anhydrous methanol or DMSO. Once reconstituted, the stock solution is stable for 3–4 months when stored frozen at \leq -20°C, desiccated and protected from light.

Experimental Protocol

Jasplakinolide can be used for *in vitro* and *in vivo* assays using working concentrations of 50 nM to 5 μ M and incubation times ranging from a few minutes to 24 hours. ^{4,5} For use on live cells, jasplakinolide can be added to a variety of cell culture media with added serum and antibiotics. The optimal concentration and incubation time must be determined empirically.

References

1. Tetrahedron Lett 27, 2797 (1986); **2.** J Am Chem Soc 108, 3123 (1986); **3.** Antimicrob Agents Chemother 32, 1154 (1988); **4.** J Biol Chem 275, 5163 (2000); **5.** Nature Cell Biology 4, 513 (2002).

Product List Current prices may be obtained from our Web site or from our Customer Service Department.

Cat #Product NameUnit SizeJ7473jasplakinolide100 µg

MP 07473 Jasplakinolide

Contact Information

Further information on Molecular Probes' products, including product bibliographies, is available from your local distributor or directly from Molecular Probes. Customers in Europe, Africa and the Middle East should contact our office in Leiden, the Netherlands. All others should contact our Technical Assistance Department in Eugene, Oregon.

Please visit our Web site — www.probes.com — for the most up-to-date information.

Molecular Probes. Inc.

29851 Willow Creek Road, Eugene, OR 97402 Phone: (541) 465-8300 • Fax: (541) 335-0504

Customer Service: 6:00 am to 4:30 pm (Pacific Time)

Phone: (541) 335-0338 • Fax: (541) 335-0305 • order@probes.com

Toll-Free Ordering for USA and Canada:

Order Phone: (800) 438-2209 • Order Fax: (800) 438-0228

Technical Assistance: 8:00 am to 4:00 pm (Pacific Time) Phone: (541) 335-0353 • Toll-Free (800) 438-2209

Fax: (541) 335-0238 • tech@probes.com

Molecular Probes Europe BV

Poortgebouw, Rijnsburgerweg 10 2333 AA Leiden, The Netherlands

Phone: +31-71-5233378 • Fax: +31-71-5233419

Customer Service: 9:00 to 16:30 (Central European Time)

Phone: +31-71-5236850 • Fax: +31-71-5233419

eurorder@probes.nl

Technical Assistance: 9:00 to 16:30 (Central European Time)

Phone: +31-71-5233431 • Fax: +31-71-5241883

eurotech@probes.nl

Molecular Probes' products are high-quality reagents and materials intended for research purposes only. These products must be used by, or directly under the supervision of, a technically qualified individual experienced in handling potentially hazardous chemicals. Please read the Material Safety Data Sheet provided for each product; other regulatory considerations may apply.

Several of Molecular Probes' products and product applications are covered by U.S. and foreign patents and patents pending. Our products are not available for resale or other commercial uses without a specific agreement from Molecular Probes, Inc. We welcome inquiries about licensing the use of our dyes, trademarks or technologies. Please submit inquiries by e-mail to busdev@probes.com. All names containing the designation ® are registered with the U.S. Patent and Trademark Office.

Copyright 2004, Molecular Probes, Inc. All rights reserved. This information is subject to change without notice.