

Pierce Premium Grade Sulfo-NHS

PG82071 PG82072 _{2535.0}

Number Description

PG82071 Premium Grade Sulfo-NHS (*N*-hydroxysulfosuccinimide), 500mg

PG82072 Premium Grade Sulfo-NHS, 10g

Molecular Weight: 217.14 CAS # 106627-54-7

AS # 100027-54-7

Storage: Upon receipt store at 4°C. Product is shipped at ambient temperature.

Introduction

Thermo ScientificTM PierceTM Premium Grade Reagents are high-quality formulations of selected chemical modification reagents, specially characterized for applications where product integrity and risk minimization are critical. Compared to standard grade equivalents, Pierce Premium Grade Reagents provide more clearly defined quality and product support by including: (a) increased analytical testing and product characterization, (b) greater batch-specific information and quality assurance review, (c) extensive lot sample retention, and (d) change control notification.

Thermo ScientificTM PierceTM Premium Grade Sulfo-NHS is used to prepare amine-reactive esters of carboxylate groups for chemical labeling, crosslinking and solid-phase immobilization applications. Carboxylates (-COOH) may be reacted with Sulfo-NHS in the presence of a carbodiimide such as Thermo ScientificTM PierceTM Premium Grade EDC (Product No. PG82079, PG82073 and PG82074), resulting in a semi-stable Sulfo-NHS ester, which may then be reacted with primary amines (-NH₂) to form amide crosslinks (Figure 1). Although Sulfo-NHS is not required for carbodiimide reactions, its use greatly enhances coupling efficiency. Furthermore, Sulfo-NHS makes it possible to perform a two-step reaction.

Sulfo-NHS is soluble in aqueous and organic solvents. Activation with Sulfo-NHS preserves or increases water-solubility of the modified molecule, by virtue of the charged sulfonate group. Although prepared Sulfo-NHS esters are sufficiently stable to process in a two-step reaction scheme, they will hydrolyze within hours or minutes, depending on water content and pH of the reaction solution (NHS esters have a half-life of 4-5 hours at pH 7, 1 hour at pH 8 and only 10 minutes at pH 8.6.). ¹⁻³

The activation reaction with EDC and Sulfo-NHS is most efficient at pH 4.5-7.2, and EDC reactions are often performed in MES buffer (Thermo Scientific, Product No. 28390) at pH 4.7-6.0. Reaction of Sulfo-NHS-activated molecules with primary amines is most efficient at pH 7-8, and Sulfo-NHS-ester reactions are usually performed in phosphate-buffered saline (PBS) at pH 7.2-7.5. For best results in two-step reactions, perform the first reaction in MES buffer (or other non-amine, non-carboxylate buffer) at pH 5-6, then raise the pH to 7.2-7.5 with phosphate buffer (or other non-amine buffer) immediately before reaction to the amine-containing molecule. EDC reactions can be quenched with 2-mercaptoethanol (2-ME), or the excess reagent can simply be removed (as well as the reaction pH adjusted) by buffer-exchange with a desalting column (see Related Thermo Scientific Products).



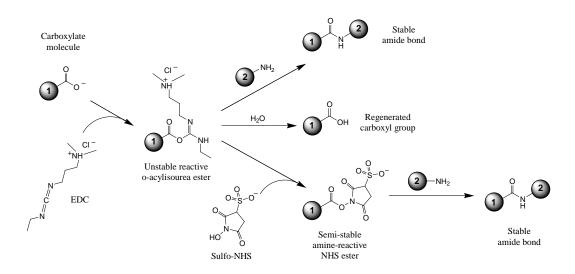


Figure 1. Reactions involving EDC, including activation as an NHS ester.

Procedure for EDC/Sulfo-NHS Crosslinking of Carboxylates with Primary Amines

A. Additional Materials Required

- Activation Buffer: 0.1M MES (2-[morpholino]ethanesulfonic acid), 0.5M NaCl, pH 6.0. Alternatively, use Thermo ScientificTM BupHTM MES Buffered Saline (Product No. 28390)
- Phosphate-buffered Saline (PBS): 0.1M sodium phosphate, 0.15M NaCl, pH 7.2-7.5 (e.g., Thermo Scientific, Product No. 28372)
- Protein #1: Prepare 1mL of Protein #1 in activation buffer at ~10mg/mL
- Protein #2: Lyophilized or dissolved at 1-10mg/mL in PBS or other amine-free buffer, pH 7-8
- Pierce Premium Grade EDC (1-ethyl-3-[3-dimethylaminopropyl]carbodiimide) (Product No. PG82079, PG82073, PG82074) for best results, use a 10-fold molar excess of EDC (MW = 191.7) to Protein #1
- (Optional) 2-Mercaptoethanol (e.g., Thermo Scientific, Product No. 35600) for quenching EDC activation reaction
- (Optional) Desalting column of appropriate size for the volume of final activation reaction (e.g., Thermo ScientificTM ZebaTM Spin Desalting Columns). If intending to use this method for clean-up and buffer exchange of the activation reaction, be sure to equilibrate the desalting column so that it is ready for use when needed in Section C.
- (Optional) Hydroxylamine (e.g., Thermo Scientific, Product No. 26103) for quenching the amine reaction

B. Sulfo-NHS-ester Activation (for larger-scale activations, scale accordingly)

- 1. Add 0.4mg of Pierce Premium Grade EDC (final concentration 2mM) directly to 1mL of Protein #1, which, based on a 50kDa protein, results in a 10-fold molar excess of EDC to Protein #1.
- 2. Add 1.1mg of Pierce Premium Grade Sulfo-NHS to the reaction (final concentration 5mM).
- 3. Mix reaction components well and react for 15 minutes at room temperature.
- 4. (Optional): Add 1.4μL of 2-mercaptoethanol (final concentration of 20mM) to inactivate the EDC.
- 5. (Optional): Separate activated Protein #1 from excess EDC, EDC by-products, Sulfo-NHS and, if used, 2-mercaptoethanol using an appropriate size desalting column that has been equilibrated with PBS. Follow desalting column instructions and recover the fraction containing the activated protein. If using absorbance at 280nm to identify fractions containing protein, be aware that Sulfo-NHS absorbs strongly at 260-280nm.



C. Amine Reaction

- 1. If Step B.5 was not performed (i.e., buffer not exchanged using a desalting column), then increase buffer pH above 7.0 using concentrated PBS or other non-amine buffer such as sodium bicarbonate.
- 2. Add Protein #2 to the solution containing activated Protein #1.
- 3. Mix the solution well and then allow reaction to proceed for 2 hours at room temperature.
- 4. (Optional): Quench reaction by adding hydroxylamine to a final concentration of 10mM. The excess hydroxylamine reacts with all Sulfo-NHS esters remaining on the surface of Protein #1, resulting in conversion of the original carboxyl groups to a hydroxamic acid. Alternative quenching reagents include 20-50mM Tris, lysine, glycine and ethanolamine. Addition of base to raise the pH > 8 will promote hydrolysis of the Sulfo-NHS esters, thereby regenerating the original carboxyl groups.

Related Thermo Scientific Products

PG82079	Pierce Premium Grade EDC (1-ethyl-3-[3-dimethylaminopropyl]carbodiimide), 1g
PG82073	Pierce Premium Grade EDC, 25g
PG82074	Pierce Premium Grade EDC, 500g
77149	Pierce EDC, 10mg
22980	Pierce EDC, 5g
22981	Pierce EDC, 25g
24510	Pierce Sulfo-NHS (N-hydroxysulfosuccinimide), 500mg
24525	Pierce Sulfo-NHS, 5g
24520	Pierce Sulfo-NHS, No-Weigh TM Format, 8 × 2mg microtubes
28390	BupH TM MES Buffered Saline, 10 packs, each pack results in 0.1M MES, 0.9% NaCl, pH 4.7 when dissolved in 500mL water
28372	BupH Phosphate Buffered Saline Packs, 40 packs, each pack results in 0.1M sodium phosphate, 0.15M NaCl, pH 7.2 when dissolved in 500mL water
89891	Zeba Spin Desalting Columns, 7K MWCO, 5mL, 5/pkg
89892	Zeba Spin Desalting Columns, 7K MWCO, 5mL, 25/pkg
89893	Zeba Spin Desalting Columns, 7K MWCO, 10mL, 5/pkg
89894	Zeba Spin Desalting Columns, 7K MWCO, 10mL, 25/pkg

Cited References

- Lomant, A.J. and Fairbanks, G. (1976). Chemical probes of extended biological structures: synthesis and properties of the cleavable cross-linking reagent [35S] dithiobis(succinimidyl propionate). J Mol Biol 104:243-61.
- Staros, J.V., et al. (1986). Enhancement by N-hydroxysulfosuccinimide of water-soluble carbodiimide-mediated coupling reactions. Anal Biochem 156:220-2.
- 3. Cuatrecaseas, P. and Parikh, I. (1972). Adsorbents for affinity chromatography. Use of *N*-hydroxysulfosuccinimide esters of agarose. *Biochemistry* 11:291-9.
- Grabarek, Z. and Gergely, J. (1990). Zero-length crosslinking procedure with the use of active esters. Anal Biochem 185:131-5.

Product References

Boels, K., et al. (2001). The neuropeptide head activator induces activation and translocation of the growth-factor-regulated Ca²⁺-permeable channel GRC. *J Cell Sci* **114**:3599-606.

Byun, C.H., et al. (2001). Identification of the peptides that inhibit the function of human monoclonal thyroid-stimulating antibodies from phage-displayed peptide library. J Clin Endocrinol Metab 86:3311-18.

Epstein, A. L. (2003). Identification of a protein fragment of interleukin 2 responsible for vasopermeability. J Natl Cancer Inst 95:741-9.

Nogami, K., et al. (2004). Mechanisms of interactions of Factor X and Factor Xa with the acidic region in the Factor VIII A1 domain. J Biol Chem 279:3104-13.

Nyman, T., *et al.* (2002). A cross-linked profilin-actin heterodimer interferes with elongation at the fast-growing end of F-actin. *J Biol Chem* **277:**15828-33. Remmert, K., *et al.* (2004). CARMIL is a bona fide capping protein interactant. *J Biol Chem* **279:**3068-77.



Romaniouk, A.V., et al. (2004). Synthesis of a novel photoaffinity derivative of 1-deoxynojirimycin for active site-directed labeling of glucosidase I. Glycobiology 14:301-10.

Van Troys, M., et al. (2004). TetraThymosinbeta is required for actin dynamics in *Caenorhabditis elegans* and acts via functionally different actin-binding repeats. *Mol Biol Cell* **15**:4735-48.

Witt, S., et al. (2004). Comparative biochemical analysis suggests that vinculin and metavinculin cooperate in muscular adhesion sites. J Biol Chem 279:31533-43.

Products are warranted to operate or perform substantially in conformance with published Product specifications in effect at the time of sale, as set forth in the Product documentation, specifications and/or accompanying package inserts ("Documentation"). No claim of suitability for use in applications regulated by FDA is made. The warranty provided herein is valid only when used by properly trained individuals. Unless otherwise stated in the Documentation, this warranty is limited to one year from date of shipment when the Product is subjected to normal, proper and intended usage. This warranty does not extend to anyone other than Buyer. Any model or sample furnished to Buyer is merely illustrative of the general type and quality of goods and does not represent that any Product will conform to such model or sample.

NO OTHER WARRANTIES, EXPRESS OR IMPLIED, ARE GRANTED, INCLUDING WITHOUT LIMITATION, IMPLIED WARRANTIES OF MERCHANTABILITY, FITNESS FOR ANY PARTICULAR PURPOSE, OR NON INFRINGEMENT. BUYER'S EXCLUSIVE REMEDY FOR NON-CONFORMING PRODUCTS DURING THE WARRANTY PERIOD IS LIMITED TO REPAIR, REPLACEMENT OF OR REFUND FOR THE NON-CONFORMING PRODUCT(S) AT SELLER'S SOLE OPTION. THERE IS NO OBLIGATION TO REPAIR, REPLACE OR REFUND FOR PRODUCTS AS THE RESULT OF (I) ACCIDENT, DISASTER OR EVENT OF FORCE MAJEURE, (II) MISUSE, FAULT OR NEGLIGENCE OF OR BY BUYER, (III) USE OF THE PRODUCTS IN A MANNER FOR WHICH THEY WERE NOT DESIGNED, OR (IV) IMPROPER STORAGE AND HANDLING OF THE PRODUCTS.

Unless otherwise expressly stated on the Product or in the documentation accompanying the Product, the Product is intended for research only and is not to be used for any other purpose, including without limitation, unauthorized commercial uses, in vitro diagnostic uses, ex vivo or in vivo therapeutic uses, or any type of consumption by or application to humans or animals.

Current product instructions are available at www.thermoscientific.com/pierce. For a faxed copy, call 800-874-3723 or contact your local distributor. © 2013 Thermo Fisher Scientific Inc. All rights reserved. Unless otherwise indicated, all trademarks are property of Thermo Fisher Scientific Inc. and its subsidiaries. Printed in the USA.