Eutech EcoTestr pH 2 pocket tester

Instruction Manual

ECOPHTEST2
Revision Date: 13-11-2015

Language Index	
English	Page 2
Deutsch	Page 3
Français	Page 4
Nederlands	Page 5
Portuguese	Page 6
Español	Page 7

Waterproof EcoTestr pH 2 Microprocessor-based Pocket pH Tester

Note: DO NOT BE ALARMED if white crystals form around the cap, this is normal.

CONDITIONING before first use. Fill the cap with water, close the cap and soak the sensor for 1 hour to activate the sensor and dissolve the crystals. *Keep the tester upright during soaking, as the cap is not leak proof.*

CALIBRATION: This tester allows up to 3-point calibration. Start with pH 7 buffer. To cancel calibration, press **cal** to escape. *If no button is pressed for 15 seconds, tester will automatically confirm calibration.*

1. Press on/off button to switch the tester on	5. Press CAL, display will flash
2. Rinse the sensor with De-ionized water	6. Press HOLD/ENT to confirm
3. Put the sensor in pH7 buffer (2cm deep)	7. Repeat steps 2 to 6 for pH4 and/or pH10
4. Wait for the value to stabilize	

MEASUREMENT: Remove cap and press **on/off** to turn on the tester • Dip the sensor about 20 mm into the test solution • Stir once and let the display stabilize. Note the pH value • Press **hold/ent** to hold the reading. Press **hold/ent** again to release the held reading • Press **on/off** to shut off • Tester automatically shuts off after 8.5 minutes of non-use.

MAINTENANCE: Rinse the sensor in tap water after use. In harsh samples, take readings quickly, rinse sensor immediately afterwards with de-ionized water to remove residues and prevent sensor contamination. To maximize sensor life, fill the cap with 20 mm clean tap water (NOT DE-IONIZED WATER), and replace cap.

CHANGING BATTERIES: Lift up front battery cover and two sides of the pocket clip to remove battery cover. Remove old batteries and replace with fresh ones noting polarity as shown in battery compartment. You do not need to re-calibrate.

ERROR MESSAGES: **Lb** - Weak batteries. Need replacement. **Er** - Wrong or bad buffer value (out of range) or the sensor is failing. **Or / Ur** – Over and Under range signal or sensor is contaminated / damaged.

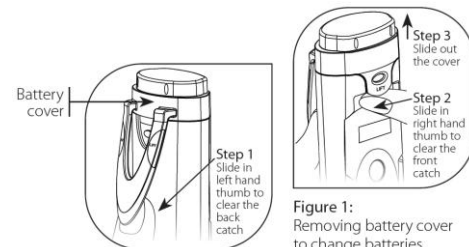
Figure 1:
Removing battery cover to change batteries

RESET option restores the calibration back to factory default settings • Turn off the tester • Press and hold **cal** button and then switch on. The display shows a blinking 'rS' (reset) • Press **hold/ent** to reset or **cal** to cancel reset.

WARRANTY: The tester has a warranty against manufacturing defects for ONE year from date of purchase.

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