

GeneArt™ Gibson Assembly® EX Cloning Kits

USER GUIDE

For highly-efficient, simultaneous, and seamless *in vitro* assembly of up to 15 DNA fragments plus a vector in a pre-determined order

for use with any of these products:

- GeneArt™ Gibson Assembly® EX Cloning Kit, Chemically Competent Cells (Cat. No. A46633)
- GeneArt™ Gibson Assembly® EX Cloning Kit, Electrocompetent Cells (Cat. No. A46634)
- GeneArt™ Gibson Assembly® EX Master Mix (Cat No. A46635, A46636, and A46637)

Catalog Numbers A46633, A46634, A46635, A46636, and A46637

Publication Number MAN0019063

Revision C.0



Life Technologies Corporation | 29851 Willow Creek | Eugene, OR 97402

For descriptions of symbols on product labels or product documents, go to [thermofisher.com/symbols-definition](https://www.thermofisher.com/symbols-definition).

The information in this guide is subject to change without notice.

DISCLAIMER: TO THE EXTENT ALLOWED BY LAW, THERMO FISHER SCIENTIFIC INC. AND/OR ITS AFFILIATE(S) WILL NOT BE LIABLE FOR SPECIAL, INCIDENTAL, INDIRECT, PUNITIVE, MULTIPLE, OR CONSEQUENTIAL DAMAGES IN CONNECTION WITH OR ARISING FROM THIS DOCUMENT, INCLUDING YOUR USE OF IT.

Revision history: Pub. No. MAN0019063

Revision	Date	Description
C.0	28 September 2020	Corrected the storage temp for Master mix A and B in contents and storage topic.
B.0	06 April 2020	Adding several edits.
A.0	13 February 2020	New document.

Important Licensing Information: This product may be covered by one or more Limited Use Label Licenses. By use of this product, you accept the terms and conditions of all applicable Limited Use Label Licenses.

TRADEMARKS: All trademarks are the property of Thermo Fisher Scientific and its subsidiaries unless otherwise specified. Gibson Assembly® is a registered trademark of SGI-DNA, Inc.

©2020 Thermo Fisher Scientific Inc. All rights reserved.

Contents

■ CHAPTER 1 Product information	5
Product description	5
Contents and storage	5
Required materials not supplied	6
Experimental outline	7
Workflow	8
■ CHAPTER 2 Methods	9
Prepare DNA inserts	9
DNA fragment design	9
Guidelines for designing PCR primers	9
PCR primers for attaching the DNA fragment to the vector	11
PCR primers for attaching adjacent DNA fragments	11
Guidelines for PCR	12
Generate a linearized <i>E. coli</i> cloning vector	12
Vector considerations	12
Guidelines for generating a linearized <i>E. coli</i> cloning vector	12
GeneArt™ Gibson Assembly® EX reaction	13
Transform One Shot™ TOP10 Chemically Competent <i>E. coli</i> cells	15
Materials required	15
Prepare reaction	15
Transform One Shot™ TOP10 Chemically Competent <i>E. coli</i> cells	16
Transform ElectroMAX™ DH10B electrocompetent <i>E. coli</i> cells	16
Materials required	16
Electroporate ElectroMAX™ DH10B cells	17
Analyze transformants	18
Analyze positive clones	18
Analyze transformants by colony PCR	19
Materials required	19
Long-term storage of positive clones	19
References	19

■	CHAPTER 3	Troubleshooting	20
■	APPENDIX A	Safety	22
		Chemical safety	23
		Biological hazard safety	24
		Documentation and support	25
		Customer and technical support	25
		Limited product warranty	25



Product information

Product description

The Gibson Assembly[®] method is an established DNA assembly reaction that allows multiple overlapping DNA fragments to be seamlessly linked in a single-tube reaction. DNA fragments of different lengths are uniformly assembled using complementary overlaps between fragments. The inherent flexibility of this approach is suitable for small and large DNA constructs and includes both single and multiple inserts. The resulting products can be used for a variety of downstream applications including transformation, PCR and rolling circle amplification (RCA).

Engineering DNA molecules by homologous *in vitro* recombination presents a seamless alternative to traditional methods using restriction endonucleases and ligases. GeneArt[™] Gibson Assembly[®] EX Master Mix contains a proprietary mixture of enzymes and reagents optimized to facilitate assembly of overlapping DNA fragments. Products assembled with GeneArt[™] Gibson Assembly[®] demonstrate low rates of junction error and high sequence fidelity.

Key features of the GeneArt[™] Gibson Assembly[®] EX Cloning Kits are:

- Accurate
- Seamless
- Ideal for complex assembly reactions
- Assemble up to 15 fragments

To seamlessly assemble up to 15 DNA fragments plus a vector into a single recombinant DNA molecule using the GeneArt[™] Gibson Assembly[®] EX Cloning Kits:

1. Combine your own linearized *E. coli* vector and the DNA fragments to assemble in a microcentrifuge tube.
2. Add appropriate amounts of the GeneArt[™] Gibson Assembly[®] EX Master Mix A and incubate the tube for in a thermocycler (chew back creates single-strand DNA 5' overhangs).
3. Slowly cool (anneal complementary overhangs).
4. Add appropriate amounts of the GeneArt[™] Gibson Assembly[®] EX Master Mix B and incubate the tube in a thermocycler (fill in gaps and seal nicks).
5. Transform the assembled DNA molecule into chemically or electrocompetent *E. coli* cells.

Contents and storage

Contents shown are for the individual kits including:

- GeneArt[™] Gibson Assembly[®] EX Cloning Kit, Chemically Competent Cells (Cat. No. A46633)
- GeneArt[™] Gibson Assembly[®] EX Cloning Kit, Electrocompetent Cells (Cat. No. A46634)
- GeneArt[™] Gibson Assembly[®] EX Master Mix (Cat. No. A46635, A46636, A46637)

Component	A46633 10 reactions	A46634 10 reactions	A46635 10 reactions	A46636 50 reactions	A46637 200 reactions	Storage Temp.
GeneArt™ Gibson Assembly® EX Master Mix A	50 µL	50 µL	50 µL	250 µL	1 mL	-70°C (or -20°C for up to 2 months)
GeneArt™ Gibson Assembly® EX Master Mix B	100 µL	100 µL	100 µL	500 µL	2 mL	
Positive Control	50 µL	50 µL	50 µL	50 µL	50 µL	-20°C
One Shot™ TOP10 Chemically Competent <i>E. coli</i>	11 × 50 µL	—	—	—	—	-80°C
ElectroMAX™ DH10B Cells	—	5 × 100 µL	—	—	—	-80°C
S.O.C. Medium	6 mL	2 × 6 mL	—	—	—	4°C

The Positive Control consists of a mixture of 10 ng of a 1.5 kb insert (kanamycin cassette) and 30 ng of a 2.7 kb vector containing an ampicillin resistance gene. Select for the 4.2 kb assembled construct on LB agar plates with 100 µg/mL ampicillin or 50 µg/mL kanamycin.

Required materials not supplied

- DNA fragments for the GeneArt™ Gibson Assembly® EX cloning reaction
- Thermocycler
- Luria-Bert (LB) plates with appropriate antibiotic

If a Master Mix kit is purchased, the following materials are also required:

- S.O.C. medium
- High efficiency competent cells. We recommend One Shot™ TOP10 Chemically Competent *E. coli* or ElectroMAX™ DH10B Cells from Thermo Fisher Scientific.

Additional optional materials

- DNA polymerase (e.g., SuperFi™ II DNA Polymerase, Thermo Fisher Scientific) for producing fragments to be assembled with the GeneArt™ Gibson Assembly® EX cloning kit
- Restriction enzymes for producing linearized vector
- PureLink™ PCR Purification Kit (Thermo Fisher Scientific) or GeneJET™ Gel Extraction Kit (Thermo Fisher Scientific)
- Thermoblock or electroporator plus electroporation cuvettes


Experimental outline

DNA fragments can be obtained from a variety of methods. Building a fragment de novo such as a GeneArt™ String™ or High-Q String™ DNA fragment allows direct design of the homologous region for Gibson Assembly® cloning. It is also possible to obtain a DNA fragment via amplification.

The table below describes the major steps required to assemble your recombinant DNA molecule using the GeneArt™ Gibson Assembly® EX cloning kit. Refer to the specified pages for details to perform each step.

Step	Action	Page
1	Design your primers to amplify your fragments	“DNA fragment design” on page 9
2	Amplify your DNA fragments with DNA Polymerase (SuperFi™ II mix)	“Guidelines for PCR” on page 12
3	Prepare a linearized vector (by using DNA Polymerase or restriction digestion)	“Guidelines for generating a linearized E. coli cloning vector” on page 12
4	Perform the GeneArt™ Gibson Assembly® EX cloning reaction	“Guidelines for GeneArt™ Gibson Assembly® EX cloning reactions” on page 13
5	Transform chemically competent One Shot™ TOP10 cells - or - electrocompetent ElectroMAX™ DH10B <i>E. coli</i> cells	“Transform One Shot™ TOP10 Chemically Competent E. coli cells” on page 15 or “Transform ElectroMAX™ DH10B electrocompetent E. coli cells” on page 16
6	Analyze colonies by colony PCR, restriction analysis and/or sequencing	“Analyze transformants” on page 18

Workflow

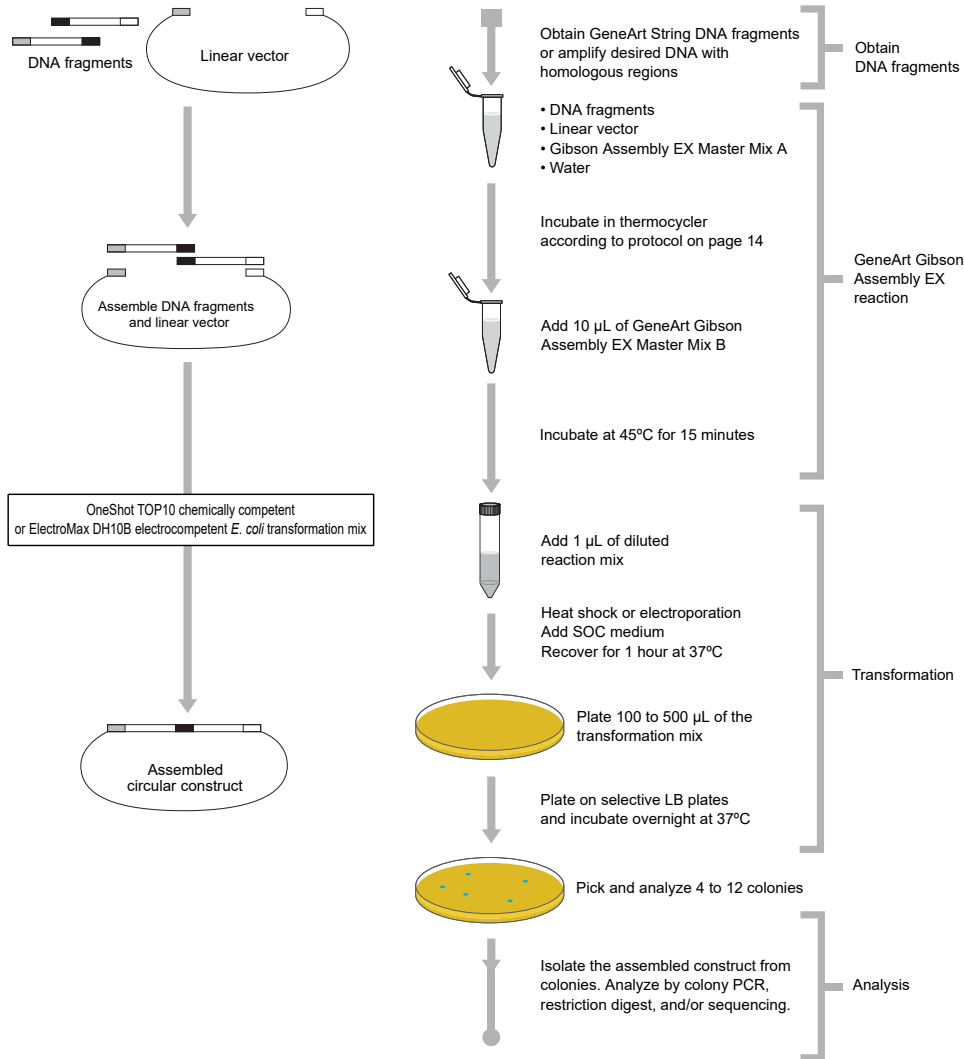

Prepare DNA inserts

DNA fragment design

GeneArt™ Gibson Assembly® technology relies on homologous recombination to assemble adjacent DNA fragments sharing end-terminal homology. The optimal length of the homologous fragment ends region depends on the number and length of the fragments in the assembly reaction. The following table lists the suggested length of the end-terminal homology.

Number of Fragments	Fragment Size	Length of Overlap Regions
1-15	≤ 8 kb	20-40 bp
	> 8 kb	40-100 bp

- For higher order assembly, longer overlap regions will result in higher efficiency.
- You may need to optimize PCR amplification reactions when using PCR primers with long homologous overlap regions.

The GeneArt™ Gibson Assembly® EX reaction requires that each DNA fragment (including the cloning vector) shares a 20-40 bp terminal homology with the adjacent fragment. When obtaining de novo DNA fragments such as GeneArt Strings™ DNA fragments, simply include homologies directly to the sequence design. If fragments are generated from existing DNA templates via PCR, primers used for generating your inserts must contain the 20-40 nt overhangs on their 5' ends to provide this homology with the adjacent fragments; however, this homology may be split between the primers used for adjacent PCR-amplified DNA fragments (see below).

Guidelines for designing PCR primers

- Design your PCR primers such that each DNA fragment to be assembled is between 0.5-5 kb in length.

Note: Large fragments (>5 kb) are more susceptible to damage in a gel extraction procedure. Furthermore, many PCR enzymes are not processive enough to amplify fragments >5 kb.

Therefore, we recommend that you assemble multiple fragments of ≤ 5 kb in one reaction rather than a single large fragment.

-
- The 5' ends of each primer pair (forward and reverse) must contain a 20-40 nt sequence (i.e., an overhang) that is homologous to the 20-40 bp at one end of the adjacent DNA fragment (i.e., the vector or another insert) while the 3' end of each primer must be specific to your DNA element you want to assemble.
 - If you are recombining the insert to the linearized vector, all 20-40 nt providing the requisite homology must be on the 5' end of the primer. To recombine two adjacent inserts, you may split the 20-40 bp homology between the fragments (e.g., 20 bp on the reverse primer of fragment 1 and 20 bp on the forward primer of fragment 2).

Note: You can split the 20-40 bp homology between adjacent fragments in any combination (e.g., 20+20 as in the example above or 10+30, 35+5, etc.).


-
- PCR primers should be up to 65 nt in length (20-40 nt to provide the requisite homology at the 5' end and 18-25 nt specific to your DNA element).

Note: If you are splitting the required 20-40 bp homology between adjacent fragments, one PCR primer may have as few as 18 nt and no overhang whatsoever. For example, in a 0+40 homology split, reverse primer of fragment 1 would not contribute to the homology and will consist of only 18-25 gene-specific nucleotides for fragment 1, while the forward primer of fragment 2 would contribute the entire 40 bp for the homology in addition to having 18-25 gene-specific nucleotides for fragment 2 (for a total length of 58-65 nt).

-
- We recommend splitting the homology between the primers to avoid long primers however.
 - You may also design your PCR primers to provide specific restriction enzyme sites or to reconstruct the restriction sites used for linearizing your cloning vector. Note that these sites will not be a part of the 20-40 nt homology.
 - Prepare each primer pair at a concentration of 100 μ M in DNase- and RNase- free water.


PCR primers for attaching the DNA fragment to the vector

The figure below shows examples of PCR primers used for preparing the DNA inserts for assembly in linear vectors with 5' overhangs, 3' overhangs, and blunt ends. The primer sequences providing the homology (in this example 25 nt) are shown in bold. Even though 5' and 3' overhangs produce slightly different lengths of homology to the amplified PCR product, we recommend to disregard the overhangs to the overall homology. The remaining sequences of each primer, i.e., the 3' end, are specific to the DNA element to assemble.

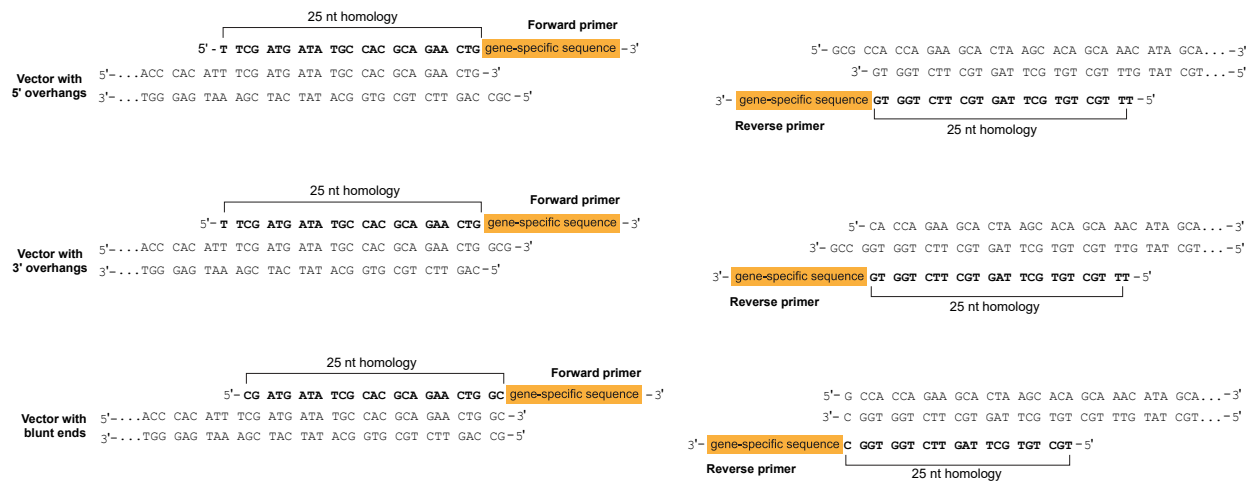

PCR primers for attaching adjacent DNA fragments

The figure below shows examples of PCR primers used for preparing adjacent DNA inserts for their assembly into a larger construct. When connecting two adjacent inserts, you may split the homology between the fragments in any combination. In the example below, the reverse primer of fragment 1 provide 12 nt and the forward primer of fragment 2 contribute 18 nt of a 30 bp homology for recombination. The primer sequences adding to the homology are shown in bold; the remaining sequences of each primer, i.e., the 3' end, are specific to the DNA element to assemble.

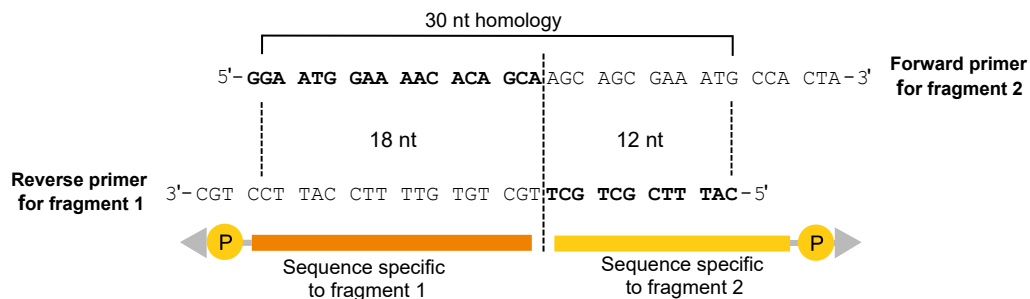

Note: While primers can be easily designed by hand, we recommend using appropriate software for primer design to avoid human error. Today, most molecular biology cloning programs support design of Gibson Assembly[®] experiments.

Guidelines for PCR

When using plasmid DNA, 5-20 ng of plasmid DNA is usually sufficient as a PCR template in a reaction volume of 100 μ L. When using *E. coli* or human genomic DNA, you can increase the amount of template DNA to 20-200 ng.

- We recommend using Platinum™ SuperFi II PCR Master Mix (Cat. No. 12368050) for achieving the best results during PCR amplification of your DNA fragments of interest.

Note: The processivity of the polymerase is crucial for obtaining full-length PCR products

- Use PCR cycling parameters that are appropriate for your template. Make sure to optimize PCR conditions to produce a single, discrete PCR product.
- PCR products may be used directly in assembly reactions without additional purification, although results may be improved by purifying the DNA fragments through a PCR cleanup kit. We recommend that you use the PureLink™ PCR Purification Kit (Cat. No. K310001).
- After preparing your DNA fragments by PCR, verify the PCR products by gel electrophoresis. If you obtain multiple bands, we recommend gel purification of your DNA fragments. We recommend using the GeneJET™ Gel Extraction Kit (Cat. No. K0691) for best results.

Generate a linearized *E. coli* cloning vector

The GeneArt™ Gibson Assembly® technology relies on homologous recombination to assemble adjacent DNA fragments sharing end-terminal homology. Because of the mechanism of action of the enzyme mix, the cloning vector and DNA fragments used with the GeneArt™ Gibson Assembly® Kit must be linear.

This section provides guidelines for generating a linearized *E. coli* cloning vector for use in the GeneArt™ Gibson Assembly® reaction.

Vector considerations

- For small inserts (<10 kb) or non-toxic gene(s), use a high copy number vector such as pUC19.
- For large insert(s) or toxic gene(s), use a low copy number or inducible vector.

Guidelines for generating a linearized *E. coli* cloning vector

- You can prepare the linearized *E. coli* cloning vector using restriction enzymes (single or double digest) or using PCR amplification.
- When generating the linearized vector by PCR amplification, we recommend treating the PCR amplification product with DpnI to reduce circular template carryover.
- When generating the linearized vector by restriction digest, we recommend that you digest the vector with two restriction enzymes rather than a single enzyme to reduce the amount of background. A double digest followed by PCR amplification of your linear vector virtually eliminates any background.
- You can use restriction enzymes that leave 3' protruding, 5' protruding, or blunt ends to linearize your cloning vector.

- It is very important to have a complete digest (i.e., very low background of uncut vector). Therefore, we recommend you increase the enzyme digestion time (2-3 hours to overnight) and the reaction volume.
- Analyze your restriction digestion products using agarose gel electrophoresis to verify that the digest is complete and then purify the digested vector using the PureLink™ PCR Purification Kit or equivalent.
- For a small scale digest, we recommend that you digest 2-5 µg of vector using 30-50 units of enzyme in a reaction volume of 100-200 µL.
- If you are planning on using the same linearized vector in multiple reactions, we recommend that you digest 20-50 µg of vector using 150-300 units of each restriction enzyme in a reaction volume of 400-800 µL. After purification, aliquot linearized vector and store at -20°C.

GeneArt™ Gibson Assembly® EX reaction

Before you set up your Gibson cloning and assembly reaction, be sure you have:

- Devised your DNA assembly strategy and possibly verified it by performing *in silico* cloning using an appropriate software tool.
- Prepared your DNA fragments (i.e., inserts) according to the guidelines in “Prepare DNA inserts” on page 9.
- Generated your linear cloning vector according to the guidelines in “Generate a linearized *E. coli* cloning vector” on page 12.

Materials required:

- DNA fragments to assemble
- Your own linearized *E. coli* cloning vector
- Positive Control
- GeneArt™ Gibson Assembly® EX Master Mix A
- GeneArt™ Gibson Assembly® EX Master Mix B
- Sterile deionized water

Guidelines for GeneArt™ Gibson Assembly® EX cloning reactions

- For maximum cloning efficiency we recommend using all fragments (vector and insert(s)) at equimolar ratio.
- Determine the concentration of your DNA insert solutions by OD₂₆₀ or fluorescence and use the concentrations to calculate the volume required to achieve the required molar ratio of insert to vector.
- Use the following formula to calculate molarities:

$$pmols \approx \frac{(weight\ in\ ng) \times 1000}{(fragment\ length\ in\ bp) \times 660}$$

- Keep GeneArt™ Gibson Assembly® EX Master Mix A and B on ice at all times and promptly return to freezer after use.
- When assembling multiple fragments, create a master mix of fragments in the proper ratios to minimize pipetting error.

GeneArt™ Gibson Assembly® EX cloning reaction

1. Thaw GeneArt™ Gibson Assembly® EX Master Mix A on ice.
2. Vortex GeneArt™ Gibson Assembly® EX Master Mix A immediately before use.
3. In a microcentrifuge tube on ice, set up the Gibson Assembly® EX cloning reaction as described in the table below:

	1-5 Inserts Assembly	Positive Control [1]
Recommended DNA molar ratio	vector:insert = 1:1	—
Amount of each fragment	0.04 pmol vector 0.04 pmol (each insert) X μ L	5 μ L
GeneArt™ Gibson Assembly® EX Master Mix A	5 μ L	5 μ L
Deionized water volume	(5 – X) μ L	—
Total volume	10 μ L	10 μ L

[1] The positive control reagents contain all necessary fragments.

(Optional) For the positive control, combine 5 μ L of the Positive Control and 5 μ L of GeneArt™ Gibson Assembly® EX Master Mix A in a tube on ice. Mix the reaction by vortexing.

Note: To assemble more than 5 inserts in a reaction, work to maintain a 1:1 molar ratio of vector to insert. Molar ratios of vector and insert can range from 0.01-0.04 pmol for each, but ideally the value is as close as possible to 0.04 pmol.

Note: For complex assemblies (e.g., > 5-15 fragments) we recommend using ElectroMAX™ DH10B electrocompetent cells for transformation.

4. Mix the reactions by vortexing and spin down.
5. Incubate in a thermocycler with the following conditions:

Temperature	For < 80bp overlaps	For \geq 80 bp overlaps
37°C	5 minutes	15 minutes
75°C	20 minutes	
Cool to 60°C at 0.1°C/second		
60°C	30 minutes	
Cool to 4°C at 0.1°C/second		

6. Thaw GeneArt™ Gibson Assembly® EX Master Mix B on ice. Vortex immediately before use and spin down.

7. Transfer the reactions from step 4 to ice and add 10 µL of GeneArt™ Gibson Assembly® EX Master Mix B. Mix by vortexing.

Note: If you increased the volume of the assembly reaction, also increase Master Mix B accordingly.

8. Incubate at 45°C for 15 minutes.
9. After incubation, place the reaction mix on ice and immediately proceed to the transformation step.

Note: Reactions can be stored at –20°C for later use.

Note: When using the GeneArt™ Gibson Assembly® EX Cloning Kits for the first time, we strongly recommend that you perform the GeneArt™ Gibson Assembly® positive control reactions in parallel with your samples to verify that the kit components are performing properly.

Transform One Shot™ TOP10 Chemically Competent *E. coli* cells

Materials required

- GeneArt™ Gibson Assembly® reaction mix
- One Shot™ TOP10 Chemically Competent *E. coli* cells
- S.O.C. medium
- 42°C heat block
- LB plates containing the appropriate selection antibiotic
- 37°C shaking and non-shaking incubator
- (Optional) pUC19 Control DNA for the transformation control reaction

Prepare reaction

1. Warm the heat block to 42°C.
2. Warm the vial of S.O.C. medium to room temperature.
3. Warm selective plates at 37°C for 30 minutes.
4. Thaw on ice 1 vial of One Shot™ Chemically Competent *E. coli* cells for each transformation.

Transform One Shot™ TOP10 Chemically Competent *E. coli* cells

1. Dilute the GeneArt™ Gibson Assembly® EX cloning reaction (“GeneArt™ Gibson Assembly® EX cloning reaction” on page 14) 1:5 in nuclease-free water (e.g., 12 µL nuclease-free water and 3 µL GeneArt™ Gibson Assembly® EX cloning reaction). Mix the reaction by vortexing, then spin down and keep the diluted reaction on ice.
2. Add 1 µL of the dilution into a vial of One Shot™ TOP10 Chemically Competent *E. coli* cells and mix gently.

IMPORTANT! Do not mix by pipetting up and down or vortexing.

(Optional) If you are performing a transformation control, add 2.5 µL of pUC19 Control DNA into a separate vial of One Shot™ TOP10 Chemically Competent *E. coli* cells and follow the transformation procedure.

3. Incubate the transformation mix on ice for 25 minutes.
4. Heat-shock the cells for 30 seconds at 42°C without shaking.
5. Immediately transfer the tubes to ice.
6. Add 450 µL of room temperature S.O.C. medium to the transformation mix.
7. Cap the tube tightly and shake at 300 rpm at 37°C for 1 hour to allow the cells to recover.
8. After incubation, spread 100 µL from each transformation on a pre-warmed selective plate. Before spreading on selective plates, mix suspension by pipetting up and down. If you have performed an assembly with ≥ 4 fragments, plate 100 µL on one selective plate and the remaining transformation on a second selective plate.
(Optional) For the positive control, plate 100 µL of the transformed reaction onto LB plates containing 100 µg/mL ampicillin or 50 µg/mL kanamycin.
9. Incubate the plates overnight at 37°C.
10. The next day, pick individual colonies and isolate the plasmid DNA or screen for the presence of the insert(s) by colony PCR, miniprep and restriction analysis or direct sequencing.

Transform ElectroMAX™ DH10B electrocompetent *E. coli* cells

Materials required

- Gibson Assembly® cloning reaction mix
- ElectroMAX™ DH10B electrocompetent *E. coli* cells
- S.O.C. medium
- Electroporation cuvettes
- Electroporator

- LB plates containing the appropriate selection antibiotic
- 37°C shaking and non-shaking incubator
- (Optional) pUC19 Control DNA for the transformation control reaction

Electroporate ElectroMAX™ DH10B cells

1. Chill microcentrifuge tubes and electroporation cuvettes on ice.
2. Thaw ElectroMAX™ DH10B cells on ice.
3. Mix the reactions by vortexing and spin down before proceeding with transformation.
4. Dilute the GeneArt™ Gibson Assembly® EX cloning reaction 1:5 in nuclease-free water (e.g., 12 µL nuclease-free water and 3 µL GeneArt™ Gibson Assembly® cloning reaction). Keep diluted reactions on ice.

Note: For assembly of large constructs, transformation efficiency could be reduced. In this case, we recommend to desalt the GeneArt™ Gibson Assembly® cloning reaction (e.g., by ethanol precipitation) and use ≥ 25% of the total reaction volume for electroporation.

5. When cells are thawed, mix cells by tapping gently. Pipet 20 µL of cells to each chilled microcentrifuge tube
6. Add 1 µL of the dilution, into the microcentrifuge tube containing ElectroMAX™ DH 10B, mix gently and return the tube to ice.

Note: If you are performing transformation control, add 1 µL of pUC19 Control DNA into a separate microcentrifuge tube containing ElectroMAX™ DH10B and follow the transformation procedure.

7. Incubate the cell/reaction mix on ice for one minute.
8. After the incubation, carefully pipette the cell/reaction mixture into a chilled 0.1 cm electroporation cuvette. Gently tap the cuvette to ensure that the cell/reaction mixture makes contact all the way across the bottom of the cuvette chamber. Avoid formation of bubbles.
9. Insert the cuvette into the electroporator and press the pulse button. If you are using the BTX™ ECM™ 630 or Bio-Rad™ Gene Pulser™ II electroporator, we recommend using the following electroporator conditions: 2.0 kV, 200 Ω, 25 µF. In case of arcing, use higher dilutions (1:10 to 1:50) or desalt the GeneArt™ Gibson Assembly® EX cloning reaction and proceed with step 6.
10. Add 800 µL of S.O.C. medium to the cuvette immediately after the end of the pulse and thoroughly pipet the mixture up and down. Transfer the solution back to the microcentrifuge tube. Repeat steps 7-8 for the remaining tubes.
11. Cap the tube tightly and shake at 300 rpm at 37°C for 1 hour to allow the cells to recover.

12. After incubation, spread 50-100 μL from each transformation on a pre-warmed selective plate. Before spreading on selective plates, mix suspension by pipetting up and down.
(Optional) For the positive control, plate 50-100 μL of the transformed reaction onto LB plates containing 100 $\mu\text{g}/\text{mL}$ ampicillin or 50 $\mu\text{g}/\text{mL}$ kanamycin.
13. Incubate the plates overnight at 37°C.
14. The next day, pick individual colonies and isolate the plasmid DNA or screen for the presence of the insert(s) by colony PCR, miniprep and restriction analysis, or direct sequencing.

Analyze transformants

Once you have performed the GeneArt™ Gibson Assembly® cloning reaction and the transformation procedure, screen for "positive" colonies containing your assembled recombinant DNA molecule by isolating the plasmid DNA and sequencing or by performing restriction analysis followed by agarose gel electrophoresis. You can also screen for the presence of the insert(s) by colony PCR. Calculate cloning efficiency using the following formula:

$$\text{Cloning efficiency (\%)} = \frac{(\text{number of positive colonies})}{(\text{number of total colonies})} \times 100$$

Typically, positive control yields cloning efficiencies >90%. Colony output is dependent on several factors, including transformation efficiency. Note that low colony output is not necessarily indicative of low cloning efficiency.

Analyze positive clones

1. Pick 4-10 colonies and culture them overnight in LB medium containing the appropriate selection antibiotic for your cloning vector.

Note: Pick at least 10 colonies for an assembly reaction involving ≥ 4 insert fragments plus the vector. For assembly reactions involving less than 4 insert fragments, 4-8 colonies should suffice.

2. Isolate plasmid DNA using your method of choice. If you need ultra-pure plasmid DNA for automated or manual sequencing, we recommend using PureLink™ HQ Mini Plasmid Purification or PureLink™ HiPure Plasmid Miniprep kits.
3. Analyze the plasmids by restriction analysis and/or by sequencing. We highly recommend that you perform sequence analysis when assembling fragments that were PCR amplified to rule out any errors made by the DNA polymerase or introduced by PCR primers during amplification.

Analyze transformants by colony PCR

1. For each sample, aliquot 19.2 μL PCR master mix into a PCR tube. Add 0.4 μL of each forward and reverse PCR primer.
2. Pick 4-10 colonies with a pipet tip and resuspend them individually in 20 μL of the PCR cocktail from step 1.
3. Using the same pipet tip, streak each colony on an LB plate containing the appropriate selection antibiotic to save for preparing glycerol stocks (“Long-term storage of positive clones” on page 19).
4. Incubate the reaction for 2-3 minutes at 94°C to lyse the cells and to inactivate the nucleases.
5. Amplify your samples for 20-30 cycles using the amplification conditions you have determined.
6. Store the reactions at 4°C.
7. Visualize the results by agarose gel electrophoresis.

Materials required

- Platinum™ Direct PCR Universal Master Mix
- Appropriate forward and reverse PCR primers (10 μM each)
- LB plates containing the appropriate selection antibiotic

Long-term storage of positive clones

After you have identified the correct clone, purify the colony and make a glycerol stock for long-term storage. Keep a DNA stock of your plasmid at -20°C .

1. Streak the original colony out on an LB agar plate containing the appropriate selection antibiotic for your cloning vector. Incubate the plate at 37°C overnight.
2. Isolate a single colony and inoculate with 1-2 mL of LB containing the appropriate selection antibiotic for your cloning vector.
3. Grow the cells until the culture reaches stationary phase ($\text{OD}_{600} = 1-2$).
4. Mix 0.85 mL of the culture with 0.15 mL of sterile glycerol and transfer the mix to a cryovial.
5. Store the glycerol stocks at -80°C .

References

- Gibson, D.G., *et al.* 2009. *Nat Methods*. 6:343-345.
Gibson, D.G., *et al.* 2010. *Science*. 329:52-56.

3

Troubleshooting

Observation	Possible cause	Recommended action
No colonies after transformation with DNA inserts and the transformation control did not work.	Low transformation efficiency.	Perform the transformation procedure exactly as described.
	Competent <i>E. coli</i> cells were handled incorrectly.	Competent <i>E. coli</i> cells are very fragile. Handle the cells gently and resuspend them by pipetting up and down gently.
		Do not vortex the competent <i>E. coli</i> cells.
		Do not freeze/thaw competent <i>E. coli</i> cells. They can only be thawed once without dramatic loss in competency.
	Store competent <i>E. coli</i> cells at -80°C .	
Insufficient amount of <i>E. coli</i> cells plated	Increase the amount of <i>E. coli</i> cells plated.	
Transformants plated on selective plates containing the wrong antibiotic.	Use the appropriate antibiotic for selection.	
No colonies after transformation with DNA inserts, but transformation with the control assembly reaction is successful.	PCR products were not pure enough.	Repeat PCR amplification and purify product using a different method of purification.
	DNA fragments do not share the required end-terminal homology.	Make sure that your DNA fragments and the linearized cloning vector share the required 20-40 bp end-terminal homology. Refer to "Guidelines for designing PCR primers" on page 9 for the requirements on PCR primer design.
	DNA fragment ends generated by PCR were damaged.	Employ extra caution to minimize any potential damage to the ends of your DNA fragments by leaving the gel on the gel tray when exposing to UV light, using low UV power, and minimizing the time the gel is exposed to UV light. Note that an additional isopropanol precipitation after gel purification might be required to obtain the best results.
	Incorrect amounts of DNA fragments and/or vector were used.	Make sure that you use the correct amounts of DNA fragments and/or vectors for cloning. For maximum cloning efficiency, use a 1:1 molar ratio of vector:insert(s).

Observation	Possible cause	Recommended action
No colonies after transformation with DNA inserts, but transformation with the control assembly reaction is successful. <i>(continued)</i>	The GeneArt™ Gibson Assembly® EX Master Mixes A and B were not handled correctly.	Quickly thaw the GeneArt™ Gibson Assembly® EX Master Mixes A and B on ice, and immediately return to –20°C after use. The enzyme mixes may be subjected to 20 freeze/thaw cycles without a loss in activity.
		Do not leave the GeneArt™ Gibson Assembly® EX Master Mixes A and B at room temperature or on ice for extended periods of time.
Large number of the transformants contain no insert.	Cloning vector was incompletely linearized.	It is crucial that your cloning vector is fully linearized and any uncut vector is removed prior to the cloning and assembly reaction. If necessary, recut your vector and gel purify.
	Plates were too old or contained incorrect antibiotic.	Make sure to use freshly prepared LB plates containing the selection antibiotic appropriate for your cloning vector.
	The GeneArt™ Gibson Assembly® EX cloning reaction was not set up correctly.	Make sure to set up the reaction according to the guidelines in “Guidelines for GeneArt™ Gibson Assembly® EX cloning reactions” on page 13.
	Incubation time was too short or too long.	Make sure you incubate the cloning and assembly reaction mix for 15 or 60 minutes, depending on the number of fragments according to the table in “Guidelines for GeneArt™ Gibson Assembly® EX cloning reactions” on page 13. After incubation, immediately proceed to transformation.
	Cloning and assembly reaction was not performed at the correct temperature.	Make sure to perform the seamless cloning and assembly reaction according to the procedure described in “Guidelines for GeneArt™ Gibson Assembly® EX cloning reactions” on page 13.
Large number of the transformants contain incorrect insert.	PCR products were not pure enough.	If your PCR product is not a single distinct band, then it may be necessary to gel purify the PCR product to ensure cloning of the correct insert.



Safety



WARNING! GENERAL SAFETY. Using this product in a manner not specified in the user documentation may result in personal injury or damage to the instrument or device. Ensure that anyone using this product has received instructions in general safety practices for laboratories and the safety information provided in this document.

- Before using an instrument or device, read and understand the safety information provided in the user documentation provided by the manufacturer of the instrument or device.
- Before handling chemicals, read and understand all applicable Safety Data Sheets (SDSs) and use appropriate personal protective equipment (gloves, gowns, eye protection, and so on). To obtain SDSs, see the “Documentation and Support” section in this document.

Chemical safety



WARNING! GENERAL CHEMICAL HANDLING. To minimize hazards, ensure laboratory personnel read and practice the general safety guidelines for chemical usage, storage, and waste provided below. Consult the relevant SDS for specific precautions and instructions:

- Read and understand the Safety Data Sheets (SDSs) provided by the chemical manufacturer before you store, handle, or work with any chemicals or hazardous materials. To obtain SDSs, see the "Documentation and Support" section in this document.
- Minimize contact with chemicals. Wear appropriate personal protective equipment when handling chemicals (for example, safety glasses, gloves, or protective clothing).
- Minimize the inhalation of chemicals. Do not leave chemical containers open. Use only with sufficient ventilation (for example, fume hood).
- Check regularly for chemical leaks or spills. If a leak or spill occurs, follow the manufacturer cleanup procedures as recommended in the SDS.
- Handle chemical wastes in a fume hood.
- Ensure use of primary and secondary waste containers. (A primary waste container holds the immediate waste. A secondary container contains spills or leaks from the primary container. Both containers must be compatible with the waste material and meet federal, state, and local requirements for container storage.)
- After emptying a waste container, seal it with the cap provided.
- Characterize (by analysis if needed) the waste generated by the particular applications, reagents, and substrates used in your laboratory.
- Ensure that the waste is stored, transferred, transported, and disposed of according to all local, state/provincial, and/or national regulations.
- **IMPORTANT!** Radioactive or biohazardous materials may require special handling, and disposal limitations may apply.



WARNING! HAZARDOUS WASTE (from instruments). Waste produced by the instrument is potentially hazardous. Follow the guidelines noted in the preceding General Chemical Handling warning.



WARNING! 4L Reagent and Waste Bottle Safety. Four-liter reagent and waste bottles can crack and leak. Each 4-liter bottle should be secured in a low-density polyethylene safety container with the cover fastened and the handles locked in the upright position.

Biological hazard safety



WARNING! Potential Biohazard. Depending on the samples used on this instrument, the surface may be considered a biohazard. Use appropriate decontamination methods when working with biohazards.



WARNING! BIOHAZARD. Biological samples such as tissues, body fluids, infectious agents, and blood of humans and other animals have the potential to transmit infectious diseases. Conduct all work in properly equipped facilities with the appropriate safety equipment (for example, physical containment devices). Safety equipment can also include items for personal protection, such as gloves, coats, gowns, shoe covers, boots, respirators, face shields, safety glasses, or goggles. Individuals should be trained according to applicable regulatory and company/ institution requirements before working with potentially biohazardous materials. Follow all applicable local, state/provincial, and/or national regulations. The following references provide general guidelines when handling biological samples in laboratory environment.

- U.S. Department of Health and Human Services, *Biosafety in Microbiological and Biomedical Laboratories (BMBL)*, 5th Edition, HHS Publication No. (CDC) 21-1112, Revised December 2009; found at:
<https://www.cdc.gov/labs/pdf/CDC-BiosafetymicrobiologicalBiomedicalLaboratories-2009-P.pdf>
- World Health Organization, *Laboratory Biosafety Manual*, 3rd Edition, WHO/CDS/CSR/LYO/2004.11; found at:
www.who.int/csr/resources/publications/biosafety/Biosafety7.pdf

Documentation and support

Customer and technical support

Visit [thermofisher.com/support](https://www.thermofisher.com/support) for the latest service and support information.

- Worldwide contact telephone numbers
- Product support information
 - Product FAQs
 - Software, patches, and updates
 - Training for many applications and instruments
- Order and web support
- Product documentation
 - User guides, manuals, and protocols
 - Certificates of Analysis
 - Safety Data Sheets (SDSs; also known as MSDSs)

Note: For SDSs for reagents and chemicals from other manufacturers, contact the manufacturer.

Limited product warranty

Life Technologies Corporation and/or its affiliate(s) warrant their products as set forth in the Life Technologies' General Terms and Conditions of Sale at www.thermofisher.com/us/en/home/global/terms-and-conditions.html. If you have any questions, please contact Life Technologies at www.thermofisher.com/support.

