



PRODUCT INFORMATION

Klenow Fragment

Pub. No. MAN0012882

Rev. Date 30 September 2016 (Rev. B.00)

#_

Lot: _ Expiry Date: _

Store at -20 °C

Components	#EP0051	#EP0052	#EP0054
Klenow Fragment	300 U 10 U/μL	1500 U 10 U/μL	300 U 2 U/μL
10X Reaction Buffer	1 mL	5 × 1 mL	1 mL

www.thermofisher.com

For Research Use Only. Not for use in diagnostic procedures.

Description

Klenow Fragment is the Large Fragment of DNA Polymerase I, *E.coli*. It exhibits 5'→3' polymerase activity and 3'→5' exonuclease (proofreading) activity, but lacks 5'→3' exonuclease activity of DNA Polymerase I.

Applications

- DNA blunting by fill-in of 5'-overhangs or removal of 3'-overhangs. (1), see protocols on back page.
- Random-primed DNA labeling (2-4).
- Labeling by fill-in 5'-overhangs of dsDNA.
- DNA sequencing by the Sanger method (5).
- Site-specific mutagenesis of DNA with synthetic oligonucleotides (6).
- Second strand synthesis of cDNA (7).

Source

E.coli cells with a cloned fragment of the *polA* gene.

Molecular Weight

68 kDa monomer.

Definition of Activity Unit

One unit of the enzyme catalyzes the incorporation of 10 nmol of deoxyribonucleotides into a polynucleotide fraction in 30 min at 37 °C.

Storage Buffer

The enzyme is supplied in: 25 mM Tris-HCl (pH 7.5), 0.1 mM EDTA, 1 mM DTT and 50% (v/v) glycerol.

10X Reaction Buffer

500 mM Tris-HCl (pH 8.0 at 25 °C), 50 mM MgCl₂, 10 mM DTT.

Inhibition and Inactivation

- Inhibitors: metal chelators, PP_i, P_i (at high concentrations) (8).
- Inactivated by heating at 75 °C for 10 min or by addition of EDTA.

Note

- Activity of Klenow Fragment in Thermo Scientific buffers (in comparison to activity in assay buffer):

Buffers	Activity, %
for restriction enzymes: Thermo Scientific™ FastDigest™, FastDigest™ Green, O, R, 1x Thermo Scientific™ Tango™, 2x Tango™, BamHI, EcoRI Ecl136II, PaeI, SacI, KpnI B G	100 50-75 25-50 20-50
for PCR buffers: Taq buffer with KCl, Taq buffer with (NH ₄) ₂ SO ₄ , Pfu buffer	100
RT buffers	100

CERTIFICATE OF ANALYSIS

Endodeoxyribonuclease Assay

No detectable degradation was observed after incubation of supercoiled plasmid DNA with Klenow Fragment.

Quality authorized by:

 Jurgita Zilinskiene

(continued on back page)

Protocol for DNA 3'-end labeling by fill-in of 5'-overhangs

1. Prepare the following reaction mixture:

Linear DNA	0.1-4 µg
10x reaction buffer for Klenow Fragment	2 µL
[α-³²P]-dNTP, ~15-30 TBq/mmol (400-800 Ci/mmol) or [α-³²P]-dNTP, ~110 TBq/mmol (3000 Ci/mmol)	0.74 MBq (20 µCi) 2.96 MBq (80 µCi)
3 dNTP Mix, 2 mM each (without a labeled dNTP)	2.5 µL (0.25 mM final concentration)
Klenow Fragment	0.1 µL (1 U)
Water, nuclease-free (#R0581)	to 20 µL
Total volume	20 µL

2. Incubate at 37 °C for 15 min.

3. Stop the reaction by heating at 75 °C for 10 min.

Note

This protocol is suitable for labeling of the following DNA markers, composed of DNA fragments with 5'-overhangs:

Lambda DNA EcoRI Marker, #SM0281

Lambda DNA HindIII Marker, #SM0101

Lambda DNA EcoRI/HindIII Marker, #SM0191

- The modified version of this protocol can be used for nonradioactive labeling of DNA markers. Substitute a part of dTTP with a modified nucleotide (e.g. Biotin-11-dUTP or Fluorescein-12-dUTP) at a molar ratio of 1:2.

Protocol for DNA Blunting by fill-in of 5'-overhangs or removal of 3'-overhangs

1. Prepare the following reaction mixture:

Linear DNA	10-15 µL (0.1-4 µg)
10X reaction buffer for Klenow Fragment	2 µL
dNTP Mix, 2mM each (#R0241)	0.5 µL (0.05 mM final concentration)
Klenow Fragment	0.1-0.5 µL (1-5 U)
Water, nuclease-free (#R0581)	to 20 µL
Total volume	20 µL

2. Mix thoroughly, spin briefly and incubate at 37 °C for 10 min.

3. Stop the reaction by heating at 75 °C for 10 min.

Note

The enzyme incorporates modified nucleotides (e.g. biotin-, digoxigenin-, fluorescently-labeled nucleotides).

References

1. Sambrook, J., Fritsch, E.F. and Maniatis, T. (1989). *Molecular Cloning: A Laboratory Manual*. (2nd ed.), 5.40-5.43. Cold Spring Harbor: Cold Spring Harbor Laboratory Press.
2. Feinberg, A.P., Vogelstein, B., A technique for radiolabeling DNA restriction endonucleases fragments to high specific activity, *Anal. Biochem.*, 132, 6-13, 1983.
3. Feinberg, A.P., Vogelstein, B., Addendum to: A technique for radiolabeling DNA restriction endonuclease fragments to high specific activity, *Anal. Biochem.*, 137, 266-267, 1984.
4. Yu, H., et al., Cyanine dye dUTP analogs for enzymatic labeling of DNA probes, *Nucleic Acids Res.*, 22, 3226-3232, 1994.
5. Sanger, F., et al., DNA sequencing with chain-terminating inhibitors, *Proc. Natl. Acad. Sci. USA*, 74, 5463-5467, 1977.
6. Wallace, R.B., et al., Directed deletion of a yeast transfer RNA intervening sequence, *Science*, 209, 1396-1400, 1980.
7. Rougeon, F., et al., Insertion of rabbit β -globin gene sequence into an *E.coli* plasmid, *Nucleic Acids Res.*, 2, 2365-2378, 1975.
8. Eun, H-M., *Enzymology Primer for Recombinant DNA Technology*, Academic Press, Inc., 1996.

LIMITED USE LABEL LICENSE: Internal Research and Development Use Only.

The purchase of this product conveys to the buyer the limited, non-exclusive, non-transferable right (without the right to resell, repackage, or further sublicense) to use this product for internal research and development purposes. No other license is granted to the buyer whether expressly, by implication, by estoppel or otherwise. In particular, the purchase of the product does not include or carry any right or license to use, develop, or otherwise exploit this product commercially and no rights are conveyed to the buyer to use the product or components of the product for purposes including but not limited to provision of services to a third party, generation of commercial databases or clinical diagnostics. This product is sold pursuant to authorization from Thermo Fisher Scientific and Thermo Fisher Scientific reserves all other rights. For information on purchasing a license for uses other than internal research and development purposes, please contact outlicensing@lifetech.com or Out Licensing, Life Technologies Inc., 5791 Van Allen Way, Carlsbad, California 92008.

PRODUCT USE LIMITATION

This product is developed, designed and sold exclusively *for research purposes and in vitro use only*. The product was not tested for use in diagnostics or for drug development, nor is it suitable for administration to humans or animals.

Please refer to www.thermofisher.com for Material Safety Data Sheet of the product.

© 2016 Thermo Fisher Scientific Inc. All rights reserved. All trademarks are the property of Thermo Fisher Scientific Inc. and its subsidiaries.