RNase Inhibitor (Cloned)

Catalog Number AM2682, AM2684

Pub. No. 4393874 Rev. B

Contents	Quantity	Storage conditions
RNase-Inhibitor (Cloned), 40 U/µL	Cat. no. AM2692: 2,500 Units	Store at –20°C. <i>Do not store in a frost-free freezer.</i>
	Cat. no. AM2694: 10,000 Units	

WARNING! Read the Safety Data Sheets (SDSs) and follow the handling instructions. Wear appropriate protective eyewear, clothing, and gloves. Safety Data Sheets (SDSs) are available from **www.lifetechnologies.com**/**support**.

Product description

RNase Inhibitor, a recombinant human protein produced in *E. coli* (Blackburn, 1979), is a potent inhibitor of neutral pancreatic ribonucleases, including RNases A, B, and C. The mode of inhibition is noncompetitive; the inhibitor tightly binds RNases in a 1:1 ratio. The enzyme has been shown to inactivate a variety of RNases that are present in many tissues and cell types. RNase Inhibitor does not inhibit RNase T1, RNase 1, RNase H, S1 Nuclease, or RNase from *Aspergillus*.

Source: Recombinant E. coli strain.

Unit (U) definition: One unit is the amount of protein required to inhibit the activity of 5 ng of RNase A by 50%. Unit assay conditions: 100 mM Tris-acetate (pH 6.5), 1 mM EDTA, 1 mM cyclic 2',3'-CMP, and RNase inhibitor. Addition of RNase A initiates the reaction. RNase inhibitor activity is measured by the inhibition of hydrolysis of cyclic 2',3'-CMP by RNase A.

Storage buffer (*not included***):** 20 mM HEPES-KOH (pH 7.6). 50 mM KCl, 5 mM DTT and 50% (v/v) glycerol. If opened frequently during storage, we recommend adding DTT to maintain the optimal 5 mM level.

Using RNase Inhibitor

Addition of RNase Inhibitor has been shown to be useful whenever the integrity of RNA must be maintained, such as in the preparation of cDNA by reverse transcription (deMartynoff, et. al., 1980), *in vitro* RNA transcription (Melton, et. al.1984), and *in vitro* protein synthesis (Scheele and Blackburn, 1979).

RNase Inhibitor requires a minimum of 1 mM DTT to maintain activity and requires a pH of 5–8, with maximal activity between

pH 7 and 8. Since the mode of inhibition is the formation of a 1:1 complex with RNases, avoid denaturation or oxidation of RNase Inhibitor, which would result in the release of active RNase. The half-life of RNase Inhibitor:RNase A binding is approximately 8 hr.

Guidelines for RNA transciption, protein translation, and cDNA synthesis

Add RNase Inhibitor to transcription, translation, and cDNA synthesis reactions at a final concentration of 1 U/ μ L. RNase Inhibitor requires a minimum of 1 mM DTT and is active over a broad pH range. Avoid denaturation of RNase Inhibitor by SDS, urea, etc., which could result in a release of active RNases.

Limited product warranty

Life Technologies Corporation and/or its affiliate(s) warrant their products as set forth in the Life Technologies' General Terms and Conditions of Sale found on Life Technologies' website at **www.lifetechnologies.com/termsandconditions**. If you have any questions, please contact Life Technologies at **www.lifetechnologies.com/support**.

References

Blackburn, P. 1979. Ribonuclease inhibitor from human placenta: rapid purification and assay. *J. Biol. Chem.* 254: 12484–12487.

deMartynoff, G., Pays, E. and Vassart, G. 1980. Synthesis of a full length DNA complementary to thyroglobulin 33 S messenger RNA. *B.B.R.C.* 93: 645–653.

Melton, D.A., Krieg, P.A., Rebagliati, M.R., Maniatis, T., Zinn, K. and Green, M.R. 1984. Efficient in vitro synthesis of biologically active RNA and RNA hybridization probes from plasmids containing a bacteriophage SP6 promoter. *Nucl. Acids Res.* 12, 70335–70356.

Scheele, G. and Blackburn, P. 1979. Role of mammalian RNase inhibitor in cell-free protein synthesis. *P.N.A.S. USA* 76: 4898–4902.

For Research Use Only. Not for use in diagnostic procedures.

The information in this guide is subject to change without notice.

DISCLAIMER: LIFE TECHNOLOGIES CORPORATION AND/OR ITS AFFILIATE(S) DISCLAIM ALL WARRANTIES WITH RESPECT TO THIS DOCUMENT, EXPRESSED OR IMPLIED, INCLUDING BUT NOT LIMITED TO THOSE OF MERCHANTABILITY, FITNESS FOR A PARTICULAR PURPOSE, OR NON-INFRINGEMENT. TO THE EXTENT ALLOWED BY LAW, IN NO EVENT SHALL LIFE TECHNOLOGIES AND/OR ITS AFFILIATE(S) BE LIABLE, WHETHER IN CONTRACT, TORT, WARRANTY, OR UNDER ANY STATUTE OR ON ANY OTHER BASIS FOR SPECIAL, INCIDENTAL, INDIRECT, PUNITIVE, MULTIPLE OR CONSEQUENTIAL DAMAGES IN CONNECTION WITH OR ARISING FROM THIS DOCUMENT, INCLUDING BUT NOT LIMITED TO THE USE THEREOF.

© 2013 Life Technologies Corporation. All rights reserved. The trademarks mentioned herein are the property of Life Technologies Corporation and/or its affiliate(s) or their respective owners.

For support visit lifetechnologies.com/support or email techsupport@lifetech.com



