# thermo scientific

## PRODUCT INFORMATION Thermo Scientific Phusion Blood Direct PCR Kit

Pub. No. MAN0016315

Rev. Date 25 November 2016 (Rev. A.00)

#\_

Lot:

## Store at -20 °C

#### Ordering information

Component	#F-547S 100 rxns	#F-547L 500 rxns
Phusion Blood II DNA Polymerase	40 µL	200 µL
2X Phusion Blood PCR Buffer (includes dNTPs and MgCl <sub>2</sub> )	1 .25 mL	5 × 1.25 ml
Universal control primer mix	40 µL	40 µL
50 mM EDTA (pH 8.0)	0.5 mL	0.5 mL
50 mM MgCl <sub>2</sub> solution	1.5 mL	1.5 mL
100 % DMSO	0.5 mL	0.5 mL

Expiry Date: \_\_\_\_

## Storage

Upon arrival, store the components at -20 °C.

## **CERTIFICATE OF ANALYSIS**

Performance in PCR is tested by the amplification of 7.5 kb fragment from human genomic DNA and 20 kb fragment from  $\lambda$  DNA.

Quality authorized by: Jurgita Zilinskiene

www.thermofisher.com

For Research Use Only. Not for use in diagnostic procedures.

## 1. Introduction

Thermo Scientific<sup>™</sup> Phusion<sup>™</sup> Blood Direct Kit is designed to perform PCR directly from whole blood with no prior DNA extraction or sample preparation. Blood stored at +4 °C or frozen, and preserved with EDTA, citrate or heparin are all suitable for this kit, as is blood dried onto commercially available cards such as Whatman 903<sup>™</sup> and FTA<sup>™</sup> Cards. Both genomic and exogenous target DNA can be amplified. The Phusion Blood Direct PCR Kit employs a modified Thermo Scientific<sup>™</sup> Phusion<sup>™</sup> Hot Start II High-Fidelity DNA Polymerase that exhibits extremely high resistance to inhibitors found in blood. The kit has been optimized to give excellent results from mammalian blood with amplicons up to 7.5 kb. The recommended blood concentration is 1-20%, although robust amplification with up to 40% blood can often be achieved with some optimization. This kit is recommended for end-point PCR protocols.

The Phusion Blood Direct PCR Kit includes a pair of universal control primers that are compatible with a number of mammalian and other vertebrate species (see Section 7).

## 2. Important Notes

- The recommended starting amount is 5% blood added directly to the reaction without further modification.
- Use 98 °C for denaturation.
- The annealing rules are different from many common DNA polymerases (such as *Taq* DNA polymerases).
   Read Section 5.2 carefully.
- For extension, use 15 s for amplicons ≤1 kb or 30 s/kb for amplicons >1 kb.
- After PCR, spin the reactions at 1 000 × g (about 4 000 rpm) for 1–3 minutes to pellet debris from blood.
- For blood concentrations greater than 10%, optimization of MgCl<sub>2</sub> concentration may be required (see 4.3).
- Phusion DNA Polymerases produce blunt end PCR products.

## 3. Guidelines for PCR

Carefully mix and spin down all tubes before opening to ensure homogeneity and improve recovery. The PCR setup can be performed at room temperature.

Table 1. Pipetting in	nstructions		
Component	20 µL rxn	50 µL rxn	Final conc.
H <sub>2</sub> O	Add to 20 $\mu L$	add to 50 µL	
2X Phusion Blood PCR Buffer	10 µL	25 µL	1X
Primer A	XμL	XμL	0.5 µM
Primer B	XμL	XμL	0.5 µM
Phusion Blood II DNA Polymerase	0.4 µL	1 µL	
Whole Blood	1μL	2.5 µL	(See 4.1)
Optional components for reactions optimization			
50 mM MgCl <sub>2</sub>	0.6 µL	1.5 µL	(See 4.3)
50 mM EDTA	0.5-1.0 µL	1.25-2.5 µL	(See 4.3)

Table 2.Recommended cycling protocol

1.0 uL

Cycle step	2-step protocol		3-step protocol		Cycles
Cycle step	Temp.	Time	Temp.	Time	Uycies
Lysis of cells (see 5.1)	98 °C	5 min	98 °C	5 min	1
Denaturation Annealing (see 5.2) Extension	98 °C -	1 s -	98 °C X °C	1 s 5 s	35-40
(see 5.3)	72 °C	15-30 s/kb	72 °C	15-30 s/kb	
Final Extension	72 °C 4 °C	1 min hold	72 °C 4 °C	1 min hold	1

2.5 uL

5% (See 4.4)

## 4. Notes about reaction components

## 4.1. Blood sample

DMSO

With the Phusion Blood Direct PCR Kit, it is possible to use a wide range of blood concentrations in the reaction (from 1% to 20%, or in some cases even up to 40%). The recommended starting point is 5%.

In general, if higher blood percentage (>10%) is used, higher reaction volume (up to 50  $\mu$ L) is recommended, and some optimization may be required (in particular the MgCl<sub>2</sub> concentration may need to be increased up to 4.5 mM final concentration). For small amplicons, whole blood up to 40% can be used (please note that with mouse blood 20% should not be exceeded). More information about blood volumes for specific species available at <u>www.thermofisher.com/directpcr</u>.

If blood dried onto a Whatman card is used, a 1 mm punch can be added directly to the reaction without pretreatment. The optimal reaction volume varies depending on the card type used:

- Whatman 903 Card: 10-50 μL
- Whatman FTA Elute Card: 25-50 μL
- Whatman FTA Gene Card: 50 μL

If larger punches or smaller volumes are desired for the FTA cards, the punches may need to be pretreated by washing with 20  $\mu$ L H<sub>2</sub>O at 50 °C for 3 minutes. After washing, the water can simply be removed and the PCR mix added directly to the rinsed punch.

## Non-mammalian blood

The Phusion Blood Direct PCR Kit has been successfully tested with a variety of mammalian species. In addition, good results have been obtained with several different bird species. For birds and other species with nucleated blood cells, it may be necessary to reduce the amount of blood in the PCR reaction.

## Blood storage

For long term storage, it is recommended to store blood at -20 °C or dried on Whatman FTA/903 Cards. For short term storage (less than 3 months), blood can be stored at 4 °C. EDTA, sodium citrate and heparin all work well with the Phusion Blood Direct PCR Kit.

## 4.2. Enzyme

Physion Blood II DNA Polymerase is a proofreading polymerase that contains a unique processivityenhancing domain, making this polymerase accurate and rapid. Phusion Blood II DNA Polymerase is a hot start polymerase utilizing a reversibly binding Affibody<sup>™</sup> protein<sup>1,2</sup>. Phusion Blood II DNA Polymerase possesses 5' $\rightarrow$ 3' DNA polymerase activity and 3' $\rightarrow$ 5' exonuclease activity. The error rate of Phusion Blood II DNA Polymerase is approximately 25-fold lower than that of Thermus aquaticus DNA polymerase when determined with a modified lacl-based method<sup>3</sup>. The recommended amount of Physion Blood II DNA Polymerase is optimized to work with most amplicons. If high blood volume is used, increasing the enzyme amount may improve the results. The volume per reaction can be increased up to 0.6 µL per 20 µL reaction, or up to 1.5 µL per 50 µL reaction. When cloning fragments amplified with Phusion Blood II DNA Polymerase, blunt end cloning is recommended. If TA cloning is required, it can be performed by adding A overhangs to the blunt PCR product with Tag DNA Polymerase, for example. However, before adding the overhangs it is very important to remove all Phusion Blood II DNA Polymerase by purifying the PCR product carefully. Any remaining Phusion Blood II DNA Polymerase will degrade the A overhangs, creating blunt ends again. A detailed protocol for TA cloning of PCR fragments amplified with any of the Phusion DNA Polymerases can be found on our website: www.thermofisher.com.

(continued on reverse page)

#### 4.3. MgCl<sub>2</sub> concentration

The 2X Phusion Blood PCR buffer included in the kit provides 3.0 mM MgCl<sub>2</sub> in the final reaction. Extensive testing has shown 3.0 mM MgCl<sub>2</sub> to be optimal for most PCR reactions using whole blood. There is a separate tube of 50 mM MgCl<sub>2</sub> (F-510MG) included in the kit for adjusting the MgCl<sub>2</sub> concentration (up to 4.5 mM if necessary) for reactions containing very high amounts of blood. It should be noted, though, that excess MgCl<sub>2</sub> may result in spurious PCR products.

If unspecific products are created with the default reaction buffer (providing 3.0 mM MgCl<sub>2</sub>), the effective MgCl<sub>2</sub> concentration can be decreased by adding the chelating agent EDTA, which is included in the kit. Typically, adding 0.5 to 1.0  $\mu$ L of 50 mM EDTA to a 20  $\mu$ L reaction is sufficient to eliminate non-specific products. Note that the optimal conditions will depend on the primers, the percentage of blood in the reaction and/or the type of card used, since anticoagulants and other chemicals impregnating the cards can alter the available Mg<sup>2+</sup> concentration.

#### 4.4. DMSO

DMSO has been found to improve PCR results for some amplicons (particularly GC-rich templates). With these amplicons it is recommended to add 1–5% DMSO to the reaction. Note that if high DMSO concentration is used, the annealing temperature must be decreased, as DMSO alters the melting point of the primers. It has been reported that 10% DMSO decreases the annealing temperature by 5.5-6.0 °C<sup>4</sup>.

Other PCR additives such as formamide, glycerol, and betaine are also compatible with Phusion Blood II DNA Polymerase.

#### 5. Notes about cycling conditions

#### 5.1. Lysis of cells and DNA denaturation

The initial 5-minute incubation step at 98 °C allows the lysis of leukocytes, making genomic DNA available for PCR. The DNA denaturation step can be very short. It is sufficient that the reaction mixture reaches the required 98 °C.

#### 5.2. Primer annealing

Note that the optimal annealing temperature for Phusion Blood II DNA Polymerase may differ significantly from that of *Taq*-based polymerases. Always use the Tm calculator and instructions on our website: <u>www.thermofisher.com/tmcalculator</u> to determine the Tm values of primers and optimal annealing temperature. As a basic rule, for primers >20 nt, anneal for 5 seconds at a Tm +3 °C of the lower Tm primer. For primers ≤20 nt, use an annealing temperature equal to the Tm of the lower Tm primer. If necessary, use a temperature gradient to find the optimal annealing temperature for each template-primer pair combination. The annealing gradient should extend up to the extension temperature (two-step PCR). Two-step cycling without an annealing step is recommended for high-Tm primer pairs (Tm at least 69–72 °C).

#### 5.3. Extension

The extension is performed at 72 °C. The optimal extension time varies depending on the amplicon length and the blood concentration in the reaction. The recommended extension time is 15 seconds for amplicons  $\leq$ 1 kb, and 30 s/kb for amplicons >1 kb. When higher blood concentrations are used, longer extension times may improve the results.

#### 6. Post-amplification analysis

After amplification, centrifuge the PCR reaction at 1 000 × g for 1–3 minutes to collect the supernatant for analysis. This step separates the various components of blood that might interfere with subsequent assays, e.g. gel electrophoresis. This is especially important when high blood concentrations are used, as there can be a substantial amount of cell debris, etc. in the tube after the PCR reaction.

If post-PCR enzyme treatment is performed (e.g. PCR-RFLP), it may be necessary to dilute the PCR product 2–4-fold in order to dilute the salts and other inhibitors from the PCR reaction.

## 7. Control reactions

### 7.1. Universal control primer mix

This component is a mix of degenerate primers in H<sub>2</sub>O that amplify a 237 bp fragment of mammalian genomic DNA. The amplified region is a highly conserved non-coding region upstream of the SOX21 gene,<sup>5</sup> and the primers are designed to amplify this region from a wide range of vertebrate species. Each primer concentration is  $25 \ \mu M$ .

#### Primer #1 (24-mer)

5'- AGCCCTTGGGGASTTGAATTGCTG -3' Melting point: 73.5 °C

#### Primer #2 (27-mer) 5'- GCACTCCAGAGGACAGCRGTGTCAATA -3'

Melting point: 72.2°C (R=A), 75.3 °C (R=G)

#### Table 3. Pipetting instructions for control reactions

Component	20 µL rxn	50 µL rxn
H <sub>2</sub> O	Add to 20 µL	add to 50 µL
2X Phusion Blood PCR Buffer	10 µL	25 µL
Universal Control Primer Mix	0.4 µL	1 µL
Phusion Blood II DNA Polymerase	0.4 µL	1 µL
Whole Blood (see 4.1)	Up to 1 µL	Up to 10 µL

#### Table 4. Cycling instructions for control reactions

Cycle step	Temp.	Time	Cycles
Lysis of cells	98 °C	5 min	1
Denaturation Annealing/Extension	98 °C 72 °C	1 s 15 s	35
Final Extension	72 °C 4 °C	1 min hold	1

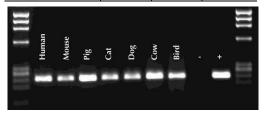


Figure 1. Amplification of the 237 bp control DNA fragment directly from whole blood (5 % in the reaction) of different species using universal control primers included in the Phusion Blood Direct PCR Kit. The blood samples used were stored at -20 °C. The + and - denote control reactions with and without purified DNA.

#### 8. Troubleshooting

o. mousicanooung	
No product at all or low yield	
<ul> <li>Repeat and make sure that there are no pipeting errors.</li> <li>Make sure that the cycling protocol was performed as recommended.</li> <li>Optimize annealing temperature.</li> <li>Titrate the amount of blood in the reaction.</li> <li>Increase the number of cycles.</li> <li>Check the purity and concentration of the primers.</li> </ul>	<ul> <li>Check primer design.</li> <li>Increase extension time.</li> <li>Increase denaturation time up to 5 seconds.</li> <li>Add 1-5 % DMSO (especially for high GC amplicons).</li> <li>Add MgCl<sub>2</sub> (1-2 µL/50 µL reaction) if high blood concentration is used.</li> <li>Add EDTA (1.25-2.5 µL/50 µL reaction).</li> </ul>
Non-specific products - High m	olecular weight smears
<ul> <li>Make sure that the extension time used was not too long. (Recommended extension time is 15–30 s/kb).</li> <li>Increase annealing temperature or perform a temperature gradient PCR.</li> </ul>	<ul> <li>Titrate the amount of blood in the reaction.</li> <li>Reduce the total number of cycles.</li> <li>Decrease primer concentration.</li> <li>Add EDTA (1.25-2.5 µL/50 µL reaction).</li> </ul>
Non-specific products - Low mol	ecular weight discrete bands
<ul> <li>Increase annealing temperature.</li> <li>Titrate the amount of blood in the reaction.</li> <li>Shorten extension time.</li> </ul>	<ul> <li>Perform a temperature gradient PCR.</li> <li>Decrease primer concentration.</li> <li>Add EDTA (1.25-2.5 µL/50 µL reaction).</li> <li>Design new primers.</li> </ul>

#### 9. References

- 1. Nord K. et al. (1997) Nature Biotechnol. 15: 772–777.
- 2. Wikman M. et al. (2004) Protein Eng. Des. Sel. 17: 455–562.
- 3. Frey B.F. & Suppmann B. (1995) *Biochemica* 2: 34–35.
- Chester N. & Marshak D.R. (1993) Analytical Biochemistry 209: 284–290.
- 5. Woolfe A. et al. (2005) PLoS Biology 3: 116-130.

#### NOTICE TO PURCHASER:

The purchase price of this product includes a limited, non-transferable license under U.S. and foreign patents owned by BIO-RAD Laboratories, Inc., to use this product. No other license under these patents is conveyed expressly or by implication to the purchaser by the purchase of this product.

## LIMITED USE LABEL LICENSE: Internal Research and Development Use Only.

The purchase of this product conveys to the buyer the limited, nonexclusive, non-transferable right (without the right to resell, repackage, or further sublicense) to use this product for internal research and development purposes. No other license is granted to the buyer whether expressly, by implication, by estoppel or otherwise. In particular, the purchase of the product does not include or carry any right or license to use, develop, or otherwise exploit this product commercially and no rights are conveyed to the buyer to use the product or components of the product for purposes including but not limited to provision of services to a third party, generation of commercial databases or clinical diagnostics. This product is sold pursuant to authorization from Thermo Fisher Scientific and Thermo Fisher Scientific reserves all other rights. For information on purchasing a license for uses other than internal research and development purposes, please contact outlicensing@lifetech.com or Out Licensing, Life Technologies Inc., 5781 Van Allen Way, Carlsbad, California 92008.

#### PRODUCT USE LIMITATION

This product is developed, designed and sold exclusively for research purposes and in vitro use only. The product was not tested for use in diagnostics or for drug development, nor is it suitable for administration to humans or animals. Please refer to <u>www.thermofisher.com</u> for Material Safety Data Sheet of the product.

© 2016 Thermo Fisher Scientific, Inc. All rights reserved. Affibody is a registered trademark of Affibody AB, Sweden. Whatman, FTA and 903 are trademarks of GE Healthcare and its subsidiaries. All other trademarks are the property of Thermo Fisher Scientific, Inc. and its subsidiaries.