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## Transferpipetten

<b>DE</b>	Transferpipetten – Gebrauchshinweis	2 - 5
<b>EN</b>	Transfer pipettes – Instructions for Use	6 - 9
<b>BG</b>	Пастърор пипети – Инструкции за употреба	10 - 13
<b>CS</b>	Transferové pipety – Návod k obsluze	14 - 17
<b>DA</b>	Transfer pipette – Brugsanvisning	18 - 21
<b>EL</b>	Πινέτες μεταφοράς – Οδηγός χρήσεως	22 - 25
<b>ES</b>	Pipetas de transferencia – Instrucciones de uso	26 - 29
<b>ET</b>	Ülekandepipett – Kasutusjuhend	30 - 33
<b>FI</b>	Pasteur pipetti – Käyttöohje	34 - 37
<b>FR</b>	Pipettes de transfert – Mode d'emploi	38 - 41
<b>HR</b>	Transfer pipete – Uputa za upotrebu	42 - 45
<b>HU</b>	Transzferpipetták – Használati utasítás	46 - 49
<b>IT</b>	Pipette di trasferimento – Istruzioni d'uso	50 - 53
<b>KO</b>	이송 피펫 – 사용 설명서	54 - 57
<b>LT</b>	Perkėlimo pipetės – Naudojimo instrukcijos	58 - 61
<b>LV</b>	Pastēra pipetes – Lietošanas norādes	62 - 65
<b>NL</b>	Transferpipet – Gebruiksaanwijzing	66 - 69
<b>NO</b>	Transferpipette – Bruksanvisning	70 - 73
<b>PL</b>	Pipety transferowe – Instrukcja obsługi	74 - 77
<b>PT</b>	Pipetas de transferência – Instruções de utilização	78 - 81
<b>RO</b>	Pipete de transfer – Instrucțiuni de utilizare	82 - 85
<b>RU</b>	Пастеровские пипетки – Инструкция по применению	86 - 89
<b>SK</b>	Prenosové pipety – Návod na Použitie	90 - 93
<b>SL</b>	Pipete za prenos – Navodila za uporabo	94 - 97
<b>SV</b>	Transferpipetter – Bruksanvisning	98 - 101
<b>TH</b>	ปีเปตที่แบบปริมาตร – คำแนะนำในการใช้งาน	102 - 105
<b>TR</b>	Transfer pipetleri – Kullanım için talimatlar	106 - 109
<b>ZH</b>	移液器 – 使用说明	110 - 113

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EN: Instructions for Use

## Intended use

Transfer pipettes are used for transferring diagnostic sample material from one vessel to another. This allows fluids such as urine, blood, serum and plasma to be transferred for subsequent analysis.

Transfer pipettes are used when it is not necessary for the sample volume to be exact. If exact volumes are needed, piston pipettes with suitable pipette tips should be used.

No special ambient conditions are required.

Transfer pipettes should be used by medical professionals.

## Product description

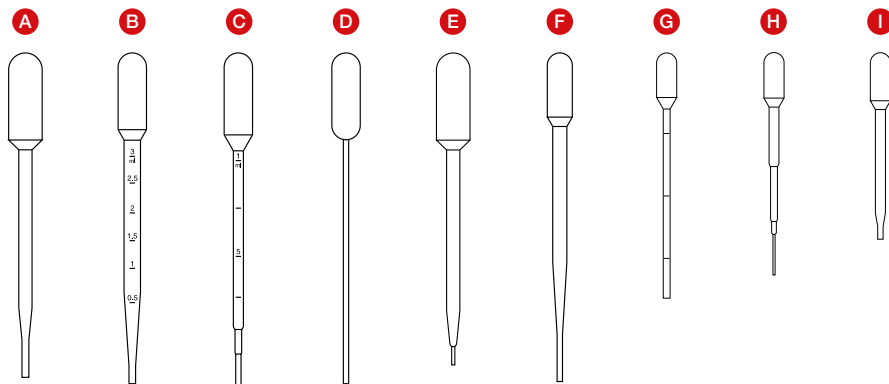
Transfer pipettes with an integrated suction bulb are available in nine different versions (see figure 1 on the next page).

The different versions (A-I) vary in terms of suction volume, droplet size, length and graduation. In addition, sterile and non-sterile variants are available, as well as different packaging options.

Sterile transfer pipettes are used, for example, for transferring sample material for microbial analyses. Here, there is a choice between individual sterile packs and sterile packs of 10 transfer pipettes.

SARSTEDT

Type	Art. no.	Name
<b>A</b>	86.1174.XXX	6 ml LD-PE transfer pipette, without graduation, Length: 152 mm, droplet size: 40 - 55 $\mu$ l
<b>B</b>	86.1171.XXX	3.5 ml LD-PE transfer pipette, with graduation 3: 0.50 ml (0.5 - 3.0 ml), Length: 155 mm, droplet size: 35 - 55 $\mu$ l
<b>C</b>	86.1172.XXX	3.5 ml LD-PE transfer pipette, with graduation 1: 0.25 ml (0.25 - 1.0 ml), Length: 156 mm, droplet size: 30 - 45 $\mu$ l
<b>D</b>	86.1173.XXX	3.5 ml LD-PE transfer pipette, without graduation, Length: 155 mm, droplet size: 30 - 45 $\mu$ l
<b>E</b>	86.1175.XXX	6 ml LD-PE transfer pipette, without graduation, Length: 146 mm, droplet size: 20 - 30 $\mu$ l
<b>F</b>	86.1176.XXX	2 ml LD-PE transfer pipette, without graduation, Length: 154 mm, droplet size: 35 - 45 $\mu$ l
<b>G</b>	86.1170.XXX	1 ml LD-PE transfer pipette, with graduation 1: 0.10 ml (0.1 - 0.3 ml), Length: 115 mm, droplet size: 45 - 55 $\mu$ l
<b>H</b>	86.1180.XXX	1 ml LD-PE transfer pipette, without graduation, Length: 104 mm, droplet size: 17 - 22 $\mu$ l
<b>I</b>	86.1179.XXX	1 ml LD-PE transfer pipette, without graduation, Length: 87 mm, droplet size: 30 - 45 $\mu$ l

LD-PE = low density polyethylene

## Handling

1. For blister pack variants: Open the blister at the opening tab marked with corrugations. This tab is located on the side of the suction bulb of the transfer pipette.  
The transfer pipette can be grasped by the suction bulb and removed from the blister pack.
2. Press in the suction bulb of the transfer pipette.
3. Dip the tip of the transfer pipette into the fluid you intend to transfer.
4. Release the suction bulb, allowing the vacuum created to fill the transfer pipette with fluid.
5. Transfer the tip of the transfer pipette into the target vessel and press the suction bulb again to release the fluid.
6. Dispose of the transfer pipette in accordance with the regulations of your facility.

## Restriction

- The accuracy of the droplet size and the graduation is  $\pm 15\%$ . Transfer pipettes are therefore intended for transferring undefined fluid quantities. For pipetting exact volumes, a piston pipette with a suitable pipette tip should be used.

## Disposal

- Wearing disposable gloves and other personal protective equipment reduces the risk of infection. Follow the regulations of your facility.
- General hygiene guidelines and legal regulations for the proper disposal of infectious material should be observed and complied with.
- Contaminated transfer pipettes must be disposed of in appropriate biohazard disposal containers. Follow the regulations of your facility.

## Safety information/warnings

- Personal protective equipment must be used when handling diagnostic samples and hazardous materials.
- Do not use transfer pipettes after the expiry date has passed
- The handling of biological samples must take place in accordance with the applicable hygiene and safety regulations.
- Transfer pipettes are intended for single use. After use they should be disposed of according to the applicable regulations. There is a risk of contamination if reused.

All serious incidents relating to the product shall be notified to the manufacturer and the competent national authority.

 SARSTEDT AG & Co. KG  
Sarstedtstr. 1  
D-51588 Nümbrecht  
[www.sarstedt.com](http://www.sarstedt.com)

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