# RiboLock RNase Inhibitor

Catalog Number EO0381, EO0382, EO0384

Pub. No. MAN0012010 Rev. C.00

 $\dot{\mathbb{N}}$ 

**WARNING!** Read the Safety Data Sheets (SDSs) and follow the handling instructions. Wear appropriate protective eyewear, clothing, and gloves. Safety Data Sheets (SDSs) are available from **thermofisher.com/support**.

# **Product description**

Thermo Scientific RiboLock RNase Inhibitor inhibits the activity of RNases A, B and C by binding them in a noncompetitive mode at a 1:1 ratio. It does not inhibit eukaryotic RNases: T1, T2, U1, U2, CL3 as well as prokaryotic RNases I and H.

#### Contents and storage

Cat. No.	Contents	Source	Molecular Weight	Amount	Storage
EO0381	RiboLock RNase Inhibitor	E.coli cells with a cloned gene encoding a mammalian	49.6 kDa	2500 U, 40 U/μL	-25 °C to -15 °C
EO0382				4 x 2500 U, 40 U/µL	
EO0384		ribonuclease inhibitor.	monomer	24 x 2500 U, 40 U/µL	

# **Applications**

- Inhibition of RNA degradation in the following:
  - in vitro transcription (1);
  - cDNA synthesis (2);
  - in vitro translation (3);
  - isolation of mammalian cell fractions that contain mRNA-protein complex (3);
  - RNA amplification (4).
- RNA purification and storage.
- Separation and identification of specific ribonuclease activities (5).
- Studies of tumor suppression (6).

## **Definition of Activity Unit**

One unit of the RiboLock™ RNase Inhibitor inhibits the activity of 5 ng of RNase A by 50 %.

Inhibitor activity is assayed in the following mixture:

100 mM Tris-HCI (pH 7.5), 1.2 mM EDTA, 0.1 mg/mL BSA, 100 ng/mL RNase A, 0.1 mg/mL [3H]-RNA, 50 mg/mL yeast RNA, 8 mM DTT.

#### Storage Buffer

The protein is supplied in: 20 mM HEPES-NaOH (pH 7.5), 50 mM NaCl, 8 mM DTT, 0.03 % (v/v) ELUGENT Detergent and 50% (v/v) glycerol.

#### Inhibition and Inactivation

- Inhibitors: common denaturants (SDS, urea and all oxidizing reagents (p-chloromercuribenzoate, dissolved oxygen, ions in their higher oxidation states) strongly inhibit RiboLock RNase Inhibitor and release the RNase bound.
- Inactivated by heating at 75 °C for 10 min. Residual activity detectable after 10 min heating at 70 °C.

### Note

- DTT provided in the Storage Buffer ensures stability during long term storage, but is not necessary for inhibitor activity.
- Recommended concentration 1 U/µL of a reaction mixture.



#### References

- 1. Nielsen, D.A., Shapiro, D.J., Preparation of capped RNA transcripts using T7 RNA polymerase, Nucleic Acids Res., 14, 5936, 1986.
- 2. Martynoff, G., et al., Synthesis of a full length DNA complementary to thyroglobulin 33S messenger RNA, Biochem. Biophys. Res. Commun., 93, 645-653, 1980.
- 3. Scheele, G., Blackburn, P., Role of mammalian RNase inhibitor in cell-free protein synthesis, Proc. Natl. Acad. Sci. USA, 76, 1898-1902. 1979.
- 4. Van Gelder., et al., Amplified RNA synthesized from limited quantities of heterogeneous cDNA. Proc. Natl. Acad. Sci. USA, 87, 1663-1667, 1990.
- 5. Eichler, D.C., et al., Effect of human placental ribonuclease inhibitor in cell-free ribosomal RNA synthesis, Biochem. Biophys. Res. Commun., 101, 396-403, 1981.
- 6. Polakowski, I.J., et al., A ribonuclease inhibitor expresses anti-angiogenic properties and leads to reduced tumor growth in mice, Amer. J. Pathol., 143, 507-517, 1993.

### Limited product warranty

Life Technologies Corporation and/or it affiliate(s) warrant their products as set forth in the Life Technologies' General Terms and Conditions of Sale at www.thermofisher.com/us/en/home/global/terms-and-conditions.html. If you have any questions, please contact Life Technologies at www.thermofisher.com/support.



Thermo Fisher Scientific Baltics UAB | V.A. Graiciuno 8, LT-02241 Vilnius, Lithuania For descriptions of symbols on product labels or product documents, go to thermofisher.com/symbols-definition.

The information in this guide is subject to change without notice.

DISCLAIMER: TO THE EXTENT ALLOWED BY LAW, THERMO FISHER SCIENTIFIC INC. AND/OR ITS AFFILIATE(S) WILL NOT BE LIABLE FOR SPECIAL, INCIDENTAL, INDIRECT, PUNITIVE, MULTIPLE, OR CONSEQUENTIAL DAMAGES IN CONNECTION WITH OR ARISING FROM THIS DOCUMENT, INCLUDING YOUR USE OF IT.

Important Licensing Information: These products may be covered by one or more Limited Use Label Licenses. By use of this product, you accept the terms and conditions of all applicable Limited Use Label Licenses.

©2019 Thermo Fisher Scientific Inc. All rights reserved. All trademarks are the property of Thermo Fisher Scientific and its subsidiaries unless otherwise specified.

