

# Ion 16S<sup>™</sup> Metagenomics Kit

### Catalog Number A26216

Pub. No. MAN0010801 Rev. A.0

Component	Part no. <sup>[1]</sup>	Cap color	Quantity	Volume	Storage conditions
2X Environmental Master Mix	4401975	Orange	2 tubes	0.8 mL per tube	Shipped at -20°C.
DNA Dilution Buffer	4405587	Clear	1 bottle	7.0 mL	After first use, store at 2–8°C if used frequently, otherwise store at –20°C.
					Protect the 2X Environmental Master Mix from light.
16S Primer Set V2–4–8 (10X)	100026495	Green	1 tube	300 µL	Store at –15 to –25°C.
16S Primer Set V3–6, 7–9 (10X)	100026496	Blue	1 tube	300 µL	
Negative Control	362250	White	1 tube	1.0 mL	
<i>E. coli</i> DNA control (30 μg/mL)	4458450	Red	1 tube	40 µL	Store at –15 to –25°C.
					Store separately from other reagents to prevent cross- contamination.

[1] Provided for identification purposes; the parts cannot be ordered separately from the kit.

**Note:** For safety and biohazard guidelines, refer to the "Safety" appendix in the *Ion* 165<sup>™</sup> *Metagenomics Kit User Guide* (Pub. no. MAN0010799). Read the Safety Data Sheets (SDSs) and follow the handling instructions. Wear appropriate protective eyewear, clothing, and gloves.

# **Product description**

The Ion 16S<sup>™</sup> Metagenomics Kit is designed for rapid analyses of polybacterial samples using Ion Torrent<sup>™</sup> sequencing technology. The kit includes two primer sets that selectively amplify the corresponding hypervariable regions of the 16S region in bacteria:

- Primer set V2-4-8
- Primer set V3-6, 7-9

Sequence the amplified fragments on the Ion PGM<sup>™</sup> System and analyze the results with the Ion Reporter<sup>™</sup> software Ion 16S<sup>™</sup> Metagenomics Kit analyses module. The combination of the two primer pools allows for sequence-based identification of a broad range of bacteria within a mixed population.

## Limited product warranty

Life Technologies Corporation and/or its affiliate(s) warrant their products as set forth in the Life Technologies' General Terms and Conditions of Sale found on Life Technologies' website at **www.lifetechnologies.com/termsandconditions**. If you have any questions, please contact Life Technologies at **www.lifetechnologies.com/support**.

The information in this guide is subject to change without notice.

#### DISCLAIMER

LIFE TECHNOLOGIES CORPORATION AND/OR ITS AFFILIATE(S) DISCLAIM ALL WARRANTIES WITH RESPECT TO THIS DOCUMENT, EXPRESSED OR IMPLIED, INCLUDING BUT NOT LIMITED TO THOSE OF MERCHANTABILITY, FITNESS FOR A PARTICULAR PURPOSE, OR NON-INFRINGEMENT. TO THE EXTENT ALLOWED BY LAW, IN NO EVENT SHALL LIFE TECHNOLOGIES AND/OR ITS AFFILIATE(S) BE LIABLE, WHETHER IN CONTRACT, TORT, WARRANTY, OR UNDER ANY STATUTE OR ON ANY OTHER BASIS FOR SPECIAL, INCIDENTAL, INDIRECT, PUNITIVE, MULTIPLE OR CONSEQUENTIAL DAMAGES IN CONNECTION WITH OR ARISING FROM THIS DOCUMENT, INCLUDING BUT NOT LIMITED TO THE USE THEREOF.

Important Licensing Information: These products may be covered by one or more Limited Use Label Licenses. By use of these products, you accept the terms and conditions of all applicable Limited Use Label Licenses.

©2014 Thermo Fisher Scientific Inc. All rights reserved. All trademarks are the property of Thermo Fisher Scientific and its subsidiaries unless otherwise specified.

For support visit lifetechnologies.com/support or email techsupport@lifetech.com

lifetechnologies.com

