

# AccuPrime™ *Pfx* DNA Polymerase

Catalog Numbers 12344-024 and 12344-032

Doc. Part No. 12344.pps Pub. No. MAN0001079 Rev. A.0



**WARNING!** Read the Safety Data Sheets (SDSs) and follow the handling instructions. Wear appropriate protective eyewear, clothing, and gloves. Safety Data Sheets (SDSs) are available from [thermofisher.com/support](https://www.thermofisher.com/support).

---

## Product description

The Invitrogen™ AccuPrime™ *Pfx* DNA Polymerase is a proprietary enzyme preparation containing recombinant DNA polymerase from *Thermococcus* species *KOD*. This polymerase possesses a proofreading 3' to 5' exonuclease activity that provides higher fidelity than *Pfu* DNA polymerase. The AccuPrime™ *Pfx* DNA Polymerase is a highly processive enzyme and possesses a fast chain extension capability. The 10X AccuPrime™ *Pfx* Reaction Mix contains MgSO<sub>4</sub>, dNTPs, and thermostable AccuPrime™ proteins; the thermostable proteins enhance specific primer-template hybridization during every cycle of PCR.

The AccuPrime™ *Pfx* DNA Polymerase is provided in an antibody-bound form that is inactive at room temperatures. The enzyme regains activity after the initial denaturation step at 94°C in PCR cycling, providing an automatic “hot start”. The hot start increases specificity, sensitivity, and yield, while allowing room-temperature assembly.

The high specificity, fidelity, and yield offered by AccuPrime™ *Pfx* DNA Polymerase make it ideal for demanding PCR applications such as site-directed mutagenesis and PCR expression cloning.

## Contents and storage

Contents	Cat. No. 12344-024 (200 reactions)	Cat. No. 12344-032 (1,000 reactions)	Storage
AccuPrime™ <i>Pfx</i> DNA Polymerase (2.5 U/μL)	100 μL	500 μL	-20°C
50 mM Magnesium sulfate	1 mL	2 × 1 mL	
10X AccuPrime™ <i>Pfx</i> Reaction Mix	1 mL	5 × 1 mL	

**Note: Unit (U) definition:** One unit of AccuPrime™ *Pfx* DNA Polymerase incorporates 10 nmol of deoxyribonucleotide into acid-insoluble material in 30 minutes at 74°C.

### Storage buffer

- 50 mM Tris-HCl (pH 8.0)
- 50 mM KCl
- 1 mM DTT
- 0.1 mM EDTA
- Stabilizers
- 50% (v/v) glycerol

### Procedural guidelines

- Take appropriate precautions to avoid cross-contamination.
- MgSO<sub>4</sub> is included in the 10X AccuPrime™ *Pfx* Reaction Mix at a final concentration of 1 mM, which is sufficient for most templates. For further optimization, add 0.1–1.0 μL of 50 mM MgSO<sub>4</sub> (included in the kit) to the reaction.
- dNTPs are included in the 10X AccuPrime™ *Pfx* Reaction Mix at a final concentration of 0.3 mM.
- Use an annealing temperature that is 5–10°C lower than the T<sub>m</sub> of the primers. If needed, gradually increase the annealing temperature by 2–3°C for higher specificity.
- For difficult primer sets, prepare titrations of KCl (not included) at final concentrations of 20–50 mM for further optimization.

## Perform the PCR

The following general procedure is suggested as a starting point when using AccuPrime™ *Pfx* DNA Polymerase in any PCR amplification.

1. Add the following components to an autoclaved microcentrifuge tube at room temperature or on ice:

Component	Volume for one 50- $\mu$ L reaction	Final concentration
10X AccuPrime™ <i>Pfx</i> Reaction Mix <sup>[1]</sup>	5 $\mu$ L	1X
Primer mix (10 $\mu$ M each) <sup>[1]</sup>	1.5 $\mu$ L	0.3 $\mu$ M each
Template DNA (10 pg to 200 ng)	$\geq$ 1 $\mu$ L	as required
AccuPrime™ <i>Pfx</i> DNA Polymerase <sup>[2]</sup>	0.4–1 $\mu$ L	1.0–2.5 U
Autoclaved, distilled water	to 50 $\mu$ L	—

<sup>[1]</sup> AccuPrime™ *Pfx* DNA Polymerase will not function in reactions that contain dUTP in the primers or in the dNTP mix.

<sup>[2]</sup> For most targets, 1 unit is optimal. Higher concentrations may be inhibitory. More enzyme may be required for longer targets (>3 kb).

2. Mix the tube contents; if needed, cover with mineral or silicone oil.

**Note:** The oil is not needed in thermal cyclers equipped with a heated lid.

3. Cap the tubes, then centrifuge briefly to collect the contents.
4. Perform thermal cycling as follows:

Step	Three-step cycling		Two-step cycling	
	Temp.	Time	Temp.	Time
Initial denaturation	95°C	2 minutes	95°C	2 minutes
<b>25–35 cycles of:</b>				
Denature	95°C	15 seconds	95°C	15 seconds
Anneal	55–64°C	30 seconds	—	—
Extend	68°C	1 minute per kb	68°C	1 minute per kb

**Note:** Two-step cycling can be used for long primers with high  $T_m$ .

5. Maintain the reaction at 4°C after cycling. The samples can be stored at –20°C until use.
6. Analyze the products by agarose gel electrophoresis and visualize by ethidium bromide staining.

## Limited product warranty

Life Technologies Corporation and/or its affiliate(s) warrant their products as set forth in the Life Technologies' General Terms and Conditions of Sale found on Life Technologies' website at [www.thermofisher.com/us/en/home/global/terms-and-conditions.html](http://www.thermofisher.com/us/en/home/global/terms-and-conditions.html). If you have any questions, please contact Life Technologies at [www.thermofisher.com/support](http://www.thermofisher.com/support).

---

The information in this guide is subject to change without notice.

### DISCLAIMER

TO THE EXTENT ALLOWED BY LAW, LIFE TECHNOLOGIES AND/OR ITS AFFILIATE(S) WILL NOT BE LIABLE FOR SPECIAL, INCIDENTAL, INDIRECT, PUNITIVE, MULTIPLE, OR CONSEQUENTIAL DAMAGES IN CONNECTION WITH OR ARISING FROM THIS DOCUMENT, INCLUDING YOUR USE OF IT.

**Revision history:** Pub. No. MAN0001079

Revision	Date	Description
A.0	5 May 2016	Format, style, and legal updates
—	1 June 2010	Baseline for this revision history

**Important Licensing Information:** These products may be covered by one or more Limited Use Label Licenses. By use of these products, you accept the terms and conditions of all applicable Limited Use Label Licenses.

Corporate entity: Life Technologies Corporation | Carlsbad, CA 92008 USA | Toll Free in USA  
1 800 955 6288

©2016 Thermo Fisher Scientific Inc. All rights reserved. All trademarks are the property of Thermo Fisher Scientific and its subsidiaries unless otherwise specified.

For support visit [thermofisher.com/support](http://thermofisher.com/support) or email [techsupport@lifetech.com](mailto:techsupport@lifetech.com)

[thermofisher.com](http://thermofisher.com)

5 May 2016

**ThermoFisher**  
SCIENTIFIC