

# GeneRuler 1 kb DNA Ladder, ready-to-use

Catalog Number SM0313, SM0314

Pub. No. MAN0013007 Rev. D.00



**WARNING!** Read the Safety Data Sheets (SDSs) and follow the handling instructions. Wear appropriate protective eyewear, clothing, and gloves. Safety Data Sheets (SDSs) are available from [thermofisher.com/support](https://www.thermofisher.com/support).

## Contents and storage

Cat. No.	Contents	Amount	Storage
SM0313	GeneRuler 1 kb DNA Ladder, ready-to-use	250 (5 x 50) µg (for 500 applications), 0.1 µg/µL	at room temperature or at 4 °C for periods up to 6 months. For longer periods store at -20 °C.
	6X TriTrack DNA Loading Dye	2 x 1 mL	
SM0314	GeneRuler 1 kb DNA Ladder, ready-to-use	50 µg (for 100 applications), 0.1 µg/µL	
	6X TriTrack DNA Loading Dye	1 mL	

## Description

Thermo Scientific™ GeneRuler™ 1kb DNA Ladder, ready-to-use, is designed for sizing and approximate quantification of wide range double-stranded DNA fragments on agarose gel. The ladder is composed of fourteen chromatography-purified individual DNA (in base pairs): 10000, 8000, **6000**, 5000, 4000, 3500, **3000**, 2500, 2000, 1500, **1000**, 750, 500, 250. It contains three reference bands (6000, 3000 and 1000 bp) for easy orientation. The ladder is ready to use – it is premixed with 6X TriTrack DNA Loading Dye for direct loading on gel.

## Storage and Loading Buffer

10 mM Tris-HCl (pH 7.6), 10 mM EDTA, 0.005 % bromophenol blue, 0.005 % xylene cyanol FF, 0.025 % orange G and 10 % glycerol.

## 6X TriTrack DNA Loading Dye

10 mM Tris-HCl (pH 7.6), 0.03 % bromophenol blue, 0.03 % xylene cyanol FF, 0.15 % orange G, 60 % glycerol and 60 mM EDTA.

## Protocol for Loading


**Step 1:** Mix gently

**Step 2:** Load 1 µL per 1 mm gel lane.

## Recommendations

- Do not heat before loading.
- Dilute your DNA sample with the 6X TriTrack DNA Loading Dye (#R1161, supplied with the ladder): mix 1 volume of the dye solution with 5 volumes of the DNA sample;
- Load the same volumes of the DNA sample and the DNA ladder;
- For quantification, adjust the concentration of the sample to equalize it approximately with the amount of DNA in the nearest band of the ladder.
- For DNA band visualization with SYBR™ Green and other intercalating dyes, do not add the dyes into the sample, use gel staining after electrophoresis or include dyes into agarose gel to avoid aberrant DNA migration.
- **Important note:** For DNA bands visualization with GelRed™ use gel staining after electrophoresis to avoid aberrant DNA migration.

## GeneRuler 1 kb DNA Ladder, ready-to-use

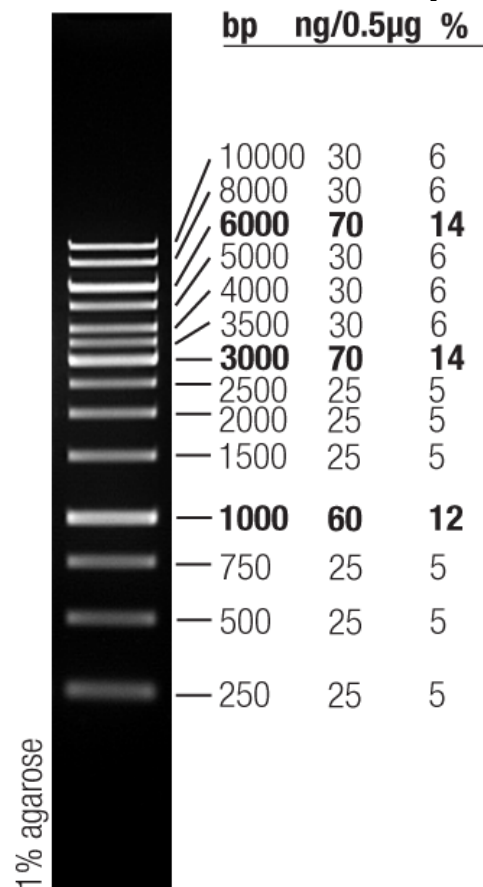

0.5 µg/lane, 8 cm length gel,  
1X TAE, 7 V/cm, 45 min

## Limited product warranty

Life Technologies Corporation and/or its affiliate(s) warrant their products as set forth in the Life Technologies' General Terms and Conditions of Sale at [www.thermofisher.com/us/en/home/global/terms-and-conditions.html](http://www.thermofisher.com/us/en/home/global/terms-and-conditions.html). If you have any questions, please contact Life Technologies at [www.thermofisher.com/support](http://www.thermofisher.com/support).



Thermo Fisher Scientific Baltics UAB | V.A. Graiciuno 8, LT-02241 Vilnius, Lithuania  
For descriptions of symbols on product labels or product documents, go to [thermofisher.com/symbols-definition](http://thermofisher.com/symbols-definition).

The information in this guide is subject to change without notice.

**DISCLAIMER:** TO THE EXTENT ALLOWED BY LAW, THERMO FISHER SCIENTIFIC INC. AND/OR ITS AFFILIATE(S) WILL NOT BE LIABLE FOR SPECIAL, INCIDENTAL, INDIRECT, PUNITIVE, MULTIPLE, OR CONSEQUENTIAL DAMAGES IN CONNECTION WITH OR ARISING FROM THIS DOCUMENT, INCLUDING YOUR USE OF IT.

**Important Licensing Information:** These products may be covered by one or more Limited Use Label Licenses. By use of this product, you accept the terms and conditions of all applicable Limited Use Label Licenses.

©2019 Thermo Fisher Scientific Inc. All rights reserved. All trademarks are the property of Thermo Fisher Scientific and its subsidiaries unless otherwise specified. GelRed is a registered trademark of Biotium Inc.

[thermofisher.com/support](http://thermofisher.com/support) | [thermofisher.com/askaquestion](http://thermofisher.com/askaquestion)

[thermofisher.com](http://thermofisher.com)

26 July 2019

**ThermoFisher**  
SCIENTIFIC