Invitrogen[™] Platinum[™] II Hot-Start PCR Master Mix (2X)

| Pub. No. MAN0017535 Rev. A.0 | |
|--|--|
| Catalog No.Size14000-01250 reactions14000-013200 reactions14000-0141000 reactions | Enzyme char Hot-start: Length: |
| All the components of the kit can be stored at 4° C for periods up to 3 months. For longer storage, keep all components at -20° C. | Fidelity vs. <i>Ta</i> Format: |
| Template: cDNA, genomic DNA, plasmid DNA, phage DNA Forward and reverse primers Invitrogen[™] E-Gel[™] EX Agarose Gels, 1% (Cat. No. G4010-01) | PCR setup Use the measured parameters in the |
| Invitrogen[™] E-Gel[™] 1 kb Plus Express DNA Ladder | Component |
| | Water, nuclease |
| Gel loading buffer | Platinum [™] II Ho Master Mix (2X) |
| Varies depending on amplicon length. | 10 µM forward |
| • Invitrogen [™] Platinum [™] II Hot-Start PCR Master Mix (2X) contains | 10 µM reverse p |
| optimized Platinum ^{\square} II PCR buffer with dNTPs. | Template DNA ² |
| Platinum[™] II <i>Taq</i> Hot-Start DNA Polymerase is an engineered <i>Taq</i> DNA polymerase that shows increased resistance to reaction inhibitors originating from sample material or DNA purification | Platinum [™] GC E (optional) ³ |
| steps. The polymerase activity is blocked at ambient temperatures and restored after the initial denaturation step at 94°C. This automatic "hot start" provides increased sensitivity, specificity, and yield, | ¹ Provides 1.5 mM ² 0.5–500 ng geno synthesis reaction ³ Recommended |
| Due to unique composition of the Platinum[™] II PCR buffer, the annealing temperature is 60°C for most primer pairs designed | PCR protocol Go to page 2 |
| Platinum[™] II <i>Taq</i> DNA polymerase extends 1 kb in 15 seconds. The extension step can be prolonged without a negative effect on | Click here for |
| specificity. The enzyme has a template independent terminal transferase activity that adds a single deoxyadenosine (A) to the 3' ends of | Optimization Click here for |
| PCR products. Like standard <i>Taq</i> , it has both 5' to 3' polymerase and 5' to 3' exonuclease activities, but lacks 3' to 5' exonuclease activity. | Troubleshoot |
| Visit our product page for additional information. | Limited w |
| | Catalog No. Size 14000-012 50 reactions 14000-013 200 reactions 14000-014 1000 reactions All the components of the kit can be stored at 4°C for periods up to 3 months. For longer storage, keep all components at -20°C. • Template: cDNA, genomic DNA, plasmid DNA, phage DNA • Forward and reverse primers • Invitrogen [™] E-Gel [™] EX Agarose Gels, 1% (Cat. No. G4010-01) • Invitrogen [™] E-Gel [™] 1 kb Plus Express DNA Ladder (Cat. No. 10488-091) • 0.2 or 0.5-mL nuclease-free microcentrifuge tubes • Gel loading buffer Varies depending on amplicon length. • Invitrogen [™] Platinum [™] II Hot-Start PCR Master Mix (2X) contains Platinum [™] II Taq Hot-Start DNA Polymerase premixed in an optimized Platinum [™] II PCR buffer with dNTPs. • Platinum [™] II Taq Hot-Start DNA Polymerase is an engineered Taq DNA polymerase that shows increased resistance to reaction inhibitors originating from sample material or DNA purification steps. • The polymerase activity is blocked at ambient temperatures and restored after the initial denaturation step at 94°C. This automatic "hot start" provides increased sensitivity, specificity, and yield, while allowing reaction assembly at room temperature. • Due to unique composition of the Platinum [™] II PCR buffer, the annealing temperature is 60°C for most primer pairs designed following general primer design rules. • Platinum [™] II Taq DNA polymerase extends 1 kb in 15 seconds |

aracteristics

| Hot-start: | Antibody |
|---------------------------|------------|
| Length: | Up to 5 kb |
| Fidelity vs. <i>Taq</i> : | 1X |
| Format: | Master mix |
| | |

rements below to prepare your PCR experiment, or enter your own he column provided.

| Component | 20-µL rxn | 50-µL rxn | Custom | Final conc. |
|--|-----------|-----------|--------|-------------|
| Water, nuclease-free | to 20 µL | to 50 µL | to µI | |
| Platinum [™] II Hot-Start PCR Master Mix (2X) ¹ | 10 µL | 25 µL | μL | 1X |
| 10 µM forward primer | 0.4 µL | 1 µL | μL | 0.2 µM |
| 10 µM reverse primer | 0.4 µL | 1 µL | μL | 0.2 µM |
| Template DNA ² | varies | varies | μL | <500 ng/rxn |
| Platinum [™] GC Enhancer (<i>optional</i>) ³ | 4 µL | 10 µL | μL | 1X |

nM MgCl₂ in final reaction concentration.

nomic DNA, 1 pg–50 ng plasmid or viral DNA, or 1–5 µL of cDNA on per 50-µL PCR reaction.

d for targets with >65% GC sequences.

ЪL

2 for instructions to prepare and run your PCR experiment.

uidelines

for important PCR guidelines.

n strategies

for guidelines to optimize your PCR experiment.

oting

to troubleshoot your PCR experiment.

warranty, disclaimer, and licensing information



For Research Use Only. Not for use in diagnostic procedures.

PCR protocol

The example procedure below shows the appropriate volumes for a single **50-µL** reaction. For multiple reactions, prepare a master mix of components common to all reactions to minimize pipetting error, then dispense the appropriate volumes into each 0.2-mL or 0.5-mL PCR tube before adding template DNA and primers.

| | Steps | Action | | Pro | cedure det | tails | | |
|---|-------|------------------------|---|----------------------|----------------|---------------------|--------------------------------|----|
| 1 | | Thaw reagents | Thaw, mix, and briefly centrifuge | each component b | efore use. | | | |
| | | | a. Add the following component | s to each reaction t | rube. | | | |
| | | | Note: Consider the volumes for required to reach your final read | | isted in steps | s 2 and 3 to determ | ine the correct amount of wate | er |
| | | | Component | | Volume | for 50-µL rxn | Final concentration | |
| 2 | | Prepare PCR master mix | Water, nuclease-free | | t | ο 50 μL | _ | |
| | | | Platinum [™] II Hot-Start PCR Ma | aster Mix (2X) | | 25 µL | 1X | |
| | | | Platinum [™] GC Enhancer (opti | ional)1 | | 10 µL | 1X | |
| | | | ¹ Recommended for targets with | a >65% GC sequence | es. | | | |
| | | | b. Mix, then briefly centrifuge the | e components. | | | | |
| | | | a. Add your template DNA and j | primers to each tub | pe for a final | reaction volume o | f 50 µL. | |
| | 191 | | Component | Volume for 50 |)-μL rxn | Final concentra | ation | |
| 2 | R | Add template DNA and | 10 μM forward primer | 1 µL | | 0.2 µM | | |
| 3 | | primers | 10 μM reverse primer | 1 µL | | 0.2 μM | | |
| | | | Template DNA | varies | ; | <500 ng/rxi | 1 | |
| | | | b. Cap each tube, mix, then brief | y centrifuge the co | ontents. | | | |

| Steps | Action | Procedure details | | | | | |
|---------|---|--|---------------------|-----------------|---------------|------------------------------|------------|
| | | Step | | 3-step protocol | | 2-step protocol ¹ | |
| | | | | Temperature | Time | Temperature | Time |
| | | Initial denaturation | | 94°C | 2 minutes | 94°C | 2 minutes |
| | | | Denature | 94°C | 15 seconds | 98°C | 5 seconds |
| - | Incubate reactions in a | 25–35 PCR cycles | Anneal ² | 60°C | 15 seconds | 60°C | 1E accordo |
| 4 (• ^ | | Cycles | Extend | 68°C | 15 seconds/kb | | 15 seconds |
| 7 | thermal cycler | Но | ld | 4°C | hold | 4°C | hold |
| | ¹ Recommended for simple amplicons up to 1 kb with 45–65% GC sequences. For longer, GC-rich, and complex amplicons, or cDNA targets, use the 3-step cycling protocol. ² 60°C annealing temperature works for most primers. In cases when annealing temperature requires additional optimization, we recommend performing gradient PCR or rdesigning the primers. Note: Refer to "Optimization strategies", page 1, for guidelines to optimize cycling conditions. | | | | | | |
| 5 | Add gel loading buffer and analyze with gel electrophoresis | a. Add gel loading buffer to 10 µL of PCR sample, mix, and briefly centrifuge the contents. Note: Dilute the PCR sample 2- to 20-fold for optimal separation on E-Gel[™] agarose gels. b. Analyze the sample using agarose gel electrophoresis. c. Use your PCR product immediately in down-stream applications, or store it at -20°C. | | | | | |

