

TaqMan® GMO Maize 35S Detection Kit

Catalog Number 4327690

Pub. No. 4327599 Rev. D

| Contents | Quantity | Storage conditions |
|-------------------------------|----------|--------------------------------------|
| TaqMan® GMO Maize 35S PCR Mix | 3 tubes | Upon receipt, store at -25 to -15°C. |
| AmpliTaq Gold® DNA Polymerase | 2 tubes | |
| Positive Control DNA | 1 tube | |
| Negative Control | 1 tube | |

Note: For safety and biohazard guidelines, refer to the “Safety” appendix in the *TaqMan® Genetically Modified Organism (GMO) Detection Kits User Guide* (Pub. no. 4327687). Read the Safety Data Sheets (SDSs) and follow the handling instructions. Wear appropriate protective eyewear, clothing, and gloves.

Product description

The TaqMan® GMO Maize 35S Detection Kit (Cat. no. 4327690) contains sufficient reagents to perform 100 reactions of 50- μ L reaction volume, including the recommended controls.

Additional information and procedures can be found in the *TaqMan® Genetically Modified Organism (GMO) Detection Kits User Guide* (Pub. no. 4327687).

Limited product warranty

Life Technologies Corporation and/or its affiliate(s) warrant their products as set forth in the Life Technologies' General Terms and Conditions of Sale found on Life Technologies' website at www.lifetechnologies.com/termsandconditions. If you have any questions, please contact Life Technologies at www.lifetechnologies.com/support.

For testing of Food and Environmental samples only.

For testing of Food and Environmental samples only.

The information in this guide is subject to change without notice.

DISCLAIMER

LIFE TECHNOLOGIES CORPORATION AND/OR ITS AFFILIATE(S) DISCLAIM ALL WARRANTIES WITH RESPECT TO THIS DOCUMENT, EXPRESSED OR IMPLIED, INCLUDING BUT NOT LIMITED TO THOSE OF MERCHANTABILITY, FITNESS FOR A PARTICULAR PURPOSE, OR NON-INFRINGEMENT. TO THE EXTENT ALLOWED BY LAW, IN NO EVENT SHALL LIFE TECHNOLOGIES AND/OR ITS AFFILIATE(S) BE LIABLE, WHETHER IN CONTRACT, TORT, WARRANTY, OR UNDER ANY STATUTE OR ON ANY OTHER BASIS FOR SPECIAL, INCIDENTAL, INDIRECT, PUNITIVE, MULTIPLE OR CONSEQUENTIAL DAMAGES IN CONNECTION WITH OR ARISING FROM THIS DOCUMENT, INCLUDING BUT NOT LIMITED TO THE USE THEREOF.

LIMITED USE LABEL LICENSE No. 492: Environmental Testing, Quality Control/Quality Assurance Testing, Food and Agricultural Testing :

Notice to Purchaser: The purchase of this product conveys to the purchaser the limited, non-transferable right to use the purchased amount of the product (a) to perform internal research for the sole benefit of the purchaser; and (b) for environmental testing, quality control/quality assurance testing, food and agricultural testing, including reporting results of purchaser's activities in environmental testing, quality control/quality assurance testing, food and agricultural testing for a fee or other commercial consideration. No other right is hereby granted expressly, by implication, or by estoppel. This product is for environmental testing, quality control/ quality assurance testing, food and agricultural testing and research purposes only.

The purchase of this product does not grant the purchaser any additional rights, including (without limitation) the right to transfer or resell the product in any form or the right to use the product as a therapeutic agent or diagnostics test component. For information on obtaining additional rights, please contact outlicensing@lifetech.com or Out Licensing, Life Technologies, 5791 Van Allen Way, Carlsbad, California 92008.

Life Technologies is a Thermo Fisher Scientific brand. ©2014 Thermo Fisher Scientific Inc. All rights reserved. All trademarks are the property of Thermo Fisher Scientific and its subsidiaries. AmpliTaq Gold® is a registered trademark of Roche Molecular Systems, Inc. TaqMan® is a registered trademark of Roche Molecular Systems, Inc., used under permission and license.

For support visit lifetechnologies.com/support or email techsupport@lifetech.com

lifetechnologies.com

25 February 2014

