VetMAX[™] Xeno[™] Internal Positive Controls and Assays for PCR and RT-PCR USER GUIDE

Catalog Numbers A29761, A29762, A29763, A29764, A29765, A29766, A29767, and A29768 Publication Number MAN0014500 Revision C.0



The information in this guide is subject to change without notice.

DISCLAIMER

TO THE EXTENT ALLOWED BY LAW, LIFE TECHNOLOGIES AND/OR ITS AFFILIATE(S) WILL NOT BE LIABLE FOR SPECIAL, INCIDENTAL, INDIRECT, PUNITIVE, MULTIPLE, OR CONSEQUENTIAL DAMAGES IN CONNECTION WITH OR ARISING FROM THIS DOCUMENT, INCLUDING YOUR USE OF IT.

Revision history

Table 1 Pub. No. MAN0014500

Revision	Date	Description
C.0	14 April 2016	Correction to troubleshooting.
B.0	January 2016	Changed regulatory label
A.0	October 2015	New document

Limited Use Label License No. 569: Veterinary and Research Use

Notice to Purchaser : The purchase of this product conveys to the purchaser the limited, non-transferable right to use the purchased amount of the product for research and veterinary use only. No other right is hereby granted expressly, by implication, or by estoppel. The purchase of this product does not grant the purchaser any additional rights, including (without limitation) the right to transfer or resell the product in any form or the right to use the product as a therapeutic agent or human diagnostics test component. For information on obtaining additional rights, please contact **outlicensing@lifetech.com** or Out Licensing, Life Technologies, 5791 Van Allen Way, Carlsbad, California 92008.

Corporate entity

Life Technologies Corporation | Carlsbad, CA 92008 USA | Toll Free in USA 1 800 955 6288

TRADEMARKS

All trademarks are the property of Thermo Fisher Scientific and its subsidiaries unless otherwise specified.

©2016 Thermo Fisher Scientific Inc. All rights reserved.

Contents

Product information	4
Product description	4
Methods	7
Workflow A (add Xeno [™] IPC to samples)	7
Lyse the samples	7
Perform the PCR or RT-PCR reaction	7
Workflow B (add Xeno [™] Assav to PCR/RT-PCR mix)	8
Perform the PCR or RT-PCR reaction	8
Workflow C (add Xeno ^{M} IPC to samples and Xeno ^{M} Assay to PCR/RT-PCR mix)	
I vee the comples	0
Perform the PCP or PT_PCP reaction	/
Analyze results	9
APPENDIX A Troubleshooting	10
APPENDIX B Safety	. 11
Chemical safety	. 12
Biological hazard safety	. 13
Documentation and support	. 14
Customer and technical support	. 14
Limited product warranty	. 14



Product information

IMPORTANT! Before using this product, read and understand the information in the "Safety" appendix in this document.

Product description

Applied Biosystems[™] VetMAX[™] Xeno[™] Internal Positive Controls (IPC) and Assays provide a complete solution to qualify the effectiveness of sample preparation and real-time PCR and RT-PCR workflows.

- Add VetMAX[™] Xeno[™] IPC to the lysis solution used for nucleic acid isolation, to serve as a positive control for recovery of nucleic acid.
- Additionally, use diluted VetMAX[™] Xeno[™] IPC as a positive control for PCR and RT-PCR.
- Multiplex VetMAX[™] Xeno[™] IPC Assay with the assay of interest, to detect Xeno[™] IPC in test samples, to enable monitoring of nucleic acid recovery and PCR inhibition.

VetMAX[™] Xeno[™] IPC and VetMAX[™] Xeno[™] IPC Assays enable the user to distinguish true target negatives from samples affected by PCR inhibition.

VetMAX[™] Xeno[™] IPC is available in DNA and RNA formulations. VetMAX[™] Xeno[™] IPC Assays are available with VIC[™] or LIZ[™] reporter dyes, for flexibility in multiplex reactions. The assay formulations are optimized for use with challenging animal samples to detect RNA or DNA pathogens.

For these reagents	You need	And follow
 Applied Biosystems[™] VetMAX[™] Reagents that do not include Xeno[™] IPC (RNA or DNA), such as: VetMAX[™] NA and EU PRRSV Reagents (Cat. no. 4468465) VetMAX[™] M. hyopneumoniae Reagents (Cat. no. 4415218) 	VetMAX [™] Xeno [™] IPC (RNA or DNA)	Workflow A (page 7)



For these reagents	You need	And follow
 Applied Biosystems[™] VetMAX[™] Plus reagents that do not include Xeno[™] Assay: Master mix with Xeno[™] DNA control: VetMAX[™]-Plus qPCR Master Mix (Cat. no. 4415327) Master mixes with Xeno[™] RNA control: VetMAX[™]-Plus Multiplex One-Step RT-PCR Kit (Cat. no. 4415320) VetMAX[™]-Plus One-Step RT-PCR Kit (Cat. no. 4415328) 	VetMAX [™] Xeno [™] IPC Assay (VIC [™] or LIZ [™] dyes)	Workflow B (page 8)
 Master mix reagents and assays that do not include Xeno[™] IPC (RNA or DNA) and Xeno[™] Assay: Applied Biosystems[™] PCR Master Mix: Path-ID[™] qPCR Master Mix (Cat. no. 4388643) Applied Biosystems[™] RT-PCR Master Mixes, such as: 	VetMAX [™] Xeno [™] IPC (RNA or DNA) and VetMAX [™] Xeno [™] IPC Assay (VIC [™] or LIZ [™] dyes)	Workflow C (page 8)

Kit contents and storage

Iable 2 VetMAX Xeno Internal Positive Control RN	able 2 V	VetMAX [™] Xer	io [™] Internal	Positive	Control	RNA
--	----------	-------------------------	--------------------------	----------	---------	------------

Component	Cat. no. A29763 (100 rxns)	Cat. no. A29761 (500 rxns)	Storage
Xeno [™] RNA Control (10,000 copies/µL)	250 µL	1250 µL	−25°C to −15°C
Nucleic Acid Dilution Solution	250 μL	1250 µL	



Component	Cat. no. A29764 (100 rxns)	Cat. no. A29762 (500 rxns)	Storage
Xeno [™] DNA Control (10,000 copies/µL)	250 µL	1250 µL	-25°C to -15°C
Nucleic Acid Dilution Solution	250 μL	1250 µL	

Table 3 VetMAX[™] Xeno[™] Internal Positive Control DNA

 Table 4
 VetMAX[™] Xeno[™] Internal Positive Control – VIC[™] Assay

Component	Cat. no. A29765 (100 rxns)	Cat. no. A29767 (500 rxns)	Storage
Xeno [™] VIC [™] Primer Probe Mix	110 µL	550 μL	–25°C to –15°C

Table 5 VetMAX[™] Xeno[™] Internal Positive Control – LIZ[™] Assay

Component	Cat. no. A29766 (100 rxns)	Cat. no. A29768 (500 rxns)	Storage
Xeno [™] LIZ [™] Primer Probe Mix	110 µL	550 μL	–25°C to –15°C

Methods



reaction

Workflow A (add Xeno ^{TM} IPC to samples)	7
Workflow B (add Xeno [™] Assay to PCR/RT-PCR mix)	8
Workflow C (add Xeno [™] IPC to samples and Xeno [™] Assay to PCR/RT- PCR mix)	8
Analyze results	9

Workflow A (add Xeno[™] IPC to samples)

Follow this procedure when:

- The PCR/RT-PCR reagent mix contains Xeno[™] Assay
- The samples do not contain Xeno[™] IPC (DNA or RNA)
- Lyse the samples
 1. Add 2 μL (20,000 copies) per reaction of Xeno[™] IPC (DNA or RNA depending on the nucleic acid isolated) to the lysis solution.
 2. Perform the nucleic acid isolation according to the standard procedure.
- Perform the PCR1. Program the real-time PCR instrument according to the manufacturer's
recommendation.
 - Dilute the Xeno[™] IPC (DNA or RNA depending on the nucleic acid isolated) 10fold (1,000 copies per µL) to use as a PCR control.
 - **3.** Assemble the PCR or RT-PCR reaction mix according to the master mix used, reaction size, and number of reactions.
 - **4.** Distribute the PCR or RT-PCR reaction mix to the wells of a PCR plate or to PCR tubes.
 - 5. Add purified nucleic acid and controls according to the following table.

Reaction type	Component
Test sample	Purified nucleic acid
Positive control ^[1]	Diluted Xeno [™] IPC
Negative control ^[1]	Nuclease-Free Water

^[1] Use the same volume of control as of test sample.

6. Load the plate in the instrument and run the thermal cycling program.

Workflow B (add Xeno[™] Assay to PCR/RT-PCR mix)

Follow this procedure when:

- The PCR/RT-PCR reagent mix does not contain Xeno[™] Assay
- The samples contain Xeno[™] IPC

Note: We recommend that you perform appropriate testing to evaluate the performance of your target assay in presence of the Xeno[™] Assay (i.e. the performance of your target remains the same after addition of the Xeno[™] Assay to the master mix).

Perform the PCR or RT-PCR reaction Program the real-time PCR instrument according to the manufacturer's recommendation.
 IMPORTANTI Make sure to set up the reporter corresponding to the Xero

IMPORTANT! Make sure to set up the reporter corresponding to the XenoTM Assay used (VICTM or LIZTM dye).

- 2. Dilute the Xeno[™] IPC (DNA or RNA depending on the nucleic acid isolated) 10fold (1,000 copies per μL) to use as a PCR control.
- **3.** Assemble the PCR or RT-PCR reaction mix according to the master mix used, reaction size, and number of reactions.
- 4. Add 1 µL of Xeno[™] Assay per 25 µL of PCR or RT-PCR mix.
- **5.** Distribute the PCR or RT-PCR reaction mix to the wells of a PCR plate or to PCR tubes.
- **6.** Add purified nucleic acid and controls according to the following table.

Reaction type	Component
Test sample	Purified nucleic acid
Positive control ^[1]	Diluted Xeno [™] IPC
Negative control ^[1]	Nuclease-Free Water

^[1] Use the same volume of control as of test sample.

7. Load the plate in the instrument and run the thermal cycling program.

Workflow C (add Xeno[™] IPC to samples and Xeno[™] Assay to PCR/RT-PCR mix)

Follow this procedure when:

- The PCR/RT-PCR reagent mix does not contain Xeno[™] Assay
- The samples do not contain Xeno[™] IPC

Note: We recommend that you perform appropriate testing to evaluate the performance of your target assay in presence of the Xeno[™] Assay (i.e. the performance of your target remains the same after addition of the Xeno[™] Assay to the master mix).



- **Lyse the samples** 1. Add 2 μL (20,000 copies) per reaction of Xeno[™] IPC (DNA or RNA depending on the nucleic acid isolated) to the lysis solution.
 - 2. Perform the nucleic acid isolation according to the standard procedure.

Perform the PCR or RT-PCR reaction

1. Program the real-time PCR instrument according to the manufacturer's recommendation.

IMPORTANT! Make sure to set up the reporter corresponding to the XenoTM Assay used (VICTM or LIZTM dye).

- 2. Dilute the Xeno[™] IPC (DNA or RNA depending on the nucleic acid isolated) 10fold (1,000 copies per μL) to use as a PCR control.
- **3.** Assemble the PCR or RT-PCR reaction mix according to the master mix used, reaction size, and number of reactions.
- **4.** Add 1 μ L of XenoTM Assay per 25 μ L of PCR or RT-PCR mix.
- **5.** Distribute the PCR or RT-PCR reaction mix to the wells of a PCR plate or to PCR tubes.
- 6. Add purified nucleic acid and controls according to the following table.

Reaction type	Component
Test sample	Purified nucleic acid
Positive control ^[1]	Diluted Xeno [™] IPC
Negative control ^[1]	Nuclease-Free Water

^[1] Use the same volume of control as of test sample.

7. Load the plate in the instrument and run the thermal cycling program.

Analyze results

The general process for analyzing results from real-time PCR and RT-PCR experiments is:

- 1. View the amplification plots for all reactions to make sure they appear normal.
- 2. Set the baseline and threshold values.
- **3.** Use the relative standard curve or the comparative C_T method to analyze your data.

The details of data analysis depend on the real-time PCR instrument that you use; refer to the appropriate user guide for instructions on how to analyze your data.



Troubleshooting

Observation	Possible cause	Recommended action
Test samples	The Xeno [™] Assay components	No action required.
Xeno [™] IPC—no or low signal Target—high signal	the PCR. High levels of target in a sample can reduce	No or low signal from Xeno [™] IPC is expected in a reaction that has a strong signal for the
	amplification of Xeno [™] IPC.	larget.
Test samples	Poor nucleic acid recovery.	Check the C _T values of Xeno [™] IPC in the
Xeno™ IPC—no or low signal		negative extraction control.
Target—no signal or inconclusive-range signal		A high C _T value for Xeno [™] IPC indicates that the nucleic acid recovery was poor.
		Repeat the purification of the original sample.
	The RNA samples contain inhibitors of RT-PCR.	Dilute purified nucleic acid two- and four-fold and check if the signal strength increases due to the dilution of inhibitors.

Safety





WARNING! GENERAL SAFETY. Using this product in a manner not specified in the user documentation may result in personal injury or damage to the instrument or device. Ensure that anyone using this product has received instructions in general safety practices for laboratories and the safety information provided in this document.

- Before using an instrument or device, read and understand the safety information provided in the user documentation provided by the manufacturer of the instrument or device.
- Before handling chemicals, read and understand all applicable Safety Data • Sheets (SDSs) and use appropriate personal protective equipment (gloves, gowns, eye protection, etc). To obtain SDSs, see the "Documentation and Support" section in this document.

Chemical safety



WARNING! GENERAL CHEMICAL HANDLING. To minimize hazards, ensure laboratory personnel read and practice the general safety guidelines for chemical usage, storage, and waste provided below, and consult the relevant SDS for specific precautions and instructions:

- Read and understand the Safety Data Sheets (SDSs) provided by the chemical manufacturer before you store, handle, or work with any chemicals or hazardous materials. To obtain SDSs, see the "Documentation and Support" section in this document.
- Minimize contact with chemicals. Wear appropriate personal protective equipment when handling chemicals (for example, safety glasses, gloves, or protective clothing).
- Minimize the inhalation of chemicals. Do not leave chemical containers open. Use only with adequate ventilation (for example, fume hood).
- Check regularly for chemical leaks or spills. If a leak or spill occurs, follow the manufacturer's cleanup procedures as recommended in the SDS.
- · Handle chemical wastes in a fume hood.
- Ensure use of primary and secondary waste containers. (A primary waste container holds the immediate waste. A secondary container contains spills or leaks from the primary container. Both containers must be compatible with the waste material and meet federal, state, and local requirements for container storage.)
- After emptying a waste container, seal it with the cap provided.
- Characterize (by analysis if necessary) the waste generated by the particular applications, reagents, and substrates used in your laboratory.
- Ensure that the waste is stored, transferred, transported, and disposed of according to all local, state/provincial, and/or national regulations.
- **IMPORTANT!** Radioactive or biohazardous materials may require special handling, and disposal limitations may apply.

Biological hazard safety



 U.S. Department of Health and Human Services, *Biosafety in Microbiological* and *Biomedical Laboratories (BMBL)*, 5th Edition, HHS Publication No. (CDC) 21-1112, Revised December 2009; found at:

www.cdc.gov/biosafety/publications/bmbl5/BMBL.pdf

• World Health Organization, *Laboratory Biosafety Manual*, 3rd Edition, WHO/CDS/CSR/LYO/2004.11; found at:

www.who.int/csr/resources/publications/biosafety/Biosafety7.pdf

Documentation and support

Customer and technical support

Visit **thermofisher.com/support** for the latest in services and support, including:

- Worldwide contact telephone numbers
- Product support, including:
 - Product FAQs
 - Software, patches, and updates
- Order and web support
 - Product documentation, including:
 - User guides, manuals, and protocols
 - Certificates of Analysis
 - Safety Data Sheets (SDSs; also known as MSDSs)

Note: For SDSs for reagents and chemicals from other manufacturers, contact the manufacturer.

Limited product warranty

Life Technologies Corporation and/or its affiliate(s) warrant their products as set forth in the Life Technologies' General Terms and Conditions of Sale found on Life Technologies' website at **www.thermofisher.com/us/en/home/global/ terms-and-conditions.html**. If you have any questions, please contact Life Technologies at **www.thermofisher.com/support**.



