

RNaseZap[®] Wipes

RNase Decontamination Wipes

Part Number AM9786, AM9788



The Ambion[®] popular RNaseZap[®] solution is now provided in a convenient wipe format. RNaseZap Wipes are designed to eliminate RNase contamination on work surfaces, pipettors and equipment that must be RNase-free. RNaseZap solution is a formulation of three ingredients known to be active against RNase. It effectively removes even high levels of RNase contamination that similar products cannot. With thorough rinsing, RNaseZap solution leaves no residues that are inhibitory to enzymatic reactions. Always wear gloves while using RNaseZap Wipes because prolonged contact with skin may cause irritation.

A. Instructions for Use

Preparing RNaseZap Wipes for use

1. Place the wipe roll into the empty canister.
2. Pour RNaseZap solution over the wipes in the canister.
3. Wait 5 minutes to allow the RNaseZap solution to saturate the wipes.
4. Pull the innermost wipe up from the center of the roll and thread it through the opening in the lid of the canister. Secure the lid tightly on the canister.

Decontaminating work surfaces

Wipe the surface to be cleaned thoroughly with RNaseZap Wipes, rinse with distilled water and then dry with a clean paper towel.

Decontaminating lab equipment

Thoroughly wipe all exposed surfaces with RNaseZap Wipes, rinse with distilled water and then dry with a clean paper towel.



CAUTION

Do not expose corrodible metal surfaces to RNaseZap solution for extended periods of time.

Decontaminating plastic and glass vessels

Large vessels may be cleaned by wiping the entire inner surface with RNaseZap Wipes. Then rinse the vessel thoroughly with distilled water and dry.

Decontaminating pipettors

Clean the shaft of the pipettor thoroughly with a RNaseZap Wipe, then rinse with distilled water and dry.

B. RNaseZap® Wipes Specifications

Contents, Storage & Stability

Store at room temperature.

P/N AM9786	P/N AM9788	Component
1	–	RNaseZap Wipes canister (empty)
1	3	Silver pouch containing RNaseZap Wipes
1	3	Bottle containing RNaseZap

C. Safety Information

This product can be harmful if swallowed or inhaled. It causes irritation to skin, eyes and respiratory tract. It may cause allergic skin reactions or respiratory reactions. Please see the RNaseZap Wipes MSDS for more information.

Chemical safety guidelines

To minimize the hazards of chemicals:

- Read and understand the Material Safety Data Sheets (MSDS) provided by the chemical manufacturer before you store, handle, or work with any chemicals or hazardous materials.

- Minimize contact with chemicals. Wear appropriate personal protective equipment when handling chemicals (for example, safety glasses, gloves, or protective clothing). For additional safety guidelines, consult the MSDS.
- Minimize the inhalation of chemicals. Do not leave chemical containers open. Use only with adequate ventilation (for example, fume hood). For additional safety guidelines, consult the MSDS.
- Check regularly for chemical leaks or spills. If a leak or spill occurs, follow the manufacturer's cleanup procedures as recommended on the MSDS.
- Comply with all local, state/provincial, or national laws and regulations related to chemical storage, handling, and disposal.

Obtaining the MSDS

To obtain Material Safety Data Sheets (MSDSs) for any chemical product supplied by Applied Biosystems or Ambion:

- At www.appliedbiosystems.com, select **Support**, then **MSDS**. Search by chemical name, product name, product part number, or MSDS part number. Right-click to print or download the MSDS of interest.
- At www.ambion.com, go to the web catalog page for the product of interest. Click **MSDS**, then right-click to print or download.
- E-mail (MSDS_Inquiry_CCRM@appliedbiosystems.com) or telephone (650-554-2756; USA) your request, specifying the catalog or part number(s) and the name of the product(s). We will e-mail the associated MSDSs unless you request fax or postal delivery. Requests for postal delivery require 1–2 weeks for processing.

For the MSDSs of chemicals not distributed by Applied Biosystems or Ambion, contact the chemical manufacturer.



Manual 9786M Revision C

Revision Date: September 30, 2008

For research use only. Not for use in diagnostic procedures.

Information in this document is subject to change without notice. Applied Biosystems assumes no responsibility for any errors that may appear in this document.

Applied Biosystems disclaims all warranties with respect to this document, expressed or implied, including but not limited to those of merchantability or fitness for a particular purpose. In no event shall Applied Biosystems be liable, whether in contract, tort, warranty, or under any statute or on any other basis for special, incidental, indirect, punitive, multiple or consequential damages in connection with or arising from this document, including but not limited to the use thereof.

Literature Citation: When you are describing a procedure utilizing this product in a Materials and Methods Section for publication, we would appreciate that you refer to it as RNaseZap® Wipes.

Warranty and Liability: Applied Biosystems is committed to delivering superior product quality and performance, supported by industry-leading global service and technical support teams. Warranty information for the accompanying consumable product is available at www.ambion.com/info/warranty in “Limited Warranty for Consumables,” which is subject to the exclusions, conditions, exceptions, and limitations set forth under the caption “EXCLUSIONS, CONDITIONS, EXCEPTIONS, AND LIMITATIONS” in the full warranty statement. Please contact Applied Biosystems if you have any questions about our warranties or would like information about post-warranty support.

Trademarks: Applied Biosystems, AB (Design), Ambion, and RNaseZap are registered trademarks of Applied Biosystems, Inc. or its subsidiaries in the US and/or certain other countries.

All other trademarks are the sole property of their respective owners.

© 2008 Ambion, Inc. All Rights Reserved.

AB Applied Biosystems

Ambion®

U.S./Canada
(English) 800-327-3002
(French) 800-668-6913
Fax +1-512-651-0201

Europe
tel +44 (0)1480-373-020
fax +44 (0)1480-373-010

Japan
tel 81 (0) 3 5566 6230
fax 81 (0) 3 5566 6507

Direct free phone numbers, distributors:
www.appliedbiosystems.com