

3730/3730xl DNA Analyzer Sequencing Standards, BigDye™ Terminator v3.1

Catalog Number 4336943

Pub. No. 4336884 Rev. B

WARNING! Read the Safety Data Sheets (SDSs) and follow the handling instructions. Wear appropriate protective eyewear, clothing, and gloves. Safety Data Sheets (SDSs) are available from thermofisher.com/support.

Contents and storage

Contents	Amount	Storage
3730/3730xl DNA Analyzer Sequencing Standards, BigDye™ Terminator v3.1	4 tubes ^[1]	-25°C to -15°C

^[1] Contents are lyophilized.

Resuspend the standard

- Resuspend one tube of sequencing standard with 1 mL of Hi-Di™ Formamide (Cat. No. 4311320).
- Vortex for 1 minute, then centrifuge briefly in a microcentrifuge to collect contents at the bottom of the tube.
- Denature the DNA fragments.
 - Incubate the microcentrifuge tube at 95°C for 2 minutes.
 - Incubate the microcentrifuge tube at 4°C, or on ice, for 2 minutes. Immediately proceed to the next step.

Prepare the plates

- Dispense the standard.

For	Dispense
96-well plates	10 µL of standard into every well.
384-well plates	5 µL of standard into the wells corresponding to a single 96 injection (that is, A1, C1, E1, G1, I1, K1, M1, O1, A3, C3, and so on)

- Cover the plate with a 96-well septa (Cat. No. 4315933) or heat-seal film (Cat. No. 4337570).
- Briefly centrifuge the plate to bring the samples to the bottom of the wells and eliminate air bubbles.
- Run the plate on the instrument.

Instrument	Run type	Array size	Polymer	Dye set	Module	
3730/3730xl	Spectral	36 cm	POP-7™	Z	Spect36_SeqStd_POP7	
		50 cm			Spect50_SeqStd_POP7	
		36 cm	POP-6™		Spect36_SeqStd_POP6	
		50 cm			Spect50_SeqStd_POP6	
	Sequencing	36 cm	POP-7™		StdSeq36_POP7	
					RapidSeq36_POP7	
		TargetSeq36_POP7				
		XLRSq50_POP7				
		LongSeq50_POP7				
		FastSeq50_POP7				
		36 cm			POP-6™	StdSeq36_POP6
						LongSeq50_POP6

For more information, see your instrument user guide.

Control sequence

```
1 AATTCCTGC AGGCGTGGCT GCAGCCTGGT TATGATTACT GTTAATGTTG CTACTIONACTGC TGACAATGCT GCTGCTGCTT
81 CTCCTCACTG TCTCCACTTC CTTGAACAAT GCGCCGTCAT GCTTCTTTTG CCTCCCGCTG CTCCAGAAAG CTAGGCCGCA
161 GATCAGAACC ACCACAGTCA ATATCACCAC CTTCTCTTTA TAGATTGCGA ATCTCATGAT AGGGGCTCAG CCTCTGTGCG
241 AGTGGAGAGA AGTTTGCAGG CGAGCTGAGG AGCAATTGCA GGTGATATGA TGTGCTCGGC TCAAGAAGCG GGCCCGGAGA
321 GGAAGAAGTC GTGCCGGGGC TAATTATTGG CAAAACGAGC TCTTGTGTGA AACATTGATC CAACTGGAAT GTCACAAATG
401 GCGAATCAAT ATTCCATAAG GCATGATGGT TGCTCAGAGG CAGGAGAAGA GCAACGAATA CGATCCTATA AAAGATAAAA
481 CATAAATAAA CAGTCTTGAT TATATTCTGG GTATTAAAGC CACAATCAGA ACAAATATAT GCTTTGTATC TTTTCTTGCC
561 TTCTTCATTA CCAACTGCTT CCGCGGCCAC ATTAAGAGAA CTTGTGGTAA GATAAGAAGA TATTTTATTC GTTCTGCTGA
641 CTTGCTGGAT GTCGGGAAAT ATTCTGCATT TGATAAGAGG CGGTTAATTG CAGATATAAT TGGTAGTGAA AAGGGTCGTT
721 GCTATGGTCA CCGTGAAGCG AGTACAGCAG CACAAGAATG TGTGCCGTTT TCAGTTAATA TTGTTTGAAT ATGGTAACTT
801 GTTTTAGTCG GTTTAAAGGT AAGAAGATCT AACCAAAAAC AACACTGCAG TGACTGATTG TAGTATTTAT TTTTCTACTT
881 AATCTTAATT TTGGTGTAAT CATCAACGGC GCACCTCAAC CAATACTCCA ATGTTTATC CATCGACATG ACGTTCGAGA
961 TAGGGTTGAG TGTTGTCCA GTTTGGAACA AGAGTCCACT ATTAAGAAGC GTGGACTCCA ACGTCAAAGG GCGAAAACC
1041 GTCTATCAGG GCGATGGCC ACTACGTGAA CCATCACCCA AATCAAGTTT TTTGGGGTCG AGGTGCCGTA AAGCACTAAA
1121 TCGGAACCT AAAGGGAGCC CCCGATTTAG AGCTTGACGG GGAAAGCCGG CGAACGTGGC GAGAAAGGAA GGAAGAAAG
1201
```

Limited product warranty

Life Technologies Corporation and/or its affiliate(s) warrant their products as set forth in the Life Technologies' General Terms and Conditions of Sale at www.thermofisher.com/us/en/home/global/terms-and-conditions.html. If you have any questions, please contact Life Technologies at www.thermofisher.com/support.



Life Technologies Corporation | 35 Wiggins Avenue | Bedford, MA 01730

For descriptions of symbols on product labels or product documents, go to thermofisher.com/symbols-definition.

The information in this guide is subject to change without notice.

DISCLAIMER: TO THE EXTENT ALLOWED BY LAW, THERMO FISHER SCIENTIFIC INC. AND/OR ITS AFFILIATE(S) WILL NOT BE LIABLE FOR SPECIAL, INCIDENTAL, INDIRECT, PUNITIVE, MULTIPLE, OR CONSEQUENTIAL DAMAGES IN CONNECTION WITH OR ARISING FROM THIS DOCUMENT, INCLUDING YOUR USE OF IT.

Revision history: Pub. No.

Revision	Date	Description
B	14 November 2018	Rebranded and updated.
A	13 October 2002	New document.

Important Licensing Information: These products may be covered by one or more Limited Use Label Licenses. By use of these products, you accept the terms and conditions of all applicable Limited Use Label Licenses.

©2018 Thermo Fisher Scientific Inc. All rights reserved. All trademarks are the property of Thermo Fisher Scientific and its subsidiaries unless otherwise specified.