



pOG44

Flp-recombinase expression vector designed for use with the Flp-In[™] System

Catalog number V6005-20

Revision date 1 August 2012 **Publication Part number** 25-0352

MAN0000170



For Research Use Only. Not for human or animal therapeutic or diagnostic use.

Contents

Kit Contents and Storage	iv
Introduction	1
About the kit	1
Methods	
Using pOG44	3
Appendix	5
pOG44 Vector	
Accessory Products	7
Technical Support	9
Purchaser Notification	10
References	12

Kit Contents and Storage

Contents	40 μL of 0.5 $\mu g/\mu L$ pOG44 in TE, pH 8.0.
Shipping and Storage	pOG44 is shipped at room temperature. Store at -30° C to -10° C.

Introduction

About the kit

Introduction	pOG44 is a 5.8 kb Flp recombinase expression vector designed for use with the Flp-In [™] System (Catalog nos. K6010-01 and K6010-02). When cotransfected with the pcDNA5/FRT plasmid into a Flp-In [™] mammalian host cell line, the Flp recombinase expressed from pOG44 mediates integration of the pcDNA5/FRT vector containing the gene of interest into the genome via Flp Recombination Target (FRT) sites. The vector contains the following elements:
	• The human cytomegalovirus (CMV) immediate-early enhancer/promoter for high-level constitutive expression of the Flp recombinase in a wide range of mammalian cells (Andersson <i>et al.</i> , 1989; Boshart <i>et al.</i> , 1985; Nelson <i>et al.</i> , 1987).
	• Synthetic intron to enhance expression of the <i>FLP</i> gene (Huang and Gorman, 1990; O'Gorman <i>et al.</i> , 1991).
	• <i>FLP</i> gene encoding the Flp recombinase (Buchholz <i>et al.</i> , 1996) to mediate integration of the pcDNA5/FRT expression plasmid into the genome.
	For more information about the Flp-In [™] System, the pcDNA5/FRT plasmid, and generation of the Flp-In [™] host cell line, refer to the Flp-In [™] System manual. The Flp-In [™] System manual is supplied with the Flp-In [™] Complete or Core Systems, but is also available at www.lifetechnologies.com/manuals or by contacting Technical Support (see page 9).
<i>FLP</i> Gene	The <i>FLP</i> gene was originally isolated from the <i>Saccharomyces cerevisiae</i> 2µ plasmid (Broach et al., 1982; Broach and Hicks, 1980), and encodes a site-specific recombinase that is a member of the integrase family of recombinases (Argos et al., 1986). The Flp recombinase mediates a site-specific recombination reaction between interacting DNA molecules via the pairing of interacting FRT sites. For more information about site-specific recombination, refer to the next page and published reviews (Craig, 1988; Sauer, 1994).
	The native <i>FLP</i> gene encodes a protein of 423 amino acids with a calculated molecular weight of 49 kDa. The <i>FLP</i> gene expressed from pOG44 encodes a temperature-sensitive Flp recombinase which carries a point mutation (flp-F70L) that results in a change in amino acid 70 from phenylalanine to leucine (Buchholz <i>et al.</i> , 1996). For more information about the properties of the flp-F70L protein, see below and Buchholz <i>et al.</i> , 1996.
Activity of the Flp Recombinase	When tested in mammalian cells, the native Flp recombinase has been shown to possess optimum recombination activity near 30°C and relatively low activity at 37°C, a result consistent with its physiological role in yeast (Buchholz et al., 1996).
	The flp-F70L protein expressed from pOG44 exhibits increased thermolability at 37°C in mammalian cells when compared to the native Flp recombinase (Buchholz et al., 1996). Studies have shown that the Flp recombinase expressed from pOG44 possesses only 10% of the activity of the native Flp recombinase at 37°C (Buchholz et al., 1996).

About the kit, Continued

Flp Recombinase- Mediated DNA Recombination	 In the Flp-In[™] System, integration of the pcDNA5/FRT expression construct containing your gene of interest into the genome occurs via Flp recombinase-mediated intermolecular DNA recombination. The hallmarks of Flp-mediated recombination are: Recombination occurs between specific FRT sites (see below) on the interacting DNA molecules. Recombination is conservative and requires no DNA synthesis; the FRT sites are preserved following recombination and there is minimal opportunity for introduction of mutations at the recombination site. Strand exchange requires only the small 34 bp minimal FRT site. For more information about the Flp recombinase and conservative site-specific recombination, refer to published reviews (Craig, 1988; Sauer, 1994).
FRT Site	The FRT site, originally isolated from <i>Saccharomyces cerevisiae</i> , serves as a binding site for Flp recombinase and has been well-characterized (Gronostajski and Sadowski, 1985; Jayaram, 1985; Sauer, 1994; Senecoff et al., 1985). The minimal FRT site consists of a 34 bp sequence containing two 13 bp imperfect inverted repeats separated by an 8 bp spacer that includes an <i>Xba</i> I restriction site (see the following figure). An additional 13 bp repeat is found in most FRT sites, but is not required for cleavage (Andrews et al., 1985). While Flp recombinase binds to all three of the 13 bp repeats, strand cleavage actually occurs at the boundaries of the 8 bp spacer region (Andrews et al., 1985; Senecoff et al., 1985). Minimal FRT site
	CS = cleavage site In the Flp-In [™] System, the pFRT/ <i>lac</i> Zeo and pcDNA5/FRT vectors each contain a single FRT site. The pFRT/ <i>lac</i> Zeo plasmid is used to generate the Flp-In [™] host cell line and the pcDNA5/FRT plasmid is used to express the gene of interest in the Flp-In [™] host cell line. For more information about pFRT/ <i>lac</i> Zeo, pcDNA5/FRT, and the Flp-In [™] System, refer to the Flp-In [™] System manual.
Generating Stable Expression Cell Lines	You will cotransfect the pOG44 plasmid and your pcDNA5/FRT construct into your Flp-In [™] host cell line(s) to generate stable cell lines that express your protein of interest. Cotransfection of pOG44 and pcDNA5/FRT allows expression of Flp recombinase resulting in integration of the pcDNA5/FRT plasmid into the genome via the FRT sites. Once the pcDNA5/FRT construct has integrated into the genome, the Flp recombinase is no longer required. The continued presence of Flp recombinase would actually be detrimental to the cells because it could mediate excision of the pcDNA5/FRT construct. For this reason, the pOG44 plasmid lacks an antibiotic resistance marker for selection in mammalian cells. When generating stable expression cell lines, the pOG44 plasmid and, therefore, Flp recombinase expression, will gradually be lost from transfected cells as they are cultured and selected.

Methods

Using pOG44

Introduction	General guidelines to transform pOG44 into <i>E. coli</i> are provided in this section.
General Molecular Biology Techniques	For help with <i>E. coli</i> transformation, restriction enzyme analysis, and DNA biochemistry, refer to <i>Molecular Cloning: A Laboratory Manual</i> (Sambrook <i>et al.,</i> 1989) or <i>Current Protocols in Molecular Biology</i> (Ausubel <i>et al.,</i> 1994).
<i>E. coli</i> Strain	Many <i>E. coli</i> strains are suitable for the propagation and maintenance of the pOG44 vector including TOP10, DH5 α , and JM109. We recommend that you propagate the vector in <i>E. coli</i> strains that are recombination deficient (<i>rec</i> A) and end nuclease A deficient (<i>end</i> A).
	For your convenience, TOP10 is available as chemically competent or electrocompetent cells (see page 7 for ordering information).
Transformation Method	You may use any method of your choice for transformation. Chemical transformation is the most convenient method for many researchers. Electroporation is the most efficient and the method of choice for large plasmids.
Maintenance of Plasmid	To propagate and maintain the pOG44 vector, transform a <i>recA</i> , <i>endA E</i> . <i>coli</i> strain like TOP10, DH5 α , JM109, or equivalent. Select transformants on LB agar plates containing 50 to 100 µg/mL ampicillin. Be sure to prepare a glycerol stock of the plasmid for long-term storage (see the following section).
Preparing a Glycerol Stock	After identifying the correct clone, purify the colony and make a glycerol stock for long-term storage. You should keep a DNA stock of your plasmid at –20°C.
•	 Streak the original colony out on an LB plate containing 50 μg/mL ampicillin. Incubate the plate at 37°C overnight.
	 Isolate a single colony and inoculate into 1–2 mL of LB containing 50 μg/mL ampicillin.
	• Grow the culture to mid-log phase ($OD_{600} = 0.5-0.7$).
	• Mix 0.85 mL of culture with 0.15 mL of sterile glycerol and transfer to a cryovial.
	• Store at -80° C.
Plasmid Preparation	Plasmid DNA for transfection into eukaryotic cells must be very clean and free from phenol and sodium chloride. Contaminants will kill the cells and salt will interfere with lipids decreasing transfection efficiency. We recommend isolating plasmid DNA using the PureLink [®] HiPure Miniprep Kit or the PureLink [®] HiPure Midiprep Kit (see page 7 for ordering information).

Using pOG44, Continued


Note	Several Flp-In [™] host cell lines which stably express the <i>lacZ-Zeocin</i> [™] fusion gene and contain a single integrated FRT site are available (see page 7 for ordering information). To express your gene of interest in 293, CV-1, CHO, 3T3, BHK, or Jurkat cells, you may want to use one of the Flp-In [™] host cell lines to establish your expression cell line.
IMPORTANT!	We have observed down-regulation of the viral CMV promoter and subsequent loss of gene expression when pcDNA5/FRT-based expression constructs are introduced into 3T3 or BHK cells. This behavior is not observed with pEF5/FRT-based expression constructs.
	If you are generating Flp-In [™] expression cell lines using a 3T3 or BHK host cell line, we recommend that you clone your gene of interest into a pEF5/FRT-based expression plasmid (e.g. pEF5/FRT/V5-D-TOPO [®] or pEF5/FRT/V5-DEST). For more information, refer to www.lifetechnologies.com or call Technical Support (see page 9).
MMENO OUTO OUTO OUTO OUTO OUTO OUTO OUTO	Because correct integration of your pcDNA5/FRT construct into the genome is dependent upon Flp recombinase, the expression levels of Flp recombinase in the cell will determine the efficiency of the recombination reaction. Flp recombinase levels must be sufficiently high to mediate recombination at the FRT sites (single recombination event) and overcome the low intrinsic activity of the enzyme (see page 1). We have varied the ratio of pOG44 and pcDNA5/FRT expression plasmid that we cotransfect into mammalian Flp-In [™] host cells to optimize the recombination efficiency. We recommend that you cotransfect you Flp-In [™] host cell line with a ratio of <i>at least</i> 9:1 (w/w) pOG44:pcDNA5/FRT plasmid. Note that this ratio may vary depending on the nature of the cell line. You may want to determine this ratio empirically for your cell line.
IMPORTANT!	When transfecting your Flp-In [™] host cell line, be sure to use supercoiled pOG44 and pcDNA5/FRT plasmid DNA. Flp-mediated recombination between the FRT site on pcDNA5/FRT and the integrated FRT site in the Flp-In [™] host cell line will only occur if the pcDNA5/FRT plasmid is circularized. The pOG44 plasmid should be circularized to minimize the possibility of the plasmid integrating into the genome.
Cotransfection	Once you have cloned your gene of interest into pcDNA5/FRT and have prepared clean plasmid preparations of pOG44 and your pcDNA5/FRT construct, cotransfect the plasmids into your mammalian Flp-In [™] host cell line to generate your stable Flp-In [™] expression cell line. We recommend that you include the appropriate positive and negative controls to help you evaluate your results. Specific guidelines and protocols for generation of the Flp-In [™] expression cell line can be found in the Flp-In [™] System manual.
	Reminder: The pcDNA5/FRT plasmid contains the hygromycin resistance gene to allow selection of transfectants using hygromycin. The pOG44 plasmid does not contain an antibiotic resistance gene for selection in mammalian cells (see pages 5–6).

Appendix

pOG44 Vector

Map of pOG44

pOG44 is a 5785 bp vector that expresses a temperature-sensitive Flp recombinase (flp-F70L) under the control of the human CMV promoter as previously described (O'Gorman *et al.*, 1991). The vector contains a synthetic intron to enhance expression of the *FLP* gene. Note that the vector does not contain an antibiotic resistance marker to allow stable selection in mammalian cells. The following figure summarizes the features of the pOG44 vector. **The sequence for pOG44 is available from www.lifetechnologies.com or by contacting Technical Support (see page 9).**

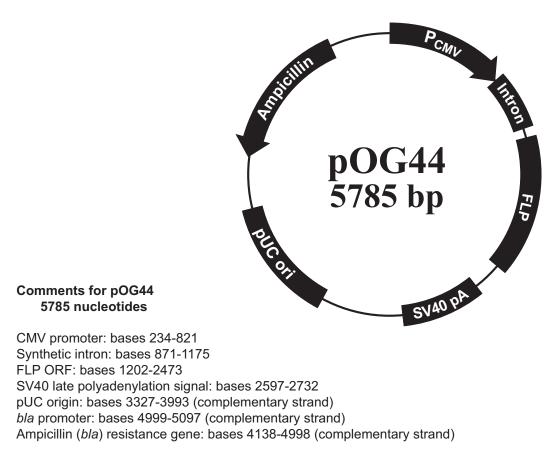

pOG44 Vector, Continued

Features of pOG44 The following table describes the relevant features of pOG44. All features have been functionally tested.

Feature	Benefit
Human cytomegalovirus (CMV) immediate early promoter	Allows high-level expression of the <i>FLP</i> gene (Andersson <i>et al.</i> , 1989; Boshart <i>et al.</i> , 1985; Nelson <i>et al.</i> , 1987).
Synthetic intron	Hybrid fragment which contains sequences derived from the adenovirus major late region and an IgG variable region (Huang and Gorman, 1990; O'Gorman <i>et al.</i> , 1991) and functions to enhance expression of the <i>FLP</i> gene.
<i>FLP</i> ORF (flp-F70L)	Encodes a temperature-sensitive Flp recombinase (Buchholz <i>et al.</i> , 1996) that mediates conservative recombination via FRT sites (O'Gorman <i>et al.</i> , 1991).
SV40 late polyadenylation signal	Allows polyadenylation of mRNA.
pUC origin	Allows high-copy number replication and growth in <i>E. coli.</i>
<i>bla</i> promoter	Allows expression of the ampicillin (<i>bla</i>) resistance gene.
Ampicillin (<i>bla</i>) resistance gene (ß-lactamase)	Allows selection of transformants in <i>E. coli.</i>

Accessory Products

Additional products

Many of the reagents used with the Flp-In[™] System are available separately. See the following table for ordering information.

Product	Quantity	Catalog no.
pFRT/ <i>lac</i> Zeo	20 µg	V6015-20
pFRT/ <i>lac</i> Zeo2	20 µg	V6022-20
pcDNA5/FRT	20 µg	V6010-20
T7 Promoter Primer	2 µg	N560-02
Zeocin™ Selection Reagent	8 × 1.25 mL	R250-01
	50 mL	R250-05
Hygromycin B	20 mL	10687-010
PureLink [®] HiPure Plasmid Miniprep Kit	100 preps	K2100-03
PureLink [®] HiPure Plasmid Midiprep Kit	25 preps	K2100-04
One Shot [®] TOP10F´ (chemically competent cells)	21 x 50 µL	C4040-03
One Shot® TOP10 Electrocomp™ (electrocompetent cells)	21 x 50 µL	C4040-52
Electrocomp [™] TOP10 (electrocompetent cells)	5 x 80 µL	C664-55

Flp-In[™] Expression Vectors

Additional Flp-In[™] expression vectors are available. For more information about the features of each vector or to download a manual for a vector, refer to **www.lifetechnologies.com/manuals** or call Technical Support (see page 9).

Product	Quantity	Catalog no.
pcDNA5/FRT/V5-His TOPO® TA Expression Kit	1 kit	K6020-01
pSecTag/FRT/V5-His TOPO® TA Expression Kit	1 kit	K6025-01
pEF5/FRT/V5 Directional TOPO® Expression Kit	1 kit	K6035-01
pEF5/FRT/V5-DEST Gateway™Vector Pack	6 µg	V6020-20

Accessory Products, Continued

Flp-In[™] Host Cell Lines

For your convenience, Life Technologies has available several mammalian Flp-In[™] host cell lines that stably express the *lacZ-Zeocin*[™] fusion gene from pFRT/*lac*Zeo or pFRT/*lac*Zeo2. Each cell line contains a single integrated FRT site as confirmed by Southern blot analysis. The cell lines should be maintained in medium containing Zeocin[™]. For more information about the Flp-In[™] Cell Lines, see www.lifetechnologies.com or call Technical Support (see page 9).

Cell Line	Quantity	Catalog no.
Flp-In™-293	3 x 10º cells, frozen	R750-07
Flp-In [™] -CV-1	3 x 10⁴ cells, frozen	R752-07
Flp-In [™] -CHO	3 x 10⁴ cells, frozen	R758-07
Flp-In [™] -BHK	3 x 10⁴ cells, frozen	R760-07
Flp-In [™] -3T3	3 x 10⁴ cells, frozen	R761-07
Flp-In [™] -Jurkat	3 x 10⁴ cells, frozen	R762-07

Technical Support

Obtaining support	For the latest services and support information for all locations, go to www.lifetechnologies.com/support .
	At the website, you can:
	 Access worldwide telephone and fax numbers to contact Technical Support and Sales facilities
	 Search through frequently asked questions (FAQs)
	 Submit a question directly to Technical Support (techsupport@lifetech.com)
	 Search for user documents, SDSs, vector maps and sequences, application notes, formulations, handbooks, certificates of analysis, citations, and other product support documents
	Obtain information about customer training
	 Download software updates and patches
Safety Data Sheets (SDS)	Safety Data Sheets (SDSs) are available at www.lifetechnologies.com/support .
Certificate of Analysis	The Certificate of Analysis provides detailed quality control and product qualification information for each product. Certificates of Analysis are available on our website. Go to www.lifetechnologies.com/support and search for the Certificate of Analysis by product lot number, which is printed on the box.
Limited product warranty	Life Technologies Corporation and/or its affiliate(s) warrant their products as set forth in the Life Technologies' General Terms and Conditions of Sale found on Life Technologies' website at www.lifetechnologies.com/termsandconditions . If you have any questions, please contact Life Technologies at www.lifetechnologies.com/support .

Purchaser Notification

Limited Use Label License No. 64: Flp-In™ System

Life Technologies Corporation ("Life Technologies") has a license to sell the Flp-In[™] System and its components ("System") to scientists for research purposes only, under the terms described below. Use of the System for any Commercial Purpose (as defined below) requires the user to obtain commercial licenses as detailed below. Before using the System, please read the terms and conditions set forth below. Your use of the System shall constitute acknowledgment and acceptance of these terms and conditions. If you do not wish to use the System pursuant to these terms and conditions, please contact Life Technologies' Technical Services within 10 days to return the unused and unopened System for a full refund. Otherwise, please complete the User Registration Card and return it to Life Technologies.

Life Technologies grants you a non-exclusive license to use the enclosed System for research purposes only. The System is being transferred to you in furtherance of, and reliance on, such license. You may not use the System, or the materials contained therein, for any Commercial Purpose without licenses for such purpose. Commercial Purpose includes: any use of the System or Expression Products in a Commercial Product; any use of the System or Expression Products in the manufacture of a Commercial Product; any sale of the System or Expression Products; any use of the System or Expression Products to facilitate or advance research or development of a Commercial Product; and any use of the System or Expression Products to facilitate or advance any research or development program the results of which will be applied to the development of a Commercial Product. "Expression Products" means products expressed with the System, or with the use of any vectors or host strains in the System. "Commercial Product" means any product intended for sale or commercial use.

Access to the System must be limited solely to those officers, employees and students of your entity who need access to perform the aforementioned research. Each such officer, employee and student must be informed of these terms and conditions and agree, in writing, to be bound by same. You may not distribute the System or the vectors or host strains contained in it to others. You may not transfer modified, altered, or original material from the System to a third party without written notification to, and written approval from Life Technologies. You may not assign, sub-license, rent, lease or otherwise transfer any of the rights or obligations set forth herein, except as expressly permitted by Life Technologies. This product is licensed under U.S. Patent Nos. 5,654,182 and 5,677,177 and is for research purposes only. Inquiries about licensing for commercial or other uses should be directed to: The Salk Institute for Biological Studies, 10010 North Torrey Pines Road, La Jolla, CA 92037, Attn.: Department of Intellectual Property and Technology Transfer. Phone: 858-453-4100 ext 1703; Fax: 858-450-0509; Email: mwhite@salk.edu .

Purchaser Notification, Continued

Limited Use Label License No. 284: Flp-In Services This product is the subject of intellectual property owned by The Salk Institute for Biological Studies, and licensed to Life Technologies Corporation. The purchase of this product conveys to the buyer the non-transferable right to use the purchased amount of the product and components of the product in research conducted by the buyer (whether the buyer is an academic or for-profit entity). The buyer cannot sell or otherwise transfer (a) this product (b) its components or (c) materials made using this product or its components to a third party or otherwise use this product or its components or materials made using this product or its components for Commercial Purposes. The buyer may transfer information or materials made through the use of this product to a scientific collaborator, provided that such transfer is not for any Commercial Purpose, and that such collaborator agrees in writing (a) not to transfer such materials to any third party, and (b) to use such transferred materials and/or information solely for research and not for Commercial Purposes. Commercial Purposes means any activity by a party for consideration and may include, but is not limited to: (1) use of the product or its components in manufacturing; (2) use of the product or its components to provide a service, information, or data to a third party; (3) use of the product or its components for therapeutic, diagnostic or prophylactic purposes; and/or (4) resale of the product or its components, whether or not such product or its components are resold for use in research. In addition, any use of the FLP technology outside of this product or the product's intended use, requires a separate license from the Salk Institute. Life Technologies will not assert a claim against the buyer of infringement of patents owned by Life Technologies and claiming this product based upon the manufacture, use or sale of a therapeutic, clinical diagnostic, vaccine or prophylactic product developed in research by the buyer in which this product or its components was employed, provided that neither this product nor any of its components was used in the manufacture of such product or for a Commercial Purpose. If the purchaser is not willing to accept the limitations of this limited use statement, Life Technologies is willing to accept return of the product with a full refund. For information on purchasing a license to this product for purposes other than research, contact Licensing Department, Life Technologies Corporation, 5791 Van Allen Way, Carlsbad, California 92008, Phone (760) 603-7200. Fax (760) 602-6500, or The Salk Institute for Biological Studies, 10010 North Torrey Pines Road, La Jolla, CA 92037, Attn.: Office of Technology Management, Phone: (858) 453-4100 extension 1275, Fax: (858) 546-8093.

Limited Use Label License: Research Use Only

The purchase of this product conveys to the purchaser the limited, nontransferable right to use the purchased amount of the product only to perform internal research for the sole benefit of the purchaser. No right to resell this product or any of its components is conveyed expressly, by implication, or by estoppel. This product is for internal research purposes only and is not for use in commercial services of any kind, including, without limitation, reporting the results of purchaser's activities for a fee or other form of consideration. For information on obtaining additional rights, please contact **outlicensing@lifetech.com** or Out Licensing, Life Technologies, 5791 Van Allen Way, Carlsbad, California 92008.

- Andersson, S., Davis, D. L., Dahlbäck, H., Jörnvall, H., and Russell, D. W. (1989). Cloning, Structure, and Expression of the Mitochondrial Cytochrome P-450 Sterol 26-Hydroxylase, a Bile Acid Biosynthetic Enzyme. J. Biol. Chem. 264, 8222-8229.
- Andrews, B. J., Proteau, G. A., Beatty, L. G., and Sadowski, P. D. (1985). The FLP Recombinase of the 2 Micron Circle DNA of Yeast: Interaction with its Target Sequences. Cell 40, 795-803.
- Argos, P., Landy, A., Abremski, K., Egan, J. B., Ljungquist, E. H., Hoess, R. H., Kahn, M. L., Kalionis, B., Narayana, S. V. L., and Pierson, L. S. (1986). The Integrase Family of Site-Specific Recombinases: Regional Similarities and Global Diversity. EMBO J. 5, 433-440.
- Ausubel, F. M., Brent, R., Kingston, R. E., Moore, D. D., Seidman, J. G., Smith, J. A., and Struhl, K. (1994). Current Protocols in Molecular Biology (New York: Greene Publishing Associates and Wiley-Interscience).
- Boshart, M., Weber, F., Jahn, G., Dorsch-Häsler, K., Fleckenstein, B., and Schaffner, W. (1985). A Very Strong Enhancer is Located Upstream of an Immediate Early Gene of Human Cytomegalovirus. Cell 41, 521-530.
- Broach, J. R., Guarascio, V. R., and Jayaram, M. (1982). Recombination Within the Yeast Plasmid 2mu Circle is Site-specific. Cell 29, 227-234.
- Broach, J. R., and Hicks, J. B. (1980). Replication and Recombination Functions Associated with the Yeast Plasmid, 2 mu Circle. Cell 21, 501-508.
- Buchholz, F., Ringrose, L., Angrand, P. O., Rossi, F., and Stewart, A. F. (1996). Different Thermostabilities of FLP and Cre Recombinases: Implications for Applied Site-specific Recombination. Nuc. Acids Res. 24, 4256-4262.
- Craig, N. L. (1988). The Mechanism of Conservative Site-Specific Recombination. Ann. Rev. Genet. 22, 77-105.
- Gronostajski, R. M., and Sadowski, P. D. (1985). Determination of DNA Sequences Essential for FLPmediated Recombination by a Novel Method. J. Biol. Chem. 260, 12320-12327.
- Huang, M. T. F., and Gorman, C. M. (1990). Intervening Sequences Increase Efficiency of RNA 3' Processing and Accumulation of Cytoplasmic RNA. Nuc. Acids Res. 18, 937-947.
- Jayaram, M. (1985). Two-micrometer Circle Site-specific Recombination: The Minimal Substrate and the Possible Role of Flanking Sequences. Proc. Natl. Acad. Sci. USA *82*, 5875-5879.
- Nelson, J. A., Reynolds-Kohler, C., and Smith, B. A. (1987). Negative and Positive Regulation by a Short Segment in the 5'-Flanking Region of the Human Cytomegalovirus Major Immediate-Early Gene. Molec. Cell. Biol. 7, 4125-4129.
- O'Gorman, S., Fox, D. T., and Wahl, G. M. (1991). Recombinase-Mediated Gene Activation and Site-Specific Integration in Mammalian Cells. Science 251, 1351-1355.
- Sambrook, J., Fritsch, E. F., and Maniatis, T. (1989). Molecular Cloning: A Laboratory Manual, Second Edition (Plainview, New York: Cold Spring Harbor Laboratory Press).

References, Continued

- Sauer, B. (1994). Site-Specific Recombination: Developments and Applications. Curr. Opin. Biotechnol. 5, 521-527.
- Senecoff, J. F., Bruckner, R. C., and Cox, M. M. (1985). The FLP Recombinase of the Yeast 2-micron Plasmid: Characterization of its Recombination Site. Proc. Natl. Acad. Sci. USA *82*, 7270-7274.

©2012 Life Technologies Corporation. All rights reserved.

The trademarks mentioned herein are the property of Life Technologies Corporation or their respective owners.

LIFE TECHNOLOGIES CORPORATION AND/OR ITS AFFILIATE(S) DISCLAIM ALL WARRANTIES WITH RESPECT TO THIS DOCUMENT, EXPRESSED OR IMPLIED, INCLUDING BUT NOT LIMITED TO THOSE OF MERCHANTABILITY, FITNESS FOR A PARTICULAR PURPOSE, OR NON-INFRINGEMENT. TO THE EXTENT ALLOWED BY LAW, IN NO EVENT SHALL LIFE TECHNOLOGIES AND/OR ITS AFFILIATE(S) BE LIABLE, WHETHER IN CONTRACT, TORT, WARRANTY, OR UNDER ANY STATUTE OR ON ANY OTHER BASIS FOR SPECIAL, INCIDENTAL, INDIRECT, PUNITIVE, MULTIPLE OR CONSEQUENTIAL DAMAGES IN CONNECTION WITH OR ARISING FROM THIS DOCUMENT, INCLUDING BUT NOT LIMITED TO THE USE THEREOF.

Notes

 Headquarters

 5791 Van Allen Way | Carlsbad, CA 92008 USA | Phone +1 760 603 7200 | Toll Free in USA 800 955 6288

 For support visit www.lifetechnologies.com/support or email techsupport@lifetech.com



www.lifetechnologies.com